



March 4, 2022

Christopher Dere, EPA Region II
UWS EPA Project Manager
290 Broadway
New York, NY 10007
Dere.Christopher@epa.gov

Supriya Rao, P.E., C.M., EPA Quality Assurance Officer
Air & Water Quality Assurance Team
Monitoring & Assessment Branch
2890 Woodbridge Avenue, Edison, NJ 08837
732-321-4461
Rao.Supriya@epa.gov

Dear Mr. Dere and Ms. Rao:

Circumstances surrounding Covid-19 are improving but the Unified Water Study (UWS) will continue to offer trainings and field audits via virtual platforms for the 2022 UWS Season. Specifically, annual trainings are being offered remotely via a virtual online platform with modules and annual field audits will commence in June utilizing video conferencing technology.

I have consulted with the UWS Science Advisors and the UWS Laboratory Coordinator and we agree that moving the remaining 2019, 2020, and 2021 Tier II nutrient sample analyses to the University of Connecticut Center for Environmental Sciences and Engineering Laboratory from the currently listed laboratory in the approved 2019, 2020, and 2021 UWS QAPPs is a prudent course of action and necessary to get these samples processed and results out in a timely manner. The UWS Science Advisors, UWS Laboratory Coordinator, and I agree that the certifications of the new laboratory meet or exceed certifications of the previous project laboratory as required by the 2019, 2020, and 2021 UWS QAPPs. This letter will be added to each of those QAPPs and distributed to the QAPP distribution list. Save the Sound has secured funding, and is currently working towards, outfitting our laboratory with capacity to analyze Tier II nutrient parameters in the UWS and are in the process of acquiring Environmental Laboratory Approval Program certification for these parameters.

Please accept this letter as an addendum to the 2022 version of the Quality Assurance Project Plan titled: *Long Island Sound Embayments Water Quality Monitoring QAPP for Monitoring Activities Conducted in the Unified Water Study: Long Island Sound Embayment Research*.

Best Regards,

A handwritten signature in black ink, appearing to read 'Peter Linderoth', is written over a light blue horizontal line.

Peter H. Linderoth
Save the Sound
UWS Monitoring Program Project Manager
plinderoth@savethesound.org

U.S. EPA REGION 2 - QUALITY ASSURANCE PROJECT PLAN APPROVAL FORM

PROJECT INFORMATION

Quality Assurance Officer: Supriya Rao

Project Officer: Christopher Dere

Title of Quality Assurance Project Plan: Long Island Sound Embayments Water Quality Monitoring QAPP

Assistance Agreement or Contract #: 96244321-0

QA File Number: 2022-31

REGIONAL QA MANAGER OR DELEGATED APPROVER

Approved ☒ Conditionally Approved* ☐

* conditional approval may be provided when there are unresolved comments that do not impact the data collection or the quality of the data and where the project has a small window of opportunity to collect such data. Conditional Approval expires 90 days from the signature date. If updated quality documentation (qdp) is not provided by the expiration date or another due date is not agreed upon by EPA, then the QD will be considered delinquent.

Comments:

Signature EPA QA Officer:

SUPRIYA RAO

Digitally signed by SUPRIYA RAO
Date: 2022.03.15 15:22:13 -04'00'

Signature EPA PO or Project lead

Christopher E. Dere

Digitally signed by Christopher E.
Dere
Date: 2022.03.15 15:50:00 -04'00'

REVIEW/ SUMMARY:

A review was conducted on the above referenced Quality Assurance Project Plan. The subject QAPP was reviewed for conformance with the [EPA Requirements for Quality Assurance Project Plans](#) (EPA QA/R-5), EPA/240/B-01/003, March 2001; USEPA Region 2 Guidance for the Development of QAPPs for Environmental Monitoring Projects, April 2004 and other EPA QAPP guidance documents as appropriate.

This approval form documents EPA's decision of approval or conditional approval* for the aforementioned QAPP. After the QAPP is approved by EPA via this approval form, obtain the required signatures from your organization on the QAPP Title/Signature page. Send the signed QAPP to the EPA Project Officer and others on the QAPP distribution list within the timeframe stipulated in the AA terms and conditions.

Long Island Sound Embayments Water Quality Monitoring QAPP

***For monitoring activities conducted in the
Unified Water Study: Long Island Sound Embayment Research.***

Monitoring Organizations

Ash Creek Conservation Association, Bronx River Alliance, Clean up Sound and Harbors (CUSH), Coalition to Save Hempstead Harbor, Connecticut River Conservancy, Cornell Cooperative Extension of Suffolk County Marine Program, Derecktor Shipyards, Friends of the Bay, Friends of the Farm River Estuary, North Fork Environmental Council, Earthplace, Inc. (Harbor Watch), Interstate Environmental Commission, New England Science & Sailing Foundation, Salonga Wetland Advocates Network, Save the River – Save the Hills, Inc., Save the Sound, Setauket Harbor Task Force, SoundWaters, The Maritime Aquarium at Norwalk, Town of Darien, Town of Fairfield – Conservation Department, Town of Stratford – Conservation Department, River Advocates of South Central Connecticut, Guardians of Flushing Bay, East River Watershed Research Institute, SUNY Maritime

Coordinating Organization

Save the Sound

Funded By

The United States Environmental Protection Agency – Long Island Sound Study

Version Date (yyyy.mm.dd)

2022.03.04

Date Approved

2022.03.15

Prior Associated Approved QAPP:

Mamaroneck Harbor and Little Neck Bay, NY, UWS Water Quality Monitoring QAPP

Approved by Kathryn Drisco, Quality Assurance Officer, EPA, Region 2, 8/3/2017

NFWF grant 53526, Connecticut Fund for the Environment, Water Quality Monitoring Initiative for Long Island Sound Embayments (NY), EPA Cooperative Agreement LI-00A00129-0 (FC.R278). Long Island Sound Embayment Water Quality Monitoring QAPP. For monitoring activities conducted as part of the Long Island Sound Tier 1 Unified Water Study (UWS). Approved by Esther Nelson, Quality Assurance Officer, EPA, Region 2, 6/5/2018. National Fish and Wildlife Foundation (NFWF), US EPA recipient via Cooperative Agreement LI 00A00382 (NFWF FC.R334). Long Island Sound Embayments Water Quality Monitoring QAPP. For monitoring activities conducted in the Unified Water Study: Long Island Sound Embayment Research Approved by Esther Nelson, Quality Assurance Officer, EPA Region 2, 5/2/2019, EPA Agreement No. LI96259818. Approved by Esther Nelson, Quality Assurance Officer, EPA Region 2, 3/16/2020, EPA Agreement No. LI96259818. Approved by Esther Nelson, Quality Assurance Officer, EPA Region 2, 4/20/2021, EPA Agreement No. LI96259818.

A. Project Management

A.1. APPROVAL PAGE



date: 4/6/2022

Peter Linderth, Monitoring Program Project
Manager Save the Sound




date: 4/6/2022

Peter Linderth, Monitoring Program Coordinator
Save the Sound



date: 4/6/2022

Jamie Vaudrey, UWS Science Advisor
Department of Marine Sciences, University of Connecticut

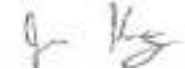

date: 4/6/2022

Jason Krumholz, UWS Science Advisor
McLaughlin Research Corporation

date: _____

Christopher E. Dere, Onondaga Lake Program Manager
US Environmental Protection Agency, Region 2

date: _____

Supriya Rao, P.E., C.M., Air & Water Quality Assurance Team
US Environmental Protection Agency, Region 2

A.2. Table of Contents

A. Project Management	2
A.1. APPROVAL PAGE.....	2
A.3. DISTRIBUTION LIST	5
A.4. PROJECT / TASK ORGANIZATION	12
A.5. PROBLEM DEFINITION / BACKGROUND	13
A.6. PROJECT / TASK DESCRIPTION	16
<i>A.6.a. Sampling Types Covered by this QAPP.....</i>	<i>20</i>
<i>A.6.b. Maps of Study Area.....</i>	<i>22</i>
<i>A.6.c. Annual Task Calendar.....</i>	<i>66</i>
A.7. DATA QUALITY OBJECTIVES.....	67
A.8. SPECIAL TRAINING / CERTIFICATION.....	76
A.9. DOCUMENTS AND RECORDS.....	78
B. Data Generation and Acquisition	80
B.1. SAMPLING PROCESS DESIGN (EXPERIMENTAL DESIGN)	80
B.2. SAMPLING METHODS	85
B.3. SAMPLE HANDLING AND CUSTODY	89
B.4. ANALYTICAL METHODS	90
B.5. QUALITY CONTROL.....	90
B.6. INSTRUMENT / EQUIPMENT TESTING, INSPECTION AND MAINTENANCE.....	91
B.7. INSTRUMENT / EQUIPMENT CALIBRATION AND FREQUENCY	92
B.8. INSPECTION / ACCEPTANCE OF SUPPLIES AND CONSUMABLES	93
B.9. NON-DIRECT MEASUREMENTS.....	93
B.10. DATA MANAGEMENT	94
C. Assessment and Oversight	97
C.1. ASSESSMENT AND RESPONSE ACTIONS.....	97

C.2. REPORTS TO MANAGEMENT.....98

D. Data Validation and Usability 99

D.1. DATA REVIEW, VERIFICATION, AND VALIDATION99

D.2. VERIFICATION AND VALIDATION METHODS100

D.3. RECONCILIATION WITH USER REQUIREMENTS.....100

E. Appendices 101

A.3. DISTRIBUTION LIST

Peter Linderoth, Monitoring Program Project Manager
Save the Sound
545 Tompkins Avenue, 3rd Floor
914-381-3140
tbrown@savethesound.org

Peter Linderoth, Monitoring Program Coordinator
Save the Sound
545 Tompkins Avenue, 3rd Floor
914-381-3140
plinderoth@savethesound.org

Elena Colón, Monitoring Program Field Coordinator
Save the Sound
545 Tompkins Avenue, 3rd Floor
914-381-3140
ecolon@savethesound.org

Peter Linderoth, Monitoring Program Quality Assurance Officer
Save the Sound
545 Tompkins Avenue, 3rd Floor
914-381-3140
plinderoth@savethesound.org

Elena Colón, Monitoring Program Lab Coordinator
Save the Sound
545 Tompkins Avenue, 3rd Floor
914-381-3140
plinderoth@savethesound.org

Elena Colón, Monitoring Program Data Management Coordinator
Save the Sound
545 Tompkins Avenue, 3rd Floor
914-381-3140
ecolon@savethesound.org

Jamie Vaudrey, UWS Science Advisor
Department of Marine Sciences, University of Connecticut
1080 Shennecossett Road, Groton, CT 06335
860-405-9149
jamie.vaudrey@uconn.edu

Jason Krumholz, UWS Science Advisor
McLaughlin Research Corporation
132 Johnnycake Hill Rd., Middletown RI 02842
401-787-0944
jkrumholz@gmail.com

Christopher E. Dere, EPA Project Officer
Onondaga Lake Program Manager
USEPA Region 2
290 Broadway, 24th Floor
New York, NY
212-637-3828
dere.christopher@epa.gov

Supriya Rao, P.E.,C.M., EPA Quality Assurance Officer Air
& Water Quality Assurance Team
Monitoring & Assessment Branch
2890 Woodbridge Avenue, Edison, NJ 08837
732-321-4461
Rao.Supriya@epa.gov

Bill Lucey
Save the Sound, Long Island Soundkeeper
900 Chapel Street, Upper Mezzanine
New Haven, CT 06510
blucey@savethesound.org

Katie Obrien-Clayton, CT DEEP Contact
CTDEEP WPLR, Environmental Analyst 2
79 Elm Street
Hartford, CT 06106
860-424-3176
Katie.obrien-clayton@ct.gov

Kelly Streich, CT DEEP Contact
Bureau of Water Management and Land Reuse
79 Elm Street
Hartford, CT 06106
860-424-3176
Kelly.streich@ct.gov

Alene Onion, NYS DEC Contact
NYSDEC, WAVE Coordinator
625 Broadway
Albany, NY 12233
518-402-8166
Alene.onion@dec.ny.gov

Ryan O'Donnell, UWS Monitoring Group Lead
Connecticut River Conservancy
15 Bank Row, Greenfield, MA 01301
860-704-0057
rodonnell@ctriver.org

Carolyn Sukowski, UWS Monitoring Group Lead
Cornell Cooperative Extension of Suffolk County Marine Program
423 Griffing Ave, Suite 100, Riverhead, NY 11901
631-293-1800 ext.21
cs424@cornell.edu

Michael Waterman, UWS Monitoring Group Lead
Derecktor Shipyards
311 E Boston Post Road, Mamaroneck, NY 10543
203-496-6210
mwight@derecktor.com

Heather Johnson, UWS Monitoring Group Lead
Friends of the Bay
111 South Street, Oyster Bay, NY 11771
516-922-6666
h.johnson@friendsofthebay.org

James Flynn, UWS Monitoring Group Lead
Friends of the Farm River Estuary
P.O. Box 2337
203-376-8120
farmriverfriends@gmail.com

Sarah Crosby, UWS Monitoring Group Lead
Earthplace, Inc. (Harbor Watch)
10 Woodside Lane, Westport, CT 06880
203-557-4403
s.crosby@earthplace.org

Evelyn Powers, UWS Monitoring Group Lead
and UWS Project Laboratory Manager
Interstate Environmental Commission
2800 Victory Blvd, Staten Island, NY 10314
718-982-3792
epowers@iec-nynjct.org

Megan Strand, UWS Monitoring Group Lead
New England Science & Sailing Foundation
72 Water Street, Stonington, CT 06378
860-595-9362
mstrand@nessf.org

Charles Muller, UWS Monitoring Group Lead
Salonga Wetland Advocates Network, Inc.
28 Marions Lane, Northport, NY 11768
631-626-8328
cjmull2003@yahoo.com

Deborah Moshier-Dunn, UWS Monitoring Group Lead
Save the River, Save the Hills, Inc.
35 Oswegatchie Road, Waterford, CT 06385
debm0727@sbcglobal.net

George Hoffman, UWS Monitoring Group Lead
Setauket Harbor Task Force, Inc.
146 Main Street, Setauket, NY 11733
631-786-6699
geohoff@optonline.net

Leigh Shemitz, UWS Monitoring Group Lead
SoundWaters
1281 Cove Road, Stamford, CT 06902
203-406-3304
lshemitz@soundwaters.org

Nicki Rosenfeld, UWS Monitoring Group Lead
The Maritime Aquarium of Norwalk
10 N Water Street, Norwalk, CT 06854
781-799-4868
nrosenfeld@maritimeaquarium.org

William Cavers, UWS Monitoring Group Lead
Town of Darien
2 Renshaw Road, Darien, CT 06820
203-940-0757
wwcavers@gmail.com

Timothy Bishop, UWS Monitoring Group Lead
Town of Fairfield – Conservation Department
725 Old Post Road, Fairfield, CT 06824
203-256-3071
tbishop@fairfieldct.org

Kelly Kerrigan, UWS Monitoring Group Lead
Town of Stratford – Conservation Department
550 Patterson Avenue, Stratford, CT 06614
203-385-4006
kkerrigan@townofstratford.com

Carol DiPaolo, UWS Monitoring Group Lead
Coalition to Save Hempstead Harbor
PO Box 159, Sea Cliff, NY 11579
516-801-6792
cshh@optonline.net

Maggie Greenfield, UWS Monitoring Group Lead
Bronx River Alliance
1 Bronx River Parkway
Bronx, NY 10462
718-430-4665
maggie.greenfield@bronxriver.org

Gail Robinson, UWS Monitoring Group Lead
Ash Creek Conservation Association
Bridgeport, CT
203-335-6690
ashcreekassoc@optonline.org

Mary Mushinsky, UWS Monitoring Group Lead
River Advocates of South Central Connecticut
1253 Whitney Avenue
Hamden, CT 06517
203-430-0921
marymushinsky@att.net

Rebecca Pryor, UWS Monitoring Group Lead
Guardians of Flushing Bay
782 Nostrand Ave. Apt. 2
Brooklyn, NY 11216
202-460-2065
rpryor@riverkeeper.org

Debra Okane, UWS Monitoring Group Lead
North Fork Environmental Council
PO Box 799
Mattituck, NY 11952.
dokane114@optonline.net

Kyle Rabin
Program Manager, Long Island Nitrogen Action Plan
Long Island Regional Planning Council
1864 Muttontown Road
Syosset, NY 11791
516-571-7613
krabin@lirpc.org

David Lipsky
Senior Policy Advisor
Bureau of Sustainability
NYC Department of Environmental Protection
59-17 Junction Blvd, 11th Floor
Flushing, NY 11363
718-595-5340
dlipsky@dep.nyc.gov

Beau Ranheim
Section Chief, Marine Sciences
New York City Department of Environmental Protection
718-595-5709
BeauR@dep.nyc.gov

Susan E. Van Patten, NYS DEC Contact
Watershed Section B, Section Chief
625 Broadway, Albany, NY 12233
518-402-8240
susan.vanpatten@dec.ny.gov

Michelle Golden, NYS DEC Contact
Environmental Program Specialist, Division of Water
625 Broadway, Albany, NY 12233
518-402-8225
michele.golden@dec.ny.gov

A.4. PROJECT / TASK ORGANIZATION

Table 1: Project Organization.

Key project personnel and their corresponding responsibilities.

Name(s)	<i>Project Title</i> - Responsibility
Peter Linderoth	Monitoring Program Project Manager – Oversees all aspects of project that incorporate the monitoring program including: fiscal management, project objectives, data uses, program changes, etc.
Peter Linderoth	Monitoring Program Coordinator – Monitoring Group recruitment and training. Develops the QAPP. Produces monitoring report. Produces or oversees outreach efforts, in coordination with project manager.
Elena Colón	Monitoring Program Field Coordinator – Responsible for assistance in training and quality assurance of monitoring groups for field work. Ensures field datasheets are properly filled out, samples and forms are transported to laboratories as needed, Standard Operations Procedures (SOPs) are being followed in entirety; and performs QA checks, including field audits, to make sure procedures are followed or corrected as needed (in collaboration QA officer and UWS Science Advisors).
Elena Colón	Monitoring Program Lab Coordinator – Makes arrangements with any lab(s) used to perform analyses according to QAPP. Ensures correct procedures are used, holding times are met, and adequate documentation is provided.
Elena Colón	Monitoring Program Data Management Coordinator – Maintains the data systems for the program. Performs/oversees data entry and checks entries for accuracy against field and lab forms.
Peter Linderoth	Monitoring Program Quality Assurance Officer – Runs Quality Assurance (QA) program.
Jamie Vaudrey and Jason Krumholz	UWS Science Advisors – Science consultants offering guidance and participating in trainings and station selection among other aspects of the project including quality assurance.
Christopher E. Dere	USEPA Project Officer – Oversees US EPA Cooperative Agreement compliance including processing recipient/subrecipient requests for QA/QC within EPA Regions
See Distribution List	UWS Monitoring Group Leads – Undertake UWS in their respective embayment(s) following all aspects of this QAPP.
Rao Supriya	USEPA Quality Assurance Officer – Reviews, comments and approves QAPP.
Changes by year. Individual names are not listed.	Monitoring Program Field Staff – Sample, perform field analyses, and assist in laboratory analyses and/or data entry.

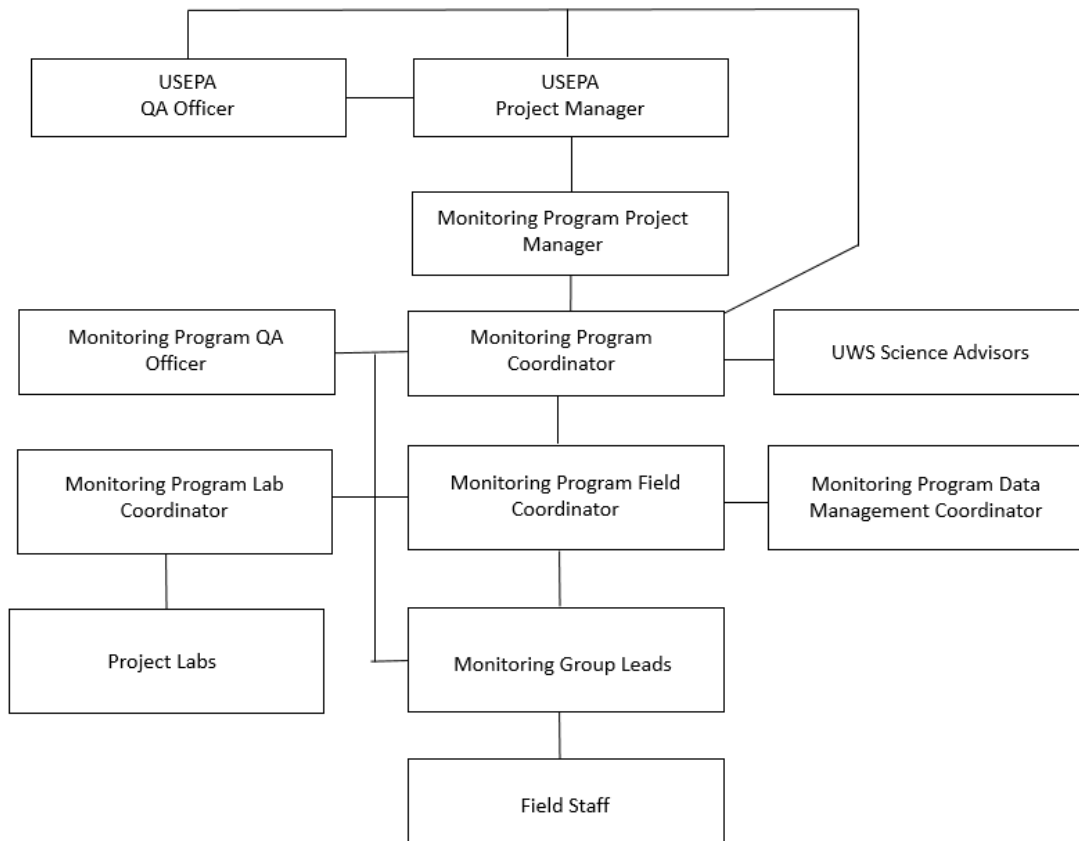


Figure 1: Organizational Chart.
Lines between boxes indicate direct communication.

A.5. PROBLEM DEFINITION / BACKGROUND

Despite three decades of effort to improve water quality, Long Island Sound (LIS) remains a severely stressed environment. In the western Sound, from Greenwich to Nassau County, dissolved oxygen concentrations—a key measure of the Sound’s health—consistently fall to levels too low to sustain aquatic wildlife. Low levels of dissolved oxygen, or hypoxia, are worsened by excess nitrogen (N) from outdated sewage collection systems, failing septic systems, contaminated stormwater runoff, and fertilizers. Moreover, there are serious

eutrophication-related impacts in embayments throughout the Sound¹. According to the EPA's Nitrogen Reduction Strategy, "Impairments linked to excess discharges of N include harmful algal blooms, low DO, poor water clarity, loss of submerged aquatic vegetation and tidal wetlands, and coastal acidification."

Main-stem Long Island Sound water quality data are abundant and readily available from Connecticut Department of Energy and Environmental Protection, New York City Department of Environmental Protection, and the Interstate Environmental Commission for analyses. However, there are limited environmental health data being collected in LIS embayments. The Unified Water Study (UWS) will fill in many of the data gaps that exist amongst LIS embayments. These data sets will have myriad of uses including comparing embayment environmental health, informing water quality management decisions, and conveying the information to the public so they can be better informed about the environmental health of LIS embayments.

The groups selected to participate in the 2018 UWS season went through an application process administered by the Monitoring Program Project Manager, Monitoring Program Coordinator, and the Monitoring Program Field Coordinator. A standardized application form was distributed and the administrators used a metric to select groups. In 2018, there were 19 monitoring groups monitoring a total of 33 embayments. In 2019, three new groups were added to the UWS. The 2019 group and embayment total is 22 groups in 37 embayments. 12 of the 37 embayments in the UWS will also receive Tier II monitoring. In 2020, one new group is being added to the UWS. The 2020 group and embayment total is 23 groups in 38 embayments. 13 of the embayments will receive Tier II monitoring. The 2021 group and embayment total is 24 groups monitoring 41 embayments. 13 of the embayments will receive Tier II monitoring. The 2022 group and embayment total is 26 groups monitoring 44 embayments. 12 of the 44 embayments will receive Tier II monitoring.

The final locations of the Tier I embayments in the UWS were dependent on the applications and respective monitoring group's interests. Tier II embayments were selected referencing priority embayment plans put forth by Connecticut Department of Energy and Environmental Protection, New York State Department of Environmental Conservation, and Long Island Sound Study. Monitoring group experience was also a factor in the decision.

¹ Vaudrey, J. M., Yarish, C., Kim, J. K., Pickerell, C., Brousseau, L., Eddings, J., & Sautkulis, M. (2016). Comparative analysis and model development for determining the susceptibility to eutrophication of Long Island Sound embayments. Connecticut Sea Grant Final Project Report, 38.

Organizational History and Mission

The mission of Save the Sound (STS) is to protect and improve the land, air and water of Connecticut and Long Island Sound. Founded in 1978, STS merged in 2004 with Save the Sound, a respected voice for the protection of Long Island Sound's shoreline, marine habitat and water quality with a track record of more than 40 years. The proposed project is in line with one of Save the Sound's strategic goals: "Our Long Island Sound, rivers and lakes are safe for people and wildlife."

Data collected under this QAPP will be collected in a manner to allow the data to be used as part of the Unified Water Study. The UWS is a coordinated effort among groups monitoring Long Island Sound with the goal of comparing water quality parameters associated with eutrophication within and among embayments. The UWS is comprised of two tiers, Tier I and Tier II. Tier I data is required for entry into the study. The parameters monitored in Tier I of the study are dissolved oxygen, conductivity (salinity), chlorophyll a, temperature, turbidity, and qualitative macrophyte assessments. Tier II parameters monitored as part of the UWS are nutrient concentrations, logged dissolved oxygen and conductivity (salinity) data, and underwater camera quantitative assessments of macrophytes. Tier I and Tier II parameters are covered in this QAPP.

Monitoring History and Status

The New York Office of Save the Sound initiated a pathogen-indicator and water quality monitoring program in 2013 and has since expanded the spatial and temporal scale of the water quality monitoring program. Save the Sound was the lead facilitator in the development of the UWS and participated in the 2016 UWS pilot season. Save the Sound continues to participate in the UWS as the coordinating organization as well as a monitoring group. Measuring the eutrophic conditions in the bays and harbors of Long Island Sound directly relates to Save the Sound's overarching goal of reducing nitrogen and other pollutants in the Sound.

The UWS conducted a pilot season for the Tier I parameters with four existing monitoring groups in 2016. The goal of the 2016 season was to develop protocols that followed standard methods for embayment monitoring. All groups involved with the 2016 season had previous experience monitoring their embayments and were involved in ongoing monitoring programs. Groups involved with the 2016 pilot season, and other advisory participants such as academics and federal and municipal agencies, assisted with developing and finalizing the UWS Tier I SOPs.

2017 marked the inaugural season of the UWS with twelve groups participating in the study. Eleven of these groups monitored in Long Island Sound and one monitored off the south shore of Long Island, NY. These groups maintained their own QAPPs based on a template. These QAPPs provided integrity to the Study but monitoring groups in the UWS now operate under a collective EPA-approved QAPP, this document. The groups are responsible for adhering to the

QAPP and Save the Sound will ensure that quality objectives are met for each embayment in the study as outlined in this document. The collective QAPP will be updated upon changes to embayments, participating groups, or procedures in the study.

Monitoring and Data Use Objectives

Data collected under this QAPP will be collected in a manner to allow the data to be used as part of the UWS. The UWS is designed to facilitate equitable water quality comparisons across Long Island Sound embayments. All data will be available to the public via request or download from Save the Sound's website. Data will also be uploaded to the Environmental Protection Agency Water Quality Portal. Save the Sound has met, and continues to meet and discuss, UWS data usage by Connecticut Department of Energy and Environmental Protection, New York State Department of Environmental Conservation, and UWS monitoring groups are welcome and encouraged to share their data on a local level to potentially elicit changes that could improve water quality in their respective embayments. There is also an overarching primary project goal to include the UWS data sets in the Long Island Sound Report Card which is issued by Save the Sound. The report card compares water quality indicators (dissolved oxygen, nutrients, chlorophyll a, and water clarity) to scientifically derived thresholds or goals. These indicators are combined into an overarching Water Quality Index, which is presented as a subregion percent score. The report card provides a geographically specific assessment of annual Long Island Sound ecosystem health.

Additionally, the activities covered under this QAPP will provide quality-controlled data that can support secondary goals such as the assessment and restoration of coastal embayments and watersheds through the implementation of programs such as but not limited to:

- EPA's 305(b) water body health assessments and 303(d) TMDL development for impaired waters
- Clean Water Act Section 319 projects
- Connecticut Department of Energy and Environmental Protection and New York Department of Environmental Conservation Watershed Management Plans
- Long Island Sound Study's Comprehensive Conservation and Management Plan
- Long Island Sound Study Environmental Indicators Project
- New York State Department of Environmental Conservation Long Island Nitrogen Action Plan

A.6. PROJECT / TASK DESCRIPTION

Five types of monitoring stations are included:

- 1) *Tier I water quality* stations are sampled within three hours of sunrise between the months of May through October using a multiparameter sonde. A minimum of four

stations per embayment are required. If multiple regions of the embayment are delineated, a minimum of three stations are required per region. Monitoring groups will plan to sample Tier I water quality stations at a minimum every two weeks from May to October. However, if unforeseeable circumstances make this plan not possible, a minimum of six sample events between May and October are required for inclusion in the UWS; including at least one sample event in each of the months of June, July, August, and September. The number and location of Tier I water quality stations for embayments in the UWS are provided in Section A.6.b.

Tier I Water Quality – Water quality parameters are selected to facilitate comparisons between embayments. Monitoring groups must collect the following data to be included in the UWS:

- for each station
 - GPS coordinates of stations, recorded each sample date
 - Date and time
 - Total water depth
 - 0.5 m below the surface, 0.5 m above bottom, mid-depth if total depth >10 m; if total depth is less than 1.5 m, only a mid-depth reading will be collected
 - Temperature
 - Conductivity (salinity)
 - Dissolved oxygen
 - Chlorophyll a
 - Turbidity
 - Once per field day undertake a replicate profile including all parameters
- Obtain from an online NOAA tide table and weather station approved by Monitoring Program Coordinator:
 - Time of high and low tide nearest time of sampling
 - High and low air temperature for 24 hours preceding field sampling
 - Precipitation out a week preceding sampling event
- Within 1 day of the field sampling day, read the GPS of a land-based reference station

Monitoring groups will collect water quality data for the Unified Water Study according to procedures provided in the UWS SOP Depth and GPS, UWS SOP Sonde Profile, and UWS SOP Filtered Chlorophyll in Appendix A.

- 2) *Qualitative Tier I Macrophyte* stations are land-based or boat-based. They are sampled only mid-summer and may be sampled on different days from the water quality stations. Sampling occurs on three separate days between July 15 and August 7. Two sample days or a date slightly outside of the date criteria may be sufficient in the event of unforeseen complications. This decision will be made by the Monitoring Program Quality Assurance Officer and UWS Science Advisor(s).

The goal of this part of the UWS is to identify potential problem areas, versus characterizing the overall condition of the embayment; field teams will look for areas with the highest macrophyte abundance they can find. Groups will also look for and note the presence of eelgrass (a beneficial condition). This is not intended to be a quantitative assessment. Macrophyte surveys will complement chlorophyll a concentrations to better understand the dominant primary producer in the system.

Qualitative Tier I Macrophytes - Monitoring groups must collect the following data to be included in the UWS:

- for each macrophyte station
 - GPS coordinates of stations, recorded each sampling date
 - Date and time
 - Photos of macrophytes
- Within 1 day of the field sampling day, read the GPS of a land-based reference station

Groups will submit data and photos to the Unified Water Study according to methods provided in the UWS SOP Qualitative Macrophytes in Appendix A.

- 3) *Quantitative Tier II Macrophyte* stations are boat-based. They are sampled only mid-summer and are typically sampled on different days from the water quality stations. Sampling occurs on one day between July 15 and August 7.

Quantitative Tier II Macrophytes - Monitoring groups must collect the following data to be included in the UWS:

- For each macrophyte station
 - GPS coordinates
 - Date and time interval
 - Video of macrophyte abundance
- Within 1 day of the field sampling day, read the GPS of a land-based reference station

Groups will submit data and photos to the Unified Water Study according to methods provided in the UWS SOP Macrophyte Percent Coverage Via Camera in Appendix A.

- 4) *Tier II nutrients* stations are sampled a minimum every two weeks from May to October. However, if unforeseeable circumstances make this plan not possible, a minimum of six sample events between May and October are required for inclusion in the UWS; including at least one sample event in each of the months of June, July, August, and September. At a minimum, there will be two stations per embayment. In embayments with multiple regions, there will be a minimum of two stations per region. The number and location of Tier II nutrients stations for embayments in the UWS are provided in Section A.6.b. Stations in tributaries are sampled for nutrients on the same day as the embayment they flow into is sampled. These stations are chosen with respect to access and salinity value <1 ppt. A Long Island Sound reference station will also be sampled for nutrients on the

same day the embayment stations are sampled.

Tier II Nutrients - Monitoring groups must collect the following data to be included in the UWS:

- For each water quality station
 - GPS coordinates of stations, recorded each sample date
 - Date and time
 - 0.5 m below the surface
 - Total Nitrogen
 - Total Dissolved Nitrogen
 - Dissolved Inorganic Nitrogen Species (nitrate, nitrite, ammonia)
 - Total Phosphorous
 - Orthophosphate (also known as dissolved inorganic phosphorus)
 - Salinity

Groups will submit data to the Unified Water Study according to methods provided in the UWS SOP Filtered Nutrients and UWS SOP Total Nitrogen and Total Phosphorous in Appendix A.

- 5) *Continuous* Dissolved Oxygen stations will be selected with consideration to representativeness and where access is granted for maintenance of the equipment. A minimum of one continuous dissolved oxygen station is required per embayment region. Logging will commence from May to October 31.

Continuous Dissolved Oxygen - Monitoring groups must collect the following data to be included in the UWS:

- for each continuous dissolved oxygen station
 - GPS coordinates of stations, recorded each sampling date
 - Date and time
 - Dissolved Oxygen
 - Conductivity (Salinity)
 - Barometric Pressure
 - Temperature

Groups will submit data to the Unified Water Study according to methods provided in the UWS SOP Continuous Dissolved Oxygen in Appendix A.

Roles of Project Participants

The Monitoring Program Coordinator and Monitoring Program Field Coordinator or designee will provide guidance and advisement to the groups participating in the UWS, conferring with the UWS Science Advisors as needed. They will conduct trainings, field audits, station selection guidance, ongoing technical support, and lab coordination among other activities. The full set of participants and their respective roles can be referenced in Table 1 of this document. Figure 1 outlines the lines of communications between project participants.

UWS Project Laboratory Manager has agreed to the UWS SOP Filtered Chlorophyll, UWS SOP Total Nitrogen and Total Phosphorous, and UWS SOP Filtered Nutrients for sample collection in the field. The project laboratory has provided their laboratory method SOPs which can be reviewed in Appendix C of this QAPP. The laboratory will adhere to both the UWS and their lab SOPs.

Participating Monitoring Groups will be responsible for conducting field work and analyses following the requirements presented in the UWS SOPs. Monitoring Group Leads or designated appointees will complete all required training. Monitoring Group members will complete all required data sheets and chain of custody forms. Any problems or deviance from this QAPP or SOPs will immediately be reported to the Monitoring Program Field Coordinator who will confer with the Monitoring Program Quality Assurance Officer on corrective course of action.

How the proposed sampling plan supports the Monitoring Program objectives

Data collected under this QAPP will be collected in a manner to allow the data to be used as part of the UWS. The UWS is a coordinated effort among groups monitoring Long Island Sound embayments with the goal of comparing water quality and macrophyte abundance within and among embayments.

Overview of data handling processes

Sampling event and field data will be collected on standardized field and instrument calibration sheets. These standardized datasheets are in Appendix B of this document.

If a field team is delivering samples to a centralized location for laboratory analysis by a member of the Monitoring Group, the field data sheet is sufficient as a chain of custody record. In this scenario a chain of custody form will not be required as sufficient information is contained on the sample event datasheet.

If a field team is delivering a sample for analysis by a lab external to the monitoring group, the UWS Chlorophyll a Chain of Custody Form or UWS Nutrient Chain of Custody Form is required. These forms are in Appendix B of this document.

A.6.a. Sampling Types Covered by this QAPP

The type of sample information that can be collected under this QAPP includes:

- GPS location to identify and track station locations
- Total water depth of the sample station; and depth of sample location
- Temperature
- Conductivity (Salinity)
- Dissolved oxygen concentration and percent saturation
- Chlorophyll a concentrations (filtered water sample)

- Chlorophyll a concentrations (*in situ* fluorescence)
- Turbidity
- Qualitative assessment of macrophytes
- Quantitative assessments of macrophytes
- Nitrogen forms to measure nutrient levels
- Phosphorous forms to measure nutrient levels

This section of the page intentionally left blank.

A.6.b. Maps of Study Area

Figure 2a-z. All UWS study sites. Reporting regions if present are clearly delineated. Red circles are water quality station locations. The maps are overlain by a hexagonal grid. The grid was used to select stations to represent the water quality of the entire embayment using a probability-based sampling design², as in the EPA National Coastal Assessment³. In some cases, hexagons have been joined to represent a local area considered similar or if a hexagon included large sections of land.

This section of the page intentionally left blank. Maps of study sites start below.

² Paul, J.F., J.L. Copeland, M. Charpentier, P.V. August, and J.W. Hollister. 2003, Overview of GIS applications in estuarine monitoring and assessment research. *Marine Geodesy Journal* 26: 63-72.

³ EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

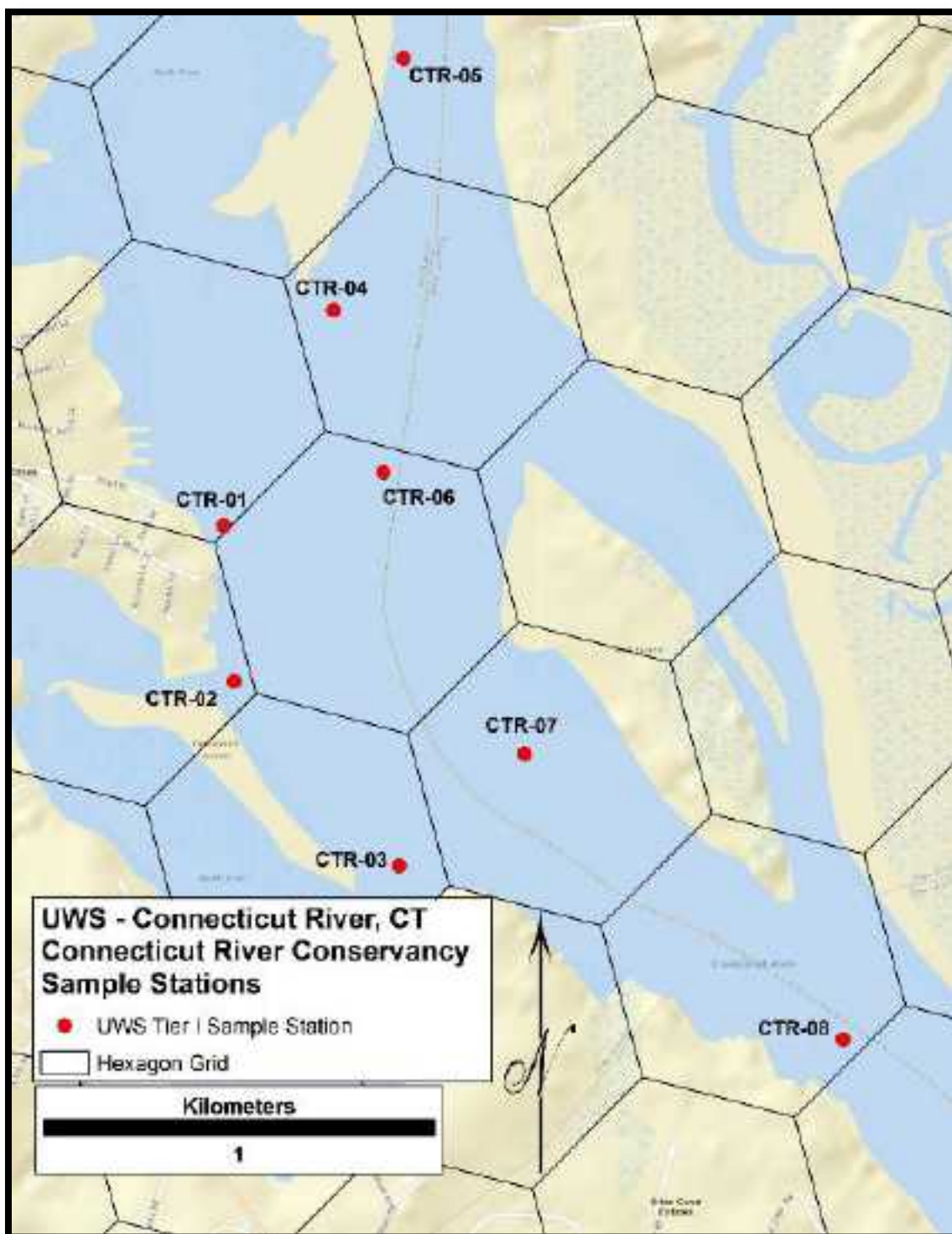


Figure 2a

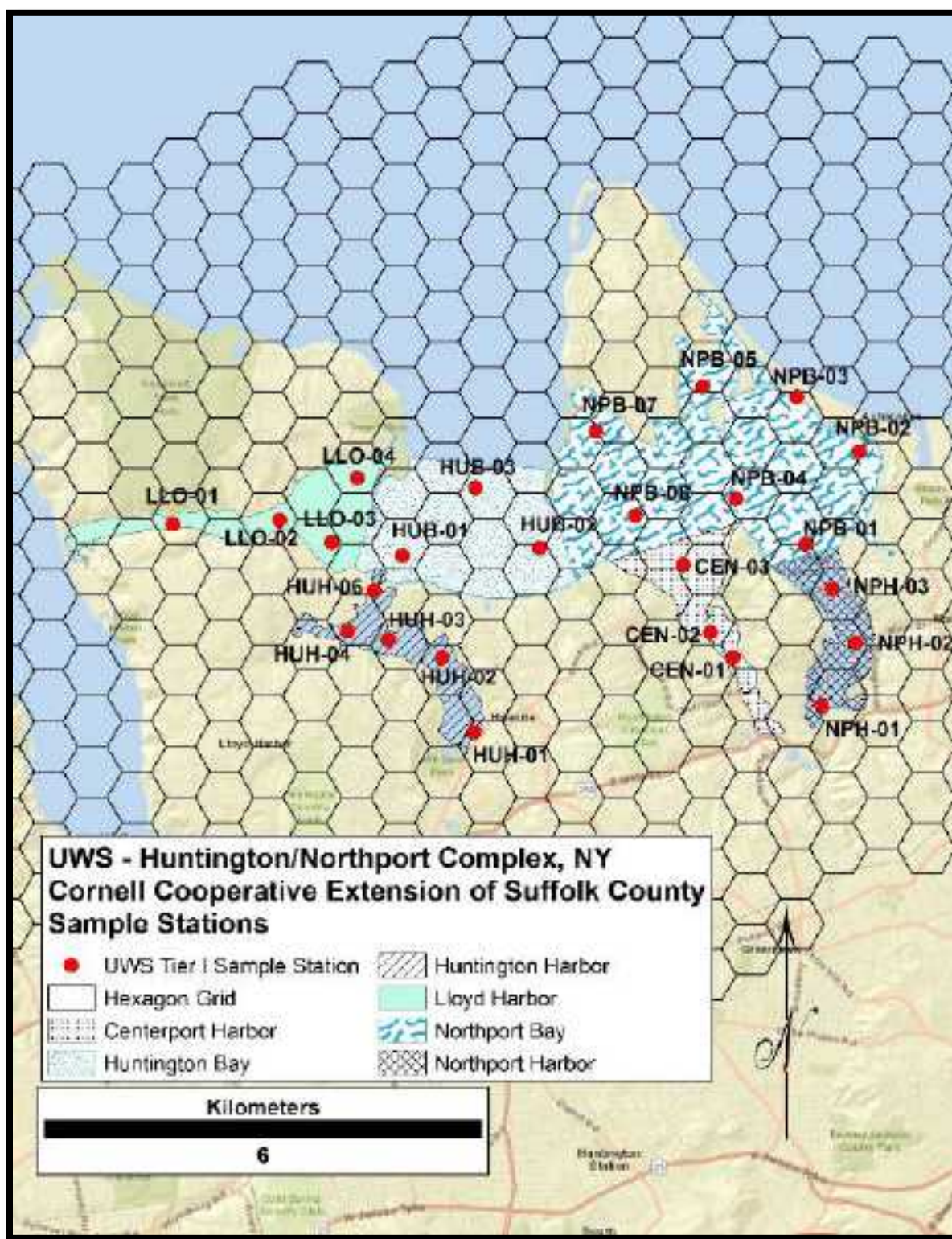


Figure 2b

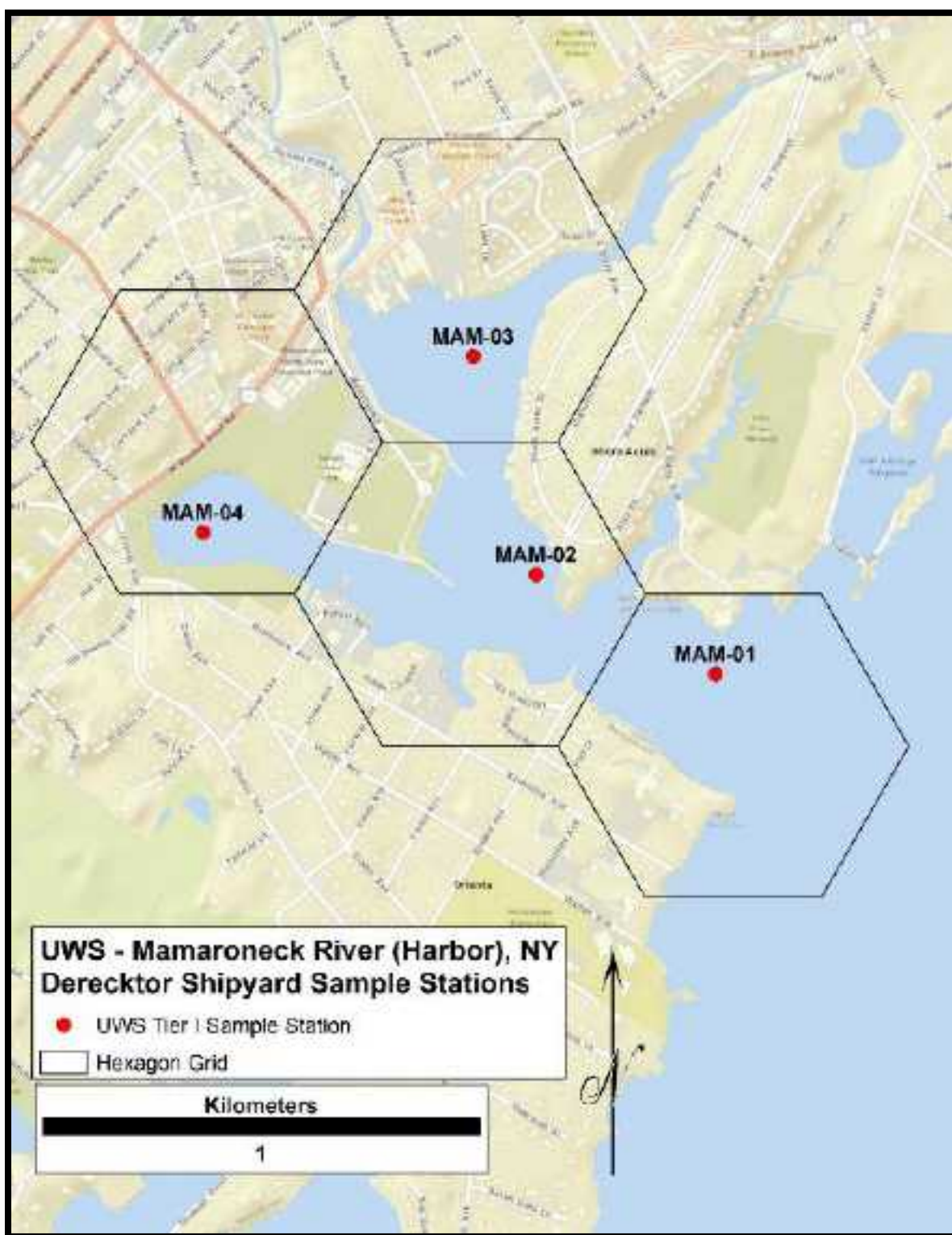


Figure 2c

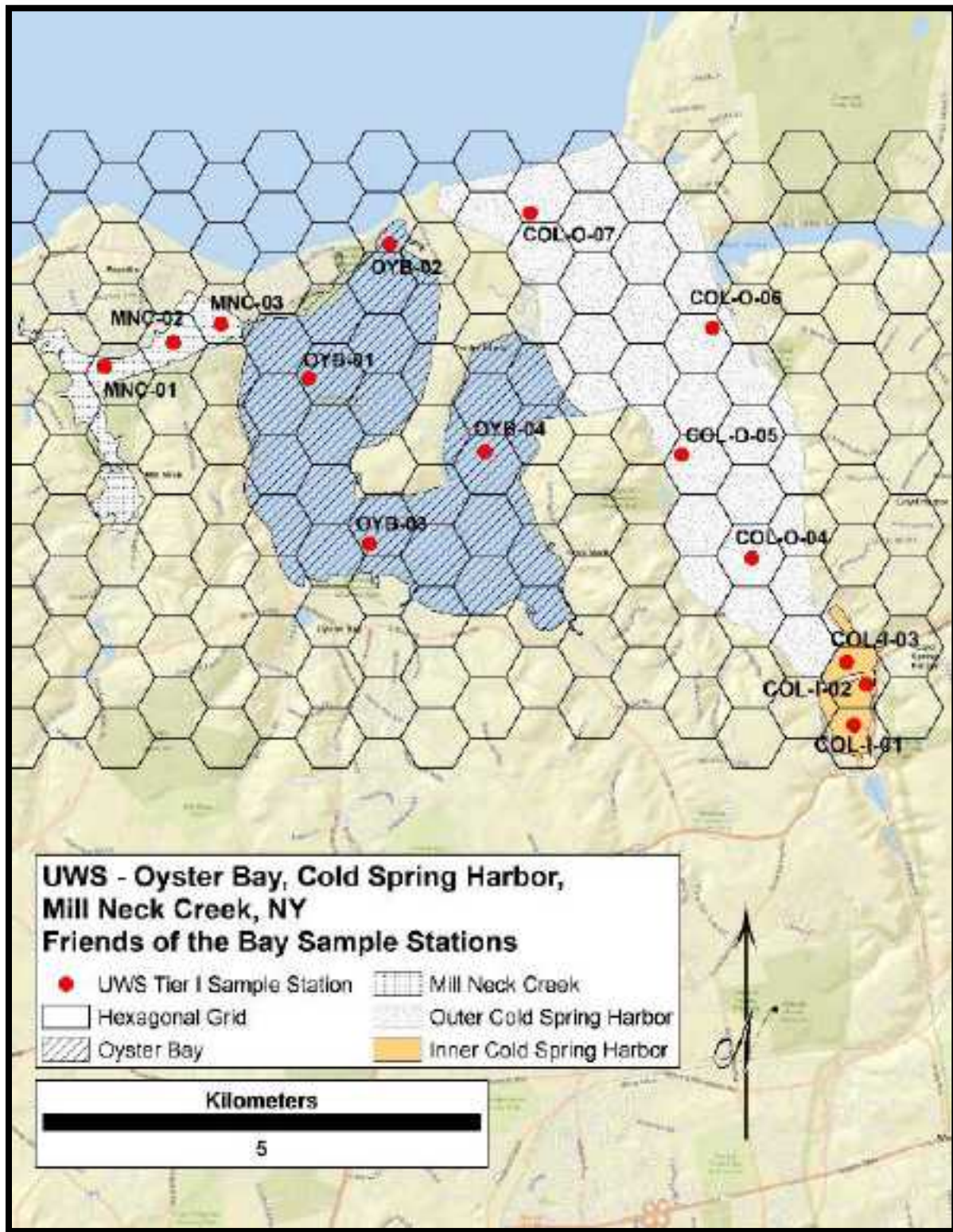


Figure 2d

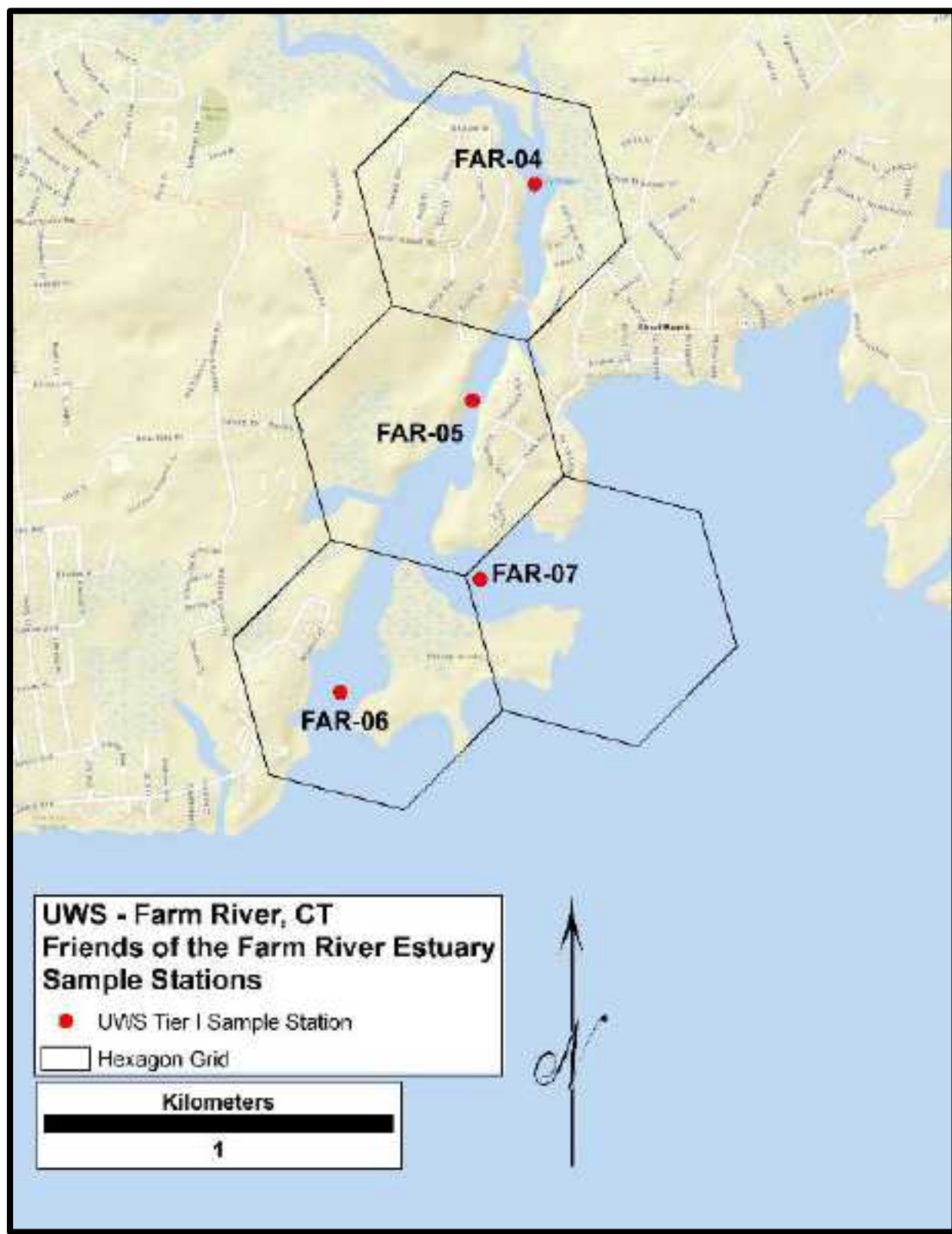


Figure 2e

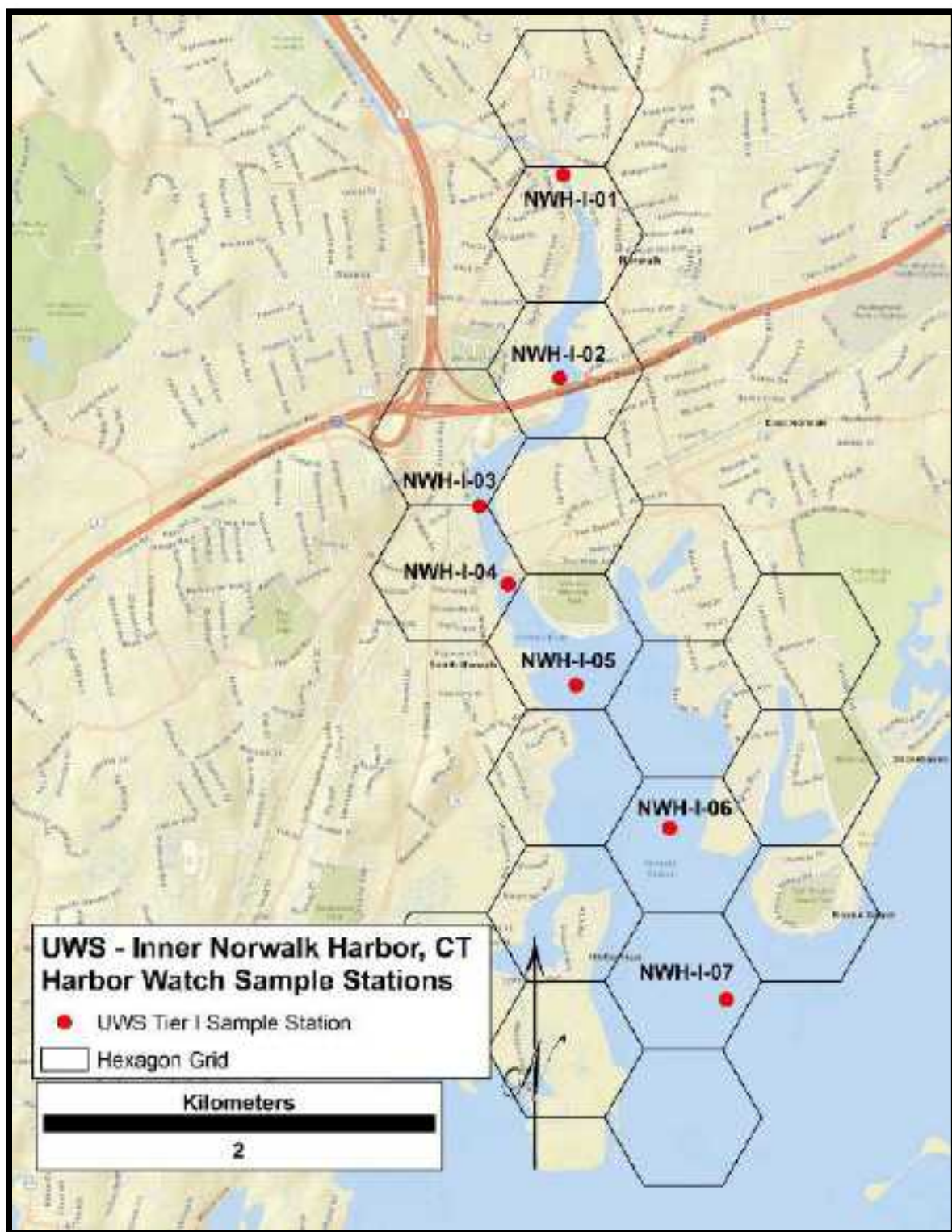


Figure 2f

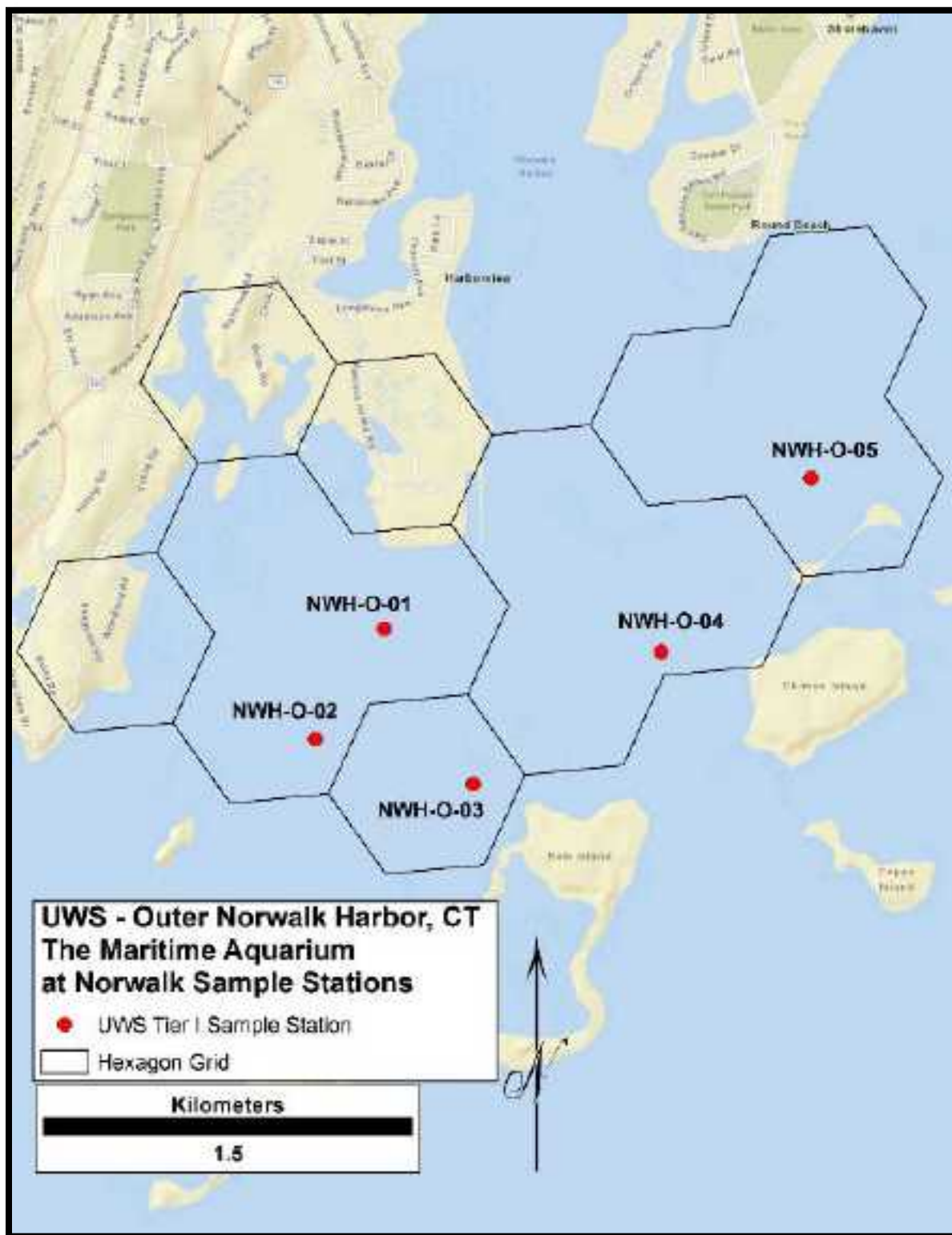


Figure 2g

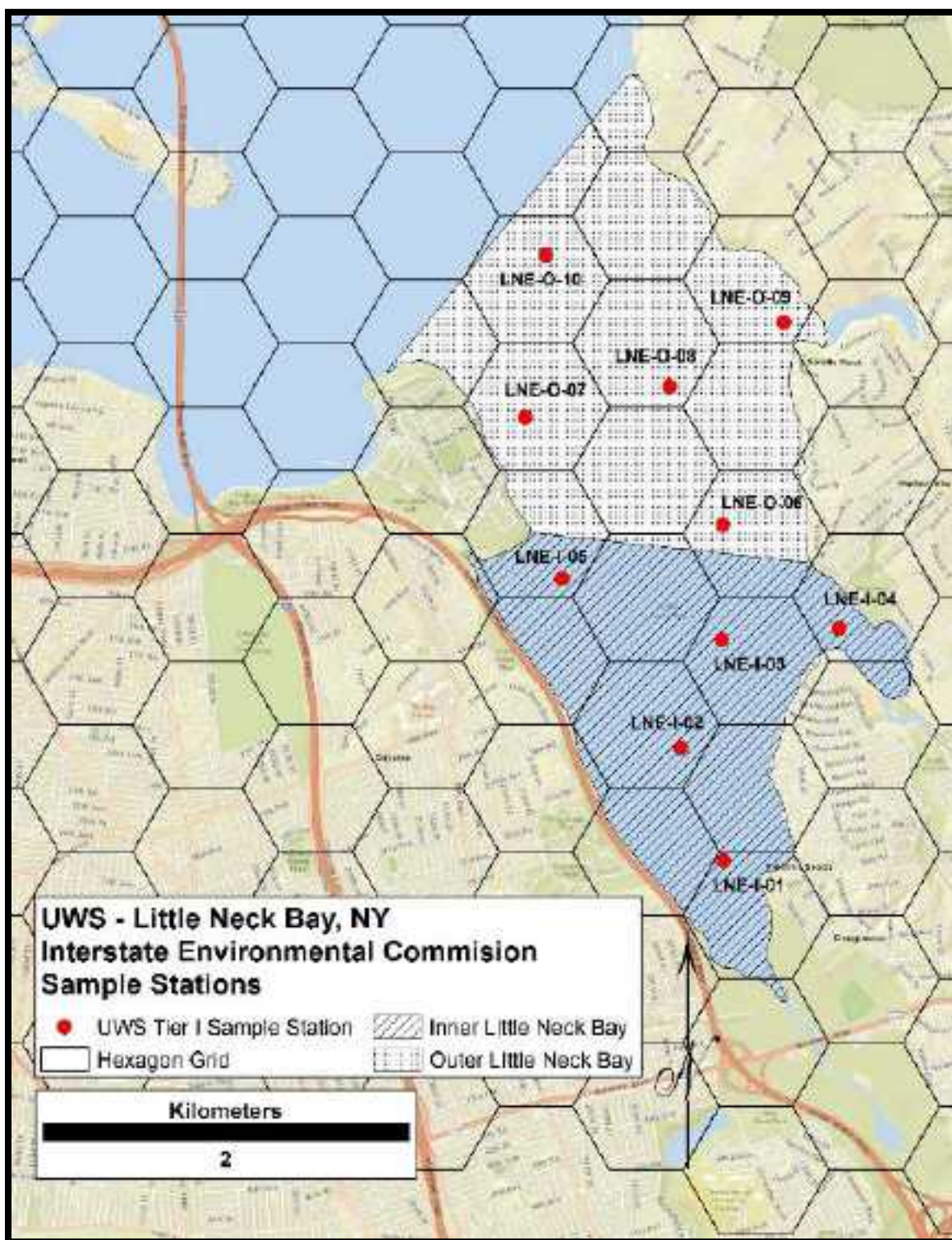


Figure 2h

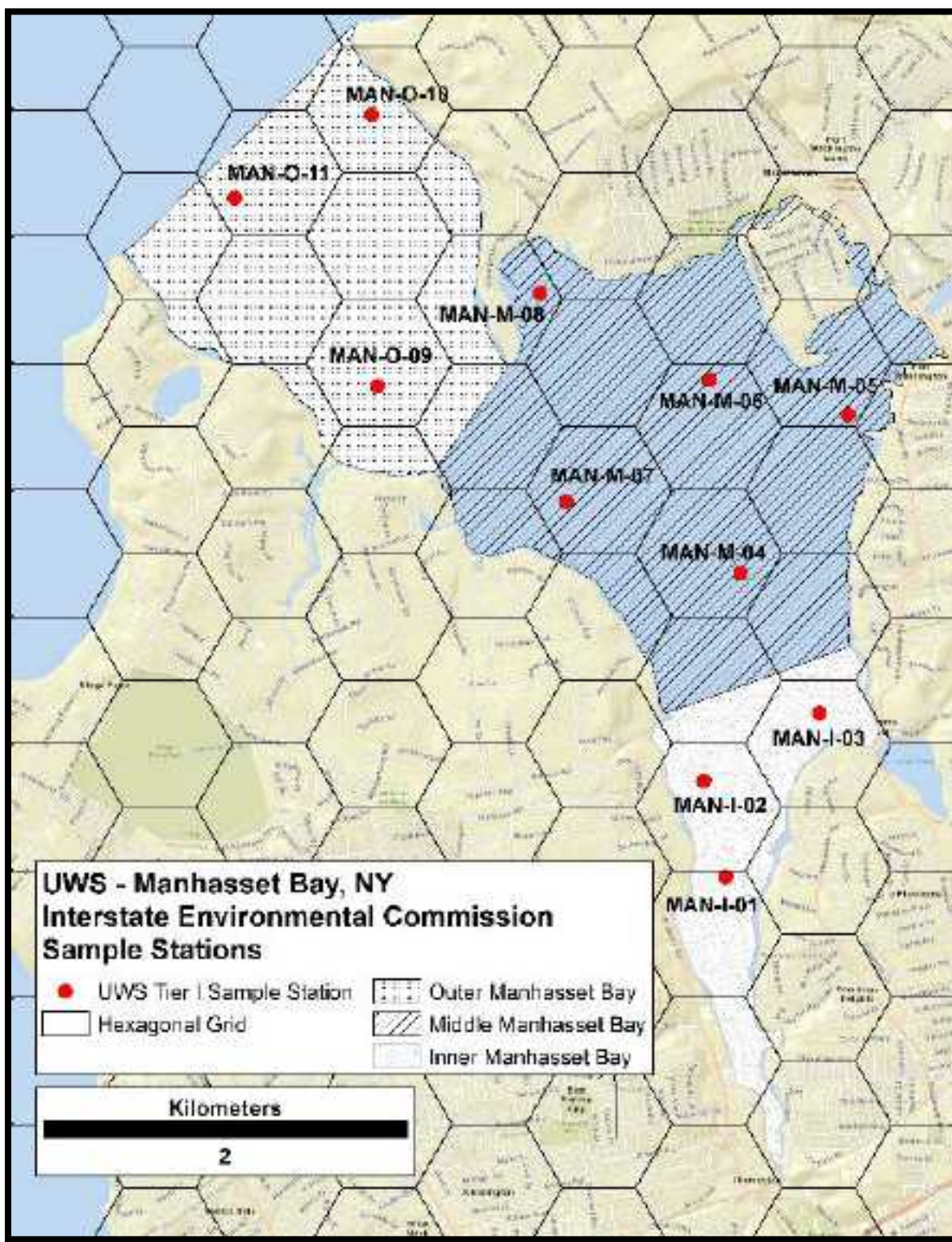


Figure 2i

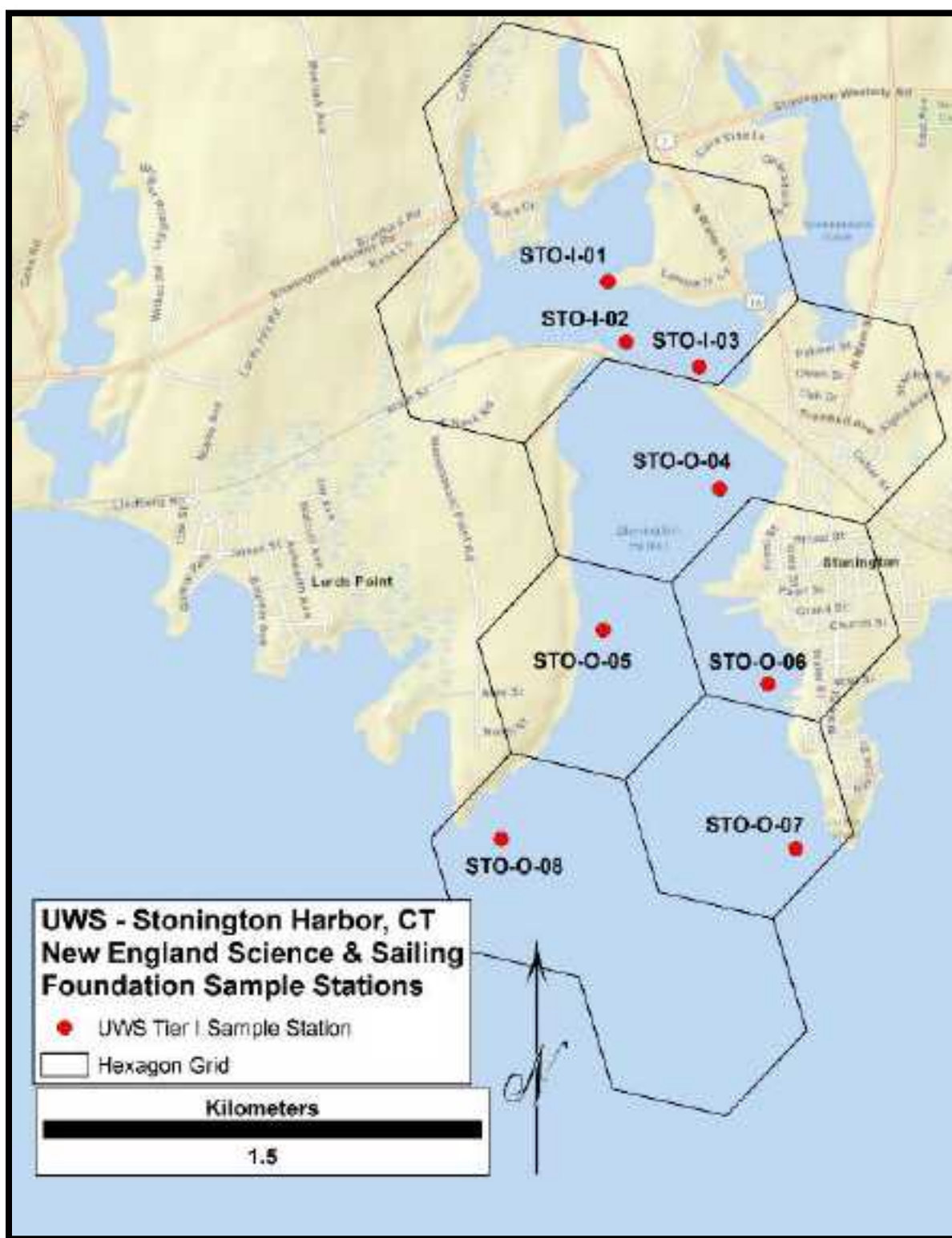


Figure 2j

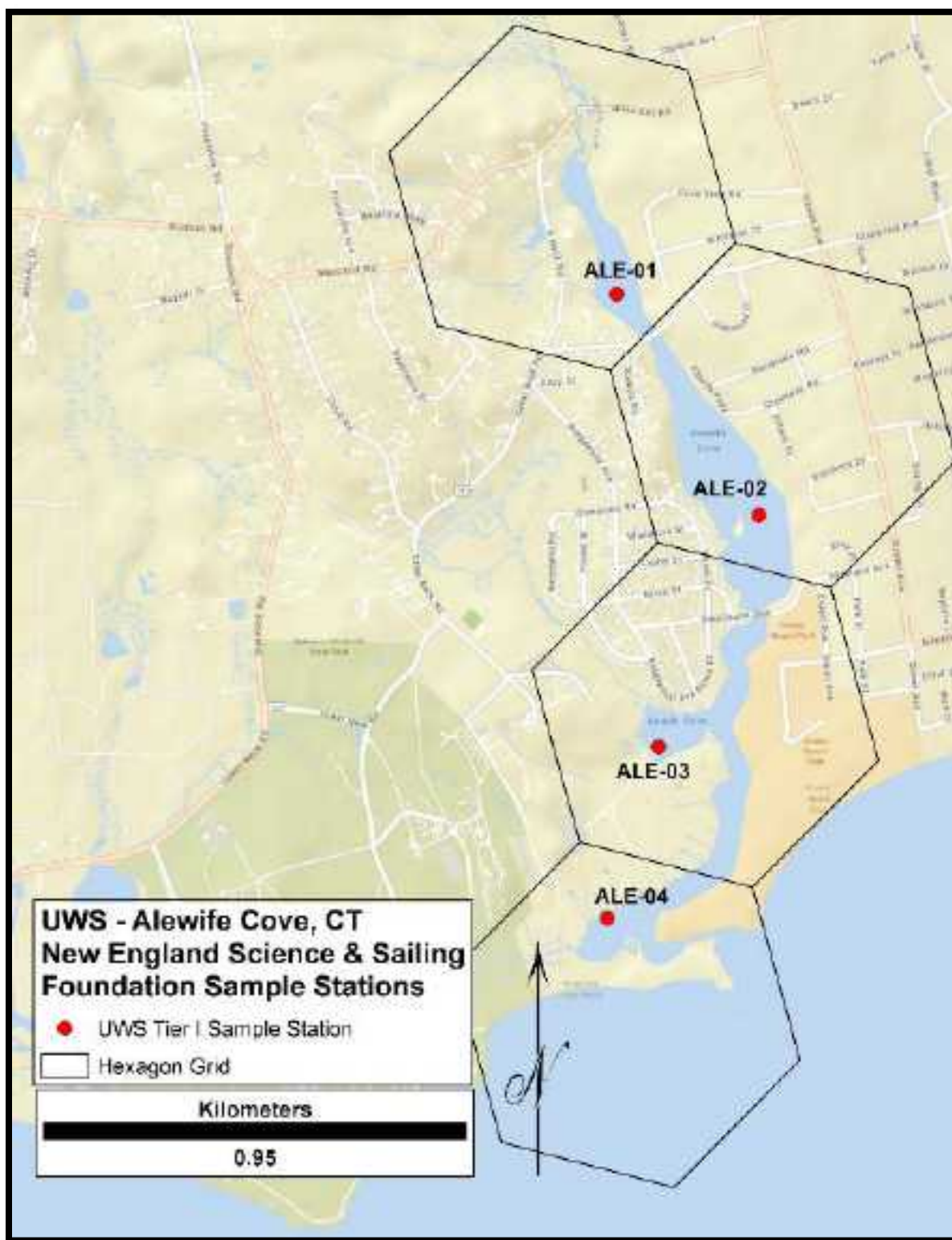


Figure 2k

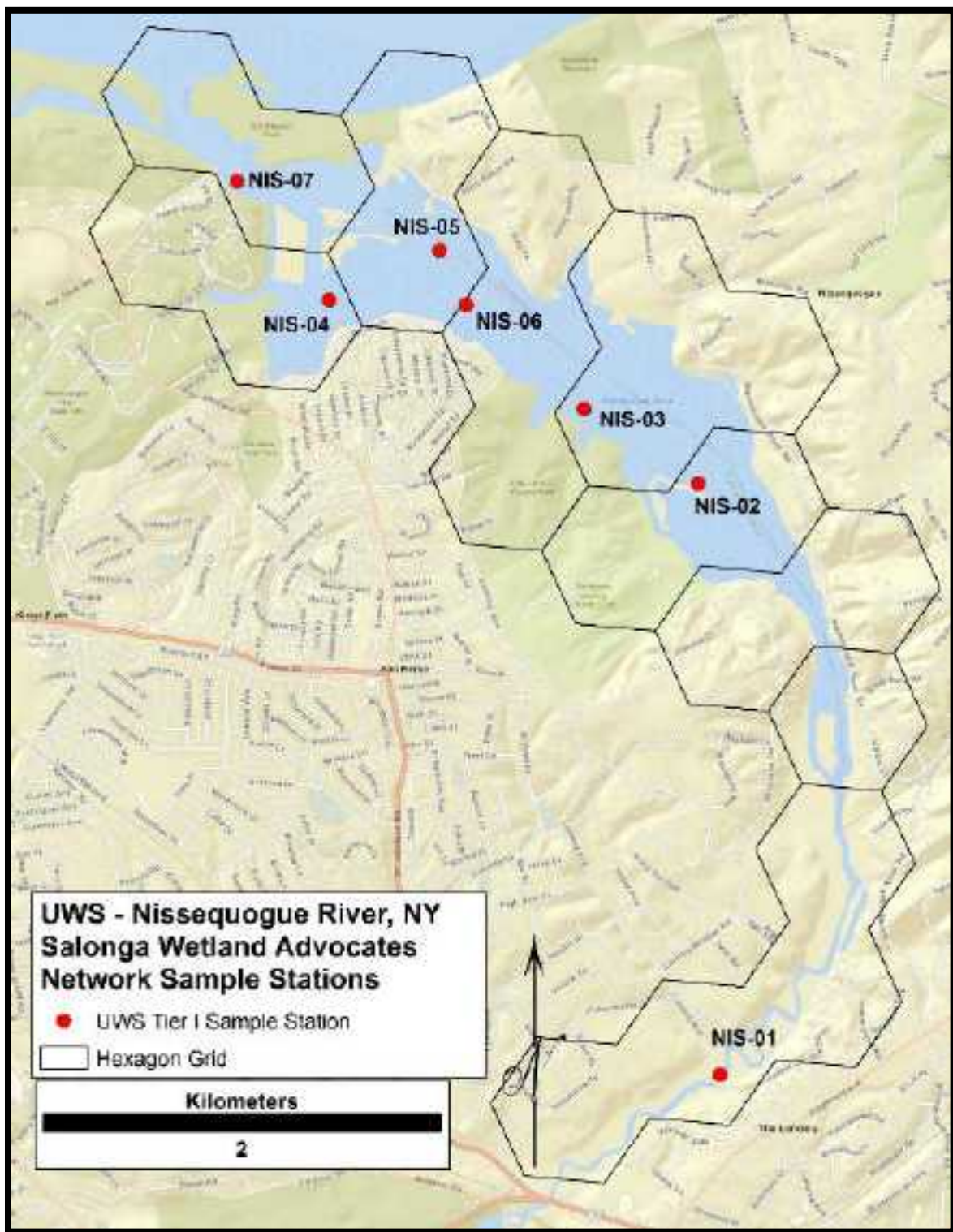


Figure 2I

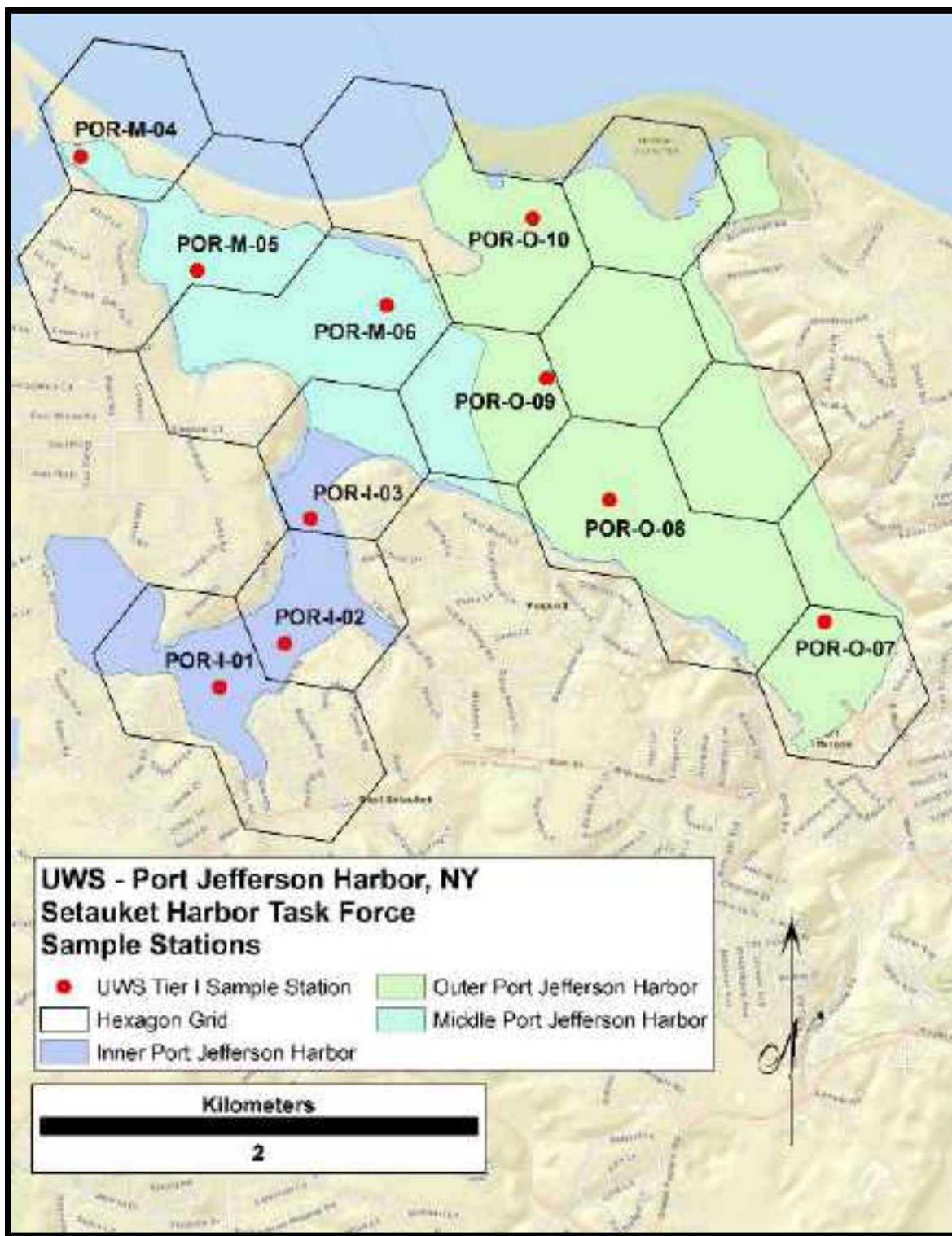


Figure 2m

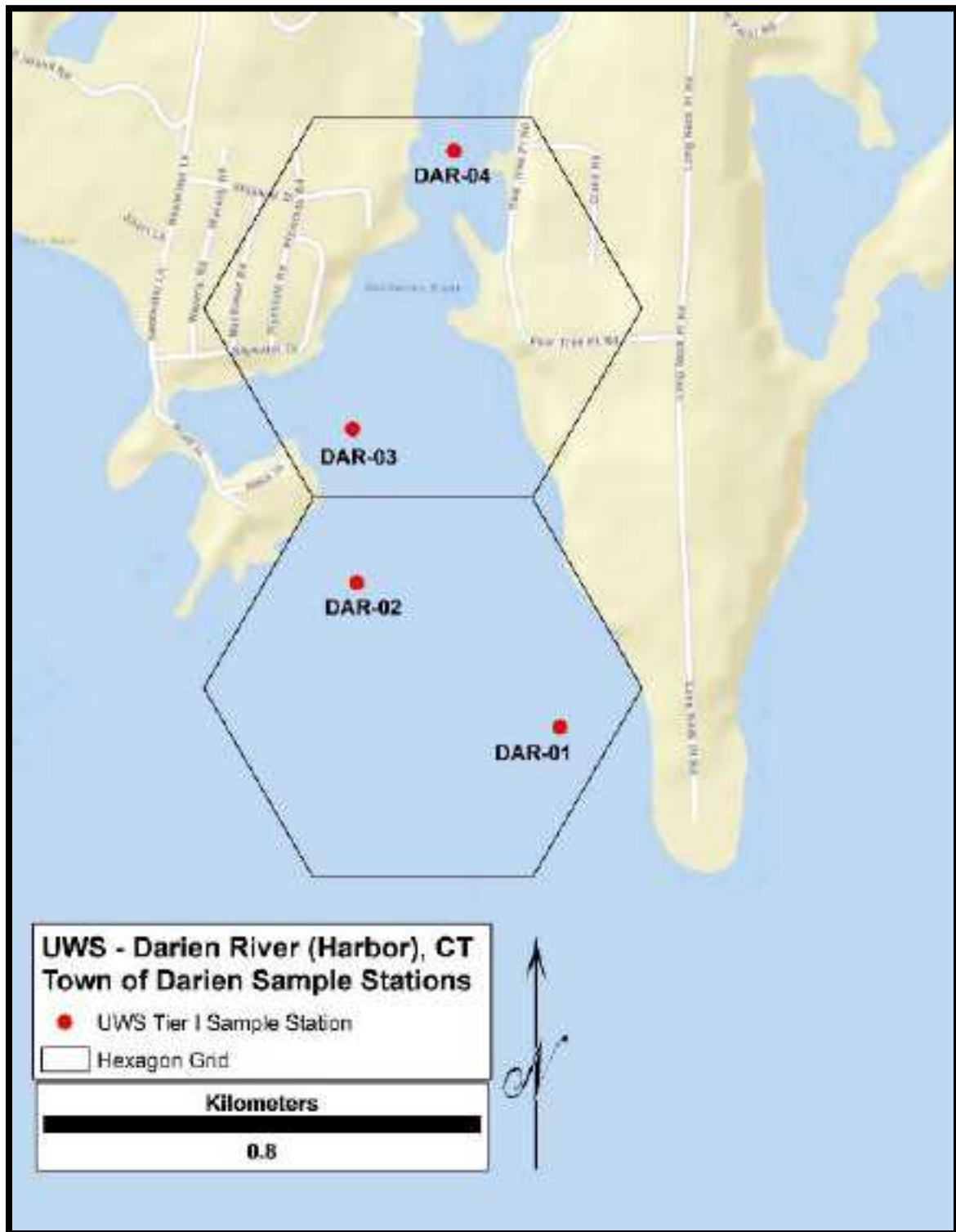


Figure 2n

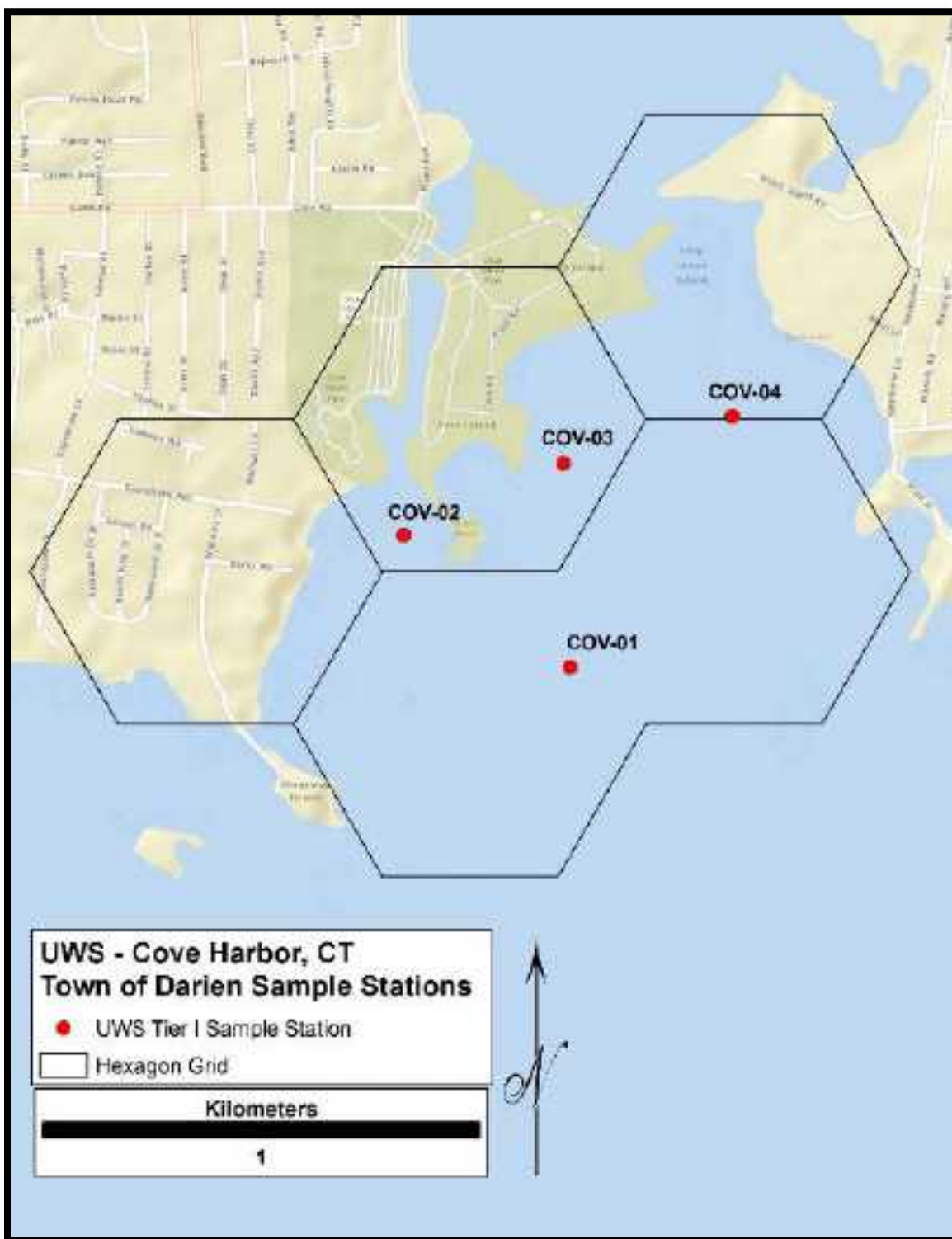


Figure 2o

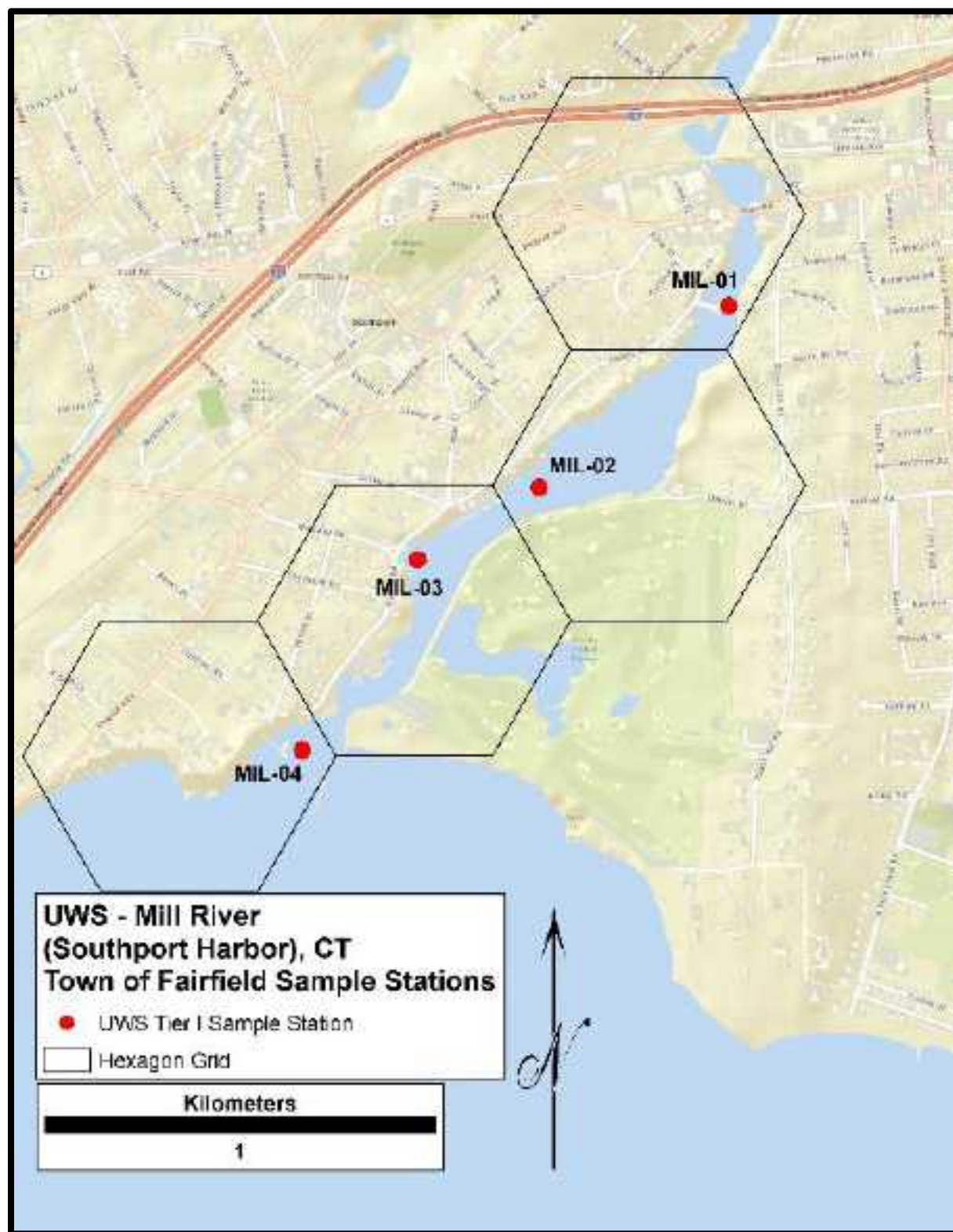


Figure 2p

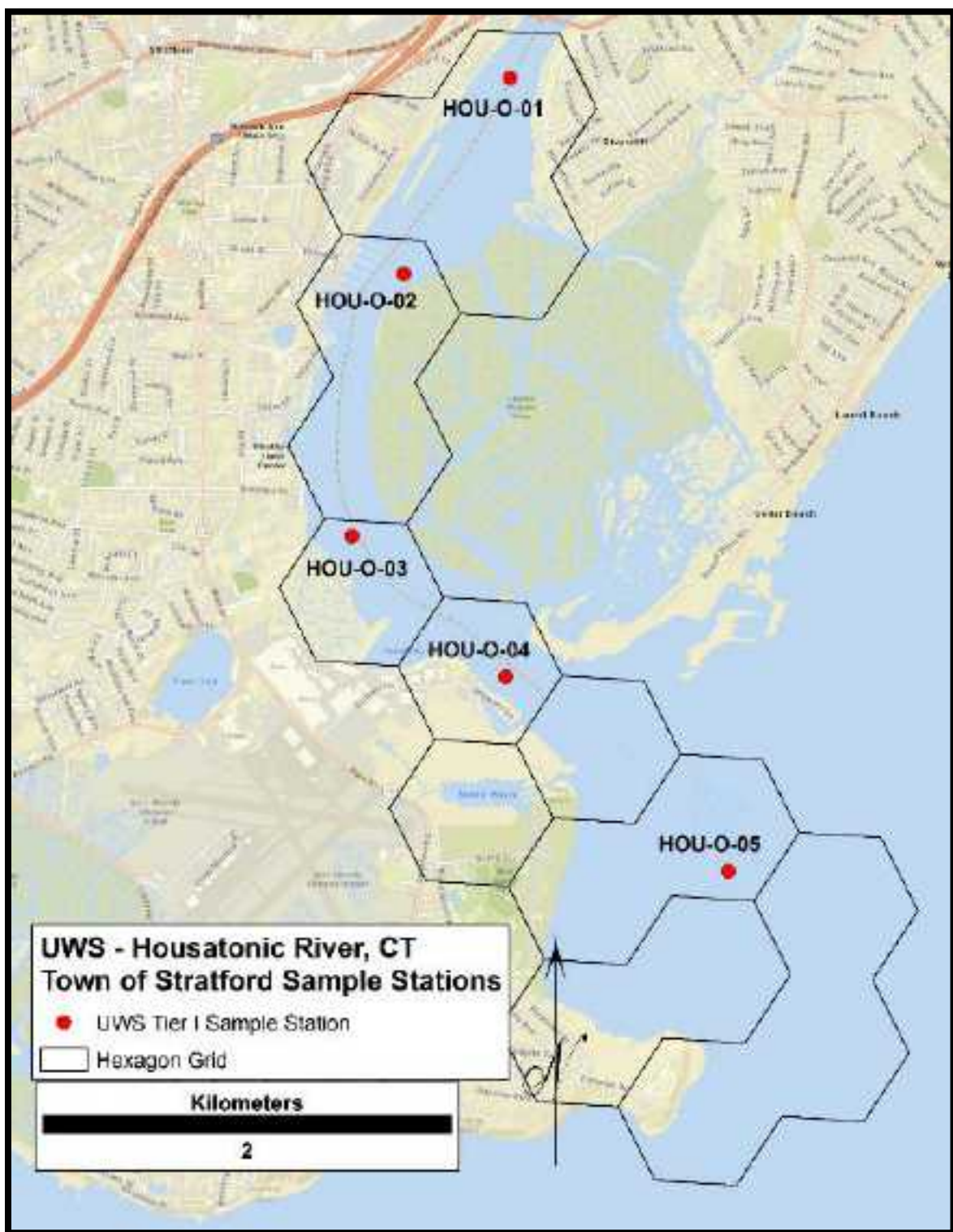


Figure 2q

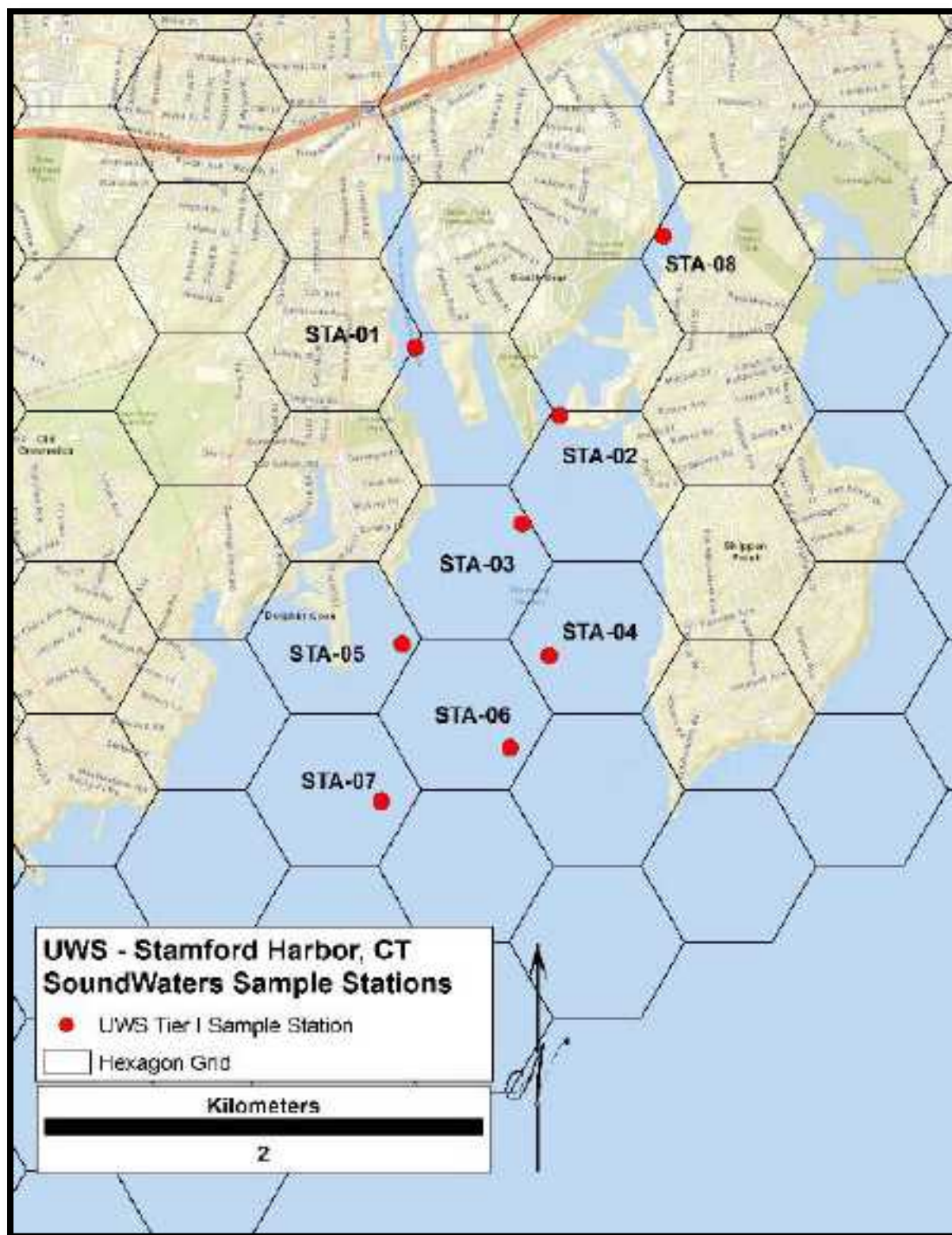


Figure 2r

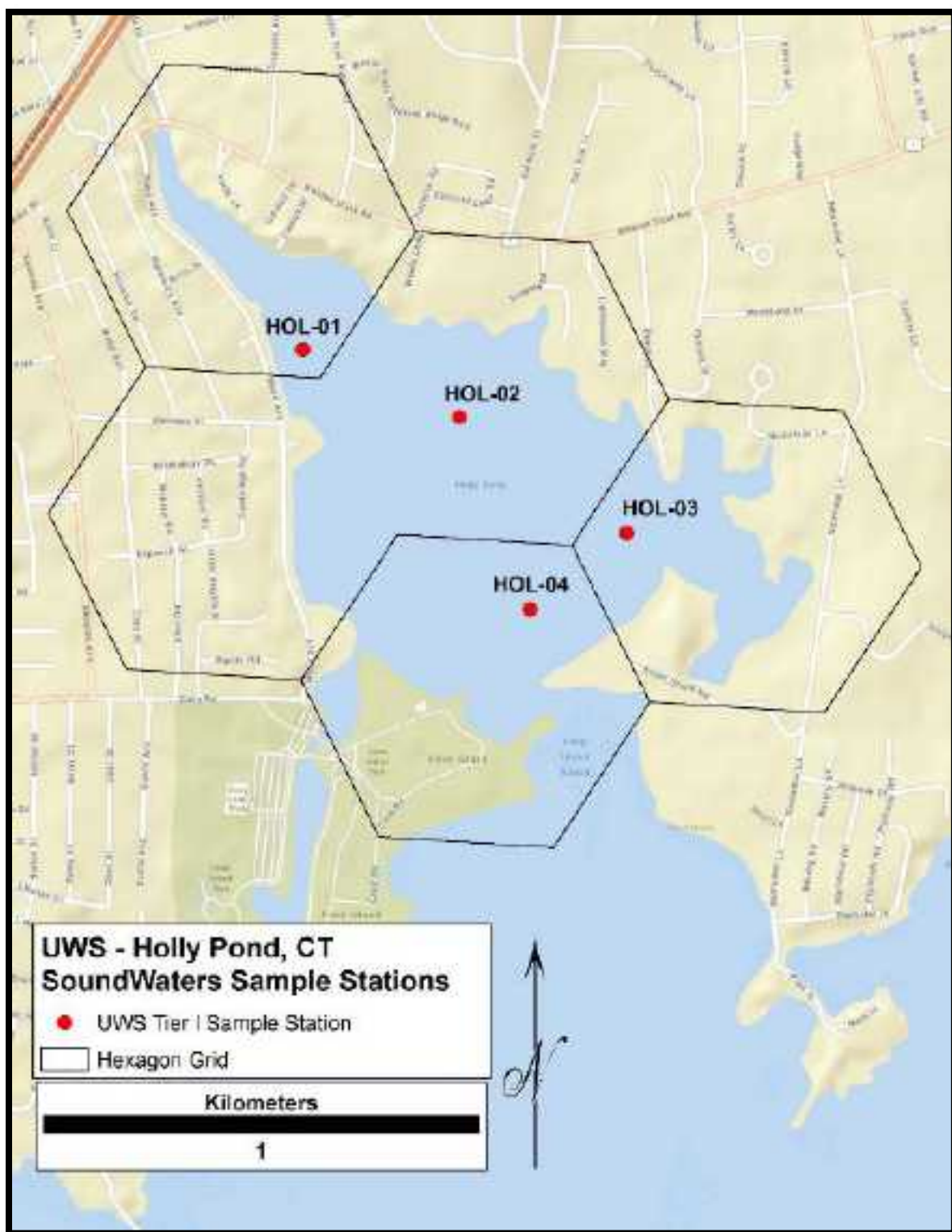


Figure 2s

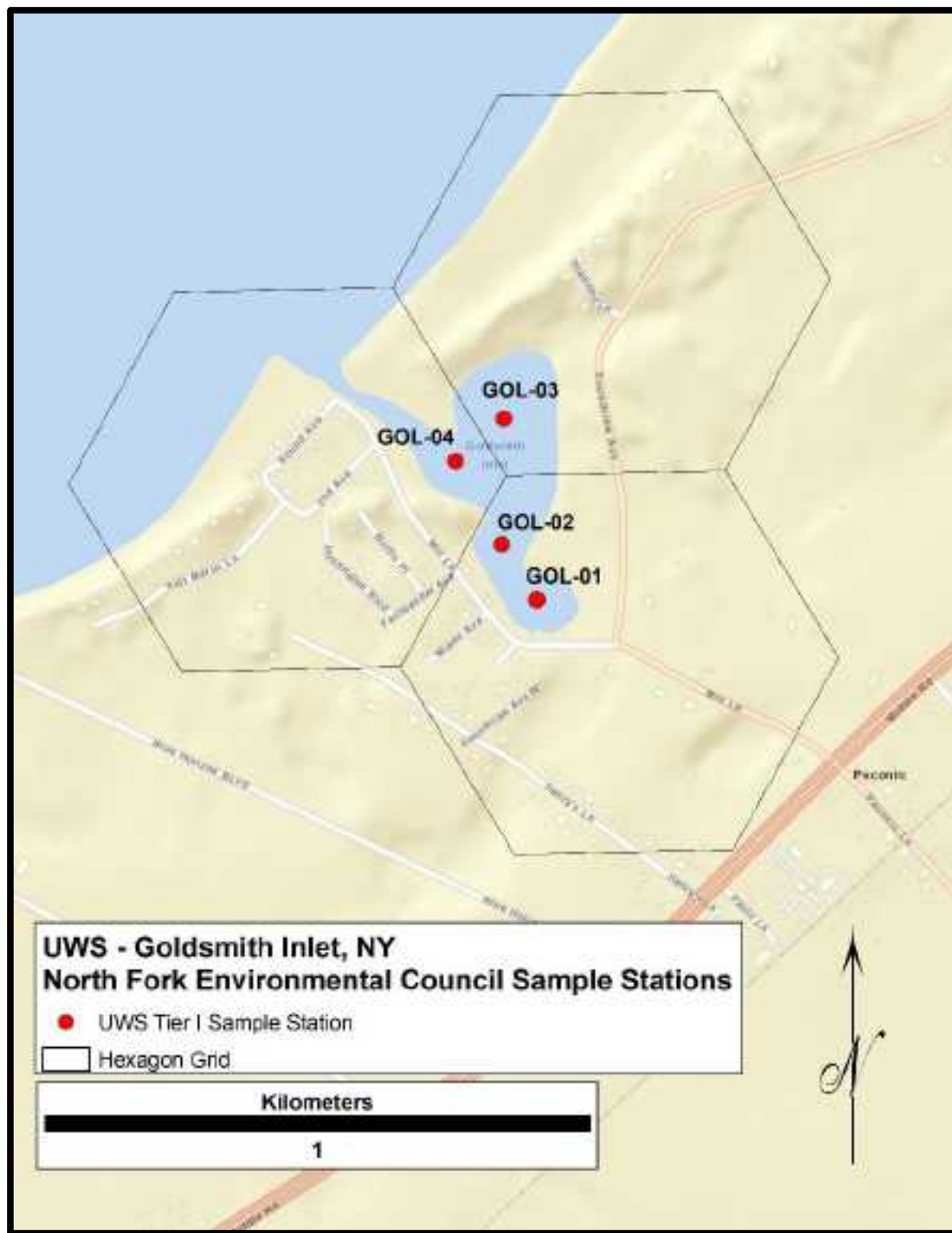


Figure 2t

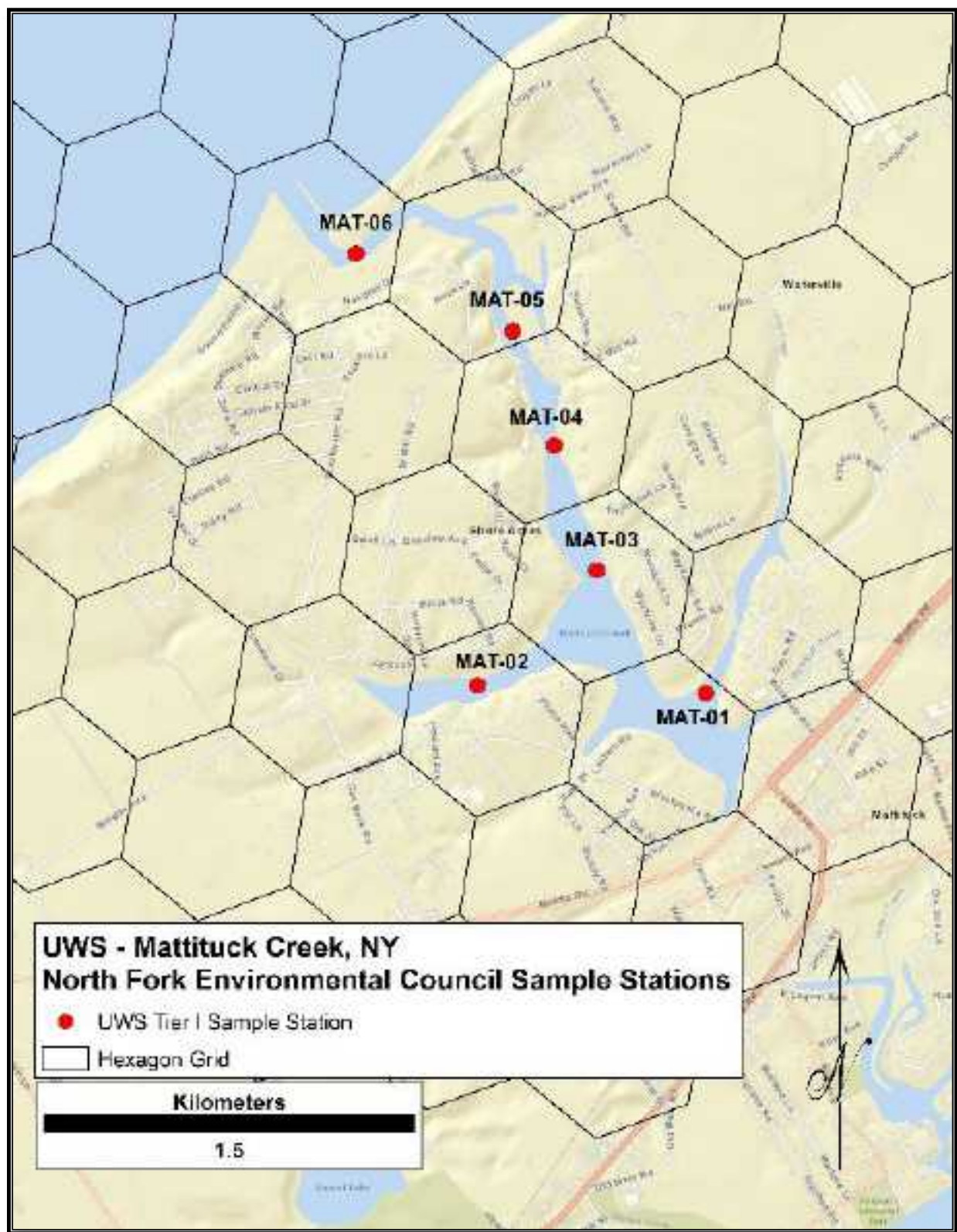


Figure 2u

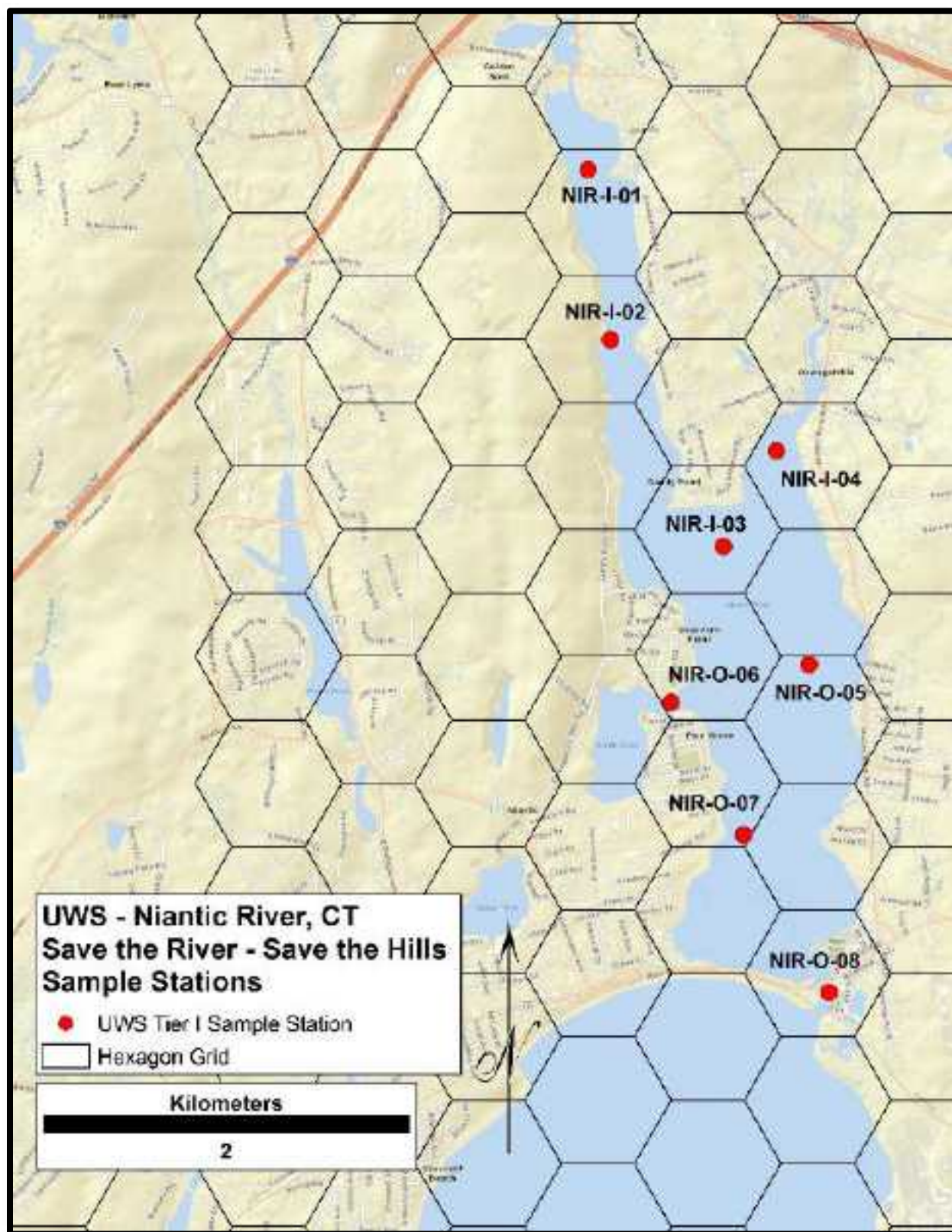


Figure 2v



Figure 2w

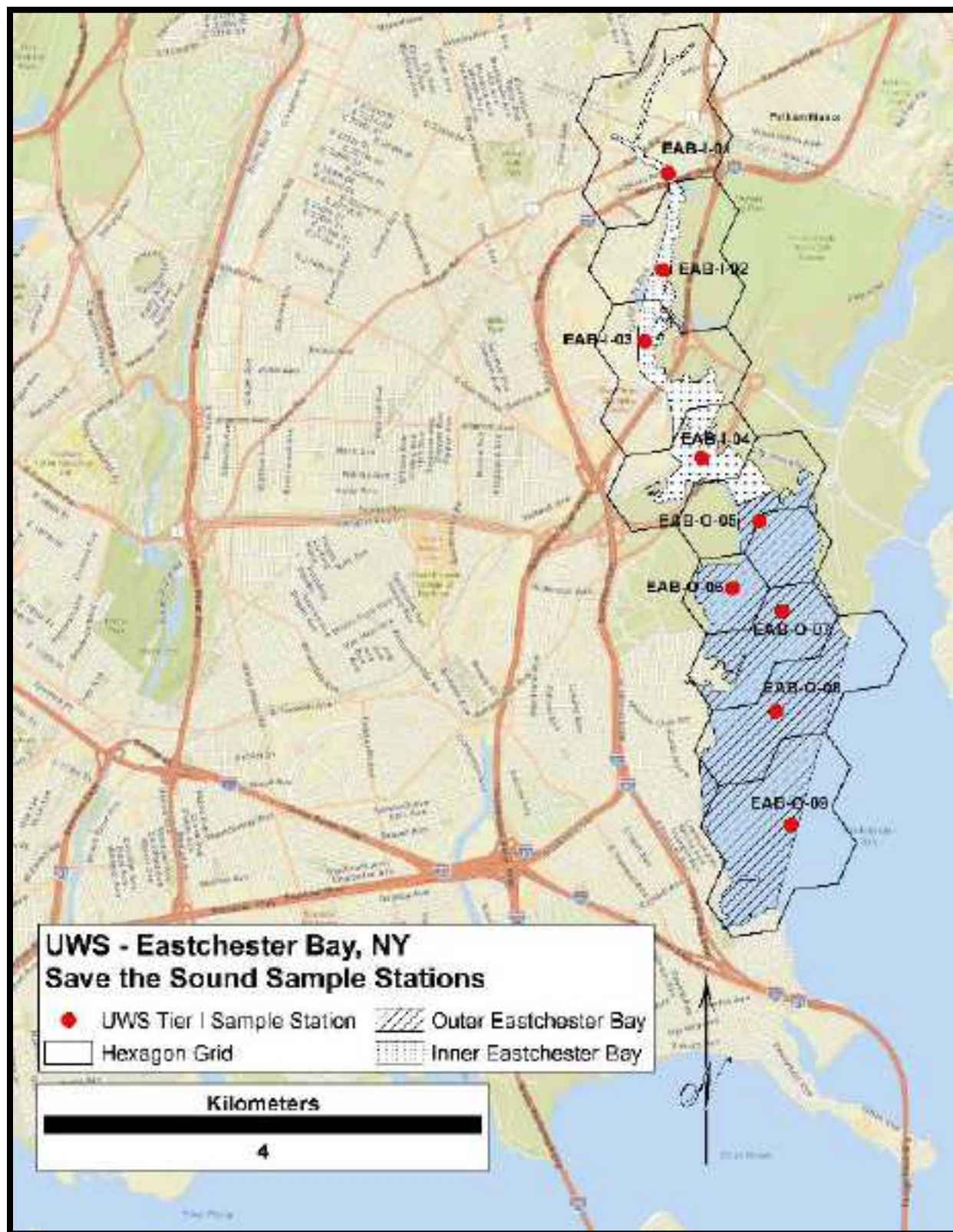


Figure 2x

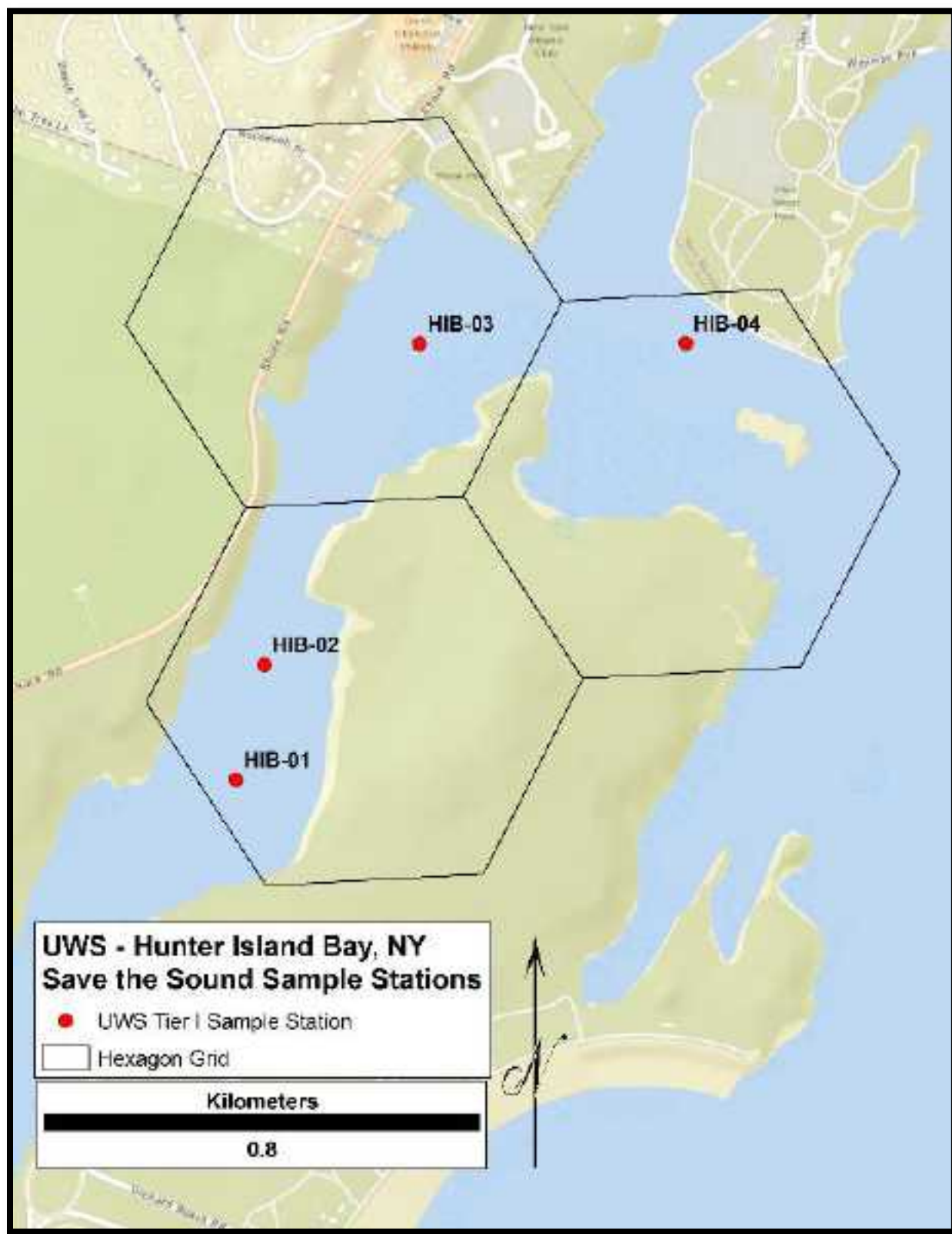


Figure 2y



Figure 2z

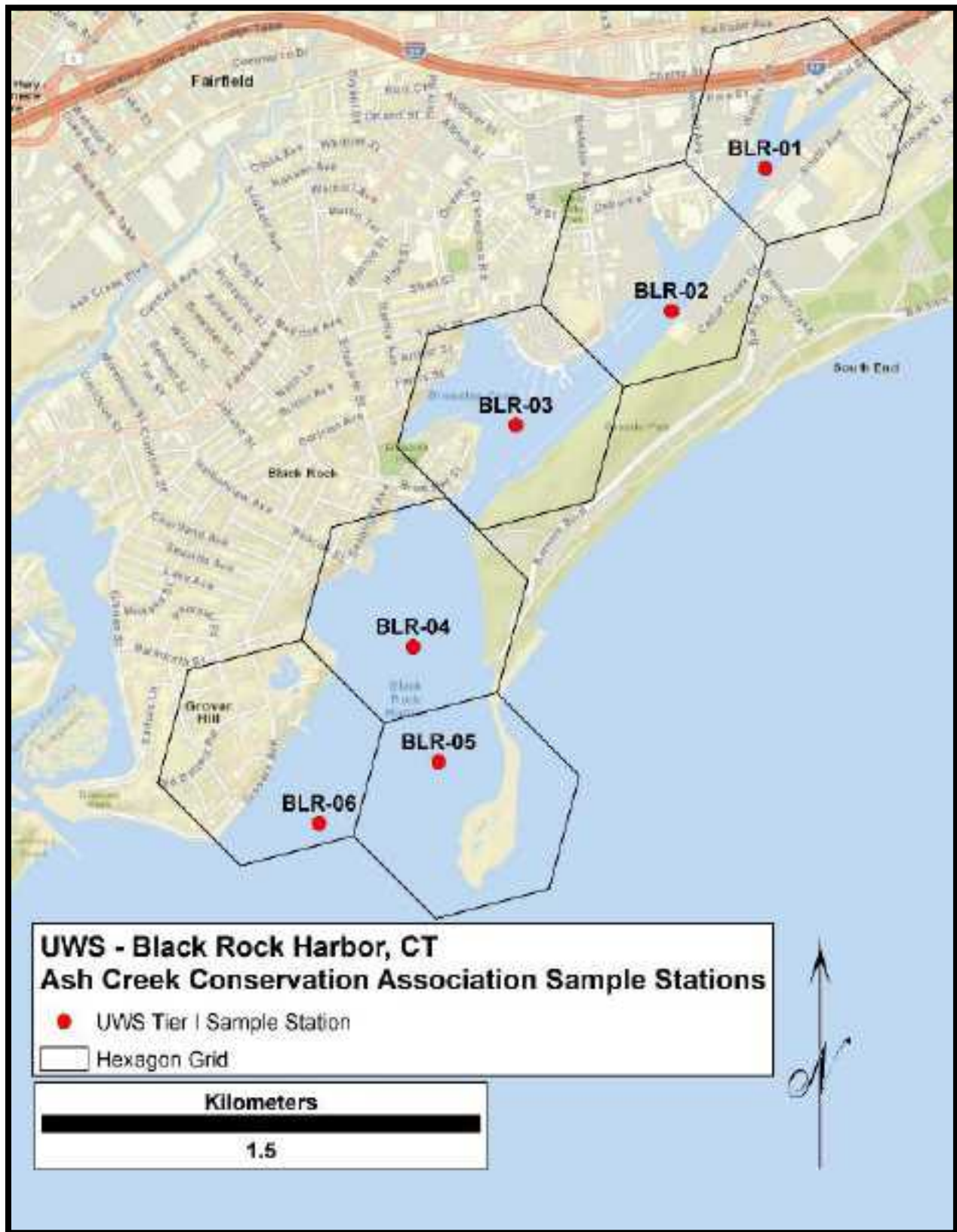


Figure 2aa



Figure 2ab

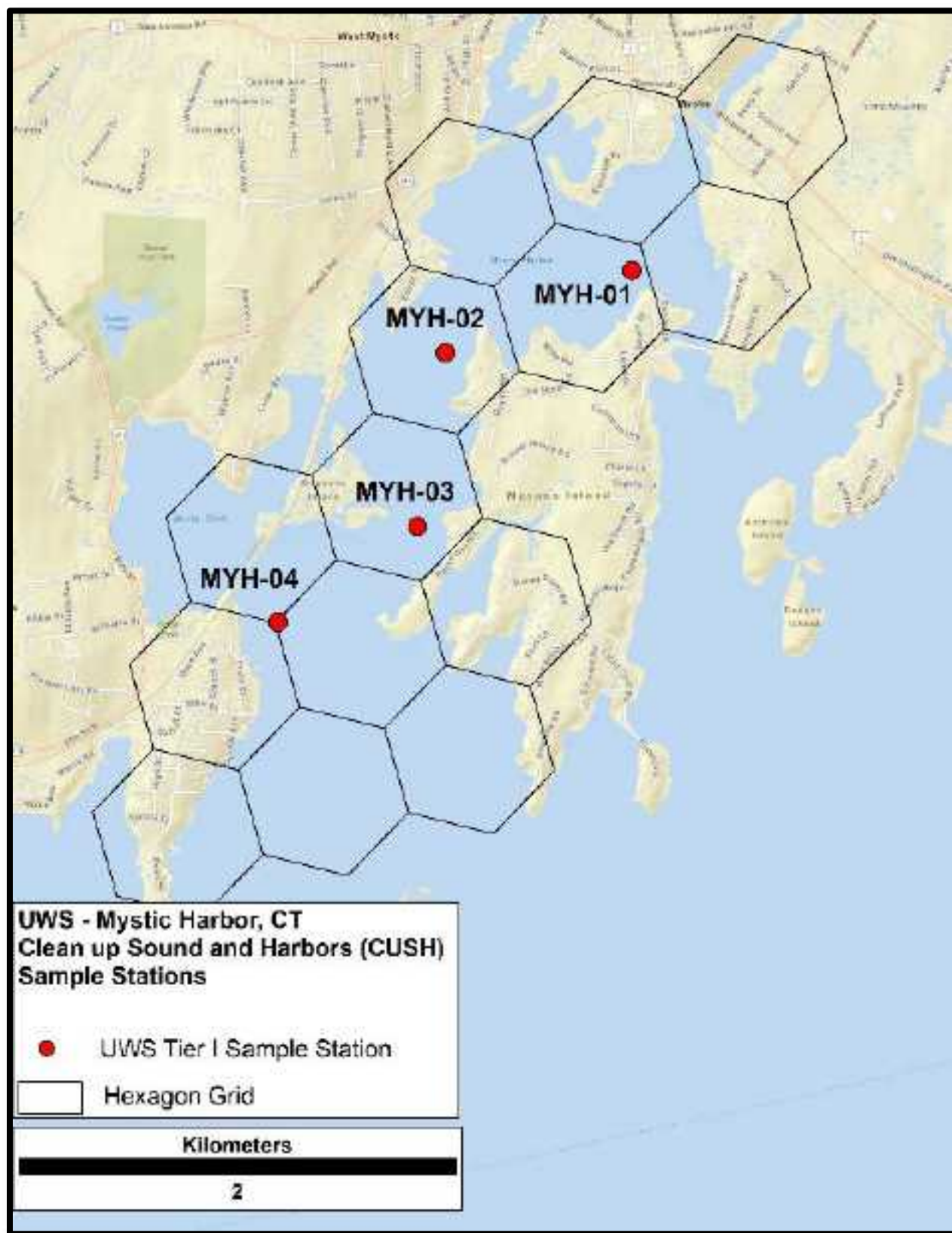


Figure 2ac

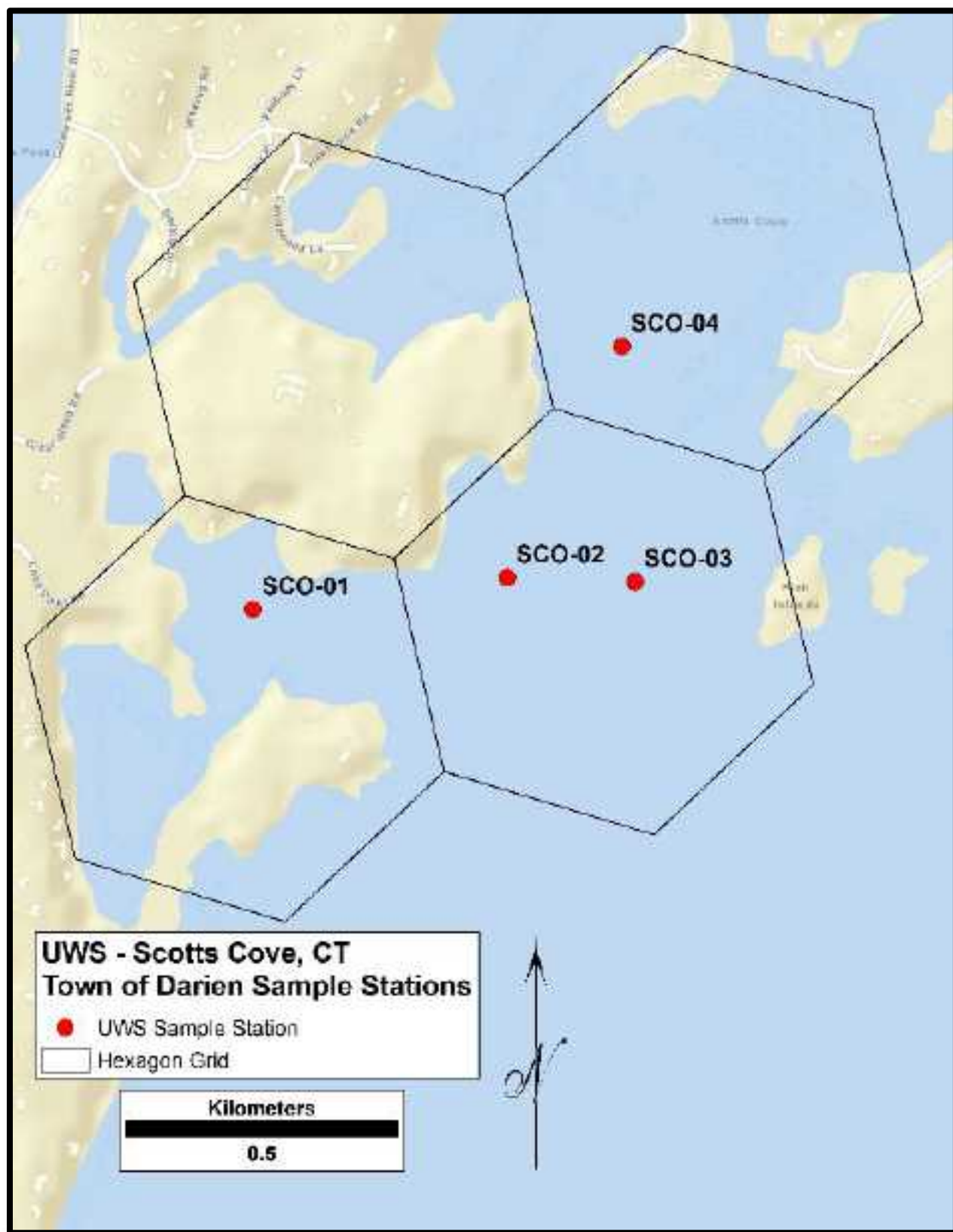


Figure 2ad

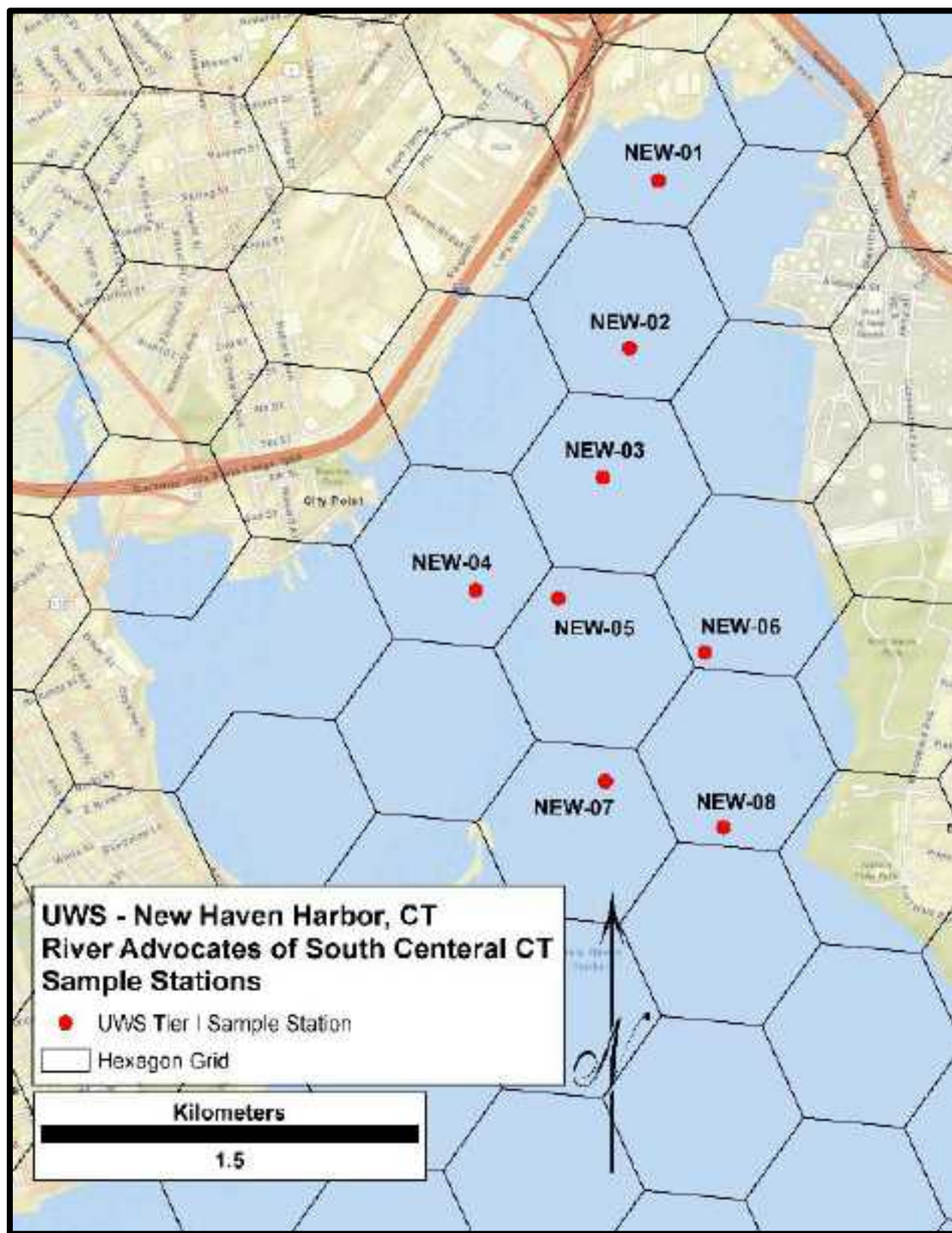


Figure 2ac

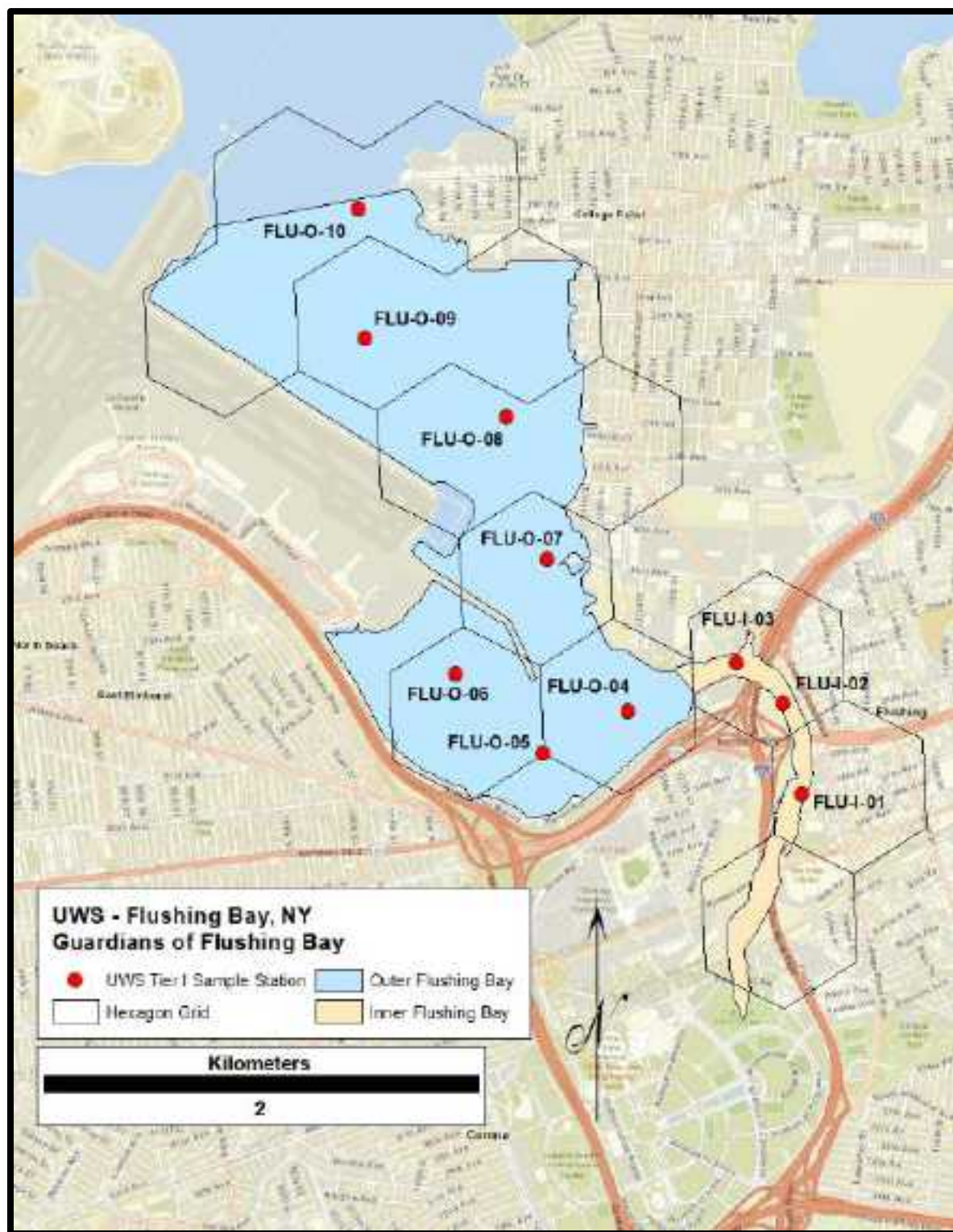


Figure 2ad

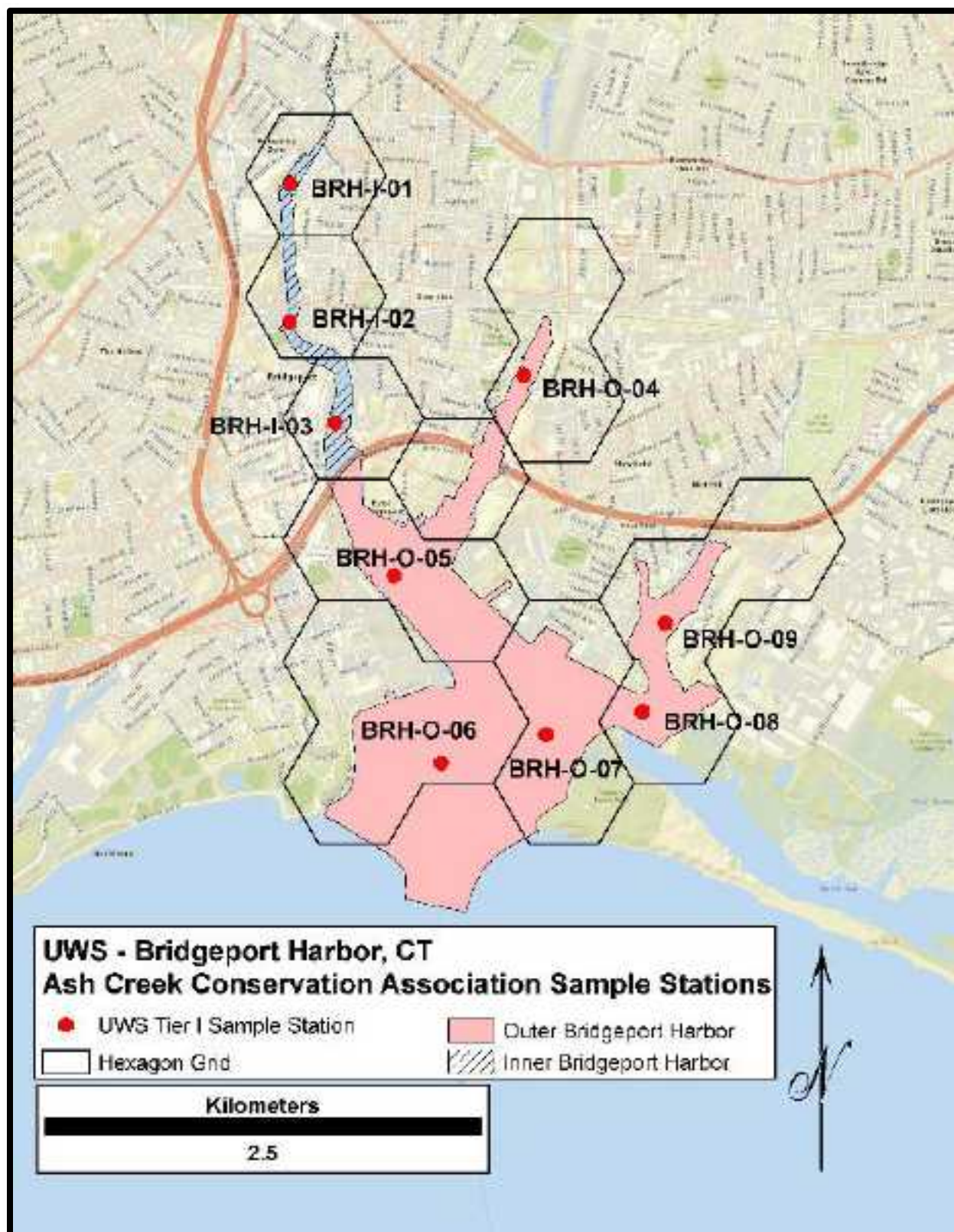


Figure 2ae

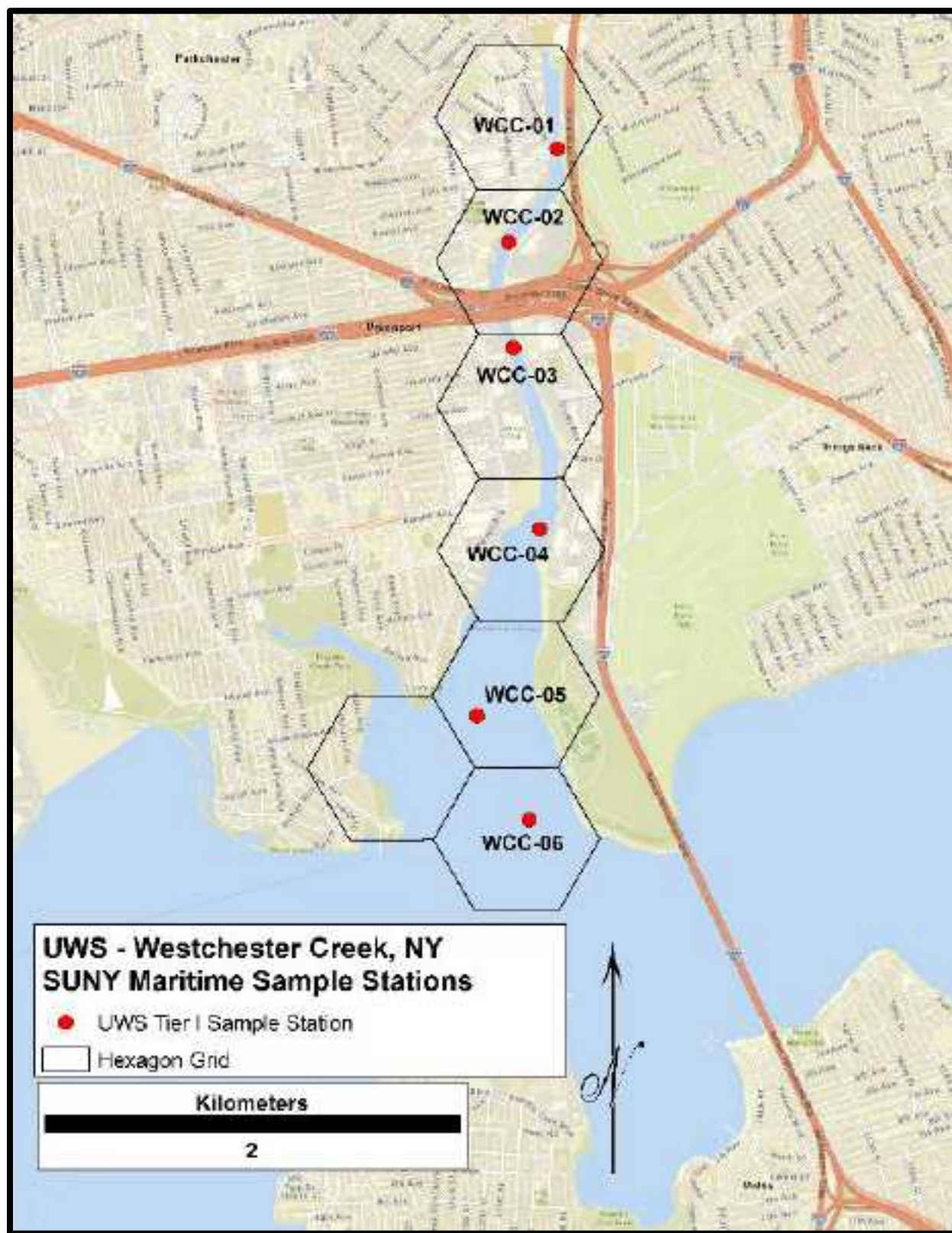


Figure 2af

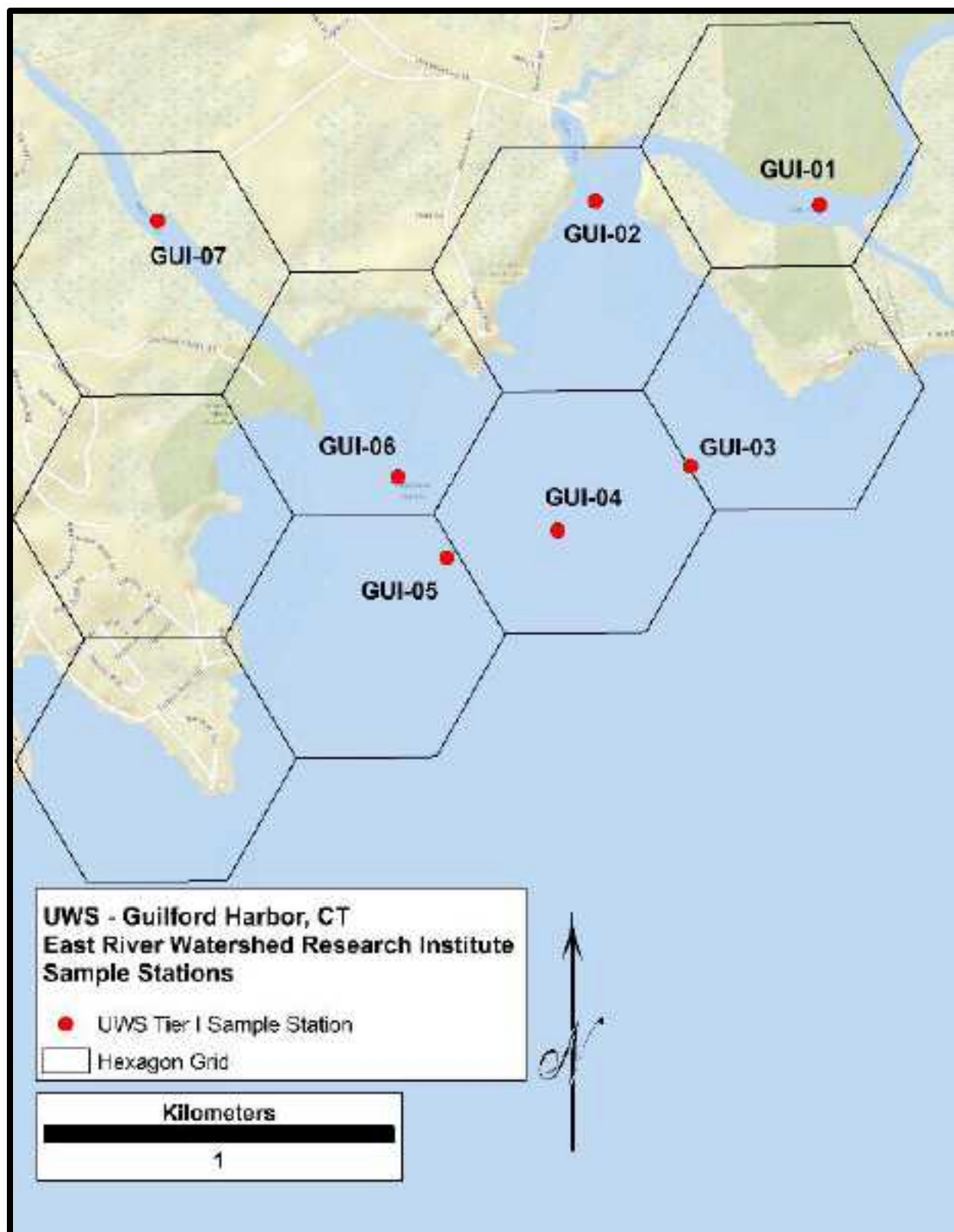


Figure 2ag

Table 2: Station coordinates in NAD_83 for Tier I water quality and Tier II nutrients stations* in the UWS.

Station ID	Embayment	Longitude	Latitude
ALE-01	Alewife Cove, CT	-72.10449	41.31814
ALE-02	Alewife Cove, CT	-72.10069	41.31364
ALE-03	Alewife Cove, CT	-72.10343	41.30898
ALE-04	Alewife Cove, CT	-72.10485	41.3055
BLR-01*	Black Rock Harbor, CT	-73.20513	41.16387
BLR-02	Black Rock Harbor, CT	-73.20895	41.15945
BLR-03*	Black Rock Harbor, CT	-73.21531	41.15589
BLR-04	Black Rock Harbor, CT	-73.21946	41.14901
BLR-05	Black Rock Harbor, CT	-73.21839	41.14545
BLR-06*	Black Rock Harbor, CT	-73.22326	41.14353
BRR-01	Bronx River, NY	-73.88403	40.82226
BRR-02	Bronx River, NY	-73.87916	40.81694
BRR-03	Bronx River, NY	-73.87346	40.81489
BRR-04	Bronx River, NY	-73.86939	40.81079
BRR-05	Bronx River, NY	-73.86751	40.80661
BRR-06	Bronx River, NY	-73.86299	40.80257
CEN-01 *	Centerport Harbor, NY	-73.37583	40.89694
CEN-02	Centerport Harbor, NY	-73.37952	40.90007
CEN-03 *	Centerport Harbor, NY	-73.38401	40.90849
COL-I-01	Cold Spring Harbor, NY	-73.46501	40.8625
COL-I-02	Cold Spring Harbor, NY	-73.46333	40.86667
COL-I-03	Cold Spring Harbor, NY	-73.46605	40.86898
COL-O-04	Cold Spring Harbor, NY	-73.47908	40.8796
COL-O-05	Cold Spring Harbor, NY	-73.48873	40.89025
COL-O-06	Cold Spring Harbor, NY	-73.48468	40.90344
COL-O-07	Cold Spring Harbor, NY	-73.50969	40.91512
CTR-01	Connecticut River, CT	-72.3842	41.352
CTR-02	Connecticut River, CT	-72.3839	41.34842

Station ID	Embayment	Longitude	Latitude
CTR-03	Connecticut River, CT	-72.37888	41.34416
CTR-04	Connecticut River, CT	-72.38082	41.35696
CTR-05	Connecticut River, CT	-72.37864	41.36275
CTR-06	Connecticut River, CT	-72.37932	41.35323
CTR-07	Connecticut River, CT	-72.37504	41.34672
CTR-08	Connecticut River, CT	-72.36533	41.34011
COV-01	Cove Harbor, CT	-73.49904	41.03958
COV-02	Cove Harbor, CT	-73.5036	41.04227
COV-03	Cove Harbor, CT	-73.49928	41.04377
COV-04	Cove Harbor, CT	-73.49468	41.04477
DAR-01	Darien River, CT	-73.4814	41.03846
DAR-02	Darien River, CT	-73.48587	41.04082
DAR-03	Darien River, CT	-73.48598	41.04336
DAR-04	Darien River, CT	-73.48382	41.04796
EAB-I-01 [*]	Eastchester Bay, NY	-73.8207	40.88621
EAB-I-02	Eastchester Bay, NY	-73.82118	40.87824
EAB-I-03	Eastchester Bay, NY	-73.82306	40.8724
EAB-I-04 [*]	Eastchester Bay, NY	-73.81672	40.8628
EAB-O-05 [*]	Eastchester Bay, NY	-73.81038	40.85766
EAB-O-06	Eastchester Bay, NY	-73.81319	40.85211
EAB-O-07	Eastchester Bay, NY	-73.80781	40.85024
EAB-O-08	Eastchester Bay, NY	-73.80829	40.84192
EAB-O-09 [*]	Eastchester Bay, NY	-73.80649	40.83259
FAR-04	Farm River, CT	-72.85192	41.26209
FAR-05	Farm River, CT	-72.85405	41.25649
FAR-06	Farm River, CT	-72.85857	41.24893
FAR-07	Farm River, CT	-72.85378	41.25186
GOL-01	Goldsmith Inlet, NY	-72.46946	41.05073
GOL-02	Goldsmith Inlet, NY	-72.47022	41.05165

Station ID	Embayment	Longitude	Latitude
GOL-03	Goldsmith Inlet, NY	-72.47017	41.05373
GOL-04	Goldsmith Inlet, NY	-72.47123	41.05303
HEM-M-01	Hempstead Harbor, NY	-73.65353	40.83189
HEM-M-02	Hempstead Harbor, NY	-73.65854	40.84172
HEM-M-03	Hempstead Harbor, NY	-73.65216	40.85365
HEM-O-04	Hempstead Harbor, NY	-73.67396	40.86077
HEM-O-05	Hempstead Harbor, NY	-73.67493	40.87349
HEM-O-06	Hempstead Harbor, NY	-73.65016	40.88365
HOL-01	Holly Pond, CT	-73.50337	41.05624
HOL-02	Holly Pond, CT	-73.49906	41.05487
HOL-03	Holly Pond, CT	-73.49446	41.0525
HOL-04	Holly Pond, CT	-73.4971	41.05092
HOU-O-01	Housatonic River, CT	-73.11245	41.1976
HOU-O-02	Housatonic River, CT	-73.11861	41.18895
HOU-O-03	Housatonic River, CT	-73.12158	41.17737
HOU-O-04	Housatonic River, CT	-73.11256	41.17121
HOU-O-05	Housatonic River, CT	-73.09952	41.16267
HIB-01	Hunter Island Bay, NY	-73.79606	40.87446
HIB-02	Hunter Island Bay, NY	-73.79547	40.87637
HIB-03	Hunter Island Bay, NY	-73.79217	40.88172
HIB-04	Hunter Island Bay, NY	-73.78636	40.88178
HUB-01 *	Huntington Bay, NY	-73.42993	40.90936
HUB-02 *	Huntington Bay, NY	-73.40746	40.91044
HUB-03	Huntington Bay, NY	-73.41805	40.91777
HUH-01 *	Huntington Harbor, NY	-73.41805	40.88749
HUH-02	Huntington Harbor, NY	-73.42333	40.89666
HUH-03	Huntington Harbor, NY	-73.43205	40.89881
HUH-04	Huntington Harbor, NY	-73.43865	40.89988
HUH-06 *	Huntington Harbor, NY	-73.43445	40.90499

Station ID	Embayment	Longitude	Latitude
LNE-I-01 *	Little Neck Bay, NY	-73.75791	40.77224
LNE-I-02	Little Neck Bay, NY	-73.7608	40.7778
LNE-I-03	Little Neck Bay, NY	-73.75823	40.78314
LNE-I-04 *	Little Neck Bay, NY	-73.75061	40.78377
LNE-I-05 *	Little Neck Bay, NY	-73.76862	40.78606
LNE-O-06 *	Little Neck Bay, NY	-73.7582	40.7888
LNE-O-07	Little Neck Bay, NY	-73.77112	40.794
LNE-O-08	Little Neck Bay, NY	-73.76179	40.79561
LNE-O-09	Little Neck Bay, NY	-73.75442	40.79884
LNE-O-10 *	Little Neck Bay, NY	-73.76992	40.80202
LLO-01 *	Lloyd Harbor, NY	-73.46734	40.91296
LLO-02	Lloyd Harbor, NY	-73.45	40.91361
LLO-03	Lloyd Harbor, NY	-73.44147	40.91093
LLO-04 *	Lloyd Harbor, NY	-73.43738	40.91889
MAM-01 *	Mamaroneck River, NY	-73.72225	40.94088
MAM-02	Mamaroneck River, NY	-73.72717	40.94288
MAM-03 *	Mamaroneck River, NY	-73.72894	40.94737
MAM-04 *	Mamaroneck River, NY	-73.73625	40.94367
MAN-I-01	Manhasset Bay, NY	-73.71316	40.80772
MAN-I-02	Manhasset Bay, NY	-73.71461	40.81244
MAN-I-03	Manhasset Bay, NY	-73.70714	40.81586
MAN-M-04	Manhasset Bay, NY	-73.71242	40.82271
MAN-M-05	Manhasset Bay, NY	-73.70551	40.83064
MAN-M-06	Manhasset Bay, NY	-73.71454	40.83228
MAN-M-07	Manhasset Bay, NY	-73.72375	40.82616
MAN-M-08	Manhasset Bay, NY	-73.72564	40.83644
MAN-O-09	Manhasset Bay, NY	-73.73613	40.83179
MAN-O-10	Manhasset Bay, NY	-73.73672	40.84517
MAN-O-11	Manhasset Bay, NY	-73.74556	40.84097

Station ID	Embayment	Longitude	Latitude
MAT-01	Mattituck Creek, NY	-72.53983	40.99671
MAT-02	Mattituck Creek, NY	-72.55082	40.99702
MAT-03	Mattituck Creek, NY	-72.54506	41.00124
MAT-04	Mattituck Creek, NY	-72.5471	41.0058
MAT-05	Mattituck Creek, NY	-72.5491	41.00997
MAT-06	Mattituck Creek, NY	-72.55664	41.01282
MNC-01	Mill Neck Creek, NY	-73.5675	40.89888
MNC-02	Mill Neck Creek, NY	-73.55809	40.90138
MNC-03	Mill Neck Creek, NY	-73.55167	40.90333
MIL-01	Mill River, CT	-73.27468	41.13761
MIL-02	Mill River, CT	-73.28045	41.13339
MIL-03	Mill River, CT	-73.28416	41.1317
MIL-04	Mill River, CT	-73.28766	41.12727
MYH-01	Mystic Harbor, CT	-71.96392	41.34344
MYH-02	Mystic Harbor, CT	-71.97418	41.34013
MYH-03	Mystic Harbor, CT	-71.97581	41.33295
MYH-04	Mystic Harbor, CT	-71.98351	41.32905
NRH-01	New Rochelle Harbor, NY	-73.77759	40.89548
NRH-02	New Rochelle Harbor, NY	-73.78096	40.89031
NRH-03	New Rochelle Harbor, NY	-73.78444	40.88806
NRH-04	New Rochelle Harbor, NY	-73.7881	40.88382
NIR-I-01	Niantic River, CT	-72.19166	41.36423
NIR-I-02	Niantic River, CT	-72.19027	41.35582
NIR-I-03	Niantic River, CT	-72.18295	41.34556
NIR-I-04	Niantic River, CT	-72.17941	41.35027
NIR-O-05	Niantic River, CT	-72.17737	41.3397
NIR-O-06	Niantic River, CT	-72.18646	41.33786
NIR-O-07	Niantic River, CT	-72.18174	41.33128
NIR-O-08	Niantic River, CT	-72.1762	41.32346

Station ID	Embayment	Longitude	Latitude
NIS-01	Nissequogue River, NY	-73.20069	40.86397
NIS-02	Nissequogue River, NY	-73.20219	40.89071
NIS-03	Nissequogue River, NY	-73.20899	40.89408
NIS-04	Nissequogue River, NY	-73.22424	40.89892
NIS-05	Nissequogue River, NY	-73.21767	40.90121
NIS-06	Nissequogue River, NY	-73.21607	40.89874
NIS-07	Nissequogue River, NY	-73.22976	40.90427
NPB-01	Northport Bay, NY	-73.36417	40.91111
NPB-02 *	Northport Bay, NY	-73.35544	40.92265
NPB-03	Northport Bay, NY	-73.36616	40.92906
NPB-04	Northport Bay, NY	-73.37555	40.91666
NPB-05 *	Northport Bay, NY	-73.38112	40.93054
NPB-06 *	Northport Bay, NY	-73.39183	40.91458
NPB-07	Northport Bay, NY	-73.39841	40.92496
NPH-01 *	Northport Harbor, NY	-73.36131	40.89117
NPH-02	Northport Harbor, NY	-73.35583	40.89888
NPH-03 *	Northport Harbor, NY	-73.35972	40.90561
NWH-I-01	Norwalk Harbor, CT	-73.41105	41.11738
NWH-I-02 *	Norwalk Harbor, CT	-73.41117	41.10799
NWH-I-03 *	Norwalk Harbor, CT	-73.416	41.10205
NWH-I-04	Norwalk Harbor, CT	-73.41419	41.09846
NWH-I-05 *	Norwalk Harbor, CT	-73.41003	41.09385
NWH-I-06 *	Norwalk Harbor, CT	-73.40425	41.08727
NWH-I-07	Norwalk Harbor, CT	-73.40073	41.07939
NWH-O-01	Norwalk Harbor, CT	-73.41195	41.06843
NWH-O-02 *	Norwalk Harbor, CT	-73.41526	41.06435
NWH-O-03	Norwalk Harbor, CT	-73.40758	41.06275
NWH-O-04	Norwalk Harbor, CT	-73.39851	41.06764
NWH-O-05 *	Norwalk Harbor, CT	-73.39131	41.07406

Station ID	Embayment	Longitude	Latitude
OYB-01	Oyster Bay, NY	-73.53963	40.89789
OYB-02	Oyster Bay, NY	-73.52878	40.91181
OYB-03	Oyster Bay, NY	-73.53113	40.88073
OYB-04	Oyster Bay, NY	-73.51553	40.89036
POR-I-01	Port Jefferson Harbor, NY	-73.10422	40.94861
POR-I-02	Port Jefferson Harbor, NY	-73.10069	40.9504
POR-I-03	Port Jefferson Harbor, NY	-73.09931	40.95557
POR-M-04	Port Jefferson Harbor, NY	-73.11192	40.97045
POR-M-05	Port Jefferson Harbor, NY	-73.10555	40.96579
POR-M-06	Port Jefferson Harbor, NY	-73.09524	40.9644
POR-O-07	Port Jefferson Harbor, NY	-73.07133	40.95141
POR-O-08	Port Jefferson Harbor, NY	-73.08307	40.95637
POR-O-09	Port Jefferson Harbor, NY	-73.08649	40.96139
POR-O-10	Port Jefferson Harbor, NY	-73.08729	40.968
STA-01	Stamford Harbor, CT	-73.54388	41.0363
STA-02	Stamford Harbor, CT	-73.53599	41.03353
STA-03	Stamford Harbor, CT	-73.53796	41.02906
STA-04	Stamford Harbor, CT	-73.53645	41.02362
STA-05	Stamford Harbor, CT	-73.54446	41.02405
STA-06	Stamford Harbor, CT	-73.53853	41.01981
STA-07	Stamford Harbor, CT	-73.54553	41.01754
STA-08	Stamford Harbor, CT	-73.53044	41.041
STO-I-01	Stonington Harbor, CT	-71.91511	41.34485
STO-I-02	Stonington Harbor, CT	-71.9144	41.34298
STO-I-03	Stonington Harbor, CT	-71.91138	41.3422
STO-O-04	Stonington Harbor, CT	-71.9106	41.33843
STO-O-05	Stonington Harbor, CT	-71.91545	41.33408
STO-O-06	Stonington Harbor, CT	-71.90871	41.33237
STO-O-07	Stonington Harbor, CT	-71.90762	41.32727

Station ID	Embayment	Longitude	Latitude
STO-O-08	Stonington Harbor, CT	-71.91973	41.32764
SCO-01	Scotts Cove, CT	-73.47318	41.04985
SCO-02	Scotts Cove, CT	-73.46762	41.05041
SCO-03	Scotts Cove, CT	-73.46483	41.05037
SCO-04	Scotts Cove, CT	-73.46516	41.05425
NEW-01	New Haven Harbor, CT	-72.91249	41.29462
NEW-02	New Haven Harbor, CT	-72.91387	41.28861
NEW-03	New Haven Harbor, CT	-72.91511	41.28396
NEW-04	New Haven Harbor, CT	-72.92178	41.28159
NEW-05	New Haven Harbor, CT	-72.91722	41.27964
NEW-06	New Haven Harbor, CT	-72.91025	41.27771
NEW-07	New Haven Harbor, CT	-72.91496	41.27309
NEW-08	New Haven Harbor, CT	-72.90936	41.27144
FLU-I-01	Flushing Bay, NY	-73.83682	40.76045
FLU-I-02	Flushing Bay, NY	-73.83789	40.76416
FLU-I-03	Flushing Bay, NY	-73.84044	40.76584
FLU-O-04	Flushing Bay, NY	-73.8463	40.76379
FLU-O-05	Flushing Bay, NY	-73.85089	40.76199
FLU-O-06	Flushing Bay, NY	-73.85573	40.76525
FLU-O-07	Flushing Bay, NY	-73.85081	40.76998
FLU-O-08	Flushing Bay, NY	-73.85314	40.77587
FLU-O-09	Flushing Bay, NY	-73.8609	40.77904
FLU-O-10	Flushing Bay, NY	-73.86131	40.78441
BRH-I-01	Bridgeport Harbor, CT	-73.18912	41.19232
BRH-I-02	Bridgeport Harbor, CT	-73.18911	41.18517
BRH-I-03	Bridgeport Harbor, CT	-73.18597	41.17997
BRH-O-04	Bridgeport Harbor, CT	-73.17312	41.1825
BRH-O-05	Bridgeport Harbor, CT	-73.1819	41.1721
BRH-O-06	Bridgeport Harbor, CT	-73.17863	41.16247

Station ID	Embayment	Longitude	Latitude
BRH-O-07	Bridgeport Harbor, CT	-73.17147	41.16398
BRH-O-08	Bridgeport Harbor, CT	-73.16489	41.16516
BRH-O-09	Bridgeport Harbor, CT	-73.16339	41.16975
WCC-01	Westchester Creek, NY	-73.83933	40.83521
WCC-02	Westchester Creek, NY	-73.84203	40.8311
WCC-03	Westchester Creek, NY	-73.84169	40.82654
WCC-04	Westchester Creek, NY	-73.84005	40.81868
WCC-05	Westchester Creek, NY	-73.8435	40.81054
WCC-06	Westchester Creek, NY	-73.84045	40.80605
GUI-01	Guilford Harbor, CT	-72.6573	41.27003
GUI-02	Guilford Harbor, CT	-72.66496	41.27013
GUI-03	Guilford Harbor, CT	-72.66174	41.26328
GUI-04	Guilford Harbor, CT	-72.66627	41.26163
GUI-05	Guilford Harbor, CT	-72.67005	41.26093
GUI-06	Guilford Harbor, CT	-72.67172	41.26301
GUI-07	Guilford Harbor, CT	-72.67993	41.26964

*These stations will be sampled for Tier II nutrients.

A.6.c. Annual Task Calendar

The annual task calendar describes when certain activities will occur.

Table 3: Annual Task Calendar

These tasks are repeated annually.

Activity	J	F	M	A	M	J	J	A	S	O	N	D
Kickoff meeting with UWS project team	x	x										
Develop draft QAPP and submit to UWS & EPA	x	x	x									
Finalize QAPP, responding to comments from EPA			x	x	x							
Application process and group admittance to UWS; includes station selection and funding	x	x	x	x			x					
Equipment inventory, purchase, inspection, and testing	x	x	x	x							x	X

Activity	J	F	M	A	M	J	J	A	S	O	N	D
Field training and database-related training session(s)			x	x								
Contact with analytical laboratory (for chlorophyll a and nutrient sampling samples)		x	x	x								
Field audits & midseason check in with Monitoring Group Leads or designated appointees					x	x		x				
Monthly check ins with Monitoring Groups					x	x	x	x	x	x	x	x
Technical support to Monitoring Groups				x	x	x	x	x	x	x	x	x
Sampling events					x	x	x	x	x	x		
Data entry					x	x	x	x	x	x	x	x
Data review and validation of data entry	o	o				x	x	x	x	x	x	x
Data uploads to STS-UWS website (must follow data review)			o									
Draft report		o										
Final annual report			o									
o indicates the year following sampling events												

A.7. DATA QUALITY OBJECTIVES

Taken together, precision, accuracy and bias, representativeness, comparability, completeness, and sensitivity comprise the major data quality indicators used to assess the quality of the program's data. A summary of criteria are provided in Table 4.

Definitions of these data quality indicator terms:

- **Precision** is the degree of agreement among repeated field measurements of the same indicator and gives information about the consistency of methods. It is typically defined as relative percent difference, or RPD.
- **Accuracy** is a measure of confidence that describes how close a measurement is to its "true" or expected value; it includes a combination of random error (precision) and systematic error (bias) components of both sampling and analytical operations.
- **Bias** is the systematic or persistent distortion of a measurement process that causes errors in one direction.
- **Representativeness** is the extent to which measurements actually represent the true environmental condition. Parameters, station selection (including location of sampling point within the water column), time, and frequency of sample collection can all play a role in determining how representative a sample is.

- **Comparability** is the extent to which data can be compared between sample locations or periods of time within a project, or between different sites.
- **Completeness** is the comparison between the amount of valid or usable data the program originally intended to collect versus how much was actually collected.
- **Sensitivity** is the capability of a method or instrument to discriminate between measurement responses representing different levels of the variable of interest.

Table 4: Measurement Performance Criteria

Data Quality Indicators	Measurement Performance Criteria	QC Sample and/or Activity Used to Assess Criteria
Precision – overall	RPD \leq value indicated in Table 5	field duplicates
Precision – analytical	RPD \leq value indicated in Table 5	analytical duplicates
Accuracy / Bias	85% \leq recovery \leq 115%	certified reference material
Comparability	standard procedures followed	NA
Completeness	data from surface, mid (if applicable) and bottom at each station meet data quality objectives	data completeness check
Sensitivity	value \geq MDL*	sample value check

* MDL = method detection limit. This is a reporting limit based on the lowest standard accurately analyzed in the analysis.

Precision - Precision objectives are listed in Table 5. Precision is evaluated in the field by participants taking replicate measurements for at least 5% of samples, where applicable.

For UWS Tier I water quality sample events, a replicate profile at one station per field day.

For water samples filtered and analyzed for extracted chlorophyll-*a*, two field replicates will be collected per sample day.

For estuary and tributary stations sampled for nutrients, a field replicate will be collected at each station.

For qualitative macrophyte stations, multiple pictures will be taken for rake toss sampling and beach sampling.

For quantitative macrophyte sampling, multiple analyses of the stills to determine percent coverage will be undertaken.

When a multiparameter sonde is used, standards will be read before and following a trip, within one day of the field day. Calibration of sondes will happen within one day of the sampling event. Multiparameter sondes can hold their calibrations for weeks. The pre and post sampling event readings can identify any potential drift outside of manufacturer recommendations for calibration. These values will be kept with all other data for review at the end of the project interval.

The Onset HOB0 loggers (dissolved oxygen) and Star-Oddi loggers (temperature, conductivity, depth) will be deployed in a common water bath before deployment and following deployment. Conductivity, temperature and oxygen will be varied in the bath, allowing for multiple values for intercomparison. The temperature, conductivity, and oxygen of the bath will be determined with the instruments being used for conducting Tier I water quality profiles. These pre- and post-baths will serve to cross-calibrate all instruments and to determine if the deployed loggers exhibited any drift over the course of the deployment. The deployed loggers will be intercalibrated by applying a multiplicative correction if initial values differ by more than 10% from the reference value (as determined from the YSI EXO1 sonde or Eureka Manta +35).

The frequency of field replicate measurements for each parameter are described in Table 7.

Relative percent difference (RPD) of replicate samples is used as one index of precision; see Table 5. This is defined as the absolute difference between the replicates divided by the average of the replicates. The allowable RPDs for each parameter are provided in Table 7. A difference greater than the designated RPD requires further investigation of the sample run. If the difference is large enough, it indicates failure (unless the average of the two samples is less than 10 times the method detection limit), and results in potential disqualification or flagging of data from that station depth, unless there is a reasonable and supported explanation for the inconsistency. Replicate precision will be analyzed by calculating the RPD using the equation:

$$\text{RPD (\%)} = |x_1 - x_2| / ((x_1 + x_2)/2) * 100$$

where x_1 is the original sample concentration and x_2 is the replicate sample concentration.

The Microsoft Excel formula for calculating the RPD is:

$$= \text{ABS}(X1-X2) / ((X1+X2) / 2) * 100$$

where $X1$ is the original sample concentration and $X2$ is the replicate sample concentration. The RPD is automatically calculated in the UWS data entry template for replicate profiles and field samples.

Accuracy and Bias - Accuracy objectives are listed in Table 4. Procedures used to test or ensure accuracy are described in Table 11. While training and audits help to ensure measurement accuracy and precision, quantitative measures of accuracy for water quality monitoring are estimated using laboratory QC data (blank results, fortified matrix results, known QC samples,

etc.). When a multiparameter sonde is used, standards will be read before and following a trip, within one day of the field day. Extracted chlorophyll a analysis will include a field replicate, laboratory blank and reference standards. Nutrient analysis will include a laboratory blank, field blank, and reference standards. Data loggers will be calibrated prior to and after deployment. Biweekly comparative readings between loggers and sonde will be recorded to keep a log of any drift occurring with loggers. These data will be evaluated with the log data in the final report and during the season.

Representativeness – Tier I water quality sample stations and quantitative Tier II macrophyte stations are selected to represent the entire embayment using a probability-based sampling design⁴, as in the EPA National Coastal Assessment⁵. In this approach, a 0.42 km² hexagonal grid is overlain on the site map. Three random stations are generated in ArcGIS in each section of the embayment within a hexagon. Large embayments will have hexagons for random station generation selected with the UWS Science Advisors. A minimum of four stations will be sampled in each embayment, with larger sites having up to twelve stations. The location of the station in each hexagon will be randomly generated, with at least two alternate locations also randomly generated, in case the original location is deemed unusable (e.g. too shallow or in the middle of a navigation channel). If none of the three random stations are accessible, a station will be determined as close as possible to a randomly generated station unless bias circumstances are identified. The Monitoring Program Coordinator will oversee station selection, providing GIS-based maps and station coordinates to groups participating in the Unified Water Study. At least one Scientific Advisor affiliated with the UWS will also advise on the choice of station locations. Stations used previously by a group are evaluated for bias before inclusion in the UWS.

Tier II nutrient stations are a subset of the Tier I water quality stations selected in consultation with the UWS Science Advisors. A pilot test of nutrient sampling was conducted in Mamaroneck River, NY and Little Neck Bay, NY in 2017. In this pilot test, all Tier I water quality stations were sampled for nutrients. The approach to selecting a subset of stations (2-3) for Tier II nutrient analysis to yield a regional average was compared to the regional average using all stations in a region (3-5). The approach described below yields the most accurate regional average when sampling a subset of Tier I water quality stations for nutrients.

⁴ Paul, J.F., J.L. Copeland, M. Charpentier, P.V. August, and J.W. Hollister. 2003, Overview of GIS applications in estuarine monitoring and assessment research. *Marine Geodesy Journal* 26: 63-72.

⁵ EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

At a minimum, there will be two stations per embayment. In embayments with multiple regions, there will be a minimum of two stations per region. These stations are located at the boundaries of region delineations or the entire embayment, when regions are not identified.

Embayment size, salinity, and best judgment of the UWS Science Advisors determine if more stations may be needed in a region. Additional stations are required when the geometry of the embayment is not linear. For example, the Y shape of Mamaroneck River, NY requires three stations as seen in this QAPP. Salinity within a region that has a range greater than 2 ppt triggers a closer inspection of available salinity data by the UWS Science Advisors, to insure selected stations are sufficient to characterize the region.

Stations that have salinity data of less than 5 ppt are not acceptable as representative embayment stations. In embayments with the minimum Tier I water quality stations (4 stations), a station where salinity is episodically below 5 ppt may be deemed acceptable if the UWS Science Advisors deem inclusion of the station as important to estimating the regional average. The next station downstream is evaluated for inclusion in cases where a station is rejected.

Tier II nutrients stations (tributary and Long Island Sound reference) are chosen based on access, location, and salinity values under 1 PPT.

Data logging stations are selected with at least one station per reporting region. Access to station is a strong consideration in the selection process. The final station locations will be conferred with UWS Science Advisors before deployment.

Qualitative macrophyte surveys are targeted qualitative assessments of areas in the embayment known to harbor macrophytes thus the random station generation does not apply to selecting these locations. Sample collection timing and frequency for water quality stations are selected to capture data that are representative of embayment conditions. While tidal stage will vary among sampling dates, the timing relative to dawn was considered of greatest importance when sampling Tier I water quality stations to evaluate hypoxia in embayments. These very shallow systems are typically dominated by benthic primary producers (macroalgae, benthic microalgae, and seagrass) versus pelagic primary producers (pelagic microalgae / phytoplankton). When the sun rises, these primary producers quickly replenish the dissolved oxygen in the water column. One of the goals of this study is to evaluate the incidence of hypoxia in embayments, thus sampling close to dawn is more important than sampling at a specific tidal stage. Time of high and low tide and precipitation volumes are recorded and will be considered in the analysis of results. Any abnormal or episodic conditions that may affect the representativeness of sample data are noted and maintained as metadata.

Comparability - The comparability of the data collected can be assured by using known protocols and documenting methods, analysis, sampling sites and stations, times and dates, sample storage and transfer, as well as laboratories and identification specialists; so that future

surveys can produce comparable data by following similar procedures. Examples of project procedures are available in the collection of Standard Operating Procedures (SOPs) provided in Appendix A of this document.

Completeness –Minimum sample events for inclusion for Tier I and Tier II monitoring are included in section A.6 of this document.

Sensitivity – Sensitivity objectives are listed in Table 5. Sensitivity is the lowest detection limit of the method or instrument for each of the measurement parameters of interest. For analytical methods, these are the method detection limits (MDLs).

Table 5: Data Quality Objectives

Parameter	Units	Accuracy	Precision (allowable RPD)	Approx. Expected Range	Sensitivity (Resolution or MDL)
Depth (calibrated line)	meters (m)	± 0.1 m	20%	0 – 50 m	0.1 m
Barometric Pressure (ONSET HOBO U20L-01)	Kilopascal (kPa)	0.62 kPa maximum error	10%	3.7 – 4.1 kPa	< 0.02 kPa
Depth (YSI EXO 1)	meters (m)	0 to 10 m ± 0.04% FS or ± 0.004 m 0 to 100 m ± 0.04% FS or ± 0.04 m 0 to 820 m ± 0.04% FS or ± 0.1 m	20%	0 – 50 m	0.001 m
Depth (Eureka Manta +35)	meters (m)	0 to 10 m ±0.02 (±0.2% of FS) 0 to 25 m ±0.05 (±0.2% of FS) 0 to 50 m ±0.1 (±0.2% of FS) 0 to 100 m ±0.2 (±0.2% of FS) 0 to 200 m ±0.4 (±0.2% of FS)	20%	0 – 50 m	0.01 m 0.01 m 0.1 m 0.1 m 0.1 m
GPS coordinates	decimal degrees (dec. deg.)	± 7.8 m http://www.gps.gov/systems/gps/performance/accuracy/	for reference point on land, within 10 m (=0.0001 dec. de g.)	NA	1.02 m

Parameter	Units	Accuracy	Precision (allowable RPD)	Approx. Expected Range	Sensitivity (Resolution or MDL)
Temperature (YSI EXO 1)	degrees Celsius (°C)	-5 to 35 °C ± 0.5 °C 35 to 50 °C ± 0.05 °C	10%	4 – 26 °C	0.001 °C
Temperature (Eureka Manta +35)	degrees Celsius (°C)	± 0.1 °C	10%	4 – 26 °C	0.01 °C
Conductivity (YSI EXO 1)	millisiemens (mS/cm)	0 to 100 mS/cm ± 5% of reading or 0.0001 mS/cm; whichever is greater 100 to 200 mS/cm ± 5% of reading	10%	0 – 50 mS/cm	0.0001 to 0.001 mS/cm, range- dependant
Conductivity (Star-Oddi DST CT)	millisiemens (mS/cm)	13-50 mS/cm ± 1.5 mS/cm	10%	13-50 mS/cm	0.01 mS/cm within range
Specific Conductance (Eureka Manta +35)	millisiemens (mS/cm)	0 to 10 mS/cm ± 1% of reading or ± 0.001 mS/cm 10 to 100 mS/cm or ± 1% of reading	10%	0 – 50 mS/cm	0.001 mS/cm 0.01 mS/cm
Dissolved oxygen (YSI EXO 1)	milligrams per liter (mg/L) = parts per million (ppm); percent saturation (% sat.)	0 to 20 mg/l ± 1% of reading or 0.1 mg/L 20 to 50 mg/l ± 5% of reading 0 to 200% ± 1% reading or 1% air saturation, whichever is greater 200 to 500% ± 5% reading	20%	0 – 14 mg/L 0 – 120 %	0.01 mg/L 0.1 % sat.
Dissolved oxygen (Eureka Manta +35)	milligrams per liter (mg/L) = parts per million (ppm); percent saturation (% sat.)	0 to 20 mg/l ± 0.2 mg/l 20 to 50 mg/l ± 10% reading 0 to 200% sat. ±1% of reading or ±0.1 % sat. 200 to 500% sat. ±10% of reading	20%	0 – 14 mg/L 0 – 120 % sat.	0.1 mg/l 0.1 % sat.

Parameter	Units	Accuracy	Precision (allowable RPD)	Approx. Expected Range	Sensitivity (Resolution or MDL)
Dissolved oxygen (ONSET HOBO U26)	milligrams per liter (mg/L) = parts per million (ppm)	0 to 8 mg/l \pm 0.2 mg/l 8 to 20 mg/l \pm 0.5 mg/l	20%	0 – 14 mg/L	0.02 mg/l
Chlorophyll a (as measured in lab)	microgram per liter (μ g/L)	75 - 125 % recovery of a lab QC sample with known μ g/L	15%	0 – 30 μ g/L; though higher concentrations may occur	0.7 μ g/L
Chlorophyll a (YSI EXO 1)	Relative Fluorescence Units (RFU), microgram per liter (μ g/L)	Chl: $R^2 > 0.999$ for serial dilution of Rhodamine WT Solution from 0 to 400 μ g/L PC equivalents	20%	0 – 30 μ g/L; though higher concentrations may occur	0.01 RFU 0.01 μ g/L
Chlorophyll a (Eureka Manta +35)	microgram per liter (μ g/L)	0.03 to 500 μ g/L \pm 3% of full scale	20%	0 – 30 μ g/L; though higher concentrations may occur	0.01 μ g/L
Turbidity (YSI EXO 1)	FNU*	0 to 999 FNU \pm 2% of reading or 0.3 FNU, whichever is greater 1000 to 4000 FNU 0.1 FNU	20%	0 – 30 FNU	0 – 999 FNU: 0.01 FNU 1000 – 9999 FNU: 0.1 FNU
Turbidity (Eureka Manta +35)	NTU	0 to 400 NTU \pm 1% of reading \pm 1 count 400 to 3000 NTU \pm 3% of reading	20%	0 – 30 NTU	4 digits 4 digits
Dissolved ammonia - NH ₃ (as measured in lab)	mg/L NH ₃ (= ppm = g/m3)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-1 mg/l	0.020 mg/l
Dissolved nitrate [†] - NO ₃ - (NO _x - NO ₂ -)	mg/l NO ₃ (= ppm = g/m3)	Value calculated from multiple N analyses	NA	0-2 mg/l	NA
Dissolved nitrite - NO ₂ - (as measured in lab)	mg/L NO ₂ (= ppm = g/m3)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-0.7 mg/l	0.004 mg/l

Parameter	Units	Accuracy	Precision (allowable RPD)	Approx. Expected Range	Sensitivity (Resolution or MDL)
Nitrate-nitrite – NO _x or NO ₃ ⁻ + NO ₂ ⁻ (as measured in lab)	mg/L NO _x (= ppm = g/m ³)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-2.5 mg/l	0.004 mg/l
Dissolved inorganic nitrogen [†] – DIN (NH ₃ +NO _x)	mg/L DIN (= ppm = g/m ³)	Value calculated from multiple N analyses	NA	0-4 mg/l	NA
Total dissolved nitrogen – TDN (as measured in lab)	Mg/l TDN (= ppm = g/m ³)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-5 mg/l	0.05 mg/l
Total Nitrogen (as measured in lab)	mg/l TDN (= ppm = g/m ³)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-8 mg/l	0.05 mg/l
Total inorganic nitrogen [†] – TIN (NH ₃ +NO _x)	mg/L TIN (= ppm = g/m ³)	value calculated from multiple N analyses	NA	0-4 mg/l	NA
Total organic nitrogen [†] – TON (TN - TIN)	mg/L TON (= ppm = g/m ³)	value calculated from multiple N analyses	NA	0-5 mg/l	NA
Dissolved organic nitrogen [†] – DON (TDN - DIN)	mg/L DON (= ppm = g/m ³)	value calculated from multiple N analyses	NA	0-4.5 mg/l	NA
Particulate nitrogen [†] – PN (TN-TDN)	mg/L PN (= ppm = g/m ³)	value calculated from multiple N analyses	NA	0-0.5 mg/l	NA
Total phosphorus – TP (as measured in lab)	mg/L TP (= ppm = g/m ³)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-0.5 mg/l	0.01 mg/l
Dissolved organic nitrogen [†] - DON (TDN - DIN)	mg/L DON (= ppm = g/m ³)	value calculated from multiple N analyses	NA	0-4.5 mg/l	NA
Dissolved orthophosphate – PO ₄ ³⁻ or DIP (as measured in lab)	mg/L PO ₄ ³⁻ mg/L DIP (= ppm = g/m ³)	85% - 115% recovery of lab fortified matrix (LFM)	Field Replicate 30% Analytical Replicate 15%	0-0.3 mg/l	0.001 mg/l

Parameter	Units	Accuracy	Precision (allowable RPD)	Approx. Expected Range	Sensitivity (Resolution or MDL)
Quantitative macrophyte amount	% coverage bare, macrophytes, and animals of bottom	Estimates from three analyses are compared. If the relative percent difference among the three estimates is greater than 5%, the Monitoring Group Lead examines the image and the three estimates, choosing the appropriate value. The three estimates will not be changed, values are retained to show the inconsistency. The Monitoring Group Lead decides on the final value for the estimate.			
Qualitative macrophyte amount	choice of: none, some, lots	This is a qualitative assessment, not quantitative. Photos are reviewed by a UWS Science Advisor or trained designee to confirm choice of amount.			

*: FNU and NTU are interchangeable in the UWS. All data reported as NTU.

†: This parameter is calculated rather than measured analytically, so MDL is not computed. RPD is also not relevant for this parameter.

A.8. SPECIAL TRAINING / CERTIFICATION

UWS trainings are hands on full day events*. They are designed for a wide range of experience in water quality monitoring; ranging from groups with an extensive existing water quality monitoring programs to groups with little to no water quality monitoring experience. The trainings help to ensure all groups are on the same understanding of project QAPP and SOPs. All Monitoring Groups are provided the SOPs and QAPP. They are required to read these documents. Monitoring Groups take notes on the project SOPs during training events to clarify any points that require extra attention. The objective of trainings is to have all Monitoring Groups, regardless of previous experience, following the project requirements in a unified manner.

The Monitoring Program Coordinator shall ensure that all UWS Monitoring Groups receive appropriate training by organizing and conducting training events. The trainings are mandatory for new and veteran groups and have hands on elements for sonde calibration and usage, filling in datasheets, macrophyte assessments, logger calibration and usage, nutrient sample collection, and filtering chlorophyll a. All topics are covered in guided step by step approach. Training will be assessed as described by checks in Section C.1 Assessment and Response Actions.

The Monitoring Program Coordinator enters training into the project database and records the following information: subject matter (i.e. what type of monitoring and procedures are covered), training course title, date and agenda, name and qualification of trainers, and names of participants trained with associated monitoring group name. The trainings and technical support offered through the Monitoring Program Coordinator, Monitoring Program Field Coordinator, and Science Advisors is in place for the duration of the project. This will ensure new and veteran groups have a reliable source for prompt answers to their inquiries. Groups are prompted to email or call the Monitoring Program Coordinator or Monitoring Program Field Coordinator with their questions. If needed, the Science Advisors will be consulted. This support

is mentioned throughout the project duration and emphasized at the trainings. Trainers remind trainees to call at any hint of a question or issue so it can be resolved.

The Monitoring Program Coordinator worked closely with the Science Advisor signatories on this QAPP to confirm procedures are appropriate. He was part of the three person team leading UWS Tier I trainings around Long Island Sound in 2017. The Coordinator oversees Save the Sound's Water Quality Program which samples water for pathogen-indicator bacteria and participates in the UWS Tier I & II monitoring.

*: UWS trainings for the 2020 season will be held remotely due to concerns and timing in regard to spread of the Coronavirus (Covid-19).

The Coordinator holds a Bachelor of Science Degree in Environmental Studies from University California Santa Barbara and a Master's of Science Degree in Environmental Science and Management from Sacred Heart University.

Project training shall take place as specified in Table 6.

Table 6: Project-Specific Training

Training: Type & Description	Trainer(s)	Training Date(s)	Trainees	Location of Training Records
<p>UWS Standard Operating Procedures and Methods, QAPP, and Data Entry Training</p> <ul style="list-style-type: none"> - General water quality parameter information - Sonde calibration and field training - Chlorophyll a field collection, filtering, preservation, & transport - Nutrient sampling collection, filtering, preservation, & transport - Qualitative macrophyte assessment procedure - Quantitative macrophyte assessment procedure - Data logger calibration and field training - QAPP review and data entry 	<p>UWS Monitoring Program Coordinator, UWS Field Coordinator, UWS Science Advisor(s), and other personnel under the supervision of the listed trainer(s)</p>	<p>Annual; Spring before sampling season commences</p>	<p>All participating UWS groups will send 1-3 representatives</p>	<p>Office of the UWS Monitoring Program Coordinator; digital record of attendees and agenda stored on computer and backed up on Save the Sound S-Drive</p>

A.9. DOCUMENTS AND RECORDS

Calibration Datasheet, Sample Event Datasheet and Field Datasheet will be completed by Monitoring Groups before, during, and after Tier I water quality station sampling event.

Qualitative Macrophyte Field Datasheet will be completed upon every qualitative macrophyte survey. Photographs must accompany and be identified in this datasheet.

UWS Nutrient Sample Event Datasheet, Calibration Datasheet, Field Datasheet will be completed upon every Tier II nutrients sampling event. The calibration datasheet will only contain the parameters being recorded in the field: conductivity (salinity).

UWS Quantitative Macrophyte Field Datasheet will be completed upon every quantitative macrophyte survey.

UWS Logger Retrieval Sample Event Datasheet, Field Datasheet, and Calibration Datasheet will be completed upon every data retrieval of data loggers in the field. The calibration datasheet will only contain the parameters being recorded in the field: dissolved oxygen and conductivity.

Sample Labels will be put on all sample containers. Labels will include the station name, organization name, date, time, sample id, and type of sample. Samples needing containers with labels are filters for extracted chlorophyll a and nutrients. Detailed instructions for chlorophyll a filters and nutrient samples are provided in the UWS SOP Chlorophyll, UWS SOP Filtered Nutrients, and UWS SOP Total Nitrogen and Total Phosphorous SOP in Appendix A.

Chain of Custody (COC) forms will accompany samples from collection sites to laboratories. COC forms will be signed by collectors and all individuals who gain custody of the samples until they arrive at a lab. Information will agree with the label information on the sample containers and field datasheet. UWS Chain of Custody forms are in Appendix B.

Training records and field audit information will be kept by the Monitoring Program Coordinator.

The electronic project database shall be organized and protected from loss and damage through proper back-up of digital data on Save the Sound's S-Drive.

No scientific collecting permits or certificates of permission are required.

The specific forms to be used for this project are provided in Appendix B.

B. Data Generation and Acquisition

B.1. SAMPLING PROCESS DESIGN (EXPERIMENTAL DESIGN)

Tier I water quality sample stations, Tier II nutrients stations (estuary), and quantitative Tier II macrophyte stations were selected to represent the water quality of the entire embayment using a probability-based sampling design⁶, as in the EPA National Coastal Assessment⁷. Monitoring Program Coordinator and a UWS Science Advisor advised on the choice of station locations. The UWS assigned unique ID codes for the embayments, reporting regions of the embayment, and stations. Water quality stations can be viewed in section A.6.b of this document.

Qualitative macrophyte stations are selected based on local knowledge and observation of the embayment during the course of the May and June sampling events. They are targeted to areas of macrophyte abundance. Monitoring Program Coordinator and a UWS Science Advisor advised on the choice of station locations. UWS SOP Qualitative Macrophytes describes the process for sampling locations for macrophytes. These stations are selected using an adaptive process that requires observations during May and June sampling events. These observations identify the best locations for high macrophyte abundance. The qualitative macrophyte sampling can be classified as judgmental design resulting in directed sampling information to complement the water quality station data. Macrophyte stations may change from year to year but records of locations are maintained by the UWS Monitoring Program Coordinator.

Data logging stations were selected with at least one station per reporting region. Access to station is a strong consideration in the selection process. The final station locations will be conferred with Monitoring Program Coordinator and the UWS Science Advisors before deployment.

Tier II nutrients stations (tributary) were selected based on access and salinity <1 ppt. Parameters, number and location of sampling sites, sampling time of day, frequency, and season are selected to meet the monitoring objectives referred to in Section A.6.a.

⁶ Paul, J.F., J.L. Copeland, M. Charpentier, P.V. August, and J.W. Hollister. 2003, Overview of GIS applications in estuarine monitoring and assessment research. *Marine Geodesy Journal* 26: 63-72.

⁷ EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

Sampling design components are described below:

Sampling Safety. Personal safety shall be a primary consideration in all activities, including selection of sampling stations, dates, and training programs. No sampling shall occur when personal safety is thought to be compromised. The Monitoring Group Lead of each participating group on this QAPP shall confer with their respective field teams before each sampling event to decide whether adverse weather or other conditions pose a threat to safety and will cancel/postpone sampling when necessary. Sampling shall take place in teams of two or more. Samplers shall wear life vests when required, in adverse conditions in boats, or wading in waters under difficult conditions. Samplers shall wear proper clothing to protect against the elements.

Design Considerations. A summary of design considerations incorporated into this project are included in Table 7. Specifics on the design approaches to the number of stations, depth of sampling, and frequency of sampling and time of day of sampling are included in the SOPs in Appendix A. A summary of general design approaches to the number of stations, depth of sampling, frequency of sampling, and time of day are included here:

There are 237 Tier I water quality stations and approximately 121 qualitative macrophyte station monitored across all the study sites. 16 data logging stations will be monitored. 75 quantitative macrophyte stations will be monitored. 49 Tier II nutrient stations will be monitored. The qualitative macrophyte stations are confirmed by the beginning of the macrophyte monitoring window described in UWS SOP Macrophytes. Sonde profiles for water quality parameters at water quality stations will be sampled 0.5 m below the surface, 0.5 m above bottom, mid-depth if total depth >10 m; if total depth is less than 1.5 m, only a mid-depth reading will be collected. Extracted chlorophyll a samples will be taken from a bucket. Two filters and a corresponding chlorophyll a sonde reading will be taken per sampling event. Land-based qualitative macrophyte stations will be photographed from land. Rake toss qualitative macrophyte stations will be photographed from land or boat. Quantitative macrophyte stations are recorded from a boat. These are the same locations as the Tier I water quality stations in the respective embayments. Tier II nutrients stations are collected 0.5 below the surface. New Tier II nutrients tributary stations are confirmed by field work prior to collecting the first batch of nutrient samples for the season. Data logging stations are 0.5 m off the bottom and record data every 15 minutes. These stations are selected and confirmed with the Monitoring Program Coordinator and a UWS Science Advisor prior to commencement of the season.

Table 7: Sampling Approaches.
Assessment Type: Tier I water quality Stations.

Indicators	Number of sample locations	Frequency, duration, special conditions	Field survey QC
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	every station	Twice a month from May - October, within 3 hours of sunrise,	repeat readings every time a station is sampled.
station depth			coordinates indicating a 100 m or greater discrepancy will be assessed and documented in final report.
sample depth	reference land site, once per sampling event		
temperature	once per field day, take readings twice at the last station sampled		
salinity			
dissolved oxygen			
turbidity			
chlorophyll a	every station , 0.5 m below surface		take readings twice for replicate at the last station sampled
	once per sampling event from bucket at reference station		calibration per SOPs
			collect filter and sonde readings at reference station

Stations are representative, defined clearly in respective SOPs

Assessment Type: Qualitative Macrophyte Surveys.

Indicators	Number of sample locations	Frequency, duration, special conditions	Field survey QC
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	Every station	Sample 3 days during the 3-week period starting July 15 and ending August 7. Try to sample once per week. If this is not possible, sample such that you maximize the days between sampling. All three days cannot be sampled in the same 7-day window. 2 sampling events or an event outside this time criteria may be accepted in unforeseeable circumstances. Consultation with Monitoring Program Coordinator and UWS Science Advisor is necessary for this decision. See UWS SOP Macrophytes for additional details.	repeat readings every time a station is sampled. coordinates indicating a 100 m or greater discrepancy will be assessed and documented in final report reference land site, once per sampling event
Macrophyte Abundance			Photos and assessment (none, some, lots) of each sample are reviewed by the Monitoring Program Coordinator and UWS Science Advisor

Stations are targeted, defined clearly in SOP

Assessment Type: Tier II nutrients Stations.

Indicators	Number of sample locations	Frequency, duration, special conditions	Field survey QC
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	Every station	Each visit to sample station; at least monthly (May – October), with 14 days separation	Repeat readings every time the station is sampled to verify coordinates. Coordinates indicating a 100 m or greater discrepancy from documented coordinates will be assessed and documented in data notes.
Salinity	Every station	Each visit to sample station; at least monthly (May – October), with 14 days separation	Probe calibration prior to survey; post sampling event readings in standard
Nutrients	Every station	Each visit to sample station; at least monthly (May – October), with 14 days separation	At minimum one field replicate per sampling event

Station are representative.

Assessment Type: Data logging stations.

Indicators	Number of sample locations	Frequency, duration, special conditions	Field survey QC
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	Every station	Each visit to sample station	Repeat readings every time the station is sampled to verify coordinates . Coordinates indicating a 100 m or greater discrepancy from documented coordinates will be assessed and documented in data notes.
Conductivity	Every station	Each visit to sample station	Probe calibration prior to survey; post sampling event readings in standard
Dissolved oxygen	Every station	Each visit to sample station	Probe calibration prior to survey; post sampling event readings in standard
Barometric pressure	Every station	Each visit to sample station	Not applicable

Stations are selected for access and other considerations addressed in this QAPP and UWS Data Logging SOP.

Assessment Type: Quantitative Macrophyte Surveys.

Indicators	Number of sample locations	Frequency, duration, special conditions	Field survey QC
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	Every station	Each visit to sample station while conducting all camera descents to bottom. Obtained from GPS track	Not applicable
Bottom coverage (% macroalgae, % bare, % eelgrass)	Every station	Each visit to sample station with specifications in SOP on image count and analysis	Not applicable

B.2. SAMPLING METHODS

Pre-coordination shall occur with the external lab to ensure that sample collection procedures meet lab needs. The project labs for this study are below:

Chlorophyll a

Save the Sound, John and Daria Barry Foundation Water Quality Laboratory, contact: Peter Linderoth & Elena Colón, plinderoth@savethesound.org; ecolon@savethesound.org, 914-381-3140, 1385 Boston Post Rd, 2nd Fl, Larchmont, NY 10538

Tier II Nutrients

Interstate Environmental Commission Lab, contact: Evelyn Powers, epowers@iec-nynjct.org, 718-982-3792, c/o College of Staten Island-CUNY, 2800 Victory Blvd., Building 6S, Room 106, Staten Island, NY 10314

A laboratory of equal or higher certification than Save the Sound or Interstate Environmental Commission can be considered if Save the Sound or Interstate Environmental Commission are unable to complete project analyses. This surrogate laboratory must adhere to analytical methods in Table 10.

To comply with UWS program guidelines, all sample collections for this project shall follow detailed methods on how samples will be collected and preserved as stated in the standard operating procedures (SOPs) contained in Appendix A of this document. The lab has reviewed the UWS SOPs and confirms they are appropriate for the select analyses.

A summary overview of sample collection methods is provided in Table 8. A summary of field sampling considerations is provided in Table 9.

Any filters or nutrient sample bottles collected for analysis will be stored in a cooler, on ice during the sampling trip. The cooler designated for these samples will not be used for the storage of macrophytes.

All deviations from the Standard Operating Procedures of this QAPP will be documented and subsequently reviewed by the Monitoring Program Coordinator and the project UWS Science Advisors. This information will be available to all signatories at the completion of this project at which time acceptability of data will be determined.

Table 8: Overview of Sample Collection Methods
Assessment Type: Tier I water quality Parameters

(NA = not applicable)

Parameter(s)	Container Type(s) and Preparation	Minimum Sample Quantity per Sample Depth (unless otherwise noted)	Sample Preservation	Maximum Holding Time
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	<i>in situ</i>	1 / station	transfer to digital format, maintain back-up copies of digital data	NA
Station depth	<i>in situ</i>	1 / station	transfer to digital format, maintain back-up copies of digital data	NA
Sample depth (metered line)	<i>in situ</i>	1, and for remainder of the sampling event if the intercomparison with project sonde and line is > 0.3 m	transfer to digital format, maintain back-up copies of digital data	NA
Multiparameter sonde: <ul style="list-style-type: none"> • depth • temperature • salinity • dissolved oxygen • turbidity • chlorophyll a 	<i>in situ</i>	1 / sample depth and a second reading for each depth at the last station of the day	transfer to digital format, maintain back-up copies of digital data	NA

Parameter(s)	Container Type(s) and Preparation	Minimum Sample Quantity per Sample Depth (unless otherwise noted)	Sample Preservation	Maximum Holding Time
chlorophyll a (extracted, fluorometric analysis)	<i>Large bucket rinsed three times with surface water. Bucket volume must accommodate sonde for reading in situ</i>	Filtered sample volume must be sufficient to provide “color” on the filter pad; 10 mL to 180 mL. 2 filters collected and one sonde reading from bucket per sampling event	GF/F filter is blown dry with a 60mL syringe and stored in the dark (foil wrapped), on ice; transferred to - 20°C freezer within 12 hours	28 days

Assessment Type: Qualitative Macrophyte Abundance

(NA = not applicable)

Parameter(s)	Container Type(s) and Preparation	Minimum Sample Quantity	Sample Preservation	Maximum Holding Time
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	<i>in situ</i>	1 / station	transfer to digital format, maintain back-up copies of digital data	NA
Macrophyte abundance	digital photos	1 / sample (a sample is a single rake toss or required distance for soft shoreline photo)	transfer to a computer, upload to online datasheet	NA

Assessment Type: Tier II nutrients Parameters

(NA = not applicable)

Parameter(s)	Container Type(s) and Preparation	Minimum Sample Quantity	Sample Preservation	Maximum Holding Time
GPS: latitude & longitude in decimal degrees; NAD83 coordinate system or record system used	<i>in situ</i>	NA	transfer to digital format; maintain back-up copies of digital data	NA
Multiparameter sonde: • Salinity	<i>in situ</i>	NA	transfer to digital format; maintain back-up copies of digital data	NA

Parameter(s)	Container Type(s) and Preparation	Minimum Sample Quantity	Sample Preservation	Maximum Holding Time
Inorganic nutrients	high density polyethylene (HDPE) polypropylene (new containers washed with ASTM Type 1 Ultrapure Water, used containers pre-acid-washed with 10% hydrochloric acid)	120 mL per station	ice or refrigerate filtered water samples at a temperature of <4 C while in the field, store at <-20 C	holding time of ~1 year once frozen
Total nutrients	high density polyethylene (HDPE) polypropylene (new containers washed with ASTM Type 1 Ultrapure Water, used containers pre-acid-washed with 10% hydrochloric acid)	120 mL per station	ice or refrigerate water samples at a temperature of <4 C while in the field, freeze at <-20 C	holding time of ~1 year once frozen

Assessment Type: Quantitative Macrophyte Abundance

Parameter(s)	Container Type(s) and Preparation	Minimum Sample Quantity	Sample Preservation	Maximum Holding Time
Macrophyte abundance (% bare, % macroalgae, % eelgrass)	Field data sheets and computer storage	20 still images per station; more if heterogeneity is observed. Detailed procedure in SOP	maintain back-up copies of digital data	NA

Table 9: Overview of Field Sampling Considerations

Sample Type	Parameter(s)	Sampling Considerations
<i>In-situ</i> sampling	Station depth	Note the tidal stage and time of day. Depth varies greatly over the tidal cycle.
<i>In-situ</i> sampling, GPS	GPS: latitude & longitude in decimal degrees; NAD83 or WGS84 coordinate system, record system used	NAD83 or WGS84 coordinate system, record system used; check GPS accuracy relative to a known, fixed location
<i>in-situ</i> Tier I water quality sampling, multiparameter sonde	Depth Temperature Salinity Dissolved oxygen Turbidity Chlorophyll a fluorescence	Sample within 3 hours of sunrise. Inspection, maintenance, pre-calibration and post-checking of probes and instruments are critical to achieving accurate and precise measurements.

Sample Type	Parameter(s)	Sampling Considerations
Data logging stations	Dissolved Oxygen, Conductivity (Salinity), Barometric Pressure	Inspection, maintenance as specified by manufacturer, and calibration of instruments are critical to achieving accurate and precise measurements, especially for DO. Loggers are rinsed and cleaned with freshwater after each retrieval and use.
Grab samples - i.e. collection of a water sample	Chlorophyll a	Keep careful and accurate track of volume of water passed through each filter pad, quantitation is impossible without this value.
Qualitative macrophyte abundance	Macrophyte abundance	Be sure to photograph all sites and samples. Record identifier for each photo on the datasheet.
Quantitative macrophyte abundance	Macrophyte abundance	Maintain low speed to minimize potential damage to camera. Monitoring Group Lead and two additional members of the sampling team, under Monitoring Group Lead supervision, will analyze the macrophyte videos as described in SOP
Grab samples - i.e. collection of a water sample in bottle	Inorganic and total nutrients	Triple-rinse sample container in ambient water immediately prior to sample collection. Care must be taken to avoid contact between fingers and inside surfaces of containers, including bottle caps. New, pre-washed bottles preferred; if not, containers for nutrient samples should be acid-washed and rinsed with deionized water. This process is overseen by the Monitoring Coordinator. These bottles will be obtained by appropriate suppliers such as Fischer Scientific. Field filtration preferred for dissolved fractions. If filtering water, triple-rinse container with <i>filtered</i> water immediately prior to sample collection, not ambient water.

B.3. SAMPLE HANDLING AND CUSTODY

Sample handling and labeling procedures shall comply with project Standard Operating Procedures (SOPs). Chlorophyll a filters and nutrient samples will be transported on ice in a cooler to the freezer on the same day as sampling occurs. Filters and nutrient samples will always be transported on ice with no more than 24 hours out of freezer to avoid thawing.

Sample labels will be associated with: station name, date, time, volume filtered, sample id, type of sample, and organization name. These details may be written on the label. Information will also be filled in the field data sheet.

Chain of Custody shall be tracked as detailed in the SOPs. The project Chain of Custody forms are provided in in Appendix B.

The following steps shall be taken to avoid sample mislabeling:

Labels will be prepared in advance and cross checked with the field datasheet before sampling event. Field team will check data sheet versus sample filter labels before storing in the cooler for transport to a freezer. A white board with name of the embayment, site and station id, and date will be filmed prior to recording every station in the macrophyte video surveys.

B.4. ANALYTICAL METHODS

To comply with the requirements of the UWS Program, all analytical methods used in the Monitoring Program, including methods used by laboratories performing analyses for the project, shall be based on standardized laboratory methods.

All analytical methods used for this project are provided in Appendix C.

Table 10 provides an overview of the analytical methods utilized in this Monitoring Program. The SOPs associated with these methods are included in Appendix C.

Table 10: Overview of Analytical Methods.

Parameter	Method #	Source of Method	MDL	Alternative Applications Special Provisions
Chlorophyll a	EPA 445.0	EPA	0.7 µg/l	
Nitrite	EPA 353.2	EPA	0.004 mg/l	
Nitrate+Nitrite	EPA 353.2	EPA	0.004 mg/l	
Ammonia	EPA 350.1	EPA	0.020 mg/l	
Total Dissolved Nitrogen	EPA 353.2	EPA	0.05 mg/l	
Orthophosphate (DIP)	EPA 365.1	EPA	0.001 mg/l	
Total Nitrogen	EPA 353.2	EPA	0.05 mg/l	
Total Phosphorous	EPA 365.1	EPA	0.01 mg/L	

B.5. QUALITY CONTROL

Lab Quality Control (QC) protocols shall be discussed with the external lab facility or contractor analyzing chlorophyll a and nutrient samples prior to sampling to ensure acceptability.

Quality control shall be discussed and defined prior to sampling (e.g., during training).

Details on quality control procedures are provided in Table 11.

Table 11: Quality Control Measures

Note that 5% of field samples equates to one station per field day sampled as replicate.

Sample Type	Instrument/ Parameter	Accuracy Checks	Precision Checks	% Field QC Samples (blanks and field duplicates)
GPS coordinates	GPS or Smart Phone app / GPS coordinates	Compare location of reference site to Google Earth coordinates	Readings at a land- based reference point and duplicate readings at one station	1 / field day
Station depth	metered line / depth	re-measure line	replicate readings at one station	5%
Multiparameter sonde and data loggers	Depth, temperature, conductivity, dissolved oxygen, turbidity, chlorophyll a	Pre-survey calibration and post-survey checks, including “zero” DO standard check	field duplicates or 3- 5 minutes stable readings recorded	5% or verify repeatability in the field
Water samples - grab	Fluorometric determination of extracted chlorophyll a	Acetone blank, standard	QC check for multiparameter sonde	100%
Qualitative Macrophyte abundance	Observation / macrophyte abundance	Photos of all assessments	Photos of all assessments	100% photos are required for inclusion of the data in the UWS
Water samples – grab	TP, P fractions TN, N fractions	Field: blanks Lab: analysis of lab-fortified matrix (spiked samples) and/or lab QC standard	Field duplicates Lab duplicates	Minimum 5%
Quantitative Macrophyte abundance	% coverage bare, macrophytes, and animals of bottom	Still images of all assessments	Still images of all assessments	100% still images are required for inclusion of the data in the UWS

B.6. INSTRUMENT / EQUIPMENT TESTING, INSPECTION AND MAINTENANCE

Maintenance of instruments and equipment shall occur as needed during the field season. Annual maintenance and intercalibration assurance will be conducted by Save the Sound.

Records of equipment inspection, maintenance, repair and replacement shall be kept in a logbook. A backup of the logbook will be kept in a separate location. If the logbook is digital,

appropriate backups of the computer files will be maintained by Monitoring Program Coordinator.

Table 12: Instrument / Equipment Inspection and Testing Procedures

Equipment Type	Inspection Frequency	Type Inspection	Maintenance, Corrective Action
GPS unit	before each sampling date	battery life	charge batteries
Depth line	Annually, or when a potential problem is noted	Check the calibrated line against a meter tape	Wipe tape after each use, if line has stretched or is damaged, replace immediately and note recent data as questionable
Multiparameter sonde	Before each sampling date	Battery life, electrical connections, sensor condition	Charge batteries, spare sensors as appropriate, batteries
Filtering apparatus (chlorophyll a)	Before each use	Proper functioning, clean storage	Spare filters and syringe
Collection rake, rope	Before each collection	Visually for integrity	Repair, replace keep spares on hand
Filtering apparatus (nutrients)	Before each use	Proper functioning, clean storage	Spare syringe, spare filters, spare pump tubing
Logging sensors	Every 7-10 days or as needed	Biofouling and battery check	Clean off fouling organisms, check battery life from data log
Underwater camera and equipment	Before each use	Battery life, test video	Recharge/replace batteries and clean lens if required

B.7. INSTRUMENT / EQUIPMENT CALIBRATION AND FREQUENCY

Calibration shall occur within a day prior to a sampling trip.

Records of calibration shall be kept in a logbook (hard copy or digital, with back-ups). Calibration records shall be maintained for a minimum of four years, ideally longer. Monitoring Groups will deliver calibration records to the Monitoring Program Coordinator. These records and digital backups will be saved on Save the Sound's S-Drive for duration of the project.

A summary of calibration procedures for instruments and equipment is provided in Table 13.

Detailed calibration procedures are described in SOPs contained in Appendices A.

Table 13: Instrument / Equipment Calibration Procedures

Instrument	Inspection and Calibration Frequency	Standard of Calibration Instrument Used	Calibration Acceptance Criteria	Corrective Action
Calibrated lines (for depth)	Annually	Tape measure	Within 0.1 m of tape measure	Recalibrate or replace with calibrated line

Instrument	Inspection and Calibration Frequency	Standard of Calibration Instrument Used	Calibration Acceptance Criteria	Corrective Action
Multiparameter sonde	Before each sampling run	Standard solutions	According to manufacturer's instruction or when not provided a maximum difference of %10 of the calibration standard value	According to UWS and manufacturer's instruction
Logging sensors	Before and after deployment	Standard solutions, according to manufacturer's recommendations	According to manufacturer's instruction or when not provided a maximum difference of %10 of the calibration standard value	According to UWS and manufacturer's instruction

B.8. INSPECTION / ACCEPTANCE OF SUPPLIES AND CONSUMABLES

The procedures for inspection and acceptance of supplies and consumables listed in Table 14 shall be followed by the Monitoring Groups.

Table 14: Supplies Inspection and Acceptance Procedures

Supplies	Inspection Frequency	Type of Inspection	Available Parts	Maintenance
Calibration standards	Before each sampling date	Visual inspection of quantity and expiration date	Spare, fresh solutions	Storage according to manufacturer's recommendations, annual replacement at beginning of sampling season
Sonde sensors, filters	Before each sampling date	Visual inspection of quantity, integrity	Spares	Storage according to manufacturer's recommendations
Field and lab sample sheets	Before each sampling date	Visual	Additional copies	
Cooler	Before each sampling date	Cleanness, ice packs		Annually or as needed
Sample bottles	Before each sampling date	Integrity, cleanness and seal for nutrient bottles, verified sterility of bacterial sample bottles	One set of spare bottles	Clean after use (note that nutrient bottles require acid washing before reuse)

B.9. NON-DIRECT MEASUREMENTS

To provide high-quality data to enhance the interpretation of data collected as part of this Monitoring Program, data may be acquired from qualified sources approved by Monitoring Program Coordinator. NOAA tide gauges will be used for tide information. Precipitation will be

acquired from local weather stations that log reasonable (in respect to northeastern USA conditions) volumes. Precipitation data out of the expected annual volumes and the observed conditions will be flagged and discussed with Monitoring Program Coordinator and shared with quality assurance personnel for review and potential disqualification. External data sources are described in Table 15.

Table 15: Non-Project Data Validity

The following data will be used as part of the Monitoring Program. This is a secondary use of data.

Title or descriptive name of data document.	Source of data.	QAPP written? Y/N	Notes on quality of data.	Planned restrictions in use of the data due to questions about data quality.
Time of low and high tide	NOAA tide gauges recorded on field data sheet	N	NOAA has internal requirements for data suitability.. High and low tide data are not generally available at the embayment. Data from NOAA tide gauges are acceptable; data are used in broad scale, to determine the potential impact of stage in tidal cycle on the day of sampling.	Data quality is acceptable. However, local tidal stage will differ from the nearest NOAA gauge even when corrected for difference in location. These data are rough predictors only.
High and low temperature and precipitation within the 24 hours prior to the field trip	Local weather station recorded on field data sheet	N	Air temperature within the last 24 hours are not generally available at the embayment. Data from nearby weather stations are acceptable; data are used in broad scale, to determine the potential impact of weather on the day of sampling.	Data will be used in comparing among embayments or among dates, as a general indication of weather during the day prior to sampling.

B.10. DATA MANAGEMENT

Field teams shall record data on field sheets, review them, and turn over to respective Monitoring Group Lead or designated appointee.

Monitoring Group Leads or designated appointees shall review sheets and confer with field teams on any needed corrective action.

The designated person shall fill out the chain-of-custody form for forwarding samples to the external laboratory. Each person who handles or transports samples shall also sign the custody form upon receipt of the samples. Chain of custody forms will follow samples to the lab and back to Monitoring Program Lab Coordinator by mail or pickup after each analysis run is completed. Alternatively, scanned copies may be emailed or faxed. These copies will be sent to Monitoring Group Leads or designated appointees.

Once laboratory analyses are complete, the laboratory personnel shall deliver (digital or hard copy) lab results to the Monitoring Program Lab Coordinator or arrange for pickup. These results will be sent to all Monitoring Group Leads or designated appointees.

The Monitoring Group Lead or other trained designee will enter raw field and lab data into the project computer system.

Computer-entered data shall be compared with field sheets for accuracy.

Original data sheets will be stored by the Monitoring Group Leads or designated appointees, following data entry into the UWS data entry template.

Digital back-ups and copies of the non-digitized data will be made and stored in a separate location designated by the Monitoring Group Lead or designated appointees and delivered to the Monitoring Program Coordinator.

Documentation of data recording and handling, including all problems and corrective actions, shall be included in all preliminary and final reports.

Table 16 in this document accurately represents the procedures utilized by the UWS for data management, review, validation, and verification.

Table 16: Data Management, Review, Validation, Verification Process Summary

The term "Field sampler" refers to the person conducting the sampling in the field.

Activity	By whom	Corrective action, if needed
Conduct field audits of Monitoring Groups performing calibrations and demonstrating field procedures.	Monitoring Program Field Coordinator or appointed designee	Correct any discrepancies with this QAPP or SOPs
Check labels just prior to sampling, to ensure correct labeling of container.	Field sampler	Correct label
At time of sampling, record data, sign field sheets.	Field sampler	Remind samplers of proper procedures; retrain if needed.
Fill out, sign chain of custody (COC) forms for any samples going to lab.	Field sampler or designated person	Remind person of proper procedures; retrain if needed.
Before turning field sheets over to Monitoring Group Lead or designated appointee, check for reasonableness to expected range, completeness.	Field sampler	Resample if feasible; otherwise, flag suspect data.
Upon receipt of field sheets, recheck for reasonableness to expected range, completeness, accuracy, and legibility.	Monitoring Group Lead or designated appointee	Confer with field sampler(s) immediately or within 24 hours. Resample if feasible; otherwise, flag suspect data.

Activity	By whom	Corrective action, if needed
Upon receipt of samples, field sheets and COC forms, check to see that sheets and forms correspond to number of samples, condition of samples as stated on COC forms. Sign COC forms. Copies of field sheets and COC forms are made, given to Monitoring Program Coordinator.	Monitoring Group Lead or designated appointee	Contact field samplers as needed to locate missing samples, data records. In case of missing/spoiled samples or data records, authorize resampling as needed and feasible. If resampling is not feasible, flag all suspect data.
Upon completion of laboratory analyses, fill out lab sheets, including data on QC tests.	External Lab	Re-analyze if possible. If not, confer with Monitoring Program Laboratory Coordinator. Flag all suspect data.
Upon receipt of lab data, review for completeness and legibility.	Monitoring Group Lead or designated appointee	Confer with Monitoring Program Laboratory Coordinator.
Upon completion of data entry, compare with field/lab sheets for accuracy.	Monitoring Group Lead or other volunteer. Data entry personnel may review their own work, but it cannot be on the same day as data entry.	Re-enter or correct data.
Translate data into preliminary data reports: run statistical analyses and/or prepare graphical summaries of data. Check for agreement with QC objectives for completeness.	Monitoring Program Coordinator	Confer with QA Officers and UWS Science Advisor(s). Flag or discard suspect data.
In-season (at least once) and end of season review of collected data sets (individual sample runs and season-total compilations); review for completeness and agreement with QC objectives and DQOs.	Monitoring Group Lead or designated appointee	Flag suspect data. Confer with Monitoring Program Quality Assurance Officer.

Data Management Systems – spreadsheets, databases, statistical or graphical software packages, location of data records (paper and electronic), are described here:

All data will be entered from field data sheets to an Excel spreadsheet for storage and retrieval by Monitoring Group Leads and appointed individuals. Digital copies of all datasheets will be kept on file on the S-Drive server in Save the Sound office for at least 4 years with a plan to keep records for duration of the project and beyond. The S-Drive is backed up weekly.

C. Assessment and Oversight

C.1. ASSESSMENT AND RESPONSE ACTIONS

The Monitoring Program Coordinator and UWS Science Advisors will identify and effectively address any issues that affect data quality, personal safety, and other important project components. The progress and quality of the monitoring program shall be assessed to ensure the objectives are being accomplished. The Monitoring Program Coordinator or appointed designees will check at the end of every month from May - October to confirm the following:

- a. Monitoring is occurring as planned.
- b. Sufficient written commentary and supporting photographs exist.
- c. Sufficient field members are available for all sampling groups.
- d. Samplers are collecting in accordance with project schedules.
- e. Datasheets and custody control sheets are being properly completed and signed.
- f. Retraining or other corrective action is implemented at the first hint of non-compliance with the QAPP or SOPs.
- g. Labs are adhering to the requirements of this QAPP in terms of work performed, accuracy, acceptable holding times, timely and understandable results and delivery process.
- h. Data management is being handled properly, i.e. data are entered on a timely basis, is properly backed up, is easily accessed, and raw data are properly stored in a safe place.
- i. Procedure for developing and reporting the results exists.

Monitoring Groups will be assessed on their ability to follow UWS procedures during field audits overseen by the Monitoring Program Quality Assurance Officer and Field Coordinator. The Monitoring Program Field Coordinator or designee will observe each monitoring group undertaking calibrations and field procedures once in May-June and follow up calls with Monitoring Groups will be scheduled after initial field audits. Field procedures will be reviewed from a set location on the water that does not need to be a UWS monitoring station. A dock or boat in a slip will be appropriate for these field audits. The CTDEEP and NYSDEC representative on this QAPP distribution list will be provided dates for field audits being held in embayments within their respective management areas. CTDEEP and NYSDEC staff have the option to attend

the field audits as observers. The Monitoring Program Quality Assurance Officer or designee will conduct a midseason check in call in August to all Monitoring Groups.

The Monitoring Program Coordinator shall confer with the UWS Science Advisors as necessary to discuss any problems that occur and what corrective actions are needed to maintain program integrity. In addition, the Monitoring Program Coordinator and UWS Science Advisors shall meet at the end of the sampling season, to review the draft report and discuss all aspects of the program and identify necessary program modifications for future sampling activities. All problems discovered and program modifications made shall be documented in the final version of the project report. If modifications require changes in the Quality Assurance Project Plan, these changes shall be submitted to the QAPP distribution list for review.

If data are found to be consistently outside the Data Quality Objectives as defined in section A.7. of this documents the Monitoring Program Coordinator shall review the program and correct problems as needed. Corrections may include retraining groups; rewriting sampling instructions; replacement of staff/Monitoring Group(s); alteration of sampling schedules, sites, stations or methods; or other actions deemed necessary. This information will be logged and maintained by the Monitoring Program Quality Assurance Officer. It will also be included in the QAPP Final Report.

C.2. REPORTS TO MANAGEMENT

Data that have passed the project quality assurance may be posted on the organization's web site, shared with the local media or at other venues (e.g. kiosks at recreation access sites), and submitted to the Long Island Sound Study, New England Interstate Water Pollution Control Commission, Interstate Environmental Commission, New York State Department of Environmental Conservation, New York City Department of Environmental Protection and/or Connecticut Department of Energy and Environmental Protection. A caveat will accompany these or any data released on a preliminary basis, explaining that they are for review purposes only and subject to correction after completion of a full data review occurring at the end of the sampling season.

The Monitoring Program Coordinator will write a final report. This will be sent to the distribution list on this QAPP. A final workbook of data from all embayments will accompany the report. The final report will also include (updated as necessary) any tables and graphs that were developed for initial data distribution efforts (i.e. the web site and media), and it will describe the program's goals, methods, quality control results, and recommendations. This report may also be used in public presentations.

All reports, preliminary or final, will include discussion of steps taken to assure data quality, findings on data quality, and decisions made on use, censorship, or flagging of questionable data. Any data that are censored in reports will be either referred to in this discussion, or presented but noted as censored.

In short, the final report will include:

- Raw data
- QC data
- Associated metadata
- Questionable data, flagged
- Identification of status as “preliminary” or “final” report

Table 17: Report Mechanisms, Responsibilities, and Distribution

Reporting Mechanism	Person Responsible for writing report	Distribution list
Monitoring Group Master Data Entry Template	Monitoring Group Lead or designated appointee	Monitoring Program Coordinator
Final Monitoring Report	Monitoring Program Coordinator	All signatories of this QAPP
Final Monitoring Data	Monitoring Program Coordinator	Signatories on this QAPP, EPA, NYS DEC, CTDEEP, NYCDEP, and other management groups

D. Data Validation and Usability

D.1. DATA REVIEW, VERIFICATION, AND VALIDATION

All project data, metadata, and quality control data shall be critically reviewed to look for problems that may compromise data usability.

Data collected before the 2018 season will be flagged as not being conducted under this QAPP when distributed. Save the Sound will be tasked with maintaining this QAPP in all aspects for the duration of the Unified Water Study.

The Monitoring Group Lead or designated appointee will review field data after each sampling run and take corrective actions as described in Table 16 of this document. At least once during the season, at the end of the season and if questions arise, the Monitoring Group Lead or designated appointee will share the data with the UWS Quality Assurance Officer to determine if the data appear to meet the objectives of the QAPP. Together, they will decide on any actions to take if problems are found.

D.2. VERIFICATION AND VALIDATION METHODS

All project data and metadata are reviewed and approved as usable data, or as un-usable data.

Data verification and validation will occur as described in Table 16, and will include checks on:

- Completion of all fields on data sheets; missing data sheets
- Completeness of sampling runs (e.g. number of stations visited / samples taken vs. number proposed, were all parameters sampled / analyzed)
- Completeness of QC checks (e.g. number and type of QC checks performed vs. number or type proposed)
- Number of samples exceeding QC limits for accuracy and precision and how far limits were exceeded.

D.3. RECONCILIATION WITH USER REQUIREMENTS

At the conclusion of the sampling season, after all in-season quality control checks, assessment actions, validation and verification checks and corrective actions have been taken, the resulting data set will be compared with the program's data quality objectives (DQOs) as defined in section A.7. This review will include, for each parameter, calculation of the following:

- Completeness goals: overall % of samples passing QC tests vs. number proposed.
- Percent of samples exceeding accuracy and precision limits.
- Average departure from accuracy and precision targets.

After reviewing these calculations, and taking into consideration such factors as clusters of unacceptable data (e.g. whether certain parameters, stations, dates, monitoring groups, etc. produced poor results), the Monitoring Program Coordinator, Quality Assurance Officer, and respective Monitor Group Lead will evaluate overall program attainment of DQOs and determine what limitations to place on the use of the data, or if a revision of the DQOs is allowable.

E. Appendices

Appendix A. Standard Operation Procedures

A-1 UWS SOP Station Selection.....	103
A-2 UWS SOP Tier I Sampling Plan.....	112
A-3 UWS SOP Depth and GPS.....	124
A-4 UWS SOP Sonde Profile.....	129
A-5 UWS SOP Filtered Chlorophyll.....	135
A-6 UWS SOP Qualitative Macrophytes.....	141
A-7 UWS SOP Filtered Nutrients.....	157
A-8 UWS SOP Total Nitrogen and Total Phosphorous.....	163
A-9 UWS SOP Continuous Dissolved Oxygen.....	169
A-10 UWS SOP Macrophytes Percentage Coverage via Camera.....	174

Appendix B. Data Forms, Checklists, and Chain of Custody Forms

B-1 Calibration Datasheet.....	181
B-2 Field Datasheet.....	183
B-3 Sample Event Datasheet.....	184
B-4 Chlorophyll Chain of Custody Form.....	185
B-5 Macrophyte Sampling Datasheet.....	186
B-6 Macrophyte Field Guide.....	187
B-7 Nutrient Chain of Custody Form.....	189
B-8 Nutrient Sample Event Datasheet.....	190
B-9 Logger Retrieval Sample Event Datasheet.....	191
B-10 Macrophyte Percent Coverage Datasheet.....	192

Appendix C. Project Laboratories - Analytical Method

C-1 Save the Sound, John and Daria Barry Foundation Water Quality Laboratory, SOP Determination of Chlorophyll a and Pheophytin a in Water.....	193
C-2 Interstate Environmental Commission Lab SOP Ammonia.....	200
C-3 Interstate Environmental Commission Lab SOP Nitrate, Nitrite, Nitrate+Nitrite and Total Nitrogen.....	213
C-4 Interstate Environmental Commission Lab SOP Total Phosphorous and Orthophosphate.....	227

Appendix A-1

UWS Standard Operating Procedure
Station Selection

January 29, 2018
Page 1 of 9

UWS Standard Operating Procedure – Station Selection

Table of Contents

1	POINT OF CONTACT	1
2	OBJECTIVE	1
3	OVERVIEW.....	2
4	SOURCES	3
5	MATERIALS AND EQUIPMENT.....	3
6	METHODS.....	3
6.1	Water Quality Station Selection.....	3
6.1.1	Assigning New Stations	4
6.1.2	Evaluating Existing Stations	5
6.2	Macrophyte Station Selection	5
6.3	GPS Coordinate Reference Station	5
7	TROUBLESHOOTING / HINTS.....	5
8	DATA PROCESSING AND STORAGE	5
9	REFERENCES	6
10	APPENDIX - LIST OF EMBAYMENT CODES.....	6

1 POINT OF CONTACT

NAME: Peter Linderoth, Save the Sound / CFE, Water Quality Program Manager
ADDRESS: 545 Tompkins Ave, 3rd Floor, Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

2 OBJECTIVE

Select sample locations to represent the entire embayment in a statistically sound manner. The Unified Water Study (UWS) will assign unique ID codes for the embayment, reporting sections of the embayment, and stations.

UWS Standard Operating Procedure Station Selection

January 29, 2018
Page 2 of 9

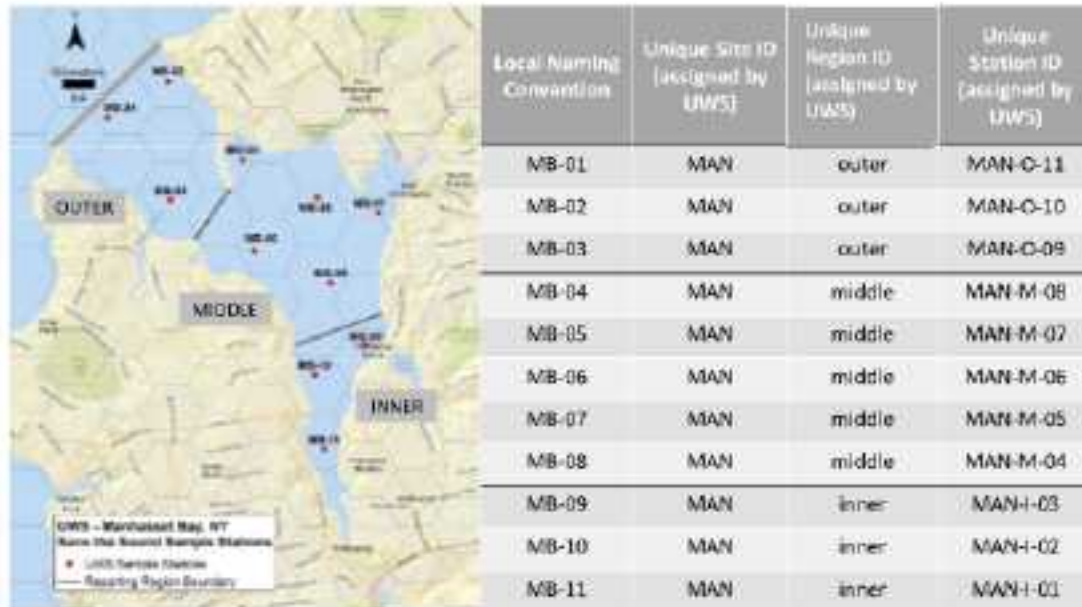


Figure 1: Visual representation of terminology. The grey line at the mouth of Manhasset Bay, NY represents the boundary of the embayment; this whole area is termed the SITE. Manhasset Bay includes three reporting regions, termed the Regions; the black lines show the divisions between the Inner, middle, and outer regions. Each region includes a minimum of three STATIONS, represented by the red dots. Monitoring groups may use their existing station IDs. When entering data into the UWS template and submitting data, the assigned UWS station ID must also be included.

3 OVERVIEW

Sample stations are selected to represent the water quality of the entire embayment using a probability-based sampling design [Paul et al., 2003], as in the EPA National Coastal Assessment (EPA, 2001). A UWS Scientific Advisor and UWS Monitoring Program Coordinator advise on the choice of station locations. Historic sites are evaluated for bias and inclusion in the UWS.

Station selection occurs only once, at the time of an embayment joining the UWS. Once stations are determined, these should be consistent in future years. Re-evaluation of stations in future years should be conducted under advisement with the UWS Monitoring Program Coordinator and the UWS Science Advisors.

Details on sampling frequency and timing during a day are covered in the UWS SOP Sampling Plan. The information below provides a general overview of the timing of sampling at the different stations.

4 SOURCES

This station selection design is based on methods used in the EPA National Coastal Assessment (EPA, 2001).

5 MATERIALS AND EQUIPMENT

- nautical charts of embayment or local knowledge of typical depth contours
- ArcGIS, access through the UWS Monitoring Program
- UWS subwatershed and embayment delineations for general reference
- hexagonal tessellation grid for use in ArcGIS

6 METHODS

6.1 Water Quality Station Selection

The location of stations will be determined or verified using a probability-based sampling design (Paul et al., 2003), as in the EPA National Coastal Assessment (EPA, 2001). This approach allows for the use of probability-based statistics in data analysis, without bias introduced by non-random station choice.

In this approach, a 0.42 km² hexagonal grid is overlain on the site map (Figure 2).

A minimum of four stations will be sampled in each embayment. If an embayment is divided into reporting regions, a minimum of three sites will be selected within each reporting region.

In cases where many hexagons are located in the embayment, sampling of some hexagons may be eliminated, if they are likely to be similar to neighboring stations. Exclusion of hexagons from the sampling plan should only be undertaken under advisement with the UWS Monitoring Program Coordinator and UWS Scientific Advisors. Nine stations per embayment are considered reasonable. Twelve stations and above are likely candidates for eliminating hexagons from the sampling plan. An example of eliminated hexagons is provided for Little Neck Bay, NY (Figure 2).

When a monitoring group has existing stations, those stations may be evaluated by the UWS Monitoring Program Coordinator and UWS Scientific Advisors for inclusion in the study.

UWS Standard Operating Procedure Station Selection

January 29, 2018
Page 4 of 9



Figure 2: Hexagons overlaying Little Neck Bay, NY. Numbers indicate station locations. Eight of the hexagons overlaying embayment water do not contain stations. This is an example of how some hexagons may be eliminated from sampling for larger embayments. Criteria for eliminating hexagons include an assessment of representability of neighboring stations, hexagons which include a large proportion of land, and stations closer to Long Island Sound likely to be well-mixed and similar to neighboring stations.

6.1.1 Assigning New Stations

Three stations are randomly generated for each hexagon targeted for sampling, identified in order of preference from low value to high. Stations will not be allowed in the center of navigational channels or in unnavigable waters. Stations will not be situated close to point sources of nutrients; this will be determined by consulting maps, local knowledge, and an on-site assessment on the first field day. The random generation of stations is conducted by the UWS Monitoring Program Coordinator in ArcGIS.

The final location of the station in each hexagon will be determined in the field. A trial sampling run is recommended to verify the stations chosen are accessible and can be sampled in the required time frame (for timeframe, see UWS SOP Sampling Plan). From the three random stations, a Monitoring Group may determine a station is unsuitable based on navigation considerations (too shallow, too much current to anchor, too rocky, etc.). If the first station is considered unsuitable, the field leader, UWS Monitoring Program Coordinator, and UWS Scientific Advisor, will evaluate the second, then third station. If all are deemed unacceptable, the collective group may choose a station within the hexagon that is close to the first random station with above considerations reviewed.

Once chosen, the station location is considered the permanent location and will not change on future sampling dates. Because the station will be permanent, the field leader should pay particular attention to the station depth in relation to the stage of the tide; the goal should be to have enough water at low tide to be able to sample the location.

UWS Standard Operating Procedure
Station Selection

January 29, 2018
Page 5 of 9

6.1.2 Evaluating Existing Stations

Existing stations in each hexagon will be evaluated for any potential biases that are present. These can include proximity to nutrient sources, disproportional sampling of beach sites, or other targeted sampling locations. Stations that are unbiased in a hexagon will be the first options for inclusion in the study. Stations deemed biased will not be included in the UWS. Monitoring

6.2 Macrophyte Station Selection

The number of stations sampled per embayment depends upon the size of the embayment, expected variability in macroalgae coverage, and access to sampling sites. An initial site visit of many locations can aid with determining the viability of sampling stations. A minimum of 3 sample stations is recommended for all embayments. More may be needed for larger embayments. The UWS Monitoring Program Coordinator, UWS Scientific Advisor, and Monitoring Group will consult to select these stations.

Station selection for macrophytes is complex. Refer to the UWS SOP Qualitative Macrophyte for this process. Macrophyte sampling occurs at different time scales and different stations than water quality parameters.

6.3 GPS Coordinate Reference Station

Select one land-based station to assess the precision of your GPS device. This station does not need to be located near the embayment being sampled. This is a land-based station that is easily accessible to the person with the GPS device.

The GPS coordinates at this station are read within 1 days of a field sampling day.

Each GPS unit or smart phone used for GPS must collect GPS coordinates form a land reference station.

7 TROUBLESHOOTING / HINTS

Station selection should be conducted in coordination with the UWS Monitoring Program Coordinator and UWS Scientific Advisors. A monitoring group is not expected to select stations without assistance.

8 DATA PROCESSING AND STORAGE

The UWS Monitoring Program Coordinator will be the custodian of the initial station selection maps and final maps determined after the first complete field day. The UWS Monitoring Program Coordinator will maintain a database which includes the unique site codes, section codes, and station codes for the embayment. Each unique station code will be affiliated with the corresponding GPS for the station.

The monitoring group is responsible for obtaining the GPS coordinates for each station during sample events.

The monitoring group is responsible for assuring that the correct unique station ID assigned by the UWS is properly matched with the local organizations station ID codes. Both codes (monitoring group's

UWS Standard Operating Procedure
Station Selection

January 29, 2018
Page 6 of 9

station code and UWS unique station ID) will be entered into the data template, along with the GPS coordinates.

9 REFERENCES

EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

Paul, J.F., J.L. Copeland, M. Charpentier, P.V. August, and J.W. Hollister. 2003, Overview of GIS applications in estuarine monitoring and assessment research. Marine Geodesy Journal 26: 63-72.

10 APPENDIX - LIST OF EMBAYMENT CODES

For record keeping purposes, each embayment is assigned a unique three letter code for the UWS. Contact the UWS Monitoring Program Coordinator to verify the names of specific waterbodies. Site names in this list may differ from the local names used for a site.

<i>sorted by RI to NY order</i>			<i>sorted by embayment name</i>		
Site Name	RI - NY order	UWS Site ID	Site Name	RI - NY order	UWS Site ID
Pawcatuck River, RI	1	PAW	Alewife Cove, CT	18	ALE
Little Narragansett Bay, CT	2	LNA	Ash Creek, CT	57	ASH
Wequetequock Cove, CT	3	WEQ	Baker Cove, CT	16	BAK
Quanaduck Cove, CT	4	QUA	Bebee Cove, CT	10	BEB
Stonington Harbor, CT	5	STO	Black Hall River, CT	28	BLH
Quiambog Cove, CT	6	QUI	Black Rock Harbor, CT	56	BLR
Wilcox Cove, CT	7	WXC	Branford Harbor, CT	45	BRA
Williams Cove, CT	8	WIL	Bride Brook, CT	25	BRB
Mystic River, CT	9	MYS	Bridgeport Harbor, CT	54	BRH
Bebee Cove, CT	10	BEB	Byram River, CT	82	BYR
West Cove, CT	11	WST	Calf Pen Meadow Creek, CT	50	CAL
Palmer Cove, CT	12	PAL	Captain Harbor, CT	81	CAP
Venetian Harbor, CT	13	VEN	Centerport Harbor, NY	106	CEN
Mumford Cove, CT	14	MUM	Clinton Harbor, CT	36	CLI
Poquonock River, CT	15	POQ	Cockenoe Harbor, CT	65	COC
Baker Cove, CT	16	BAK	Cold Spring Harbor, NY	103	COL
Thames River, CT	17	THA	Compo Cove, CT	62	COM
Alewife Cove, CT	18	ALE	Connecticut River, CT	29	CTR
Goshen Cove, CT	19	GOS	Conscience Bay, NY	112	CON
Jordan Cove, CT	20	JOR	Cove Harbor, CT	73	COV
Gardners Pond, CT	21	GAR	Darien River, CT	71	DAR
Niantic River, CT	22	NIR	Dosoris Pond, NY	99	DOS

UWS Standard Operating Procedure
Station Selection

January 29, 2018
Page 7 of 9

<i>sorted by RI to NY order</i>			<i>sorted by embayment name</i>		
Site Name	RI - NY order	UWS Site ID	Site Name	RI - NY order	UWS Site ID
Niantic Bay, CT	23	NIB	East River, NY	95	EAR
Pattagansett River, CT	24	PAT	Eastchester Bay, NY	92	EAB
Bride Brook, CT	25	BRB	Echo Bay, NY	90	ECH
Four Mile River, CT	26	FOU	Farm River, CT	47	FAR
Threemile River, CT	27	THR	Fence Creek, CT	38	FEN
Black Hall River, CT	28	BLH	Five Mile River, CT	68	FIV
Connecticut River, CT	29	CTR	Four Mile River, CT	26	FOU
South Cove, CT	30	SOU	Frost Creek, NY	100	FRO
Indiantown Harbor, CT	31	ITH	Gardners Pond, CT	21	GAR
Oyster River, Old Saybrook, CT	32	ORO	Goldsmith's Inlet, NY	116	GOL
Hagar Creek, CT	33	HAG	Gorham Pond, CT	70	GOR
Patchogue River, CT	34	PAC	Goshen Cove, CT	19	GOS
Menunkesucket River, CT	35	MEN	Greenwich Cove, CT	76	GRC
Clinton Harbor, CT	36	CLI	Greenwich Harbor, CT	80	GRH
Toms Creek, CT	37	TOM	Guilford Harbor, CT	39	GUI
Fence Creek, CT	38	FEN	Hagar Creek, CT	33	HAG
Guilford Harbor, CT	39	GUI	Hempstead Harbor, NY	98	HEM
Indian Cove, CT	40	INC	Holly Pond, CT	72	HOL
Sachem Head Harbor, CT	41	SAC	Housatonic River, CT	52	HOU
Joshua Cove, CT	42	JOS	Hunter Island Bay, NY	91	HIB
Island Bay, CT	43	ISL	Huntington Bay, NY	109	HUB
Little Harbor, CT	44	LIH	Huntington Harbor, NY	105	HUH
Branford Harbor, CT	45	BRA	Indian Cove, CT	40	INC
Pages Cove, CT	46	PAG	Indian Harbor, CT	78	INH
Farm River, CT	47	FAR	Indiantown Harbor, CT	31	ITH
New Haven Harbor, CT	48	NEW	Island Bay, CT	43	ISL
Oyster River, Milford, CT	49	ORM	Jordan Cove, CT	20	JOR
Calf Pen Meadow Creek, CT	50	CAL	Joshua Cove, CT	42	JOS
Milford Harbor, CT	51	MIF	Kirby Pond, NY	83	KIR
Housatonic River, CT	52	HOU	Larchmont Harbor, NY	88	LAR
Lewis Gut, CT	53	LEW	Lewis Gut, CT	53	LEW
Bridgeport Harbor, CT	54	BRH	Little Harbor, CT	44	LIH
Pequonnock River, CT	55	PEQ	Little Narragansett Bay, CT	2	LNA
Black Rock Harbor, CT	56	BLR	Little Neck Bay, NY	96	LNE
Ash Creek, CT	57	ASH	Lloyd Harbor, NY	104	LLO
Pine Creek, CT	58	PIN	Mamaroneck River, NY	87	MAM
Mill River, CT	59	MIL	Manhasset Bay, NY	97	MAN
Sasco Brook, CT	60	SAS	Mattituck Creek, NY	115	MAT
Sherwood Millpond, CT	61	SHE	Menunkesucket River, CT	35	MEN

UWS Standard Operating Procedure
Station Selection

January 29, 2018
Page 8 of 9

<i>sorted by RI to NY order</i>			<i>sorted by embayment name</i>		
Site Name	RI - NY order	UWS Site ID	Site Name	RI - NY order	UWS Site ID
Compo Cove, CT	62	COM	Mianus River, CT	77	MIA
Saugatuck River, CT	64	SAU	Milford Harbor, CT	51	MIF
Cockenoe Harbor, CT	65	COC	Mill Neck Creek, NY	101	MNC
Norwalk Harbor, CT	66	NWH	Mill River, CT	59	MIL
Sheffield Island Harbor, CT	67	SIH	Milton Harbor, NY	85	MIT
Five Mile River, CT	68	FIV	Mount Sinai Harbor, NY	114	MOU
Scotts Cove, CT	69	SCO	Mumford Cove, CT	14	MUM
Gorham Pond, CT	70	GOR	Mystic River, CT	9	MYS
Darien River, CT	71	DAR	New Haven Harbor, CT	48	NEW
Holly Pond, CT	72	HOL	Niantic Bay, CT	23	NIB
Cove Harbor, CT	73	COV	Niantic River, CT	22	NIR
Wescott Cove, CT	74	WES	Nissequogue River, NY	110	NIS
Stamford Harbor, CT	75	STA	Northport Bay, NY	108	NPB
Greenwich Cove, CT	76	GRC	Northport Harbor, NY	107	NPH
Mianus River, CT	77	MIA	Norwalk Harbor, CT	66	NWH
Indian Harbor, CT	78	INH	Oyster Bay, NY	102	OYB
Smith Cove, CT	79	SMI	Oyster River, Milford, CT	49	ORM
Greenwich Harbor, CT	80	GRH	Oyster River, Old Saybrook, CT	32	ORO
Captain Harbor, CT	81	CAP	Pages Cove, CT	46	PAG
Byram River, CT	82	BYR	Palmer Cove, CT	12	PAL
Kirby Pond, NY	83	KIR	Patchogue River, CT	34	PAC
Playland Lake, NY	84	PLA	Pattagansett River, CT	24	PAT
Milton Harbor, NY	85	MIT	Pawcatuck River, RI	1	PAW
Van Amringe Millpond, NY	86	VAN	Pequonnock River, CT	55	PEQ
Mamaroneck River, NY	87	MAM	Pine Creek, CT	58	PIN
Larchmont Harbor, NY	88	LAR	Playland Lake, NY	84	PLA
Premium Millpond, NY	89	PRE	Poquonock River, CT	15	POQ
Echo Bay, NY	90	ECH	Port Jefferson Harbor, NY	113	POR
Hunter Island Bay, NY	91	HIB	Premium Millpond, NY	89	PRE
Eastchester Bay, NY	92	EAB	Pugsley Creek, NY	94	PUG
Westchester Creek, NY	93	WCC	Quana duck Cove, CT	4	QUA
Pugsley Creek, NY	94	PUG	Quiambog Cove, CT	6	QUI
East River, NY	95	EAR	Sachem Head Harbor, CT	41	SAC
Little Neck Bay, NY	96	LNE	Sasco Brook, CT	60	SAS
Manhasset Bay, NY	97	MAN	Saugatuck River, CT	64	SAU
Hempstead Harbor, NY	98	HEM	Scotts Cove, CT	69	SCO
Dosoris Pond, NY	99	DOS	Sheffield Island Harbor, CT	67	SIH
Frost Creek, NY	100	FRO	Sherwood Millpond, CT	61	SHE
Mill Neck Creek, NY	101	MNC	Smith Cove, CT	79	SMI

UWS Standard Operating Procedure
Station Selection

January 29, 2018
Page 9 of 9

<i>sorted by RI to NY order</i>			<i>sorted by embayment name</i>		
Site Name	RI - NY order	UWS Site ID	Site Name	RI - NY order	UWS Site ID
Oyster Bay, NY	102	OYB	South Cove, CT	30	SOU
Cold Spring Harbor, NY	103	COL	Stamford Harbor, CT	75	STA
Lloyd Harbor, NY	104	LLO	Stonington Harbor, CT	5	STO
Huntington Harbor, NY	105	HUH	Stony Brook Creek, NY	111	SBC
Centerport Harbor, NY	106	CEN	Thames River, CT	17	THA
Northport Harbor, NY	107	NPH	Threemile River, CT	27	THR
Northport Bay, NY	108	NPB	Toms Creek, CT	37	TOM
Huntington Bay, NY	109	HUB	Van Amringe Millpond, NY	86	VAN
Nissequogue River, NY	110	NIS	Venetian Harbor, CT	13	VEN
Stony Brook Creek, NY	111	SBC	Wequetequock Cove, CT	3	WEQ
Conscience Bay, NY	112	CON	Wescott Cove, CT	74	WES
Port Jefferson Harbor, NY	113	POR	West Cove, CT	11	WST
Mount Sinai Harbor, NY	114	MOU	Westchester Creek, NY	93	WCC
Mattituck Creek, NY	115	MAT	Wilcox Cove, CT	7	WXC
Goldsmith's Inlet, NY	116	GOL	Williams Cove, CT	8	WIL

UWS Standard Operating Procedure – Tier I Sampling Plan

Table of Contents

1	POINT OF CONTACT	2
2	OBJECTIVE	2
3	DEFINITIONS AND ABBREVIATIONS	2
4	OVERVIEW	3
5	SOURCES	3
6	MATERIALS AND EQUIPMENT	3
6.1	Safety	3
6.2	Sampling Gear – All Stations	5
7	METHODS	6
7.1	Parameters to Sample	6
7.2	Timing of Sampling	6
7.2.1	Timing During the Year	6
7.2.2	Timing During a Sample Day	7
7.3	Sampling Depths	8
7.3.1	Temperature, Salinity, Dissolved Oxygen, Chlorophyll a, Turbidity	8
7.4	Required Replicates and Verification	8
7.5	Order of Events When Sampling a Water Quality Station	10
7.5.1	Prepare for Sampling Trip	10
7.5.2	Water Quality Station Sampling – within 3 hours of sunrise	10
7.5.3	End of Field Day	10
8	TROUBLESHOOTING / HINTS	10
9	DATA PROCESSING AND STORAGE	11
10	REFERENCES	11
11	Quick Sheet – Sampling Plan	12

1 POINT OF CONTACT

NAME: Peter Linderoth, Save the Sound / CFE, Water Quality Program Manager
ADDRESS: 545 Tompkins Ave, 3rd Floor, Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

2 OBJECTIVE

Sample water quality and macrophytes within an embayment following guidelines of the Unified Water Study (UWS). Frequency of sampling and daily order of events are specified.

3 DEFINITIONS AND ABBREVIATIONS

Embayment: A recess in a coastline or an indentation off a shoreline which forms a bay. In Long Island Sound, the names of embayments often include the words Harbor, River, Cove, Bay, Creek; with a few including the names Brook, Gut, Inlet, or Lake.

Field Team: Person or group of people working together to sample a station.

Macroalgae (macroalga, singular.): Commonly referred to as seaweed. This is a group of plant-like organisms. They do not have the vascular system and roots of a true plant. The “macro” prefix indicates these organisms are visible with the naked eye, no magnification is required to view the whole organism; although, magnification with a hand lens or loupe may be necessary to view the structure of the organism. In comparison, microalgae are the phytoplankton in the water which are too small to see with the naked eye.

Macrophyte: Plants and macroalgae that are viewable with the naked eye. This term includes macroalgae, seagrass (eelgrass, *Zostera marina*; widgeon grass, *Ruppia maritima*), and marsh grass.

Monitoring Group: The group conducting the field work.

Seagrass: A true plant, not an alga; they have the vascular system and roots of a land plant. These plants are fully submerged at all times (though there are a few species not found in Long Island Sound which are intertidal). Long Island Sound has two species of seagrass: *Zostera marina* (eelgrass), which is the most commonly seen seagrass in our area; and *Ruppia maritima* (widgeon grass).

Region: The reporting regions for the UWS. Each section must include a minimum of three stations. Sections will be assigned a unique name by the UWS; examples are included below.

Number of Sections	1	2	3
Name of Sections	whole	inner outer	inner middle outer
Abbreviations for Sections	W	I O	I M O

Site: The whole embayment, as defined by the UWS list. Each site has a unique three letter code assigned by the UWS; for example, Little Neck Bay, NY is “LNE”. Each site will have a minimum of four water quality stations.

Sonde: An instrument probe that automatically transmits information about its surroundings underground, under water, in the atmosphere, etc.

SOP: Standard operating procedure; this document is a SOP.

Station: The location where samples are collected, identified by a GPS location.

UWS: Unified Water Study

UWS Monitoring Program Coordinator: The person designated as the point of contact for the UWS.

UWS Scientific Advisor: Estuarine or water quality scientists designated as advisors to the UWS.

4 OVERVIEW

Sampling occurs in the months of May through October, though a shorter season is acceptable for inclusion in the UWS.

Two types of stations are included:

- 1) *Water quality stations* are sampled within three hours of sunrise between the months of May through October. A minimum of four stations per embayment are required. If multiple regions of the embayment are delineated, a minimum of three stations are required per region. Monitoring groups will plan to sample water quality stations at a minimum every two weeks. However, if unforeseeable circumstances make this plan not possible, a minimum of six sample events between May and October are required for inclusion in the UWS; including at least one sample event in each of the months of June, July, August, and September.
- 2) *Macrophyte stations* are land-based or boat-based. They are sampled only mid-summer and may be sampled on different days from the water quality stations. Sampling occurs on three separate days between July 15 and August 7. Two sample days or a date slightly outside of the date criteria may be sufficient in the event of unforeseen complications. This decision will be made by the Monitoring Program Quality Assurance Officer and UWS Science Advisor(s).

5 SOURCES

These procedures are based on the EPA Volunteer Estuary Monitoring Manual (EPA, 2007) and follows methods used in the EPA National Coastal Assessment (EPA, 2001). The EPA Volunteer Estuary Monitoring Manual (EPA, 2007) provides a wealth of specific data for monitoring groups. All groups should refer to the EPA manual for specific guidance. This SOP provides specific monitoring details relevant to the UWS.

6 MATERIALS AND EQUIPMENT

6.1 Safety

- safety plan – every volunteer should have a copy
 - Find out the location and telephone number of the nearest telephone and write it down, or have a cellular phone available.
 - Locate the nearest medical center and write down directions for guiding emergency

- personnel to your stations.
- Have each member of the sampling team complete a medical form that includes emergency contacts, and relevant health information such as allergies, diabetes, epilepsy, etc. **Please note** – this form should be kept confidential. Whoever coordinates the Monitoring Group efforts should have a copy and should review the form. A second copy should be kept with the field team in the event of an emergency; the form for each team member should be sealed in an envelope with their name.
 - Each team member should have contact information for all field team members. This list could be kept in your field box or use packing tape to affix it to the back of a clipboard or some other item you always have in the field.
 - full name
 - cell phone
 - home phone
 - email address
 - car color, make, model, and license plate
 - emergency contact information
 - Every monitoring team should have a shore-based check-in for each sampling day. This is a person who knows who is on the team, where they will be, and the time they are expected back. If the field team does not check in, the shore person should know the procedures to follow to report the team missing. The following is an example:
 - Call the cell phones and home phones of all team members.
 - Check the site for parked vehicles – if the vehicle is present, start a search.
 - If the team is considered missing at sea, contact the Coast Guard; if the team is missing on land, contact the police.
 - first aid kit – suggested items
 - telephone numbers of emergency personnel (e.g., police, ambulance service)
 - first aid manual which outlines diagnosis and treatment procedures
 - antibacterial or alcohol wipes
 - first aid cream or ointment
 - acetaminophen and ibuprofen for relieving pain and reducing fever
 - several band-aids
 - several gauze pads, 3 or 4 inches square
 - 2-inch roll of gauze bandage
 - triangular bandage
 - large compress bandage
 - 3-inch wide elastic bandage
 - needle for removing splinters
 - tick spoon for removing ticks
 - doctor-prescribed antihistamine for any participant who is allergic to bee stings
 - cell phone
 - water
 - appropriate shoes and clothes (and extras); protection from the sun
 - other items to consider:
 - flash light
 - back pack for sampling equipment, so hands are free

UWS Standard Operating Procedure
Tier I Sampling Plan

March 9, 2021
Page 5 of 12

6.2 Sampling Gear – All Stations

REQUIRED (PUT IN A SMALL TOOL BOX)

- site maps with stations on map
- list of UWS unique station IDs with GPS coordinates for the site
- clip boards
- pencils
- pencil sharpener
- permanent marker
- field data sheets
- Chlorophyll a labels

OPTIONAL, BUT VERY USEFUL (PUT IN A SMALL TOOL BOX)

- clear packing tape (can be used to cover writing on bottles or affix a label that is not sticking)
- extra batteries for any electrical sampling gear
- tailor's tape measure (seamstress measuring tape), with metric scale – comes in handy when you need to re-mark a line
- basic tools (pliers, wrench, screw drivers, etc.)
- plastic baggies
- scissors, pocket knife, nail clippers
- cable ties
- electrical tape (this tape works well underwater; comes in many colors)
- duct tape (very useful in many situations)

7 METHODS

7.1 Parameters to Sample

Monitoring groups must collect the following data to be included in the UWS:

- for each water quality station
 - GPS coordinates of stations, recorded each sample date
 - date and time
 - total water depth
 - 0.5 m below the surface, 0.5 m above bottom, mid-depth if total depth >10 m; if total depth is less than 1.5 m, only a mid-depth reading will be collected
 - temperature
 - conductivity (salinity)
 - dissolved oxygen
 - chlorophyll a
 - turbidity
 - at the last station per field day undertake a replicate profile including all parameters
- obtain from an online NOAA tide table and noted weather station:
 - time of high and low tide nearest time of sampling
 - high and low air temperature for 24 hours preceding field sampling
 - precipitation out a week preceding sampling event
- within 1 day of the field sampling day, read the GPS of a land-based reference station
- for each macrophyte station
 - GPS coordinates of stations
 - date and time
 - photos of macrophytes
- within 1 days of the field sampling day, read the GPS of a land-based reference station

7.2 Timing of Sampling

7.2.1 Timing During the Year

WATER QUALITY STATIONS

The target sampling frequency is two sampling events per month, May through October. Sampling dates should be 10 to 18 days apart.

All water quality parameters should be sampled during a sampling event. At one station per field day, conduct a replicate profile.

In the occurrence of equipment failure or other unforeseen difficulty, a minimum of six sample events between May and October are required for inclusion in the UWS; including at least one sample event in each of the months of June, July, August, and September.

UWS Standard Operating Procedure
Tier I Sampling Plan

March 9, 2021
Page 7 of 12

MACROPHYTE SURVEY SAMPLING FREQUENCY

Macrophyte surveys will occur between July 15 and August 7 of every sampling season, with 3 survey events total. One survey per week is best. If this is not possible, maximize the days between sampling. All three days cannot be sampled in the same 7-day window. 2 sampling events or one outside of the interval may be considered for inclusion in the UWS. This decision will be made by the Monitoring Program Quality Assurance Officer and UWS Science Advisor(s).

OVERVIEW OF SAMPLING FREQUENCY

The calendar below provides an example of sampling frequency throughout the season. Dates highlighted in orange are water quality sampling dates (e.g. May 11). The green text (July 15 – August 7) are potential dates for macrophyte sampling. The dates highlighted in green are the 3 planned macrophyte sampling dates (e.g. July 21).

MAY							JUNE						
S	M	T	W	T	F	S	S	M	T	W	T	F	S
1	2	3	4	5	6	7	29	30	31	1	2	3	4
8	9	10	11	12	13	14	5	6	7	8	9	10	11
15	16	17	18	19	20	21	12	13	14	15	16	17	18
22	23	24	25	26	27	28	19	20	21	22	23	24	25
29	30	31	1	2	3	4	26	27	28	29	30	1	2
5	6	7	8	9	10	11	3	4	5	6	7	8	9
JULY							AUGUST						
S	M	T	W	T	F	S	S	M	T	W	T	F	S
26	27	28	29	30	1	2	31	1	2	3	4	5	6
3	4	5	6	7	8	9	7	8	9	10	11	12	13
10	11	12	13	14	15	16	14	15	16	17	18	19	20
17	18	19	20	21	22	23	21	22	23	24	25	26	27
24	25	26	27	28	29	30	28	29	30	31	1	2	3
31	1	2	3	4	5	6	4	5	6	7	8	9	10
SEPTEMBER							OCTOBER						
S	M	T	W	T	F	S	S	M	T	W	T	F	S
28	29	30	31	1	2	3	25	26	27	28	29	30	1
4	5	6	7	8	9	10	2	3	4	5	6	7	8
11	12	13	14	15	16	17	9	10	11	12	13	14	15
18	19	20	21	22	23	24	16	17	18	19	20	21	22
25	26	27	28	29	30	1	23	24	25	26	27	28	29
2	3	4	5	6	7	8	30	31	1	2	3	4	5

7.2.2 Timing During a Sample Day

Sampling of dissolved oxygen occurs in the morning (within 3 hours of sunrise), to capture the lowest

dissolved oxygen values in the system.

Chlorophyll a and turbidity should be sampled at the same time as dissolved oxygen, to allow for linking of these parameters to their corresponding temperature, salinity, and dissolved oxygen (Tables 1).

Table 1: Text summary of required sampling times within a sample day

Parameter/Survey	Required time interval
Dissolved Oxygen (mg/l, % sat.)	Collected within 3 hours of sunrise
Temperature	Collected within 3 hours of sunrise (with dissolved oxygen)
Salinity	Collected within 3 hours of sunrise (with dissolved oxygen)
Chlorophyll a	Collected within 3 hours of sunrise (with dissolved oxygen)
Turbidity (NTU)	Collected within 3 hours of sunrise (with dissolved oxygen)
Total and Sample Station Depth	At time of sampling
Macrophyte Survey	close to low tide for wrackline survey, close to high tide for hardened shorelines, boat surveys any time of day

7.3 Sampling Depths

7.3.1 Temperature, Salinity, Dissolved Oxygen, Chlorophyll a, Turbidity

If total depth at the station is greater than 10 m, take three samples: one at 0.5 m below the surface, one at mid depth, and another at 0.5 m above the bottom.

If total depth at the station is less than 10 m, take two samples: one at 0.5 m below the surface and the other at 0.5 m above the bottom.

If total depth at the station is less than 1.5 m, take one sample halfway to the bottom and record as mid depth.

Always record the depth of the sample.

7.4 Required Replicates and Verification

During a field day, use the field data sheet as a reminder for the number of replicates required for each parameter.

Table 2: Required replicates, blanks, and verification readings:



UWS Standard Operating Procedure
Tier I Sampling Plan

March 9, 2021
Page 9 of 12

parameter & technique	replicates required	verification and/or blank
GPS coordinates	1 reading per station at 1 station per day, take a second reading before leaving the station	read a land-based reference station within 2 days of the field sampling day
sampling with multiparameter sonde	1 reading at each depth, wait for reading to stabilize before recording at 1 station per day (typically the last station), do two replicate profiles – do one complete profile, then do a second	verify depth by lowering sonde to known depth verify Chlorophyll a by reading water in a bucket and filtering 2 samples from the bucket read standards before and after a sample day (salinity, oxygen in 100% water saturated air, turbidity, chl-a); day before and after sample day is acceptable
filtered Chlorophyll a	2 filtered samples of a single water sample collected from 0.5 m below the surface at all stations	duplicate from same bucket used as a second calibration point for Chlorophyll a for a sampling event
macrophytes	3-6 stations per embayment	photos of each rake toss or the beach being sampled, reviewed by UWS Monitoring Program Coordinator and Science Advisors

7.5 Order of Events When Sampling a Water Quality Station

7.5.1 Prepare for Sampling Trip

- A. Calibrate all instruments.
- B. Gather all field supplies.
- C. Complete the pre-sampling event portions of the UWS datasheet.
- D. Arrange for a shore person. This is someone to check on you if you don't come back at the designated time.

7.5.2 Water Quality Station Sampling – within 3 hours of sunrise

- A. Record station information on the data sheet. *Be sure to complete all sections of the data sheet completely, for every data entry.*
- B. Obtain total depth of the station, determine sampling depths for water quality parameters
- C. Collect profile data using the multiparameter sonde.
 - a. Sample the surface, then mid-depth, then bottom with Sonde; or depths necessary according to total depth and respective sample depth criteria, sampling from top to bottom.
- D. At one station per day repeat measurements where only one profile is typically collected. The last station of the day is the ideal choice in terms of time management.
 - a. Repeat the sonde profile. Complete the first profile, sampling at each depth. Complete a second profile.
 - b. Record the depth and GPS coordinates a second time, just before leaving the station.
- E. Read a calibration standard just following calibration and following the field trip to verify readings, for all parameters where this applies.

7.5.3 End of Field Day

- A. If using a multiparameter sonde to sample Chlorophyll a, place the sonde in a bucket of estuarine water, read the sonde, and filter samples for Chlorophyll a analysis for verification of the sonde readings.
- B. Verify all sections of the data sheet have been completed.
- C. Store Chlorophyll a samples according to the UWS SOP Filtered Chlorophyll.
- D. Within 1 days of sampling, record the GPS of a reference station on land to verify the accuracy and precision of the GPS coordinates.

8 TROUBLESHOOTING / HINTS

- Gather field equipment the day prior to sampling. Check the field equipment in the morning, before

UWS Standard Operating Procedure
Tier I Sampling Plan

March 9, 2021
Page 11 of 12

you head out into the field. Creating a checklist is very helpful in prepping for your field day. Include personal items (sunscreen, bug spray, etc.) and safety equipment on the checklist.

- Always carry a copy of this SOP and the relevant parameter-specific SOPs.
- Print out the “quick sheets” for relevant SOPs to use as a reminder in the field. Do not laminate these as you will want to add notes. A plastic page-protector taped close can be used to keep these sheets dry.

9 DATA PROCESSING AND STORAGE

The UWS Monitoring Program Coordinator will be the custodian of the finalized data files.

The monitoring group is responsible for obtaining data, entering data into the UWS data template, and delivering the data to the UWS Monitoring Program Coordinator.

The monitoring group is responsible for assuring that the correct unique station ID assigned by the UWS is properly matched with the local organizations station ID codes. Both codes (monitoring group’s station code and UWS unique station ID) will be entered into the data template, along with the GPS coordinates.

10 REFERENCES

EPA, 2007, Volunteer Estuary Monitoring, A Methods Manual, Second Edition. Orhrel Jr., R.L., Register, K.M. (Eds.). The Ocean Conservancy & EPA. 396 p.
https://www.epa.gov/sites/production/files/2015-09/documents/2007_04_09_estuaries_monitoruments_manual.pdf

EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

11 Quick Sheet – Tier I Sampling Plan

Check your equipment list the day before a field day and again on the morning of a field day.

Calibrate instruments the day before sampling.

Read a calibration standard just following calibration and following the sampling event to verify readings, for all parameters where this applies.

Bring the relevant SOPs and quick sheets with you into the field.

Remember – rinse equipment in sample water before collecting a sample.

Water quality stations - Sample depths are shown in chart to the right.

Use the field data sheet as a reminder for the number of replicates required for each parameter. Sample 1 station / day as a replicate for sonde profiles, GPS, and depth.

For filtered Chlorophyll *a*, filter four samples per day with corresponding sonde reading.

depth < 1.5 m	1.5 m < depth < 10 m	depth > 10 m	Depth (m)
★ mid-depth	★ 0.5 m below	★ 0.5 m below	1
			2
			3
			4
			5
		★ mid-depth	6
			7
			8
	★ 0.5 m above		9
			10
		★ 0.5 m above	11
			12

MACROPHYTE STATIONS – use quick sheet from Macrophyte SOP

GPS REFERENCE CHECK

- Within 1 days of sampling, record the GPS of a reference station on land to verify the accuracy and precision of the GPS coordinates.

WATER QUALITY STATIONS (sample within 3 hours of sunrise)

- At one station per day repeat measurements where only one profile is typically collected. The last station of the day is a good choice in terms of time management. (sonde, GPS and depth)

Appendix A-3

UWS Standard Operating Procedure
Depth & GPS Coordinates

January 29, 2018
Page 1 of 5

UWS Standard Operating Procedure – Depth and GPS Coordinates

Table of Contents

1	POINT OF CONTACT	1
2	OBJECTIVE	1
3	OVERVIEW.....	1
4	SOURCES	2
5	MATERIALS AND EQUIPMENT.....	2
6	METHODS.....	2
6.1	Preparation	2
6.2	Field Collection and Processing.....	3
6.3	Sample Storage	4
6.4	Laboratory Analysis.....	4
7	TROUBLESHOOTING / HINTS.....	4
8	DATA PROCESSING AND STORAGE	4
9	REFERENCES.....	4
10	Quick Sheet – Depth & GPS Coordinates	5

1 POINT OF CONTACT

NAME: Peter Linderoth, Save the Sound / CFE, Water Quality Program Manager
ADDRESS: 545 Tompkins Ave, 3rd Floor, Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

2 OBJECTIVE

Determine the total depth and sample depth at station. Determine the GPS coordinates of the station.

3 OVERVIEW

A graduated line is weighted and used for recording total depth (m) and sample depth (m) at a station. Acceptable alternatives include multiparameter sonde depth reading.

A GPS unit is used to obtain GPS coordinates for the station on each field day. A Smart Phone app is an acceptable alternative as backup. GPS position of your actual location at each station for each sample date is required. A second reading is taken at one station per embayment per sample day to assess precision. If unforeseen circumstances make recording GPS points impossible and dead reckoning is used, your data may still be used as part of the UWS but will be flagged and reviewed in the quality assurance process.

4 SOURCES

These procedures are based on the EPA Volunteer Estuary Monitoring Manual (EPA, 2007). The EPA Volunteer Estuary Monitoring Manual (EPA, 2007) provides a wealth of specific data for monitoring groups.

5 MATERIALS AND EQUIPMENT

- Depth sampling device. Options include:
 - weighted, graduated line, with divisions every 0.1 m
 - This can be the cable attached to a multiparameter sonde
 - multiparameter sonde with depth sensor
- GPS coordinate sampling device. Options include:
 - Handheld GPS unit
 - Apps are available for your smart phone that allow you to easily log and return to GPS coordinates. The accuracy of the GPS on your phone is dependent on Wi-Fi signals and cellular signals.
 - Dead reckoning to get yourself on station should only be used if your GPS options have failed. This will need to be included in field notes for the sampling event. Data will be reviewed in the quality assurance process.

6 METHODS

6.1 Preparation

- Weighted line:
 - Examine the line for wear or damage.
 - Measure the increments to ensure line has not stretched or shrank.
- If using a multiparameter sonde:
 - Calibrate the depth sensor prior to the sampling event and the morning of on the sampling vessel according to the manufacturer's instructions.

UWS Standard Operating Procedure
Depth & GPS Coordinates

January 29, 2018
Page 3 of 5

➤ GPS device

- Check batteries
- Within 1 day of the sampling trip, read the GPS coordinates of a reference station on land (refer to "UWS Station Selection SOP" for more details on this reference station). These readings will be used to assess the precision of your GPS method.

6.2 Field Collection and Processing

METERED LINE

1. Attach weight to line.
 - a. A weighted sonde attached to a cable with 0.1 m increments is considered a weighted line.
2. Slowly lower line until the weight reaches the bottom.
3. Mark the line at the surface of the water (pinch with finger).
4. Determine the depth of the station.
5. Record max depth to the nearest 0.1 m on field data sheet.
6. Following UWS SOP Sampling Plan for number of samples and locations in the water column associated with total depth, sample all target depths using steps 2-5 replacing "weight reaches the bottom" to deliver sonde to target depth using increments on cable.
7. At least once per day per embayment (usually the last station of the day), take the reading again.

MULTIPARAMETER SONDE

1. Slowly lower until the weighted sonde reaches the bottom.
2. Determine the depth of the station by reading the display.
 - a. Add the length from the depth sensor to the weight for total depth.
3. Record max depth to the nearest 0.1 m on field data sheet.
4. Following UWS SOP Sampling Plan for number of samples and location in the water column associated with total depth, sample all target depths using steps 1-3 replacing "weight reaches the bottom" to deliver sonde to target depth using handheld display. If depth sensor is not located with sonde attachments, compensate by lowering the sonde to the depth reading required. For example, if the sonde attachments are .25 m below the depth sensor and the target sample depth is 5 m the sampler should lower sonde to a depth reading 4.75 m on the display.
5. At least once per sampling day (prior to first station), lower the sonde to a known depth and verify the reading. If the depth reading from sonde does not agree within 0.3 m, use the metered line procedures.
6. At least once per sampling day per embayment (usually the last station of the day), take the reading again.

UWS Standard Operating Procedure
Depth & GPS Coordinates

January 29, 2018
Page 4 of 5

GPS DEVICE

1. Determine the GPS coordinates of the station by reading the display.
2. At least once per sampling day per embayment (usually the last station of the day), take the reading again.
3. Within 1 day of the sampling trip, read the GPS coordinates of a reference station on land.

6.3 Sample Storage

Not applicable.

6.4 Laboratory Analysis

Not applicable.

7 TROUBLESHOOTING / HINTS

- Always carry a copy of this SOP and the relevant parameter-specific SOPs.
- Print out the “quick sheets” for relevant SOPs to use as a reminder in the field. Do not laminate these as you will want to add notes. A plastic page-protector taped close can be used to keep these sheets dry.

8 DATA PROCESSING AND STORAGE

The monitoring group is responsible for obtaining data, entering data into the UWS data template, and delivering the data to the UWS Monitoring Program Coordinator.

The monitoring group is responsible for assuring that the correct unique station ID assigned by the UWS is properly matched with the local organizations station ID codes. Both codes (monitoring group’s station code and UWS unique station ID) will be entered into the data template, along with the GPS coordinates.

9 REFERENCES

EPA, 2007, Volunteer Estuary Monitoring, A Methods Manual, Second Edition. Orhrel Jr., R.L., Register, K.M. (Eds.). The Ocean Conservancy & EPA. 396 p.
https://www.epa.gov/sites/production/files/2015-09/documents/2007_04_09_estuaries_monitoruments_manual.pdf

UWS Standard Operating Procedure
Depth & GPS Coordinates

January 29, 2018
Page 5 of 5

10 Quick Sheet – Depth & GPS Coordinates

For depth and GPS - At least once per sampling day per embayment (usually the last station of the day), take the reading again.

Preparation

- If using a weighted line:
 - Examine the line for wear or damage.
 - Measure the increments to ensure line has not stretched or shrunk.
- If using a multiparameter sonde:
 - Calibrate the depth sensor according to the manufacturer's instructions. This may be done in the field.
 - Bring a weighted line (can be sonde cable) for verification of readings (prior to sampling first station) and in case sonde depth is inaccurate.
- GPS device
 - Check batteries
 - Within 1 days of the sampling trip, read the GPS coordinates of a reference station on land. This will be used to assess the precision of your GPS method.

Field Collection and Processing

METERED LINE

1. Attach weight to line.
 - a. A weighted sonde attached to a cable with 0.1 m increments is a weighted line.
2. Slowly lower line until the weight reaches the bottom.
3. Mark the line at the surface of the water (pinch with finger).
4. Determine the depth of the station.
5. Record max depth and sample depth(s) to the nearest 0.1 m on field data sheet.

MULTIPARAMETER SONDE

1. Determine the depth of the station by reading the display.
2. Record max depth and sample depth(s) to the nearest 0.1 m on field data sheet.
3. At least once per sampling day (prior to first station), use a metered line to confirm the accuracy of the depth reading. If the depth reading from sonde does not agree within 0.3 m, use the metered line.

GPS DEVICE

1. Determine the GPS coordinates of the station by reading the display.
2. Within 1 days of the sampling trip, read the GPS coordinates of a reference station on land.

Appendix A-4

UWS Standard Operating Procedure
Sonde Water Quality Profile

January 29, 2018
Page 1 of 6

UWS Standard Operating Procedure – Sonde Profile

Table of Contents

1	POINT OF CONTACT	1
2	OBJECTIVE	1
3	OVERVIEW	2
4	SOURCES	2
5	MATERIALS AND EQUIPMENT	2
6	METHODS	2
6.1	Preparation	2
6.2	Field Collection and Processing	3
6.2.1	Sampling at a Station	3
6.2.2	Verification of Water Quality Parameters and Sonde Depth	4
6.2.3	Verification of Chlorophyll a – Bucket Sample	4
6.3	Sample Storage	4
6.4	Laboratory Analysis	4
7	TROUBLESHOOTING / HINTS	5
8	DATA PROCESSING AND STORAGE	5
9	REFERENCES	5
10	Quick Sheet – Sonde Water Quality Profile	6

1 POINT OF CONTACT

NAME: Peter Linderoth, Save the Sound / CFE, Water Quality Program Manager
ADDRESS: 545 Tompkins Ave, 3rd Floor, Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

2 OBJECTIVE

Determine the sample depth (m), temperature (°C), salinity (ppt), dissolved oxygen concentration (mg/L = ppm), and dissolved oxygen percent (% sat.) in the surface (0.5 m below surface) and bottom water (0.5 m above bottom) at the station using a multiparameter sonde. When total water depth is greater than 10 m, parameters at mid-depth will also be determined.

3 OVERVIEW

A multiparameter sonde is used to determine the sample depth (m), temperature (°C), salinity (ppt), dissolved oxygen concentration (mg/L = ppm), and dissolved oxygen percent (% sat.) in the surface (0.5 m below surface) and bottom water (0.5 m above bottom) at the station. When total water depth is greater than 10 m, parameters at mid-depth will also be determined. Multiparameter sondes may also include probes for turbidity (NTU) and chlorophyll a (RFU, µg/L). Frequency of sampling and daily order of events are specified in the "UWS Sampling Plan SOP."

At one station per day per embayment (typically the last station), do two replicate profiles – do one complete profile, then do a second. This allows for the precision of the sonde to be assessed.

4 SOURCES

These procedures are based on the EPA Volunteer Estuary Monitoring Manual (EPA, 2007) and follows methods used in the EPA National Coastal Assessment (EPA, 2001). The EPA Volunteer Estuary Monitoring Manual (EPA, 2007) provides a wealth of specific data for monitoring groups. All groups should refer to the EPA manual for specific guidance.

5 MATERIALS AND EQUIPMENT

- Eureka Manta +35 or YSI EXO1 Multiparameter Sonde, setup to record:
 - depth
 - temperature
 - conductivity (salinity)
 - dissolved oxygen
 - chlorophyll a
 - turbidity
- 5 gallon bucket
- See UWS SOP Filtered Chlorophyll. The field team will need all equipment and material listed for collecting chlorophyll a samples. Filter two samples per field day for verification of sonde readings.
- ASTM Type I reagent grade water or equivalent for 0 NTU turbidity standard solution and 0 µg/L (and 0 RFU for YSI EXO1) for chlorophyll a calibration
- YSI Turbidity Standard, 607300 (124 FNU on YSI)(100 NTU on Eureka) (YSI SKU: 607300)
- YSI Conductivity Calibrator Solution, 3169: 50,000 µS/cm (YSI SKU: 060660)

6 METHODS

6.1 Preparation

- Check the battery power. Pack extra batteries in the field tool kit.
- Calibrate the sonde the day before or morning of each field day. Follow the manufacturer's instructions for calibration steps.

UWS Standard Operating Procedure
Sonde Water Quality Profile

January 29, 2018
Page 3 of 6

- Conductivity – Conduct calibration using the 50,000 $\mu\text{S}/\text{cm}$ conductivity standard.
- Dissolved Oxygen – Conduct an air calibration, using water saturated air for calibrating 100% saturation.
- Chlorophyll a - Conduct a one-point calibration using the 0 $\mu\text{g}/\text{L}$ (ASTM Type I reagent grade water)
- Turbidity - Conduct a two-point calibration using the 0 NTU (ASTM Type I reagent grade water) and 124 FNU (100 NTU Eureka instrument) standard solutions.
- Depth – Calibrate the depth following manufacturer's guidelines.
- The instrument will be factory calibrated and serviced at the manufacturer's recommended interval.
- After each calibration, take a verification reading in each of the standard solutions. This can be done directly after the calibration with solution still in the calibration cup.
- Record calibration information in the calibration section of the UWS datasheet. The following information will be entered into the UWS data entry template:
 - dissolved oxygen - post calibration reading in 100% saturated air
 - conductivity standard for sonde - manufacturer and value ($\mu\text{S}/\text{cm}$)
 - conductivity standard for sonde - lot number
 - conductivity standard for sonde - expiration date
 - conductivity on sonde - post calibration reading in standard ($\mu\text{S}/\text{cm}$)
 - chlorophyll a fluorescence on sonde - post calibration reading ($\mu\text{g}/\text{L}$ on Eureka and $\mu\text{g}/\text{L}$, RFU on YSI)
 - turbidity standard - manufacturer and value (FNU/NTU)
 - turbidity standard - lot number
 - turbidity standard - expiration date
 - turbidity - post calibration reading in standard (FNU/NTU)

6.2 Field Collection and Processing

6.2.1 Sampling at a Station

- Turn on the sonde.
- If sampling from a boat, anchor the boat or stay on station with short engine adjustments.
- Determine sampling depths as described in the UWS SOP Sampling Plan.
- If wading in to sample, take extra precaution to not stir up sediment. Samples should be taken from an area that is not disturbed from the sampler's approach.
- Lower sonde to sample depth. Be sure that the probe ends are at the desired depth, versus other sections of the sonde. Be careful not to let the probe contact the sediments as this could foul the

UWS Standard Operating Procedure
Sonde Water Quality Profile

January 29, 2018
Page 4 of 6

probes and give inaccurate readings.

- Allow the sonde to equilibrate for at least 90 seconds. Watch the salinity first. Once salinity is steady, watch the dissolved oxygen. The value may fluctuate in response to real variability in the water column. You are waiting for the readings to hover around a value, versus consistently rising or falling. Do not wait for turbidity or chlorophyll a to stabilize, these probes respond quickly to real variability in the water column. These values should be listed as the middle point between the high and low displayed back and forth on the screen.
- Record readings on the field data sheet.
- At one station per day per embayment (typically the last station), do a replicate profile – do one complete profile, then do a second. This allows for the precision of the sonde to be assessed.

6.2.2 Verification of Water Quality Parameters and Sonde Depth

- Take and record sonde values in standards up to a day before and after a sampling day (conductivity, oxygen in 100% water saturated air, 0 µg/L chl-*a* and 124 NTU (YSI) or 100 NTU (Eureka) turbidity).
- Verify sonde depth by lowering sonde to a known depth following the procedures listed in UWS SOP Depth SOP.

6.2.3 Verification of Chlorophyll a – Bucket Sample

If a chlorophyll a probe is on the sonde, verify the chlorophyll a readings from the sonde once per day and send to an analytical lab for analysis. This may be conducted at a dock; it does not need to occur at a water quality station.

- Rinse the bucket three times with surface water.
- Fill the bucket with surface water.
- Place the sonde in the bucket and stir for 10 seconds, allow to equilibrate, record chlorophyll a reading.
- Filter two samples from the bucket following the procedures in the UWS SOP Filtered Chlorophyll. Prior to withdrawing a sample from the bucket, be sure to stir the water in the bucket, using the sonde.

6.3 Sample Storage

Not applicable for the sonde. Follow instructions in the UWS SOP Filtered Chlorophyll SOP for filtered chlorophyll a samples.

6.4 Laboratory Analysis

Not applicable for the sonde. Follow instructions in the UWS SOP Filtered Chlorophyll SOP for filtered chlorophyll a samples.

7 TROUBLESHOOTING / HINTS

- Always make sure to wait until readings have stabilized before calibrating or logging a reading.
- If the probe does not appear to be operating or calibrating properly, contact the Monitoring Group Lead to make a decision on proceeding with the sampling event. Immediately notify the UWS Monitoring Program Coordinator if a sonde continues to not operate properly.

8 DATA PROCESSING AND STORAGE

The UWS Monitoring Program Coordinator will be the custodian of the finalized data files.

The monitoring group is responsible for obtaining data, entering data into the UWS data template, and delivering the data to the UWS Monitoring Program Coordinator.

The monitoring group is responsible for assuring that the correct unique station ID assigned by the UWS is properly matched with the local organizations station ID codes. Both codes (monitoring group's station code and UWS unique station ID) will be entered into the data template, along with the GPS coordinates.

9 REFERENCES

EPA, 2007, Volunteer Estuary Monitoring, A Methods Manual, Second Edition. Orhrel Jr., R.L., Register, K.M. (Eds.). The Ocean Conservancy & EPA. 396 p.
https://www.epa.gov/sites/production/files/2015-09/documents/2007_04_09_estuaries_monitoruments_manual.pdf

EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

10 Quick Sheet – Sonde Water Quality Profile

Preparation

- Check the battery power. Pack extra batteries in the field tool kit.
- Calibrate the sonde the day before each field day. Follow the manufacturer's instructions for calibration.
- After calibrating, use the sonde to take a verification reading in each of the standard solutions.
- Record calibration information in the calibration section of the UWS datasheet.

Field Collection and Processing

- Turn on the sonde.
- If sampling from a boat, anchor the boat or stay on station with short engine adjustments.
- Determine sampling depths as described in the UWS SOP Sampling Plan.
- If wading in to sample, take extra precaution to not stir up sediment. Sample should be taken from an area that is not disturbed from the sampler's approach.
- Lower sonde to sample depth. Be sure that the probe ends are at the desired depth, versus other sections of the sonde. Be careful not to let the probe contact the sediments as this could foul the probes and give inaccurate readings.
- Allow the sonde to equilibrate for at least 90 seconds.
 - Watch the salinity first.
 - Once salinity is steady, watch the dissolved oxygen.
 - The value may fluctuate in response to real variability in the water column. You are waiting for the readings to hover around a value, versus consistently rising or falling. This will typically take about one minute, but may take longer. Do not wait for turbidity or chlorophyll a to stabilize, these probes respond quickly to real variability in the water column.
- Record readings on the field data sheet.
- At one station per day per embayment (typically the last station), do a replicate profile – do one complete profile, then do a second. This allows for the precision of the sonde to be assessed.

Verification of Water Quality Parameters and Sonde Depth

- Take and record sonde values in standards at the start and end of a sampling day (salinity, oxygen in 100% water saturated air, and turbidity).
- Verify sonde depth by lowering sonde to a known depth following the procedures listed in UWS SOP Depth GPS.

Verify chlorophyll a by sampling in a bucket.

Verify the chlorophyll a readings from the sonde once per day and send to an analytical lab for analysis. This may be conducted at a dock; it does not need to happen at a station.

- Rinse the bucket three times with surface water, fill the bucket with surface water.
- Place the sonde in the bucket, stir for 10 seconds, and then allow to equilibrate, record chlorophyll a reading.
- Filter two samples from the bucket following the procedures indicated in the UWS SOP Filtered Chlorophyll. Prior to withdrawing a sample from the bucket, be sure to stir the water in the bucket with sonde.

Appendix A-5

UWS Standard Operating Procedure
Filtered Chlorophyll a

January 29, 2018
Page 1 of 6

UWS Standard Operating Procedure – Filtered Chlorophyll a

Table of Contents

1	POINT OF CONTACT	1
2	OBJECTIVE	1
3	OVERVIEW.....	1
4	SOURCES	2
5	MATERIALS AND EQUIPMENT.....	2
6	METHODS.....	2
6.1	Preparation	2
6.2	Field Collection and Processing.....	3
6.3	Sample Storage	4
6.4	Laboratory Analysis.....	4
7	TROUBLESHOOTING / HINTS.....	4
8	DATA PROCESSING AND STORAGE	4
9	REFERENCES.....	5
10	Quick Sheet – Filtered Chlorophyll a.....	6

1 POINT OF CONTACT

NAME: Peter Linderoth, Save the Sound / CFE, Water Quality Program Manager
ADDRESS: 545 Tompkins Ave, 3rd Floor, Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

2 OBJECTIVE

Determine the concentration of chlorophyll a in the surface water, 0.5 m below the surface, following guidelines of the Unified Water Study (UWS). Frequency of sampling and daily order of events are specified in the UWS SOP Sampling Plan.

3 OVERVIEW

A water sample is collected, filtered, and analyzed at an analytical lab to determine the chlorophyll a

UWS Standard Operating Procedure
Filtered Chlorophyll a

January 29, 2018
Page 2 of 6

concentration using the fluorometric technique (90% acetone extraction). Phytoplankton, the microscopic plant-like organisms living in the water, contain chlorophyll a. Thus, chlorophyll a concentration provides a rough approximation of the amount of phytoplankton in the water column. This is considered a rough approximation because individual phytoplankton contain varying amounts of chlorophyll a based on species, size, and environmental factors.

At each station, a chlorophyll a sonde reading is collected from 0.5 m below the surface according to the UWS SOP Sonde Profile. Once per sampling event a bucket of surface water is collected. A sonde reading for chlorophyll a and two water samples from the bucket are filtered through a glass fiber filter with pore size of 0.7 μm . The filters are dried by passing air through the filter. The filters are stored in the dark on ice until it can be transferred to a freezer (within 12 hours of the initial sample collection).

4 SOURCES

These procedures are based on the EPA Volunteer Estuary Monitoring Manual (EPA, 2007) and follows methods used in the EPA National Coastal Assessment (EPA, 2001). The EPA Volunteer Estuary Monitoring Manual (EPA, 2007) provides a wealth of specific data for monitoring groups.

Analytical methods follow EPA method #445.0, "*In Vitro* Determination of Chlorophyll a and Pheophytin a in Marine and Freshwater Algae by Fluorescence" (Arar and Collins, 1997) or Standard Methods 10200 H3, "Fluorometric Determination of Chlorophyll a" (Eaton et al., 2005).

5 MATERIALS AND EQUIPMENT

- 5 gallon bucket or larger
- Glass fiber filters (Whatman GF/F with nominal pore size of 0.7 μm), 2.5 cm diameter
- Filter holders, 2.5 cm: Pall brand, 25 mm Easy Pressure Syringe Filter Holder, Delrin Plastic
- 60 mL syringe
- Forceps for handling filters
- Unused and unbleached coffee filter cut into strips
- Aluminum foil
- Airtight container
- Cooler

6 METHODS

6.1 Preparation

- Check that field equipment is prepped and operational.

UWS Standard Operating Procedure
Filtered Chlorophyll a

January 29, 2018
Page 3 of 6

- Prepare labels for filters. The pre-labeled information should include the monitoring group, UWS unique station ID, date, sample type, sample volume, and replicate ID ("a" or "b"). For example:
 - "Save the Sound LNE-I-01a 6/8/17 Chl-a Vol. 60 mL" & "Save the Sound LNE-I-01b 6/8/17 Chl-a Vol. 60 mL"
- Prep aluminum foil squares for projected number of samples (2 per reference station) plus extras.
- Load all available filter holders with filters. Have at least 4 ready for use.

6.2 Field Collection and Processing

1. Collect water in 5 gallon bucket.
 - Rinse the bucket three times with sample water from the surface. Fill bucket with sample water.
2. If not prepped, load filters into the filter holder.
 - Preloading all filter holders with filters before the sampling event begins is strongly advised.
3. Rinse the 60 mL syringe with ~5 mL of sample water. Repeat for a total of three rinses.
4. Mix the sample well by stirring for 10 seconds with sonde.
5. Record chlorophyll a data from sonde.
6. Fill the 60 mL syringe with sample water taken close to where sonde reading was recorded. Note the volume of water in the syringe – knowing the exact volume filtered is critical to analyzing the chlorophyll a concentration.
7. Connect the filter holder to the syringe.
8. Gently expel water through the filter. The pressure applied to the plunger of the syringe should be slow and steady. If you encounter more than mild resistance, the filter may be clogged (e.g. jellyfish, algae) or you may have filtered sufficient sample.
9. After filtering 60 mL, inspect the filter for color. If a visible light green or light brown color is present, there is sufficient material on the filter for analysis, skip to step 10.
 - If no color is visible, filter another 60 mL and re-inspect.
 - On some occasions, if chlorophyll a levels are very high, filtering may become difficult after as little as 30 mL. If this occurs, and the filter is colored, stop filtering and proceed to step 10.
10. Record volume recorded on field data sheet and sample label when possible. Recording the exact volume filtered is critical to analyzing the chlorophyll a concentration. If you are uncertain of the volume filtered, discard the sample and start over.
11. Use the 60 mL syringe to dry the filter by expelling air through the filter.
 - With the syringe NOT attached to the filter holder, draw air into the syringe.

UWS Standard Operating Procedure
Filtered Chlorophyll a

January 29, 2018
Page 4 of 6

- Attach syringe to holder and expel air forcefully through the filter. Do this until no “mist” is aspirated from the filter holder, a minimum of 3 times.
- DO NOT draw air backwards through the filter. Syringe should be taken off holder each time plunger is drawn up.

12. Remove filter with forceps.

- If forceps are dirty, wipe with a Kimwipe or rinse with distilled water.
- Only touch the filter on the edges. You may use clean fingers to steady the filter by touching the edge (do not touch the green or brown part).
- Fold filter in half with forceps and place it in an absorbent pad (unbleached coffee filter). Wrap in aluminum foil by folding the foil around the filter.

13. Filters should be placed in an airtight container and stored in a cooler on ice. Ice should surround the airtight container.

6.3 Sample Storage

14. Store samples in the freezer (-20°C). Frozen samples must be analyzed within 28 days.

6.4 Laboratory Analysis

15. Bring filters to partner lab for analysis or transport group; using UWS Chain of Custody. Frozen samples must be analyzed within 28 days.

7 TROUBLESHOOTING / HINTS

- Make sure filters are dry before storing.
- Gather field equipment the day prior to sampling. Check the field equipment in the morning, before you head out into the field.
- Always carry a copy of this SOP and the relevant parameter-specific SOPs.
- Print out the “quick sheets” for relevant SOPs to use as a reminder in the field. Do not laminate these as you will want to add notes. A plastic page-protector taped close can be used to keep these sheets dry.

8 DATA PROCESSING AND STORAGE

The monitoring group is responsible for obtaining data, entering data into the UWS data template, and delivering the data to the UWS Monitoring Program Coordinator.

The monitoring group is responsible for assuring that the correct unique station ID assigned by the UWS is properly matched with the local organizations station ID codes. Both codes (monitoring group's station code and UWS unique station ID) will be entered into the data template, along with the GPS coordinates.

UWS Standard Operating Procedure
Filtered Chlorophyll a

January 29, 2018
Page 5 of 6

9 REFERENCES

- Arar, E.J., and Collins, G.B. 1997. U.S. Environmental Protection Agency Method 445.0, *In vitro* determination of Chlorophyll a and pheophytin a in marine and freshwater algae by fluorescence, revision 1.2: Cincinnati, Ohio, U.S. Environmental Protection Agency National Exposure Research Laboratory, Office of Research and Development.
- Eaton, A. D., M. A. H. Franson, American Public Health Association., American Water Works Association., and Water Environment Federation. 2005. Standard methods for the examination of water and wastewater, 21st edition. New York: American Public Health Association.
- EPA, 2007, Volunteer Estuary Monitoring, A Methods Manual, Second Edition. Orhrel Jr., R.L., Register, K.M. (Eds.). The Ocean Conservancy & EPA. 396 p.
https://www.epa.gov/sites/production/files/2015-09/documents/2007_04_09_estuaries_monitoruments_manual.pdf
- EPA, U.S. 2001. National Coastal Assessment: Field Operations Manual. U. S. Environmental Protection Agency, Office of Research and Development, National Health and Environmental Effects Research Laboratory, Gulf Ecology Division, Gulf Breeze, FL. EPA 620/R-01/003. 72 p.

10 Quick Sheet – Filtered Chlorophyll a

SAMPLE from 5 gallon Bucket

Recording the exact volume filtered is critical to analyzing the chlorophyll a concentration. If you are uncertain of the volume filtered, discard the sample and start over.

The labels should include the UWS unique station ID, monitoring group, replicate ID (“a” or “b”), sample volume, and date. For example:

- station 1: Save the Sound LNE-I-01a 6/8/17 Vol. 60 mL &
Save the Sound LNE-I-01b 6/8/17 Vol. 60 mL

1. Collect surface water in a 5 gallon bucket. Rinse bucket three times with sample water prior to filling.
2. Mix the sample well by stirring for 10 seconds with sonde. Record data from sonde
3. Rinse the 60 mL syringe with ~5 mL of sample water. Repeat for a total of three rinses.
4. ***Note the volume of water in the syringe – knowing the exact volume filtered is critical to analyzing the chlorophyll a concentration.***
5. Gently expel sample water through the filter. The pressure applied to the plunger of the syringe should be slow and steady. If you encounter more than mild resistance, the filter may be clogged (e.g. jellyfish, algae) or you may have filtered sufficient sample.
 - After filtering 60 mL, inspect the filter for color. If a visible light green or light brown color is present, there is sufficient material on the filter for analysis. If no color is visible, filter another 60 mL and re-inspect. On some occasions, if chlorophyll a levels are very high, filtering may become difficult after as little as 30 mL.
6. ***Recording the exact volume filtered is critical to analyzing the chlorophyll a concentration. If you are uncertain of the volume filtered, discard the sample and start over.***
7. Use the 60 mL syringe to dry the filter by expelling air through the filter.
 - Attach syringe to holder and expel air forcefully through the filter. Do this until no “mist” is aspirated from the filter holder, a minimum of 3 times. DO NOT draw air backwards through the filter. Syringe should be taken off holder each time plunger is drawn up.
8. Remove filter with forceps.
 - Only touch the filter on the edges. You may use clean fingers to steady the filter by touching the edge (do not touch the green part).
 - Fold filter in half with forceps, place in an absorbent pad (unbleached coffee filter), wrap in aluminum foil and apply the label.
9. Labeled, foil wrapped filters should be placed in an airtight container and stored in a cooler on ice. Ice should surround the airtight container.
10. Store samples in the freezer (-20°C). Bring filters to partner lab for analysis. Frozen samples must be analyzed within 28 days.

Appendix A-6

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 1 of 16

UWS Standard Operating Procedure – Qualitative Macrophytes

Table of Contents

1	POINT OF CONTACT	1
2	OBJECTIVE	1
3	OVERVIEW	2
4	METHODS	3
4.1	Station Selection	3
4.1.1	Considerations for Choosing a Station	3
4.1.2	Station Selection Guidance for Soft Shoreline (Beaches)	4
4.1.3	Guidance for Submerged Areas (Sampling from Boats, Docks, Jetties, etc.)	4
4.1.4	Examples of Sample Locations	5
4.2	Preparation	6
4.3	Field Collection and Processing	6
4.3.1	Option 1: Soft Shoreline Station (Beaches)	7
4.3.2	Option 2: Submerged Station	9
4.4	Sample Storage	13
4.5	Laboratory Analysis	13
5	TROUBLESHOOTING / HINTS	13
6	DATA PROCESSING AND STORAGE	13
7	REFERENCES	14
8	Quick Sheet – Qualitative Macrophytes	15

1 POINT OF CONTACT

NAME: Peter Linderoth, Save the Sound / CFE, Water Quality Program Manager
ADDRESS: 545 Tompkins Ave, 3rd Floor, Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

2 OBJECTIVE

Identify problematic macroalgae blooms and the presence of eelgrass in neighboring waters by examining the wrack line on a beach or sampling submerged areas with a rake.

3 OVERVIEW

The goal is to identify areas with problematic amounts of macroalgae and areas with eelgrass (a beneficial condition). This is not intended to be a quantitative assessment. Macrophyte surveys will complement chlorophyll *a* concentrations to better understand the dominant primary producer in the system. Methods for conducting an assessment vary by the location to be assessed. If a soft shoreline is available, photos of the beach area are used to characterize the amount of macroalgae in the neighboring water. Alternatively, a rake is used to assess what is on the bottom when submerged areas are sampled.

METHOD OVERVIEW (please see text below for a full explanation)

- Sample 3 days during the 3-week period starting July 15 and ending August 7. Try to sample once per week. If this is not possible, please sample such that you maximize the days between sampling. All three days cannot be sampled in the same 7-day window. Two sample days or a date slightly outside of the date criteria may be sufficient in the event of unforeseen complications. This decision will be made by the Monitoring Program Quality Assurance Officer and UWS Science Advisor(s).
- For soft shorelines, take an overview photo of the area, focused on the beach (not the water). Take a close-up photo of each of the dominant macroalgae types on the beach. Complete the datasheet.
- For sampling submerged areas, toss the rake 6 times, 2 tosses in each of 3 directions. Photograph each rake, even empty rakes. Complete the datasheet.
- Enter the data into the online portal. Upload your photos to the online portal.

BACKGROUND

Later on in this SOP, you will see that you are characterizing macroalgae by morphology (hair-like, twig/leaf-like, sheets) and color (green, other). Macroalgae can be hard to identify, thus we are not attempting to identify the genus or species of algae present. The morphology will allow us to identify algae that are nutrient-loving, and thus may represent a water quality problem. Hair-like algae of all colors tend to be nutrient-lovers. Some branching algae in the twig/leaf-like category are also nutrient-lovers, especially if they fall into the red algae category. Sheets of green or red are problematic, while sheets of brown (e.g. kelp) can be indicative of good water quality. Based on these assessments, you may notice that color of the algae is important. Macroalgae fall into three taxonomic divisions denoted by the color of the algae: red (*Rhodophyta*), green (*Chlorophyta*), and brown (*Phaeophyta*). These colors refer to the types of photosynthetic pigments present in the organisms of each division. While a *Rhodophyta* often looks reddish, it can appear gold, green, brown, or black. The same is true of the other divisions. Green macroalgae typically appear green. Almost all green algae found in Long Island Sound is indicative of high nutrient availability. Thus, we ask that you distinguish between green algae and those of other colors. Good photos of the algae you are seeing will aid in follow-up analysis of these data.

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 3 of 16

MATERIALS AND EQUIPMENT

- GPS unit
 - GPS unit or GPS app for a smart phone
- Digital camera with resolution >5 megapixels (most smartphone cameras meet this criteria)
- Bucket or plastic bin, for putting macroalgae in if it is very muddy or falling off rake
- Bow rake with ~7 m (~20 feet) of rope attached to the tine-end of the rake and attached to the handle (cable tie, duct tape, etc.).
 - The rake should be a heavy duty bow rake with forged steel rakehead, 16 inches wide, with 15 or 16 tines. The total length (handle to tines) should be around 60 inches. A rake with an ash wood handle is appropriate; however, fiberglass or another handle material is also acceptable.
- Weight that can be attached to the rake, to help it sink to the bottom. For example, a dive weight or large fishing weights.
- Local tidal information for survey site (for soft shoreline surveys).



4 METHODS

4.1 Station Selection

The number of stations sampled per embayment depends upon the size of the embayment, expected variability in macroalgae coverage, and access to sampling areas. An initial site visit of many locations can aid with determining the viability of sampling stations. A minimum of 3 sample stations is recommended for all embayments. More may be needed for larger embayments. Consultation with the Monitoring Program Coordinator and UWS Science Advisors will aid with the determination of how many to include. **We cannot assess how much macroalgae is in an embayment using this method.** Instead, the goal is to identify embayments which have widespread macroalgae problems and those which have problem areas for macroalgae but not throughout the whole system. Additionally, eelgrass will be noted when present in the survey.

4.1.1 Considerations for Choosing a Station

Must be legally accessible. This includes areas with public access and private areas where permission has been granted by the owner or manager of the area. Roadways closed to pedestrian traffic and all train tracks ARE NOT considered areas with public access; prior permission and official escort are typically required for these areas.

Must be safe to access. The station should not endanger personnel. For example, areas to avoid include: steep inclines, roadways with insufficient buffers from passing cars, areas overrun with poison ivy, areas with excessive debris, etc.

Look for areas where macroalgae is likely to collect. The goal is to assess what is out in the system. If you

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 4 of 16

choose an area that is too deep or where the current is sweeping away all of the macroalgae, you may miss an issue that is occurring in shallower, calmer waters. However, if the overall system is deep and scoured by the current or wind driven waves, then sample these areas, as they do indeed represent the system. To assess how representative an area is, try sampling in a few spots in a general area of the embayment. If macroalgae is widespread, make a note of this. If the macroalgae you are sampling is at a spot where macroalgae tends to collect, also make a note of this.

4.1.2 Station Selection Guidance for Soft Shoreline (Beaches)

For soft shoreline assessments, the method calls for taking a photo of the beach and close-ups of the dominant macroalgae types. This is the quicker assessment when compared to a rake survey. You are looking for an area with the following characteristics:

- A beach composed of sand or gravel.
- Avoid marshes (though you may use a rake from a marsh).
- Avoid rocky intertidal zones.
- Areas with attached brown macroalgae (rockweed, knotted wrack) are fine if the area also collects free-floating macroalgae.
- An area you know is **not raked** or otherwise maintained by municipalities or local residents.

4.1.3 Guidance for Submerged Areas (Sampling from Boats, Docks, Jetties, etc.)

For sampling submerged areas, the method calls for tossing in a garden rake attached to a rope and slowly and steadily pulling in the rake. Two tosses are conducted in three directions, for a total of six tosses. A photo is taken of each rake, including when only mud is collected. Potential areas include:

- Marinas, docks, or jetties can be suitable locations, if their depth is representative of the overall area. Even deep boat slips can be acceptable as macroalgae tends to collect in the deeper locations. Avoid areas where prop scour is likely to push macroalgae away. If your rake toss looks similar to what you see in shallower areas, the areas should be suitable. *Note – when sampling from a dock, be considerate of private property: avoid areas where your actions may be misconstrued as potentially damaging to boats. In other words, choose an area of the dock where you have some room to move, not an area with little space between boats. Always check in with the office or property owner before sampling, to let them know you have permission to sample and so they know who you are and what you are doing.*
- Causeways, seawalls, and shorelines without a beach deemed safe to access are suitable sampling locations. Avoid areas scoured clean by the water flow (i.e., the mouth of a culvert). Be sure to choose areas where you will not catch brown macroalgae attached to rocks at the shoreline (rockweed, knotted wrack). We are interested in the beneficial shoreline macroalgae.
- Boat sampling is acceptable if the depth of the area is representative of the embayment. In other words, you should not sample in the deepest or the shallowest locations.




4.1.4 Examples of Sample Locations

When sampling from a dock, jetty, causeway, or other hardened shoreline, choose three locations and do two rake tosses from each location (red lines indicate direction and length of rake toss). Note that sampling is not done from the finger docks in the image below; these docks can be too skinny to safely accommodate two people and allow for proper framing of the photograph.




UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 6 of 16

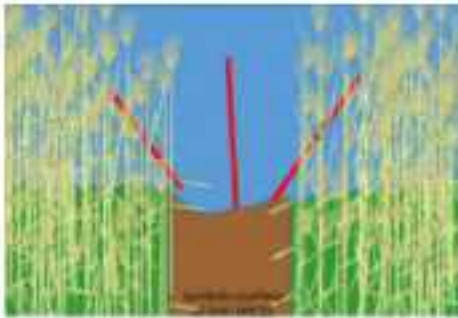

If you are in a location which inhibits moving around, do two rake tosses in each of three directions. For example, a kayak launch amidst a dense stand of *Phragmites* can provide limited access to the shoreline. In this case, two rake tosses are directed straight out into the water, two upstream, and two downstream.

When sampling from a boat at anchor, we assume the boat is shifting slightly with the wind and current. You may do all six tosses from the same location in the boat, tossing the rake in different directions. However, if you feel that you are sampling the same location six times, you will need to shift to a new location on the boat between tosses. If your boat is drifting, you may toss in the same direction each time, as you will be floating over new substrate.

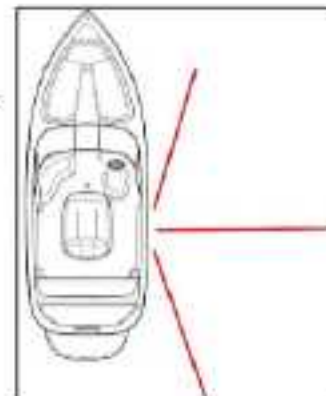

4.2 Preparation

- Ensure battery power and available storage is sufficient on camera or smartphone.
- Ensure battery power is sufficient on GPS unit or smartphone.
- Collect equipment and ensure all is in working order.
- Always carry a copy of this SOP and the relevant parameter-specific SOPs.
- Print out the "quick sheets" for relevant SOPs to use as a reminder in the field. Do not laminate these as you will want to add notes. A plastic page-protector taped close can be used to keep these sheets dry.

4.3 Field Collection and Processing

Sampling at each station should be quick. In field trials, the UWS Science Advisors completed sampling at a station in 10 to 15 minutes, though getting a rake hung up on a rock can increase the time needed. Soft shoreline (beach) areas can go even quicker. Please keep in mind this is a qualitative survey, so

spending additional time to quantify the macroalgae is not necessary. You are making a call of “not much here”, “wow, that’s a lot of macroalgae”, or “this could be a problem, but I’m not sure.” The first station may take a bit longer, as you get acquainted with the datasheet.

4.3.1 Option 1: Soft Shoreline Station (Beaches)

- Conduct the assessment within 3 hours of low tide (this gives you a 6-hour window).
- Complete the station information on the datasheet.
 - **IMPORTANT:** Record GPS points as decimal degrees (e.g. 41.56437 N, -71.87634 W). Please do not report GPS in hours and minutes. For example, a location with decimal degrees of 41.56437 N, -71.87634 W can also be reported in hours and minutes as 41° 33' 51.7314", -71° 52' 34.8234". Note that the numbers are different, decimal degrees are not just leaving out the symbols for hours (°) and minutes ('). We ask for decimal degrees because hours and minutes do not translate well in a spreadsheet. If you do not understand this distinction, please ask an advisor to explain.
- Record the amount of macroalgae on the beach as “none”, “some”, or “lots”.
 - None – essentially nothing on the beach, only attached macroalgae is present.
 - Some – ranges from a small bit of macroalgae on the beach to large wracklines of macroalgae present on the beach. This is qualitative, the Monitoring Program Coordinator and UWS Science Advisors will review when comparing stations, so do not worry too much about what truly constitutes some versus lots.
 - Lots – covers > 50% of the beach. May be a thin covering, in thick wracklines, or a thick layer covering the whole beach.
- Examine the macroalgae and use the guide to identify macroalgae morphology (hair-like, twig/leaf like, sheets) and color (green, other). You only need to record macroalgae that constitutes more than 10% of the macroalgae present on the beach.
- Take two overview photos, showing the overall appearance of the shoreline. One should include upland areas for context, the next should be slightly closer, showing macroalgae (or lack of macroalgae) on the beach (examples below).
 - Zoom in on the overview photo and confirm the beach is in focus (versus something in the background or foreground).
 - Do not submit any overview photos deemed unacceptable.
 - The two overview photos should be sufficient. However, if lighting or size of the station requires, you may take multiple photos.
 - Record an identifier for the photo on the datasheet. For a camera, this may be the filename. For a smartphone, record the time associated with the photo; this can be located by viewing the photo, the time the photo was taken should be on the screen.

UW5 Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 8 of 16

Example overview photos:



Wide scale overview shot of station X.

Closer overview shot of station X.



Wide scale overview shot of station Y.

Closer overview shot of station Y.

- For the dominant macroalgae types, take a close-up photo of each macroalgae.
 - This is a rough assessment of the dominant macroalgae types; you DO NOT need to photograph each type of macroalgae present on the beach.
 - Add something to the photo to provide a size reference. This could be a ruler, your foot, a brick, your keys, etc. It needs to be something with a known size (not a stick you pick up, which can be any size).
 - Zoom in on the macroalgae photo and confirm the macroalgae is in focus (versus something in the background or foreground).
 - Be sure that the lighting allows for realistic colors in the photo.
 - The photo should not include the shadow of the photographer.
 - Record an identifier for the photo on the datasheet. For a camera, this may be the filename. For a smartphone, record the time associated with the photo; this can be located by viewing the photo, the time the photo was taken should be on the screen.

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 9 of 16

Example close-up photo of macroalgae:



Close-up of macroalgae at station X.

Close-up of macroalgae at station Y.

4.3.2 Option 2: Submerged Station

- *Note* – a YouTube video demonstrating how to properly toss a rake is available at: <https://www.youtube.com/watch?v=yMqp5aA62VW> or find the video by searching for "macroalgae rake toss."
- The timing of sampling is not critical. The only criterion is that low tide has not exposed the sediment at the sampling station.
- Complete the station information on the datasheet. An example of a completed datasheet is included on page 14 of this SDP.
 - **IMPORTANT:** Record GPS points as decimal degrees (e.g. 41.56437 N, -71.87634 W). Please do not report GPS in hours and minutes. For example, a location with decimal degrees of 41.56437 N, -71.87634 W can also be reported in hours and minutes as 41° 33' 51.7314", -71° 52' 34.8234". Note that the numbers are different, decimal degrees are not just leaving out the symbols for hours (') and minutes ("). We ask for decimal degrees because hours and minutes do not translate well in a spreadsheet. *If you do not understand this distinction, please ask the UWS Monitoring Program Coordinator to explain.*
- Select a sampling station at your location.
 - If the station allows, conduct 2 rake tosses at each of 3 different spots. An example would be moving to three different areas on a dock. See Section 4.1.4 (page 5) for diagrams of where to sample.
 - If the station does not allow you to move around, conduct 2 tosses in each of 3 different directions. An example would be an opening in a marsh that allows access to the embayment. The rake is tossed upstream, downstream, and straight out into the water. See Section 4.1.4 for diagrams of where to sample.

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 10 of 16

- Confirm that the line is securely attached to the rake. Place your foot on the end of the rope, to keep it from going into the water.
- Face the embayment and hold the rake such that your dominant hand is at the end of the rake (ready to push the rake).
- Push the rake out into the water a distance of ~2 rake lengths. The motion is similar to an underhand slow pitch, rather than a javelin throw.
- If the toss is too short or too long, bring back and try again.
- The tines of the rake should be downward. If the rake lands with tines up, bring back and try again.
- After a successful toss, allow rake to settle on the floor then **slowly and steadily** pull the line so the teeth of the rake drag on the embayment bottom.
 - If the rake gets hooked on a rock or other object, try taking up the slack in the rope and pulling up on the rake, or moving to the left and to the right. If you feel that macroalgae was likely dislodged, try tossing again. For some rocks, freeing the rake is easy enough that macroalgae is mostly intact on the rake. For these cases, you do not need to toss again.
 - When removing the rake from the water, do so such that macroalgae does not fall off of the rake.
 - Make certain that you DO NOT pick up macroalgae removed from the rake on a previous toss. Be aware of where you discard macroalgae from the rake so that you are not capturing the same macroalgae with each toss.
- Record the amount of macroalgae on the rake as “none”, “some”, or “lots” (using sample photos shown below to assist).
 - None – no macroalgae on the rake.
 - Some – some tines are still bare.
 - Lots – all tines hold macroalgae. This includes anything from a thin coat of a slimy macroalgae to thick clumps of a wiry or leafy macroalgae.
 - If you get a lot of mud, put into a bucket or bin and rinse; mud that clumps on the rake is often being held together by fine macroalgae
- Take a photo of the rake (and macroalgae) after each toss.
 - Include all of the rake tines and macroalgae in the photo. If a portion of the rake tines is not included, take another photograph. You will want to be about 4 feet from the rake. **Do not get too close** – some cameras do not focus at shorter distances. With the resolution required, zooming in to look at macroalgae can be done on the computer. You do not need to get super close to get a good shot of the macroalgae, further back is better.
 - Frame the shot such that the light is behind the photographer (or at a slight angle), so that the light is good and the photographer’s shadow is not cast on the macroalgae.
 - Be aware of the background and change your location for the photo as needed. Light colors in the background make the macroalgae look darker; avoid light backgrounds (e.g. a white boat).
 - Zoom in on the photo and confirm the macroalgae is in focus (versus something in the background or foreground).


UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 11 of 16

- Be sure that the lighting allows for realistic colors in the photo.
- Record an identifier for the photo on the datasheet. For a camera, this may be the filename. For a smartphone, record the time associated with the photo; this can be located by viewing the photo, the time the photo was taken should be on the screen.
- Sort through the macroalgae and use the guide to identify macroalgae morphology (hair-like, twig/leaf like, sheets) and color (green, other). You only need to record macroalgae that constitutes more than 10% of the macroalgae present on the rake.

Example photos of rakes:

Station 1: Wequetequock Cove, Stonington Marina, 7/13/2016

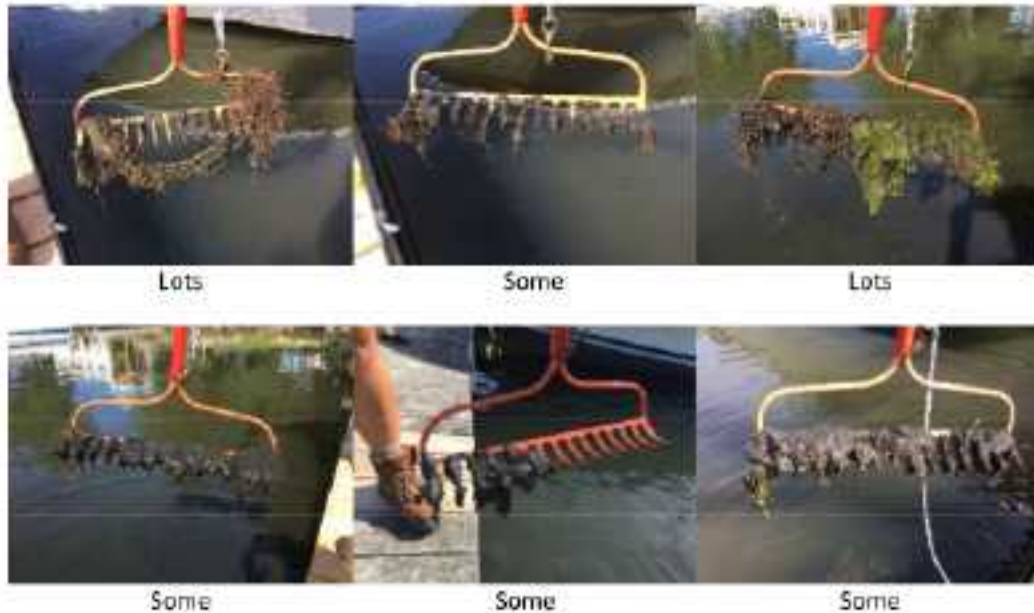

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 12 of 16

Station 2: Wequetequock Cove, kayak launch near Saltwater Farm Vineyard, 7/13/2016



Station 3: Wequetequock Cove, Elihu Island causeway, 7/13/2016



Close-up of dominant macroalgae (all beach looked like this photo, only 1 photo necessary).

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 13 of 16

The macroalgae collected on the soft shoreline at Elihu Island causeway is characterized as "lots." To verify that it represents the area, we also tossed a rake off the causeway, about midway between the mainland and the island. Each toss had "lots," confirming the beach survey was indeed representative. Eelgrass is visible on these rakes and was visible on the beach. The red branching macroalgae on the rakes was not a large component of what was on the beach. It is fine if the macroalgae composition varies, we are most interested in the amount of macroalgae.



4.4 Sample Storage

Not applicable.

4.5 Laboratory Analysis

Not applicable.

5 TROUBLESHOOTING / HINTS

- Organize all volunteers well before sampling events.
- Check tide information before initiating the surveys.
- Bring backup rake and line, if possible.
- Always carry a copy of this SOP and the relevant parameter-specific SOPs.
- Print out the "quick sheets" for relevant SOPs to use as a reminder in the field. Do not laminate these as you will want to add notes. A plastic page-protector taped close can be used to keep these sheets dry.

6 DATA PROCESSING AND STORAGE

The monitoring group is responsible for obtaining data, entering data into the UWS data template, and delivering the data to the UWS Monitoring Program Coordinator.

The monitoring group is responsible for assuring that the correct unique station ID assigned by the UWS is properly matched with the local organizations station ID codes. Both codes (monitoring group's station code and UWS unique station ID) will be entered into the data template, along with the GPS coordinates.

Macroalgae assessment data and photos will be submitted via an online portal. This will allow for consistent formatting in the data, easy transference and a storage system for the photos, and immediate access to the photos for the Monitoring Program Coordinator and UWS Scientific Advisors. We will be using a free software called KoBo Toolbox. Once uploaded by UWS participants, the data will be downloaded and stored by Save the Sound.

- Each station within your embayment will use a different form.
- Not all data on the datasheet is required in the form.
- Bookmark the form in your browser, this will make it easier to access.
- You may save an incomplete form and return to the form at a later time.
- Directions for accessing the form are included below:

How to Collect Data on a Mobile Device

Collect data on your Android device with KoBoCollect

To start collecting data with the [KoBoCollect Android app](#) you need to enter the following URL into the URL field in the KoBoCollect settings:

<https://kc.kobotoolbox.org/cfests>

Click [here](#) to read more about getting started with KoBoCollect



OR

Collect Data on Any Device in Your Browser

For iPhones or any other device that has a Web browser, use the following link to enter data - even offline:

<https://ee.kobotoolbox.org/x/11Y81e>

(Or if you have a barcode scanner on your phone, just scan the barcode on the right)



7 REFERENCES

None

8 Quick Sheet – Qualitative Macrophytes

Method Overview

- Sample 3 days during the 3-week period starting July 15 and ending August 7. Try to sample once per week. If this is not possible, please sample such that you maximize the days between sampling. All three days cannot be sampled in the same 7-day window. Two sample days or a date slightly outside of the date criteria may be sufficient in the event of unforeseen complications. This decision will be made by the Monitoring Program Quality Assurance Officer and UWS Science Advisor(s).
- For soft shorelines, take an overview photo of the area, focused on the beach (not the water). Take a close-up photo of each of the dominant macroalgae types on the beach. Complete the datasheet.
- For sampling submerged areas, toss the rake 6 times, 2 tosses in each of 3 directions (if limited by access to a single point) or 3 areas (if you have room to move to another location at the station). Photograph each rake, even empty rakes. Complete the datasheet.
- Enter the data into the online portal. Upload your photos to the online portal.

Preparation

- Gather field equipment.
- Ensure battery power and available storage is sufficient on camera or smartphone.
- Ensure battery power is sufficient on GPS unit or smartphone.
- Always carry a copy of this SOP and the relevant parameter-specific SOPs.
- Print out the “quick sheets” for relevant SOPs to use as a reminder in the field. Do not laminate these as you will want to add notes. A plastic page-protector taped close can be used to keep these sheets dry.

Field Collection and Processing

Option 1: Soft Shoreline Station (Beaches)

- Conduct the assessment within 3 hours of low tide (this gives you a 6-hour window).
- Complete the station information on the datasheet. IMPORTANT: Record GPS points as decimal degrees (e.g. 41.56437 N, -71.87634 W).
- Take two overview photos, showing the overall appearance of the shoreline. One should include upland areas for context, the next should be slightly closer, showing macroalgae (or lack of macroalgae) on the beach (examples below).
- For the dominant macroalgae types, take a close-up photo of each macroalgae.
- Record the amount of macroalgae on the beach as “none”, “some”, or “lots”.
 - None – essentially nothing on the beach, only attached macroalgae is present.
 - Some – ranges from a small bit of macroalgae on the beach to large wracklines of macroalgae present on the beach. This is qualitative, the science advisors will review when comparing stations, so do not worry too much about what truly constitutes some versus lots.
 - Lots – covers > 50% of the beach. May be a thin covering, in thick wracklines, or a thick layer covering the whole beach.
- Examine the macroalgae and use the guide to identify macroalgae morphology (hair-like, twig/leaf like, sheets) and color (green, other). You only need to record macroalgae that constitutes more than 10% of the macroalgae present on the beach.
- Remember – we are **looking for** large amounts of seaweed. If there is nothing on the beach, but you see lots of seaweed (that are not the beneficial brown seaweed) in the water, then try a rake toss in the water.

UWS Standard Operating Procedure
Qualitative Macrophytes

January 29, 2018
Page 16 of 16

Option 2: Submerged Station

- The timing of sampling is not critical. The only criterion is that low tide has not exposed the sediment at the sampling station.
- *Note* – a YouTube video demonstrating how to properly toss a rake is available at: <https://www.youtube.com/watch?v=yMqpSaA6ZWY>.
- Complete the station information on the datasheet. IMPORTANT: Record GPS points as decimal degrees (e.g. 41.56437 N, -71.87634 W).
- Confirm that the line is securely attached to the rake. Place your foot on the end of the rope, to keep it from going into the water. Face the embayment and hold the rake such that your dominant hand is at the end of the rake (ready to push the rake). Push the rake out into the water a distance of ~2 rake lengths. The motion is similar to an underhand slow pitch, rather than a javelin throw.
 - If the toss is too short or too long, bring back and try again.
 - The tines of the rake should be downward. If the rake lands with tines up, try again.
- After a successful toss, allow rake to settle on the floor then ***slowly and steadily*** pull the line so the teeth of the rake drag on the embayment bottom.
 - If the rake gets hooked on a rock or other object, try taking up the slack in the rope and pulling up on the rake, or moving to the left and to the right. If you feel that macroalgae was likely dislodged, try tossing again.
 - When removing the rake from the water, do not let macroalgae fall off of the rake.
 - Make certain that you DO NOT pick up macroalgae removed from the rake on a previous toss.
 - If you get a lot of mud, put into a bucket or bin and rinse; mud that clumps on the rake is often being held together by fine macroalgae.
- Take a photo of the rake (and macroalgae) after each toss.
 - Include all of the rake tines and macroalgae in the photo. If a portion of the rake tines is not included, take another photograph. You will want to be about 4 feet from the rake. ***Do not get too close*** – some cameras do not focus at shorter distances. With the resolution required, zooming in to look at macroalgae can be done on the computer. You do not need to get super close to get a good shot of the macroalgae, further back is better.
 - Frame the shot such that the light is behind the photographer.
 - Be aware of the background and change your location for the photo as needed. Light colors in the background make the macroalgae look darker; avoid light backgrounds (e.g. a white boat).
 - Zoom in on the photo and confirm the macroalgae is in focus (versus something in the background or foreground).
 - Be sure that the lighting allows for realistic colors in the photo.
 - Record an identifier for the photo on the datasheet. For a camera, this may be the filename. For a smartphone, record the time associated with the photo; this can be located by viewing the photo, the time the photo was taken should be on the screen.
- Record the amount of macroalgae on the rake as “none”, “some”, or “lots”.
 - None – no macroalgae on the rake.
 - Some – some tines are still bare.
 - Lots – all tines hold macroalgae. This includes anything from a thin coat of a slimy macroalgae to thick clumps of a wiry or leafy macroalgae.
 - If you get a lot of mud, put into a bucket or bin and rinse; mud that clumps on the rake is often being held together by fine macroalgae
- Examine the macroalgae and use the guide to identify macroalgae morphology (hair-like, twig/leaf like, sheets) and color (green, other). You only need to record macroalgae that constitutes more than 10% of the macroalgae present on the rake.

Appendix A-7

Standard Operating Procedure
Filtered Nutrients Sample Collection

revised January 2019
Page 1 of 6

POINT OF CONTACT

NAME: Peter Linderoth
ADDRESS: Save the Sound, 545 Tompkins Avenue, 3rd Floor
Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

I. OBJECTIVE: Determine the ammonia, nitrite, nitrate, total dissolved nitrogen, and orthophosphate concentrations in the water column.

II. OVERVIEW: Water samples are collected from the surface of the water column. This water is filtered through GF/F filters and delivered into clean HDPE bottles. The samples are analyzed on a Lachat 8500 QuikChem® Autoanalyzer, manifold reaction unit, colorimetric detector, 520 nm wavelength filter and Omnion software data system.

III. SOURCE:

IV. MATERIALS AND EQUIPMENT:

A. FIELD COLLECTION AND STORAGE OF SAMPLES

Plastic bottles, 60 mL, HDPE, (Fisher Scientific Catalog Number: 03-313-11B)

Plastic bottles, 1 L, HDPE, (Fischer Scientific Catalog Number: 02-925-3E)

Glass Fiber Filters (GF/F), 2.5 cm diameter, retains particles down to 0.7 µm, (Fisher Scientific Catalog Number: 09-874-64; Whatman Number: 1825-025)

Filter holders, 2.5 cm (*preferred*: Pall brand, 25 mm Easy Pressure Syringe Filter Holder, Delrin Plastic; *second choice*: Millipore Swinnex Filter Holder, 25 mm polypropylene with silicon gasket, Fisher Scientific Catalog Number: SX00 025 00; Millipore Number: SX0002500)

Forceps for handling filters

Acid Bath: 1 N (Certified) hydrochloric acid (Fisher Scientific Catalog Number: SA48-4)

2 L of ASTM type I (ultrapure) for field blank

Standard Operating Procedure
Filtered Nutrients Sample Collection

revised January 2019
Page 2 of 6

Equipment for getting water samples:

- option 1 – Master flex pump with silicone tubing sufficient for sample depth
- option 2 – Homemade sampling poles with 1 L Bottle

Equipment for delivering sample through filter to 60 mL bottles:

- option 1 – MasterFlex pump with silicone tubing
- option 2 – 60 mL syringe

V. METHODS

A. PREPARATION

- Check that the pump or other sampling equipment are operational
Prepare the sample bottles:
- Wear powder-free nitrile gloves
- Rinse 1 L bottles with ultrapure, ASTM type 1 (ultrapure) water. Allow to dry and store with caps on
 - If 60 mL HDPE bottles are new and un-used, rinse the vials three times with ultrapure, ASTM type I (ultrapure) water. Allow to dry and store with caps on
 - If 60 mL HDPE bottles have been used, acid wash the bottles. Acid washing entails soaking the bottles in a 1 N hydrochloric acid bath for 60 minutes followed by rinsing three times with ultrapure, ASTM type I (ultrapure) water. Allow to dry and store full of ultrapure, ASTM type 1 (ultrapure) water with caps on

B. FIELD COLLECTION AND STORAGE OF SAMPLES

A slightly different procedure is used based on the sampling method. Choose the protocols for your sampling method

MasterFlex Pump

1. Rinse the sampling equipment with sample water. This involves rinsing with three times the volume of the sampling device. Place the intake 0.5 m below the surface, allow the pump to

Standard Operating Procedure
Filtered Nutrients Sample Collection

revised January 2019
Page 3 of 6

run long enough to pass a minimum of three times the volume of the tubing through the system. Discharge this water over the boat, do not collect.

2. Assemble the filtering mechanism. Connect the filter holder (with filter loaded into the holder) to the pump. Change filter if needed during collection.
3. Rinse the sample bottles with water from target sampling depth. Cap the bottle and shake to rinse. Discard the sample used as a rinse. Repeat three times for each bottle.
4. Fill sample bottle with water from target sampling depth.
5. Sample bottle is immediately stored in cooler full of ice for transport to lab.

Note: Sample(s) can be collected in either the 1 L bottle or three 60 mL HDPE bottles. If only filling the 1 L bottle, filtering can occur when 60 mL bottles are filled as seen lower in this SOP under sample preparation. If filling 60 mL bottles in the field, only collect up to 50 mL in the bottle.

Note: About mid-way through the sampling day, perform a field blank. Follow the procedures for sampling above, but use ASTM Type I (ultrapure) water brought out on the boat in place of the field water.

Homemade Sampling Poles with 1 L Bottle

A sampling pole can be fabricated by a member of your organization. The key criteria for acceptability is the ability to sample from a specific depth without contaminating the sample. The method involves reaching into the water with the sample bottle inverted and full of air, then righting the bottle and allowing it to fill at the appropriate depth.

Homemade poles will typically work only for 0.5 m below the surface; deeper depths require alternate methods of sampling.

- Remove the cap from the bottle.
- Rinse the bottle three times with surface water. Fill bottle, cap bottle, invert, then discard rinse.

Standard Operating Procedure
Filtered Nutrients Sample Collection

revised January 2019
Page 4 of 6

- Holding the bottle with the mouth down, lower it to 0.5 m below the surface.
- Invert the bottle so that air empties out of the bottle and the bottle fills with water.
- Bring to the surface and cap.
- Sample bottle is immediately stored in cooler full of ice until being filtered.

Note – About mid-way through the sampling day, perform a field blank. Follow the procedures for sampling above, but use ASTM Type I (ultrapure) Water brought out on the boat in place of the field water.

The following steps apply to when using a MasterFlex for sample preparation:

MasterFlex Pump – after collecting sample in 1 L bottle and storing on ice

1. Rinse the sampling equipment with sample water. This involves rinsing with three times the volume of the sampling device. Place the intake into the sample bottle, allow the pump to run long enough to pass a minimum of three times the volume of the tubing through the system. Discharge this water, do not collect.
2. Assemble the filtering mechanism. Connect the filter holder (with filter loaded into the holder) to the pump. Run approximately 15 mL through the filter before next step.
3. Rinse the sample bottles. Deliver a small volume of the filtered sample into the sample bottles. Cap the vials and shake to rinse. Discard the sample used as a rinse. Repeat for a total of three small volume rinses per bottle.
4. Deliver between 45 mL and 50 mL of sample (through the filter) to the bottles. Be sure to keep the volume in the bottle under 50 mL, this head space will allow for expansion during freezing. Change the filter if needed.
5. Store the bottles *upright* in the freezer (- 20° C).

A sample, replicate, and sample in reserve are collected from each station.

The following steps apply when using a syringe for sample preparation:

Standard Operating Procedure
Filtered Nutrients Sample Collection

revised January 2019
Page 5 of 6

1. Rinse the sampling equipment with sample water. This involves rinsing with three times the volume of the sampling device.
2. Assemble the filtering mechanism. Connect the filter holder (with filter loaded into the holder). Rinse the 60 mL syringe with 5 mL of sample water. Repeat for a total of three rinses. Fill the 60 mL syringe with sample water.
3. Rinse the HDPE bottles. Deliver a small volume of the filtered sample into the HDPE bottles. Cap the bottles and shake to rinse. Discard the sample used as a rinse. Repeat for a total of three small volume rinses per bottle.
4. Deliver between 45 mL and 50 mL of sample (through the filter) to the bottles. Be sure to keep the volume below 50 mL, this head space will allow for expansion during freezing. Change the filter as needed.
5. Upon return to the lab, store the bottles in the freezer (- 20° C).

A sample, replicate, and sample in reserve are collected at each station.

C. LABORATORY ANALYSIS

- The nutrients are analyzed on a Lachat 8500 QuikChem®Autoanalyzer , manifold reaction unit, colorimetric detector, 520 nm wavelength filter and Omnion software data system.Lachat Quickchem Autoanalyzer, or similar instrument, at the Interstate Environmental Commission (IEC). Instrument operations are detailed in the IEC Laboratory SOP which is contained in the UWS QAPP.

VI. TROUBLE SHOOTING / HINTS

- Avoid contaminating the samples – do not touch: the insides of the bottle caps, the mouth of the bottles, the insides of the graduated cylinders, the filter pads
- Two people working in tandem will speed the process. One person focuses on filtering while the other handles the sample bottles and filters. The person filtering the sample can lend a hand when s/he gets ahead of the sample handler.

Standard Operating Procedure
Filtered Nutrients Sample Collection

revised January 2019
Page 6 of 6

VII. DATA PROCESSING AND STORAGE

- Enter the data on the field sheet. Be sure to fill out the data sheet completely.
- Enter data when returned from the lab into the Excel template.

VIII. QUICK SHEET

Refer to the SOP for details, this list is only a reminder!!

FIELD

Rinse everything three times before collecting sample

Store in cooler on ice while in the field

STORAGE

1 L sample bottle remains on ice immediately after collection. Filtering must commence the same day of sampling.

Store samples *upright* in the freezer

Separate replicates a and b (samples to be analyzed) from c (sample held in reserve)

ANALYSIS

Follow any potential guidelines established for analytical lab

Appendix A-8

Standard Operating Procedure
Total Nitrogen and Total Phosphorous
Sample Collection

revised January 2019
Page 1 of 6

POINT OF CONTACT

NAME: Peter Linderoth
ADDRESS: Save the Sound, 545 Tompkins Avenue, 3rd Floor
Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

I. OBJECTIVE: Determine the dissolved total nitrogen and total phosphorous concentrations in the water column.

II. OVERVIEW: Water samples are collected from the surface in the water column. The samples are analyzed on a Lachat Autoanalyzer with autosampler and dilutor, manifold reaction unit (heating unit), colorimetric detector, 880 nm wavelength filter and Omnion software data system.

III. SOURCE:

IV. MATERIALS AND EQUIPMENT:

A. FIELD COLLECTION AND STORAGE OF SAMPLES

Plastic bottles, 60 mL, HDPE, (Fisher Scientific Catalog Number: 03-313-11B)

Plastic bottles, 1 L, HDPE, (Fischer Scientific Catalog Number: 02-925-3E)

Acid Bath: 1 N (Certified) hydrochloric acid (Fisher Scientific Catalog Number: SA48-4)

2 L of ASTM type I (ultrapure) for field blank

Equipment for getting water samples:

- option 1 – MasterFlex pump with silicone tubing sufficient for sample depth
- option 2 – Homemade sampling poles with 1 L Bottle

Equipment for delivering sample to 60 mL bottles:

- option 1 – MasterFlex pump with silicone tubing
- option 2 – 60 mL syringe

Standard Operating Procedure
Total Nitrogen and Total Phosphorous
Sample Collection

revised January 2019
Page 2 of 6

B. LABORATORY ANALYSIS

Included in IEC Lab SOPs in QAPP.

V. METHODS

A. PREPARATION

- Check that the pump or other sampling equipment are operational
- Prepare the sample bottles:
- Wear powder-free nitrile gloves

Use only ultrapure, ASTM type I (ultrapure)

Rinse 1 L bottles with ASTM type I (ultrapure) Allow to dry and store with caps on.

If 60 mL HDPE bottles are new and un-used, rinse the vials three times with ultrapure, ASTM type I water. Allow to dry and store with caps on.

If 60 mL HDPE bottles have been used, acid wash the bottles. Acid washing entails soaking the bottles in a 1 N hydrochloric acid bath for 60 minutes followed by rinsing three times with ultrapure, ASTM type I water. Allow to dry and store full of ASTM type 1, ultrapure, water with caps on.

B. FIELD COLLECTION AND STORAGE OF SAMPLES

A slightly different procedure is used based on the sampling method. Choose the protocols for your sampling method.

MasterFlex Pump

1. Rinse the sampling equipment with sample water. This involves rinsing with three times the volume of the sampling device and bottle. Place the intake into the target sampling depth, allow the pump to run long enough to pass a minimum of three times the volume of the tubing through the system. Discharge this water over the boat, do not collect.
2. Rinse the sample bottle with water from target sampling depth. Cap the bottle and shake to rinse. Discard the sample used as a rinse. Repeat three times.
3. Fill sample bottle with water from target sampling depth.

Standard Operating Procedure
Total Nitrogen and Total Phosphorous
Sample Collection

revised January 2019
Page 3 of 6

4. Sample bottle is immediately stored in cooler full of ice for transport to lab.

Note: Sample(s) can be collected in either the 1 L bottle or three 60 mL HDPE bottles. If filling 60 mL bottles in the field, only collect up to 50 mL in the bottle.

Note: About mid-way through the sampling day, perform a field blank. Follow the procedures for sampling above, but use ASTM type I (ultrapure) out on the boat in place of the field water.

Homemade Sampling Poles with 1 L Bottle

A sampling pole can be fabricated by a member of your organization. The key criteria for acceptability is the ability to sample from a specific depth without contaminating the sample. The method often involves reaching into the water with the sample bottle inverted and full of air, then righting the bottle and allowing it to fill at the appropriate depth.

Homemade poles will typically work only for 0.5 m below the surface; deeper depths require alternate methods of sampling.

- Remove cap from bottle.
- Rinse the bottle three times with surface water. Fill bottle, cap bottle, invert, then discard rinse.
- Holding the bottle with the mouth down, lower it to 0.5 m below the surface.
- Invert the bottle so that air empties out of the bottle and the bottle fills with water.
- Bring to the surface and cap.
- Sample bottle is immediately stored in cooler full of ice until being filtered.

Note – About mid-way through the sampling day, perform a field blank. Follow the procedures for sampling above, but use ASTM type I (ultrapure) brought out on the boat in place of the field water.

The following steps apply when using a MasterFlex for sample preparation:

Standard Operating Procedure
Total Nitrogen and Total Phosphorous
Sample Collection

revised January 2019
Page 4 of 6

MasterFlex Pump – after collecting sample in 1 L bottle and storing on ice

1. Rinse the sampling equipment with sample water. This involves rinsing with three times the volume of the sampling device. Place the intake into the sample bottle, allow the pump to run long enough to pass a minimum of three times the volume of the tubing through the system. Discharge this water, do not collect.
2. Rinse the sample bottles. Deliver a small volume of the sample into the sample bottles. Cap the vials and shake to rinse. Discard the sample used as a rinse. Repeat for a total of three small volume rinses per bottle.
3. Deliver between 45 mL and 50 mL of sample to the bottles. Be sure to keep the volume in the bottle under 50 mL, this head space will allow for expansion during freezing. Change the filter if needed.
4. Store the bottles *upright* in the freezer (- 20° C).

A sample, replicate, and sample in reserve are collected from for station.

The following steps apply when using a syringe for sample preparation:

1. Rinse the sampling equipment with sample water. This involves rinsing with three times the volume of the sampling device.
2. Rinse the 60 mL syringe with 5 mL of sample water. Repeat for three rinses. Fill the 60 mL syringe with sample water.
3. Rinse the HDPE bottles. Deliver a small volume of the sample into the HDPE bottles. Cap the bottles and shake to rinse. Discard the sample used as a rinse. Repeat for a total of three small volume rinses per bottle.
4. Deliver between 45 mL and 50 mL of sample to the bottles. Be sure to keep the volume below 50 mL, this head space will allow for expansion during freezing.
5. Upon return to the lab, store the bottles in the freezer (- 20° C).

Standard Operating Procedure
Total Nitrogen and Total Phosphorous
Sample Collection

revised January 2019
Page 5 of 6

A sample, replicate, and sample in reserve are collected at each station.

C. LABORATORY ANALYSIS

- The nutrients are analyzed on a Lachat 8500 QuikChem®Autoanalyzer , manifold reaction unit, colorimetric detector, 520 nm wavelength filter and Omnion software data system. Lachat Quickchem Autoanalyzer, or similar instrument, at the Interstate Environmental Commission (IEC). Instrument operations are detailed in the IEC Laboratory SOP.

VI. TROUBLE SHOOTING / HINTS

- Avoid contaminating the samples – do not touch: the insides of the vial caps, the mouth of the vials, the insides of the graduated cylinders.
- Two people working in tandem will speed the process.

VII. DATA PROCESSING AND STORAGE

- Enter the data on the field sheet. Be sure to fill out the data sheet completely.
- Enter data when returned from lab into the Excel template.

Standard Operating Procedure
Total Nitrogen and Total Phosphorous
Sample Collection

revised January 2019
Page 6 of 6

VIII. QUICK SHEET

Refer to the SOP for details, this list is only a reminder!!

FIELD

Rinse everything three times before collecting sample (pump / sample collection vessel; vials)
store in cooler on ice while in the field

STORAGE

Store samples *upright* in the freezer

Separate replicates a and b (samples to be analyzed) from c (sample held in reserve)

ANALYSIS

Follow any potential guidelines established for analytical lab

Appendix A-9

Standard Operating Procedure
Continuous Dissolved Oxygen

revised February 2020
Page 1 of 5

POINT OF CONTACT

NAME: Peter Linderoth
ADDRESS: Save the Sound, 545 Tompkins Avenue, 3rd Floor
Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

I. OBJECTIVE: Determine the dissolved oxygen (DO) concentration and saturation over a period of time at select intervals.

II. OVERVIEW: Continuous dissolved oxygen sampling instruments – dissolved oxygen, conductivity, and barometric pressure loggers – are deployed in the field to log dissolved oxygen and saturation. Dissolved oxygen and conductivity are logged at approximately 0.5 meters off the bottom and barometric pressure is logged in a secure location in close proximity to the other loggers.

III. SOURCE:

IV. MATERIALS AND EQUIPMENT:

A. Continuous logging stations

Dissolved oxygen logger (HOBO U26)

Barometric pressure logger (HOBO U20L-01)

Conductivity Logger (Star-Oddi DST-CT, 13-50 mS/cm range)

Aquarium water pump for inter comparison of DO and Conductivity loggers and sonde

Aquarium air stone for inter comparison of DO and Conductivity loggers and sonde

Sodium sulfite DO calibration solution (HOBO U26-CAL-SOL)

Ice/tap water for adjusting conductivity in water bath

Multiparameter sonde (YSI EXO1 or Eureka Manta +35) with DO and conductivity attachments

Conductivity Standard (YSI: 50,000 μ S/cm) for calibrating sonde

Q tips for cleaning inside of DO loggers

Standard Operating Procedure
Continuous Dissolved Oxygen

revised February 2020
Page 2 of 5

B. LABORATORY ANALYSIS

No laboratory analysis.

V. METHODS

A. Logger Inter Comparison (Prior to Field Deployment)

- Check that all sampling equipment are operational and ready for deployment
- Calibrate conductivity and DO on multiparameter sonde to manufacturer recommendations the day before or morning of this logger comparison
- Calibrate all loggers according to manufacturer recommendations the day before or morning of this logger comparison
- Synchronize loggers and multiparameter sonde to same time. Set loggers to 15 minute interval readings
- Prepare a water bath with water similar to deployment stations; enough volume to fit all loggers for comparison and sonde
- Place DO and conductivity loggers ready for deployment in the water bath and note the time the loggers will be taking readings.
- Run the aquarium pump to keep water flowing in water bath. Do not keep loggers or sonde close enough to allow bubbles to come into contact with sensors
- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Add ice or tap water to adjust the salinity concentration in the water bath. Ensure that the salinity (conductivity) in the water bath does not go out of conductivity logger range
- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Add sodium sulfite solution to lower the DO level in the water bath. This should be done slowly
 - NOTE: Do not leave equipment in this low-oxygen solution for too long. Two readings will suffice

Standard Operating Procedure
Continuous Dissolved Oxygen

revised February 2020
Page 3 of 5

- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Logger data is retrieved from loggers and stored with multiparameter sonde readings. The salinity correction will need to be applied through the HOBOWare software for correct DO concentration

B. Logger Inter Comparison (After Field Deployment)

- Bring all loggers back to office or lab
- Prepare a water bath with water similar to deployment stations; enough volume to fit all loggers for comparison and sonde
- Place DO and conductivity loggers from the field in the water bath without cleaning any fouling and note the time the loggers will be taking readings.
- Run the aquarium pump to keep water flowing in water bath. Do not keep loggers or sonde close enough to allow bubbles to come into contact with sensors
- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Clean all sensors before proceeding to the next steps
- Place DO and conductivity loggers ready for deployment in the water bath and note the time the loggers will be taking readings.
- Run the aquarium pump to keep water flowing in water bath. Do not keep loggers or sonde close enough to allow bubbles to come into contact with sensors
- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Add ice or tap water to adjust the salinity concentration in the water bath. Ensure that the salinity (conductivity) in the water bath does not go out of conductivity logger range
- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Add sodium sulfite solution to lower the DO level in the water bath. This should be done slowly

Standard Operating Procedure
Continuous Dissolved Oxygen

revised February 2020
Page 4 of 5

- NOTE: Do not leave equipment in this low-oxygen solution for too long. Two readings will suffice
- Record DO and conductivity readings with the multiparameter sonde at the same time the loggers record data
- Logger data is retrieved from loggers and stored with multiparameter sonde readings. The salinity correction will need to be applied through the HOBOWare software for correct DO concentration

C. FIELD DATA COLLECTION

1. Logger Deployment and Data Collection

- Dissolved oxygen and conductivity loggers are deployed at approximately 0.5 m off bottom. Loggers need to be actively logging when deployed to obtain initial sonde/logger comparison in field
- Barometric pressure logger is deployed above water at a secure and close location.
 - Logger needs to be actively logging prior to deployment
- All loggers are set to record at 15 minute intervals
- Station visits are required every 7-10 days for cleaning and comparison with multiparameter sonde
 - Conductivity, dissolved oxygen, and barometric pressure are all recorded at the station at 0.5 m off bottom using sonde and handheld
 - Readings must coincide with a known time the loggers are recording data
 - Record readings on field data sheet
- Retrieve submerged loggers and clean all fouling organisms
- Couple loggers with data retrieval method of choice and collect data for storage
- Repeat the comparison sampling procedure after the loggers are cleaned and redeployed
- Bring data retrieval instrument to the laboratory/office and upload data to analysis software provided by manufacturer

Standard Operating Procedure
Continuous Dissolved Oxygen

revised February 2020
Page 5 of 5

IX. QUICK SHEET

Refer to the SOP for details, this list is only a reminder!!

INTER COMPARISON

Zero DO solution needs to be added very slowly to avoid adding too much. Obtain a full reserve of water for use in the event the DO plummets quickly and below the preferred accuracy range of the instruments.

FIELD

Secure all loggers with backup ropes and security measures. All loggers need to be set to 15 minute intervals. Inter comparison readings are taken before and after cleaning loggers of fouling organisms.

STORAGE

Data sheets are filed and digital logs of data are stored on local computer/server prior to being sent to Save the Sound.

ANALYSIS

Upload data from field onto manufacturer software for analysis.

Appendix A-10

Standard Operating Procedure
Macrophyte Percent Coverage via Camera

revised January 2019
Page 1 of 7

POINT OF CONTACT

NAME: Peter Linderoth
ADDRESS: Save the Sound, 545 Tompkins Avenue, 3rd Floor
Mamaroneck, NY 10543
EMAIL: plinderoth@savethesound.org
PHONE: 914-263-6233

I. OBJECTIVE: To quantify the percent coverage of macrophytes and determine bottom type at select locations.

II. OVERVIEW: An underwater camera (video or still) is used to capture images of the bottom in the area of a water quality station. The bottom type is identified and the percent cover of macrophytes is calculated.

III. SOURCE:

IV. MATERIALS AND EQUIPMENT:

A. FIELD COLLECTION AND STORAGE OF SAMPLES

- GoPro camera
- Setup for keeping camera a set distance from the bottom (to maintain consistent areal size in the field of vision)
- GPS unit

B. LABORATORY ANALYSIS

- not applicable

V. METHODS

A. PREPARATION

- Check the charge on the GPS battery
- Check the charge on the camera battery
- Check the camera system for proper functioning

Standard Operating Procedure
Macrophyte Percent Coverage via Camera

revised January 2019
Page 2 of 7

- Check the deployment rig for the camera system (must have a size reference in the field of vision)

B. FIELD COLLECTION AND STORAGE OF SAMPLES

- Samples stored on the SD card of the camera
- Write station, date, and time on a white board
- Approach station with camera and all equipment ready
- Start video or camera sequence with image of the white board and ensure GPS unit is tracking
- Run the boat slowly across a station starting around 15 m out and passing through it by around 15 m dropping pole to the floor waiting at least 3 seconds between drops. This can vary slightly but make sure the boat is moving. Multiple approaches from different directions to the station may be needed to collect the quantity of images required.
- Collect between 20 and 40 bottom images – in areas with greater heterogeneity, collect more samples (~40), if an area is fairly homogenous (e.g. sand, oyster, cobble) you may collect fewer (~20)
- Stop GPS track when navigating between stations
- Name GPS tracks using UWS station ID_date_time (24 hr)

C. LABORATORY ANALYSIS

- Not applicable

VI. TROUBLE SHOOTING / HINTS

- For the video system – keep the speed of the boat low, so as not to put too much tension on the pole (the pole will snap at high speeds)

VII. DATA PROCESSING AND STORAGE

- Video will be downloaded the following day and stored on computer and/or server/cloud storage.
- Download the GPS tracks the following day and store on computer as GPX file. Open GPX tracks in Garmin BaseCamp software and export as a CSV table.
- For video - A trained analyst will watch the video and do a screen-capture for all instances where the base of the pole is resting on the bottom (Figure 1).
- The bottom images will be pasted into the Excel camera work template (Figure 2).

Standard Operating Procedure
Macrophyte Percent Coverage via Camera

revised January 2019
Page 3 of 7

- The first analyst adds GPS coordinates corresponding to the image located in the CSV table exported from BaseCamp. Analyst will match the time between the video and the CSV and add to template.
- A 100 cell grid is overlaid on the image.
- The first analyst reviews the image and enters the % cover in three categories: eelgrass, macroalgae, bare sediment. The first analyst hides their columns of data, so that other analysts cannot see the first set of estimates.
- A second analyst reviews the images and makes an independent assessment of percent cover. The second analyst also checks the GPS coordinate data entry. The second analyst hides their columns of data, so that the other analysts cannot see their estimates.
- A third analyst performs the same tasks as the second analyst.
- The lab manager reviews a minimum of 10% of the images for accuracy of percent coverage estimates.
- INSERT ANY QUESTIONS INTO THE “COMMENTS” COLUMN
- In the Excel template, data are condensed into a table. Estimates from the three analysts are compared. If the relative percent difference among the three estimates is greater than 5%, the lab manager examines the image and the three estimates, choosing the appropriate value. The three estimates will not be changed, values are retained to show the inconsistency. The lab manager decides on the final value for the estimate. If a specific analyst's estimate consistently differs, the analyst will receive further training or may be removed from conducting future estimates.
- See Figure 3 for examples of coverage estimates and how to handle sparse coverage.

Macrophyte Percent Coverage via Camera

Page 4 of 7



Figure 1: Image of the bottom captured using the video camera system. Note that the black circular disc on the end of the pole keeps the pole from sinking into the sediment and provides a size reference (diameter of disc = 10 cm).

[illegible]











Figure 2: The image from Figure 1 has been entered into the data template and overlaid with the 100 cell grid.

Figure 3: Examples and tips for estimating coverage.











Standard Operating Procedure
Macrophyte Percent Coverage via Camera

revised January 2019
Page 5 of 7

		<p>eelgrass = 100 % macroalgae = 0 % bare = 0 %</p> <p>while leaves block some field of vision, this appears to be dense coverage</p>
		<p>eelgrass = 76 % macroalgae = 0 % bare = 24 %</p> <p>the turbidity of the water makes this one tricky to analyze, zoom in and look closely</p>
		<p>eelgrass = 89 % macroalgae = 0 % bare = 11 %</p> <p>harder to estimate – take care counting – you will have to estimate partial grid cells to compute bare area</p>
		<p>eelgrass = 73 % macroalgae = 0 % bare = 27 %</p> <p>harder to estimate – take care counting – you will have to estimate partial grid cells to compute bare area</p>
		<p>eelgrass = 69 % macroalgae = 0 % bare = 31 %</p> <p>harder to estimate – take care counting – you will have to estimate partial grid cells to compute bare area</p>

Standard Operating Procedure
Macrophyte Percent Coverage via Camera

revised January 2019
Page 6 of 7

		<p>eelgrass = 66 % macroalgae = 21 % bare = 13 %</p> <p>note what appears to be Ulva on the left (comment entry with any ?'s): - you will sometimes need to count more than once... (I counted macroalgae, then bare)</p>
		<p>eelgrass = 43 % macroalgae = 3 % bare = 54 %</p> <p>you can calculate the % coverage for two types, then calculate the third: $100 - 43 - 3 = 54$</p>
		<p>eelgrass = 16 % macroalgae = 1 % bare = 83 %</p> <p>in a sparse eelgrass area, do not count the leaves that extend across the frame as area – look for the base of the plants</p>
		<p>eelgrass = 0 % macroalgae = 13 % bare = 87 %</p> <p>count shells as bare area, unless colonized by macroalgae</p>

Standard Operating Procedure
Macrophyte Percent Coverage via Camera

revised January 2019
Page 7 of 7

VIII. REFERENCES:

IX. QUICK SHEET

Refer to the SOP for details, this list is only a reminder!!

FIELD

- Collect 20 – 40 images dependent on seafloor composition

STORAGE

- Download images and GPS tracks ASAP, at least by the day following field work

ANALYSIS

- Three analysts check
- Record to the nearest 1% of coverage
- Any questions go in the comments column

UWS Sonde Calibration Datasheet YSI EXO 1

- Calibrations to be completed **DAY BEFORE** or **MORNING OF** Field Sampling Date •
• Post-Readings to be completed the **AFTERNOON OF** or **DAY AFTER** Field Sampling Date •

Calibrations • Person: _____ Date: _____ Time: _____

Post-Readings • Person: _____ Date: _____ Time: _____

Handheld S/N: _____ Sonde S/N: _____

4 COMPLETE BEFORE SAMPLING 4

① Calibrate DISSOLVED OXYGEN (DO) (% sat)

Barometric Pressure (mmHg) ----

Temperature ----

Pre-Cal Value Post-Cal Value

DO Sat ----

② Calibrate CHLOROPHYLL (a) (µg/L) • 1-Point Calibration

Pre-Cal Value Post-Cal Value

③ Calibrate CHLOROPHYLL (a) (NTU) • 1-Point Calibration

Pre-Cal Value Post-Cal Value

④ Calibrate TURBIDITY • 2-Point Calibration

→ 1st Cal Value: 0 NTU (Reagent Grade Water)

Pre-Cal Value Post-Cal Value

→ 2nd Cal Value: 100 NTU (Turbidity Standard)

Pre-Cal Value Post-Cal Value

⑤ Calibrate Specific Conductance (SC) (µS/cm)

Pre-Cal Value Post-Cal Value

Cell Constant ----

⑥ Calibrate DEPTH (m)

Pre-Cal Value Post-Cal Value

5 COMPLETE AFTER SAMPLING 5

① Post-reading for DISSOLVED OXYGEN (% sat)

② Post-reading for CHLOROPHYLL (a) (µg/L)

③ Post-reading for CHLOROPHYLL (a) (NTU)

④ Post-reading for TURBIDITY (0-10 NTU)

⑤ Post-reading for TURBIDITY (1-24 NTU)

⑥ Post-reading for SP COND (50,000 µS/cm)

⑦ Post-reading for DEPTH

	Reagent Grade Water	Turbidity Standard (00 NTU)	Conductivity Standard (50,000 µS/cm)
Manufacturer			
Lot Number			
Expiration			

Accuracy Range Table	
DO (% sat)	0.1 - 100
chl a (µg/L)	0.00 - 0.50
Turbidity (0 NTU)	-1.00 - 3.00
Turbidity (100 NTU)	97.0 - 103.0
SpCond (50,000 µS/cm)	49,500 - 50,500
Depth (0 m)	-0.1 - 0.1

GPS of reference station (circle one) BND-82 WGS-84

* WITH 2 SIG OF SAMPLING DAY * IN SCIENTIFIC FORMAT *

Lat: <input type="text"/>	Long: <input type="text"/>
---------------------------	----------------------------

Appendix B-3

Sample Event Datasheet

Employment Name				GPS units (single use)			
Sample Date:				decimal degree (48.712499)			
People				degree minutes (48° 42.594')			
				degree plus sec. (48° 42' 59.56")			

Station ID									
Time									
Station Depth (m)									
GPS N									
GPS W									
	bottom (0.5 m off bottom)	mid-depth (if total depth > 10m)	surface (1m above surface)	bottom (0.5 m off bottom)	mid-depth (if total depth > 10m)	surface (1m above surface)	bottom (0.5 m off bottom)	mid-depth (if total depth > 10m)	surface (1m above surface)
Sample Depth (m)									
Temperature (°C)									
Salinity (ppt)									
Dissolved Oxygen (H)									
Dissolved Oxygen (mg/L)									
Fluorescence (R-R)									
Chl-a (µg/L)									
Turbidity (NTU)									

Enter additional field notes on back of sheet if using a different protocol than usual.

At 1 station per survey event, plot a second profile (usually at last station).

If total depth < 1.0m, use only mid-depth.

Chlorophyll Reference Check in Bucket (do once per day per outlay event)

Date & time	1st 10L	2nd 10L	3rd 10L

data entry _____ person checking _____

Chlorophyll Chain of Custody Form

Laboratory _____

Appendix B-5

Qualitative Macrophyte Sampling Datasheet

Site Name (short identifier) _____ Date _____

People (full names) _____


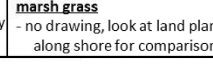

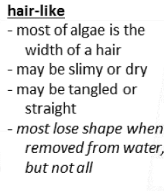
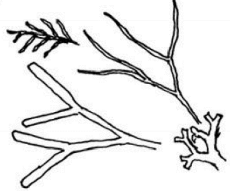

Time of Low Tide _____ Time of Sampling (now) _____

Site Description / Notes (optional) _____

GPS location (in decimal degrees) _____ N _____ W

For rake toss, complete 6 tosses. For soft shoreline, provide an overview photo and close-up photos of algae types. Remember to include something for a size reference in soft shoreline close-up photos (shoe, keys, ruler, etc.).

photo ID	NONE	SOME	LOTS	green hair-like	green twig/leaf-like	green sheets	non-green hair-like	non-green twig/leaf-like	non-green sheets	marsh grass	eelgrass

eelgrass - green when fresh - ribbon-like - up to 6 ft. long - may be brown or grey when decaying		twig/leaf like - most of algae is thicker than a hair - may have small leaf-like sections - may branch or not branch - most maintain their shape when removed from water, but not all	<i>all drawings are to scale when printed on standard paper</i>
marsh grass - no drawing, look at land plants along shore for comparison		hair-like - most of algae is the width of a hair - may be slimy or dry - may be tangled or straight - most lose shape when removed from water, but not all	sheets - may look like floppy lettuce leaves or rubbery straps (kelp) - may be very large or the size of a quarter - a few have a mid-rib (line up the middle), but most of plant does not have a "stem"
			

Appendix B-6 Macrophyte Field Guide

Long Island Sound Unified Water Study Algae Guide PAGE 1 of 2

All photos with permission from: Van Patten, Marguerite (Peg). (2009) Seaweeds of Long Island Sound, 2nd ed. Connecticut Sea Grant College Program. 104 pp.
photographer is P. Van Patten, unless otherwise indicated.

green hair-like = very thin filaments, as thick as hair; may branch



non-green hair-like = very thin filaments, as thick as hair; may branch



C. Yarnish

sheets = most of plant is sheet form, may have a small holdfast

green



non-green



Long Island Sound Unified Water Study Algae Guide PAGE 2 of 2

All photos with permission from: Van Patten, Marguerite (Peg). (2009) Seaweeds of Long Island Sound, 2nd ed. Connecticut Sea Grant College Program. 104 pp.
photographer is P. Van Patten, unless otherwise indicated.

green twig/leaf-like = thicker than a hair, may branch, typically maintains shape when removed from water



eelgrass



J. Voudrey

non-green twig/leaf-like = thicker than a hair, may branch, typically maintains shape when removed from water



"beneficial brown"

Appendix B-7
Nutrients Chain of Custody

UWS Nutrients Chain of Custody		Organization: _____	
Sample ID	Sample Date	Collection Time	Salinity (ppt)

Relinquished by: _____

Date/Time: _____

Received by: _____

Date/Time: _____

Received by: _____

Date/Time: _____

Received by: _____

Date/Time: _____

Laboratory _____

Appendix B-8

Nutrients Sampling Event Datasheet

UWS Tier II - Surface Nutrients			
Date: _____		Embayment: _____	
People: _____			
collect sample 0.5 m below surface			Notes
station (number, letter)			
GPS N			
GPS W			
time (24 hour)			
total station depth (m)			
salinity (ppt)			
60 mL bottle ID	<div style="display: flex; justify-content: space-between;"> _____ - _____ - U _____ - _____ - F </div> <small>a thru C a thru C</small>		
collect sample 0.5 m below surface			Notes
station (number, letter)			
GPS N			
GPS W			
time (24 hour)			
total station depth (m)			
salinity (ppt)			
60 mL bottle ID	<div style="display: flex; justify-content: space-between;"> _____ - _____ - U _____ - _____ - F </div> <small>a thru C a thru C</small>		
collect sample 0.5 m below surface			Notes
station (number, letter)			
GPS N			
GPS W			
time (24 hour)			
total station depth (m)			
salinity (ppt)			
60 mL bottle ID	<div style="display: flex; justify-content: space-between;"> _____ - _____ - U _____ - _____ - F </div> <small>a thru C a thru C</small>		

Appendix B-9

Logger Retrieval Sample Event Datasheet

UWS Logger Retrieval Data Sheet																				
Embayment Name: _____ Station ID: _____		GPS N <div style="border: 1px solid black; width: 60px; height: 20px;"></div>																		
Date: _____ People: _____		GPS W <div style="border: 1px solid black; width: 60px; height: 20px;"></div>																		
Barometric Pressure Unit Serial#: _____ Time of Retrieval: _____	Notes: <div style="border: 1px solid black; height: 150px; width: 100%;"></div>																			
Dissolved Oxygen Logger Serial#: _____ Time of Retrieval: _____ Time of Re-deployment: _____																				
<table border="1" style="width: 100%; border-collapse: collapse;"> <thead> <tr> <th style="text-align: left;">Sonde Readings</th> <th style="text-align: center;">Pre-retrieval</th> <th style="text-align: center;">Post-retrieval</th> </tr> </thead> <tbody> <tr><td>Time</td><td></td><td></td></tr> <tr><td>Temperature (°C)</td><td></td><td></td></tr> <tr><td>Salinity (ppt)</td><td></td><td></td></tr> <tr><td>Dissolved Oxygen (%)</td><td></td><td></td></tr> <tr><td>Dissolved Oxygen (mg/L)</td><td></td><td></td></tr> </tbody> </table>	Sonde Readings	Pre-retrieval	Post-retrieval	Time			Temperature (°C)			Salinity (ppt)			Dissolved Oxygen (%)			Dissolved Oxygen (mg/L)				
Sonde Readings	Pre-retrieval	Post-retrieval																		
Time																				
Temperature (°C)																				
Salinity (ppt)																				
Dissolved Oxygen (%)																				
Dissolved Oxygen (mg/L)																				
Embayment Name: _____ Station ID: _____		GPS N <div style="border: 1px solid black; width: 60px; height: 20px;"></div>																		
Date: _____ People: _____		GPS W <div style="border: 1px solid black; width: 60px; height: 20px;"></div>																		
Barometric Pressure Unit Serial#: _____ Time of Retrieval: _____	Notes: <div style="border: 1px solid black; height: 150px; width: 100%;"></div>																			
Dissolved Oxygen Logger Serial#: _____ Time of Retrieval: _____ Time of Re-deployment: _____																				
<table border="1" style="width: 100%; border-collapse: collapse;"> <thead> <tr> <th style="text-align: left;">Sonde Readings</th> <th style="text-align: center;">Pre-retrieval</th> <th style="text-align: center;">Post-retrieval</th> </tr> </thead> <tbody> <tr><td>Time</td><td></td><td></td></tr> <tr><td>Temperature (°C)</td><td></td><td></td></tr> <tr><td>Salinity (ppt)</td><td></td><td></td></tr> <tr><td>Dissolved Oxygen (%)</td><td></td><td></td></tr> <tr><td>Dissolved Oxygen (mg/L)</td><td></td><td></td></tr> </tbody> </table>	Sonde Readings	Pre-retrieval	Post-retrieval	Time			Temperature (°C)			Salinity (ppt)			Dissolved Oxygen (%)			Dissolved Oxygen (mg/L)				
Sonde Readings	Pre-retrieval	Post-retrieval																		
Time																				
Temperature (°C)																				
Salinity (ppt)																				
Dissolved Oxygen (%)																				
Dissolved Oxygen (mg/L)																				
Embayment Name: _____ Station ID: _____		GPS N <div style="border: 1px solid black; width: 60px; height: 20px;"></div>																		
Date: _____ People: _____		GPS W <div style="border: 1px solid black; width: 60px; height: 20px;"></div>																		
Barometric Pressure Unit Serial#: _____ Time of Retrieval: _____	Notes: <div style="border: 1px solid black; height: 150px; width: 100%;"></div>																			
Dissolved Oxygen Logger Serial#: _____ Time of Retrieval: _____ Time of Re-deployment: _____																				
<table border="1" style="width: 100%; border-collapse: collapse;"> <thead> <tr> <th style="text-align: left;">Sonde Readings</th> <th style="text-align: center;">Pre-retrieval</th> <th style="text-align: center;">Post-retrieval</th> </tr> </thead> <tbody> <tr><td>Time</td><td></td><td></td></tr> <tr><td>Temperature (°C)</td><td></td><td></td></tr> <tr><td>Salinity (ppt)</td><td></td><td></td></tr> <tr><td>Dissolved Oxygen (%)</td><td></td><td></td></tr> <tr><td>Dissolved Oxygen (mg/L)</td><td></td><td></td></tr> </tbody> </table>	Sonde Readings	Pre-retrieval	Post-retrieval	Time			Temperature (°C)			Salinity (ppt)			Dissolved Oxygen (%)			Dissolved Oxygen (mg/L)				
Sonde Readings	Pre-retrieval	Post-retrieval																		
Time																				
Temperature (°C)																				
Salinity (ppt)																				
Dissolved Oxygen (%)																				
Dissolved Oxygen (mg/L)																				

Macrophyte Percentage Coverage Datasheet

Camera Sampling		Date _____	
People		Location _____	
		Site Code _____	
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		
Sta _____	video time _____	ID on White Board _____	Notes _____ _____
	start _____		
	stop _____		

Appendix C-1

Feb 2022

Page 1 of 7

UWS Standard Operating Procedure – Determination of Chlorophyll *a* and Pheophytin *a* in Water

Table of Contents

1	POINT OF CONTACT	2
2	OBJECTIVE	2
3	OVERVIEW.....	2
4	SOURCES	2
5	MATERIALS AND EQUIPMENT	2
5.1	Field Collection	2
5.2	Lab Analysis.....	2
6	QUALITY CONTROL	3
6.1	Initial Demonstration of Performance.....	3
6.2	Calibration and Standardization	3
6.3	Continuing Calibration Checks.....	3
7	METHODS.....	4
7.1	Reagents and Standards Preparation	4
7.1.1	Aqueous Acetone Solution, 1 L.....	4
7.1.2	0.1 N Hydrochloric Acid Solution, 500 mL	4
7.1.3	Chlorophyll Stock Standard Solutions (SSS).....	4
7.1.4	Laboratory Reagent Blank (LRB)	4
7.2	Sample Collection and Storage	4
7.2.1	Sample Collection	4
7.2.2	Storage	4
7.2.3	Holding Time	5
7.3	Laboratory Analysis Procedure	5
7.3.1	Extraction.....	5
7.3.2	Sample Analysis	5
8	TROUBLESHOOTING / HINTS	6
9	DATA PROCESSING AND STORAGE	6
9.1	Sample Data Handling.....	6
9.2	Data Analysis and Calculations	6
10	REFERENCES.....	7

UWS Standard Operating Procedure
Determination of Chlorophyll *a* and Pheophytin *a* in Water

Feb 2022

Page 2 of 7

1 POINT OF CONTACT

NAME: Elena L. Colón, Save the Sound, Environmental Analyst II
ADDRESS: 1385 Boston Post Rd, 2nd Fl, Larchmont, NY 10538
EMAIL: ecolon@savethesound.org
PHONE: 914-381-3140 ext. 201

2 OBJECTIVE

Determine the concentration of chlorophyll *a* (chl *a*) and pheophytin *a* (pheo *a*) in marine or freshwater phytoplankton using fluorescence detection.

3 OVERVIEW

Fluorescence detection is used to determine the concentration of chlorophyll *a* and pheophytin *a* in the water. Samples are collected from a measured volume of water and filtered through a glass fiber filter (GF/F). The field collection method is detailed in the UWS SOP Filtered Chlorophyll *a*.

Using 90% acetone, the chlorophyll present on the GF/F is extracted for a period of no less than 18 h, but not to exceed 24 h. The acidification method will be used whereby the fluorescence is measured before and after the addition of 0.1 N HCl.

4 SOURCES

EPA Method 445.0, "In Vitro Determination of Chlorophyll *a* and Pheophytin *a* in Marine and Freshwater Algae by Fluorescence."

5 MATERIALS AND EQUIPMENT

5.1 Field Collection

See UWS SOP Filtered Chlorophyll. The field team will need all equipment and materials listed to collect chl *a* samples. Four samples (two of which will be analyzed and two saved as back-up) are to be filtered per field day for verification of sonde readings.

5.2 Lab Analysis

- Trilogy Laboratory Fluorometer
 - Chlorophyll *a* Acidification Module
- HPLC Grade Acetone, diluted to 90%
- Hydrochloric Acid, concentrated, diluted to 0.1 N
- Chlorophyll *a* Standards
- Solid Secondary Standard, Red
- 50 mL Centrifuge Tubes

UWS Standard Operating Procedure
Determination of Chlorophyll *a* and Pheophytin *a* in Water

Feb 2022

Page 3 of 7

- ASTM Type I Reagent Grade Water (for diluting acetone)
- Borosilicate Glass Tubes, 12x75mm
- Snap Caps for 12mm Glass Tubes
- Benchtop Centrifuge
- Variable Volume Pipette and Pipette Tips
- Teasing Needle Hook
- Shade Boxes (appropriately labeled with safety warning labels for acetone)
- Kimwipes

6 QUALITY CONTROL

6.1 Initial Demonstration of Performance

An initial demonstration of performance must be conducted to characterize the fluorometer's performance:

- Instrumental Detection Limit (IDL): Pure chlorophyll *a* in 90% acetone solution is serially diluted until it is no longer detected on a maximum sensitivity setting.
- Linear Dynamic Range (LDR): The LDR is determined analyzing a laboratory reagent blank and five calibration standards with concentrations ranging between 0.2 µg/L and 200 µg/L.
- Estimated Detection Limit (EDL): Pure chlorophyll *a* in 90% acetone solution is serially diluted until it yields a respond 3 times the average response of blank filters.
- Method Detection Limit (MDL): Using low-level ambient water, analyze seven replicate samples following the Laboratory Analyses Procedure in Section 7.3. After determining the concentration values, calculate the MDL as follows:

$$MDL = St_{(n-1, 1-\alpha=0.99)}$$

Where,

$t_{(n-1, 1-\alpha=0.99)}$ = the Student's t-value appropriate for a single-tailed 99th percentile t statistic and a standard deviation estimate with n-1 degrees of freedom

n = number of replicates

S = Standard Deviation of the replicate analyses

In addition to establishing the initial performance capabilities, laboratory reagent blanks and quality control samples will be used to continually check on performance. All quality control records will be maintained by the laboratory.

6.2 Calibration and Standardization

Calibration of the Turner Fluorometer should be performed on a monthly basis using Turner Designs chlorophyll *a* standards. The Turner Designs chl *a* standards package comes with two 20 mL ampoules of known concentration of chl *a* in a 90% acetone solution. Using the Turner Designs standards, a low and high concentration will be used for a two-point calibration.

6.3 Continuing Calibration Checks

A calibration standard check using Turner Designs Adjustable Solid Secondary Standard should be conducted pre- and post-readings for every batch of samples (30-40 samples) being processed.

If at any time the calibration standard check fails to meet acceptance criteria ($\pm 10\%$ of known value) the fluorometer must be re-calibrated. In the event chlorophyll *a* standards are not immediately available to perform a two-point calibration a post-correction on the RFU readings may be applied to the samples once a calibration has been able to be conducted.

7 METHODS

7.1 Reagents and Standards Preparation

7.1.1 Aqueous Acetone Solution, 1 L

Measure out 100 mL of ASTM Type I Water using a graduated cylinder or 100 mL volumetric flask. Deliver the 100 mL of water into the 90% acetone storage bottle. Measure out 900 mL of acetone in a graduated cylinder and deliver into the 90% acetone storage bottle. Cap and swirl to mix. Be sure to keep acetone solution well sealed to avoid evaporation.

Due to the resulting properties of mixing water and acetone impacting volumetric readings, it is important measure each independently and then mix the known volumes together.

7.1.2 0.1 N Hydrochloric Acid Solution, 500 mL

Add ~350 mL of ASTM Type I Water into a 500 mL volumetric flask. Add in 50 mL of 1N Hydrochloric Acid Concentrate to the water. Swirl flask gently to mix. Add ASTM Type I Water to the volumetric flask to bring the volume to 500 mL. Cap and mix the water and HCl.

7.1.3 Chlorophyll Stock Standard Solutions (SSS)

Chlorophyll *a* standard with a known concentration of chl *a* can be purchased by Turner Designs and must be kept frozen (-20°C) until ready for use. The Turner Designs standard provides two ampoules, one with a high concentration of chl *a* and the other a low concentration.

7.1.4 Laboratory Reagent Blank (LRB)

Place unused GF/F filter in glass tube and fill with 5 mL of 90% Acetone. One LRB per batch of samples should be prepared and handled in the same method as the field samples being analyzed.

7.2 Sample Collection and Storage

7.2.1 Sample Collection

Method of sample collection is detailed in the UWS SOP Filtered Chlorophyll. A measured volume of water is filtered through a 2.5 cm filter holder loaded with a circle GF/F filter (with a nominal pore size of 0.7 μm). Air is expelled through the filter several times to remove excess water. The filter is then removed from the holder, folded inward in half, placed in an unbleached coffee filter and wrapped in aluminum foil.

7.2.2 Storage

For long term storage, unprocessed filters must be kept frozen (-20°C) in the dark until extraction. Short term storage on ice for no more than 12 hours immediately after collecting the sample is acceptable.

7.2.3 Holding Time

Filters must be analyzed within 28 days from date of collection to avoid degradation of the sample material.

7.3 Laboratory Analysis Procedure

7.3.1 Extraction

IMPORTANT - Sample analyses should always be performed in a shaded and cool area of the lab with limited amount of light. Do not expose samples to direct light.

- Place filter in a 6 mL (12 x 75 mm) glass tube.
- Add 5 mL of 90% Acetone solution and seal with a snap cap.
- Swirl tube several times or vortex, mixing filter in acetone very well.
- Allow filter to steep in acetone for 18-20 hours in a dark freezer (-20°C). There must be no light exposure to the samples.

7.3.2 Sample Analysis

- Several hours (on the day of) before reading of the sample is to be conducted, it is recommended to mix the sample in the glass once more. Allow to continue to steep in the freezer.
- Remove sample from the freezer after allowing filter to steep the appropriate number of hours. Ensure light exposure is limited and wait for sample to reach room temperature before proceeding.
- Taking care to not expose the sample to light for too long, use the teasing needle to remove the filter within the tube. Place the snap cap back on the tube. The filter should be allowed to dry out in the fume hood and thereafter may be disposed of in the regular trash. Make sure to always rinse the teasing needle before using on the next sample by

UWS Standard Operating Procedure
Determination of Chlorophyll *a* and Pheophytin *a* in Water

Feb 2022

Page 6 of 7

- dipping in 90% Acetone solution.
- d. Centrifuge the sample for 5 min at 1000 g or 15 min at 675 g.
 - e. When not actively handling the sample glass tube (either to remove the filter or centrifuge), it should be covered (shade box) so as to prevent exposure to light as much as possible. Once ready to read the tube, wipe the glass tube with a kimwipe. Place in the fluorometer, close the lid and read. Record the value on your lab datasheet. This is your **F1 value** (before acid).
 - f. Once the first reading is recorded, add 0.15 µL of 0.1N HCl to the sample tube. Swirl or vortex to distribute the acid. Once acid is added to the sample tube it may be safely exposed to light without concern of compromising the sample. After adding the acid, wait for at least 90 seconds before proceeding to the next step.
 - g. Centrifuge the sample for 5 minutes at 1000 g or 15 minutes at 675 g.
 - h. Place the now acidified sample tube in the fluorometer and read. Record the value on your datasheet as your **F2 value** (after acid).

8 TROUBLESHOOTING / HINTS

- Exact volume of water filtered through the GF/F filter sample is required.
- Every attempt should be made to limit light exposure to the sample (pre-acidification) throughout the procedure.

9 DATA PROCESSING AND STORAGE

9.1 Sample Data Handling

The UWS Monitoring Program Coordinator will be the custodian of the finalized data files and final sample analysis results will be shared with the monitoring groups.

The monitoring group is responsible for recording the field sample data (date, time and filtered water volume) and providing the data to the lab via a chain-of-custody at the time the sample is transferred over.

9.2 Data Analysis and Calculations

$$F_s = \text{Calibration Response Factor} = \frac{C_a}{R_s}$$

Where:

C_a = Concentration of chl *a* Standard

R_s = Fluorometer reading of chl *a* Standard

$$F_m = \text{Acidification Ratio} = \frac{F_b}{F_a}$$

Where:

UWS Standard Operating Procedure
Determination of Chlorophyll *a* and Pheophytin *a* in Water

Feb 2022

Page 7 of 7

F_b = Fluorometer reading BEFORE acidification of chl *a* Standard
 F_a = Fluorometer reading AFTER acidification of chl *a* Standard

V_1 : Volume (mL) of acetone used to extract sample

V_2 : Volume (L) of water filtered

F_1 : Fluorescence reading of sample before acidification

F_2 : Fluorescence reading of sample after acidification

$$\frac{\mu\text{g}}{\text{L}} \text{Chlorophyll } a = \frac{F_s \cdot \frac{F_m}{(F_m - 1)} \cdot (F_1 - F_2) \cdot V_1}{V_2}$$

$$\frac{\mu\text{g}}{\text{L}} \text{Pheophytin } a = \frac{F_s \cdot \frac{F_m}{(F_m - 1)} \cdot (F_m \cdot F_2 - F_1) \cdot V_1}{V_2}$$

10 REFERENCES

EPA, U.S. 1997. Method 445.0: *In Vitro* Determination of Chlorophyll *a* and Pheophytin *a* in Marine and Freshwater Algae by Fluorescence.

Vaudrey, J. Department of Marine Sciences, UCONN. 2018. Chlorophyll – Fluorescence, Acidification and Non-acidification Methods.

Interstate Environmental Commission. 2017. Chlorophyll *a*.

Turner Designs. 2019. Trilogy Laboratory Fluorometer User's Manual, Version 1.7.

UMCES, Chesapeake Biological Laboratory. 2019. Fluorometric Determination of Chlorophyll *a* in waters and sediments of Fresh/Estuarine/Coastal Areas.

Appendix C-2

SOP XXXIX
Revision No. 3
December 2018
Page 1 of 14

Ammonia

1. TEST METHOD

Based on Lachat Method 10-107-6-1B (2001).

2. APPLICABLE MATRIX OR MATRICES

The method is applicable for non-potable water (ambient surface water and wastewater).

3. METHOD DETECTION LIMIT

The method detection limit is 0.100 mg/L as NH_3 . The reporting limit is 0.249 mg/L.

4. SCOPE AND APPLICATION

4.1 This method covers the determination of ammonia in surface waters, and domestic and industrial wastes. IEC will be using for the analysis of samples collected in western LIS (brackish samples) as well as periodic samples collected from wastewater treatment plants, as applicable.

5. SUMMARY OF METHOD

5.1 This method is based on the Berthelot reaction. Ammonia reacts with alkaline phenol, and sodium hypochlorite to form indophenol blue to form monochloramine, which, in the presence of phenol, catalytic amounts of nitroprusside (nitroferricyanide), and excess hypochlorite gives indophenol blue. The formation of monochloramine requires a pH between 8 and 11.5. At higher pH, ammonia may begin to oxidize to nitrate. At pH greater than 9.6, some precipitation of calcium and magnesium as hydroxides and carbonates occurs in seawater. EDTA added to the buffer prevents this from occurring. The indophenol blue measured at 630 nm is proportional to the original ammonia concentration.

The method is calibrated using standards prepared in deionized water. Once calibrated, samples of varying salinities (0 to 35 ppt) may be analyzed. The determination of background absorbance is necessary only for samples which have color absorbing at 630 nm.

If distillation is required, the sample is buffered at pH of 9.5 with a borate buffer to decrease hydrolysis of cyanates and organic nitrogen compounds, and is distilled into a solution of boric acid.

6. DEFINITIONS

The definitions and purposes below are specific to this method, but have been conformed to common usage as much as possible.

6.1 CALIBRATION BLANK (CB) – A volume of reagent water in the same matrix as the calibration standards, but without the analyte.

6.2 CALIBRATION STANDARD (CAL) -- A solution prepared from the primary dilution standard solution or stock standard solutions. The CAL solutions are used to calibrate the instrument response with respect to analyte concentration.

6.3 INSTRUMENT PERFORMANCE CHECK SOLUTION (ICP) -- A solution of one or more method analytes used to evaluate the performance of the instrument system with respect to a defined set of criteria.

6.4 LABORATORY SPIKED BLANK (LSB) -- An aliquot of reagent water or other blank matrices to which known quantities of the method analytes are added in the laboratory. The LSB is analyzed exactly like a sample, and its purpose is to determine whether the methodology is in control, and whether the laboratory is capable of making accurate and precise measurements

6.5 LABORATORY SPIKED SAMPLE MATRIX (LSM) -- An aliquot of an environmental sample to which known quantities of the method analytes are added in the laboratory. The LSM is analyzed exactly like sample, and its purpose is to determine whether the sample matrix contributes bias to the analytical results. The background concentrations of the analytes in the sample matrix must be determined in a separate aliquot and the measured values in the LFM correct for background concentrations.

6.6 LABORATORY REAGENT BLANK (LRB) – An aliquot of reagent water or equivalent neutral reference material treated as a sample in all aspects, except that it is not taken to the sampling site. The purpose is to determine if the analytes or interferences are present in the laboratory environment, the reagents, or the apparatus.

6.7 LINEAR CALIBRATION RANGE (LCR) – The concentration range over which the instrument response is linear.

6.8 MATERIAL SAFETY DATA SHEETS (MSDS) – Written information provided by vendors concerning a chemical's toxicity, health hazards, physical properties, fire, and reactivity data including storage, spill, and handling precautions.

6.9 METHOD DETECTION LIMIT (MDL) – The lowest level at which an analyte can be detection with 99 percent confidence that the analyte concentration is great than zero.

6.10 LIMIT OF QUANTITATION (LOQ) – The Laboratory performs a Limit of Quantitation (LOQ) study on an annual basis for analytes associated with chemistry methods. The validity of LOQ is confirmed by successful analysis of a Laboratory Fortified Blank (LFB) at approximately 2X the reporting limit. The acceptance criteria for each analyte is + 30% of the true value. After this study is completed, it is reviewed and approved by the Laboratory Management. A summary of all LOQ study performance is maintained.

6.11 QUALITY CONTROL SAMPLE (QCS) – A solution of method analytes of known concentrations that is used to spike an aliquot of LRB or sample matrix. THE QCS is obtained for a source external to the laboratory or is prepared from standards obtained from a different source than the calibration standards. The purpose is to check laboratory performance using test materials that have been prepared independently form the normal preparation process.

6.12 STOCK STANDARD SOLUTION (SSS) – A concentrated solution containing one or more method anlaytes prepared in the laboratory using assay reference materials or purchased from a reputable commercial source.

7. INTERFERENCES

7.1 Calcium and magnesium ions may precipitate if present in sufficient concentration. EDTA is added to the sample in-line to prevent this problem.

7.2 Color, turbidity, and certain organic species may interfere. Turbidity is removed by manual filtration. Sample color may be corrected for by running the samples through the manifold without color formation. See System Note 11 for specific instructions.

7.3 Sulfide may interfere at levels greater than 2 mg H₂S/L. Samples containing concentrations greater than this should be diluted.

7.4 Salinity does not normally interfere in this method. This may be verified by running the samples through the manifold with all reagents pumping, except hypochlorite, which is replaced by deionized water. The resulting concentrations are then compared to those obtained for samples determined with color formation.

7.5 The salt effect (salinity influence on absorbance) is less than 2%.

See System Note 11 for specific instructions.

8. SAFETY

8.1 The toxicity or carcinogenicity of each reagent used in this method has not been fully established. Each chemical should be regarded as a potential health hazard and exposure should be as low as reasonably achievable. Cautions are included for known extremely hazardous materials.

8.2 Each laboratory is responsible for maintaining a current awareness file of the Occupational Health and Safety Act (OSHA) regulation regarding the safe handling of the chemicals specified in this method. A reference file of Material Safety Data sheets(MSDS) is made available to all personnel involved in the chemical analysis.

8.3 The following chemicals have the potential to be highly toxic or hazardous, for detailed explanation consult the MSDS.

5.3.1. Sodium hydroxide

5.3.2. Phenol

9. EQUIPMENT AND SUPPLIES

9.1 Balance – analytical, capable of accurately weighing to the nearest 0.0001 g.

9.2 Glassware – Class A volumetric flasks and pipettes or plastic containers as required. Samples may be stored in plastic or glass.

9.3 Flow injection analysis equipment (Lachat 8500 series 2) designed to deliver and react sample and reagents in the required order and ratios.

9.3.1. Lachat Autoanalyzer

9.3.2. Multichannel proportioning pump

9.3.3. Reaction unit or manifold

9.3.4. Colorimetric detector

9.3.5. Data system

9.4 Special Apparatus

6.4.1. Seawater Accessories Kit Lachat Part No. 50969 – RAS or 50970 – ASX510
(for brackish samples)

6.4.2. PVC PUMP TUBES MUST BE USED FOR THIS METHOD.

10. REAGENTS AND STANDARDS

10.1 PREPARATION OF REAGENTS

Use deionized water (10 megohm) for all solutions.

Degassing with helium:

To prevent bubble formation, DO NOT DEGAS USING AN INVASIVE PROCEDURE SUCH AS A WAND TO AVOID CONTAMINATION. Degas by vacuum or sonication. DO NOT DEGAS PHENATE, HYPOCHLORITE, OR STANDARDS.

Reagent 1. Buffer Chelating Agent

By Volume: In a 1 L volumetric flask, dissolved 50.0 g disodium ethylenediamine tetraacetate (Na_2EDTA) and 11.0 g sodium hydroxide (NaOH) in about 900 mL DI water. Stir to mix and dilute to the mark with DI water. Degas as above. Prepare fresh monthly sample.

By Weight: To a tared 1 L container, add 966 g DI water. Add 50.0 g disodium ethylenediamine tetraacetate (Na_2EDTA) and 11.0 g sodium hydroxide (NaOH). Stir to mix. Degas as above. Prepare fresh monthly.

Preferably use Lachat Ammonia Buffer, EDTA/NaOH Solution CAT#52017.

Reagent 2. Phenate Reagent

CAUTION: Wear gloves. Phenol causes skin burns and is rapidly absorbed into the body through the skin. Do not degas this reagent.

By Volume: In a 1 L volumetric flask, dissolved 83 g crystalline phenol ($\text{C}_6\text{H}_5\text{OH}$) in approximately 500 mL DI water. While stirring, slowly add 32 g sodium hydroxide (NaOH). Cool, dilute to the mark with DI water, and invert to mix. The color of this reagent darkens with age, increasing the baseline absorbance. Prepare fresh reagent after 72 hours. Prepare fresh every 3-5 days and discard when turns dark brown.

By Weight: To a tared 1 L container, add 888 g DI water. Add 83 g crystalline phenol ($\text{C}_6\text{H}_5\text{OH}$). While stirring, slowly add 32 g sodium hydroxide (NaOH). Cool and stir to mix. The color of this reagent darkens with age, increasing the baseline absorbance. Prepare fresh every 3-5 days and discard when turns dark brown.

Preferably use Lacat Sodium Phenolate Reagent CAT#52005.

Reagent 3. Sodium Hypochlorite

By Volume: Dilute 250 mL sodium hypochlorite (SS290-1 hypochlorite solution, 4-6% NaOCl)

from Fisher) to 500 mL with DI water. Prepare fresh daily.

By Weight: Dilute 250 g sodium hypochlorite (SS290-1 hypochlorite solution, 4-6% NaOCl form Fisher) to 500 g with DI water. Prepare fresh daily.

Preferably use Lachat Hypochlorite Reagent CAT#52007.

Reagent 4. Sodium Nitroprusside

By Volume: Dissolve 1.75 g sodium nitroprusside in 500 mL DI water. Prepare fresh every 1-2 weeks.

By Weight: Dissolve 1.75 g sodium nitroprusside in 500 mL DI water. Prepare fresh every 1-2 weeks.

Preferably use Lachat Sodium Nitroprusside Reagent CAT#52006.

OR USE LACHAT AMMONIA REAGENT SET CAT # 52904 as an alternative option.

10.2 PREPARATION OF STANDARDS

Standard 1. Stock Ammonia Standard, 100 mg /L

By Volume: In a 1 L volumetric flask dissolve 0.3818 g ammonium chloride (NH₄Cl) FISHER CAT #3384-12 (Macron) in about 800 mL DI water. Dilute to the mark with DI water and invert to mix.

Standard 2. Working Stock Standard 10.0 mg /L

Dilute 10.0 mL of Stock Ammonia Standard to 100 mL with DI water.

Standards 3. Working Standards.

Dilute the following mL of working stock standard to 100 mL. The upper and lower value of the standard curve should not be modified. If project requires a different range the upper and lower standards may be modified but reporting limit must also be adjusted accordingly. Samples above the high standard (+10%) must be diluted. The mid-range standards listed below are recommended but may be modified by analyst.

mL Working Stock	mg/L Ammonia
40.0	4.00
20.0	2.00
10.0	1.00
5.00	0.50
1.00	0.10

11. SAMPLE COLLECTION, PRESERVATION AND STORAGE

For the Unified Water Study, samples are field-filtered according to UWS-approved QAPP and sampling SOP. Filters are then packaged, frozen and forwarded to IEC laboratory as per established UWS procedure and approved QAPP. See UWS QAPP and sampling SOP for details.

11.1 There is no single preservation method that may be recommended of all types of samples. The analyst must examine each situation critically and treat this information as a guide.

11.2 If samples must be chemically preserved, samples should be cooled, and adjusted to $\text{pH} < 2$ with H_2SO_4 . And store at 4°C in glass or polyethylene. Analyze within 28 days. Samples collected from western Long Island Sound sampling and stored in polyethylene bottles may be stored frozen up to 1 year. If samples are not run within 24 hours a storage stability study should be done.

11.3 Researchers have found serious errors when investigating the effects of filtration. The analyst should examine sample preparation and preservation techniques before routine testing.

12. QUALITY CONTROL

12.1 Each laboratory using this method is required to operate a formal quality control (QC) program. The minimum requirements of this program consist of an initial demonstration of laboratory capability, and the periodic analysis of laboratory reagent blanks, fortified blanks and other laboratory solutions as a continuing check on performance. The laboratory is required to maintain performance records that define the quality of the data that are generated.

12.2 INITIAL DEMONSTRATION OF PERFORMANCE

12.2.1. The initial demonstration of performance is used to characterize instrument performance (determination of LCRs and analysis of QCS) and laboratory performance (determination of MDLs) prior to performing analyses by this method.

12.2.2. Linear Calibration Range (LCR) --- The LCR must be determined initially and verified every 6 months or whenever a significant change in instrument response is observed and expected. The initial demonstration of linearity must use sufficient standards to insure that the resulting curve is linear. The verification of linearity must use a minimum of a blank and three standards. If any verification data exceeds the initial values by $\pm 10\%$, linearity must be nonlinear, sufficient standards must be used to clearly define the nonlinear portion.

12.2.3. Quality Control Sample (QCS) or Laboratory Control Sample (LCS)– When

beginning the use of this method, and as required to meet data-quality needs, verify the calibration standards are acceptable instrument performance with the preparation and analyses of a QCS. If the determined concentrations are not within +/- 10% of the stated values, performance of the determinative step of the method is unacceptable. The source of the problem must be identified and corrected before either proceeding with the initial determination of MDLs or continuing with ongoing analyses.

12.2.4. Method Detection Limit (MDL) – MDLs must be established for all analytes, using reagent water (blank) fortified at a concentration of two to three times the estimated instrument detection limit. To determine MDL values, take seven replicate aliquots of the fortified reagent water and process through the entire analytical method. Perform all calculations defined in the method and report the concentration values in the appropriate units. Calculate MDLs as follows:

$$MDL = t \times S$$

Where, t = Student's t value for a 99% confidence level and a standard deviation estimate with n-1 degrees of freedom [t= 3.14 for seven replicates, t = 2.528 for twenty-one replicates]. S= standard deviation of the replicate analyses.

MDLs should be determined every six months, when a new operator begins work, or whenever there is a significant change in the background or instrument response. The reporting limit is 5X the calculated MDL and must be verified at least annual, or as required for data-quality needs.

13. CALIBRATION AND STANDARDIZATION

13.1 Prepare a series of standards, covering the desired range, and a blank by dilution suitable volumes of standard solution (suggested range in Section 10.2).

13.2 Calibrate the instrument as described in Section.

13.3 Prepare standard curve by plotting instrument response against concentration values. A calibration curve may be fitted to the calibration solution concentration/response data using the computer. Acceptance or control limits should be established using the difference between the measured value of the calibration solution and the "true value" concentration.

13.4 After the calibration has established, it must be verified by the analysis of a suitable quality control sample (QCS). If measurements exceed +/- 10% of the established QCS value, the analysis should be terminated and the instrument recalibrated. The new calibration must be verified before continuing analysis. Periodic reanalysis of the QCS is recommended as a continuing calibration check.

14. PROCEDURE

14.1 CALIBRATION PROCEDURE

14.1.1 Prepare reagent and standards as described in section 10.

14.1.2 Set up manifold as shown in Lachat Ammonia Manual.

14.1.3 Input data system parameters as shown in Lachat Ammonia Manual.

14.1.4 Pump DI water through all reagent lines and check for leaks and smooth flow. Switch to reagents and allow the system to equilibrate until a stable baseline is achieved.

14.1.5 Place samples and/or standards. Input the information required by the data system, such as concentration, replicates, etc. (See section 14.2).

14.1.6 Calibrate the instrument by injecting the standards. The data system will then associate the concentrations with the instrument responses for each standard.

14.2 SYSTEM NOTES

14.2.1 For information on system maintenance and troubleshooting refer to the Troubleshooting Guide in the System Operation Manual. This guide is also available on request from Lachat.

14.2.2 Allow 15 minutes for heating module to warm up to 60 ° C.

14.2.3 Allow 15 minutes with reagents pumping for the system to equilibrate.

14.2.4 To check for gross carrier contamination:

- A. Monitor the baseline on the system unit screen.
- B. Equilibrate the pumping system with DI water in all reagent lines.
- C. Equilibrate the pumping system with reagents in the appropriate lines.
- D. The increase in absorbance should be 0.02 V or less.

14.2.5 A backpressure coil (100 cm x 0.5 mm (0.022") i.d. Teflon tubing) is used to prevent air bubble formation.

14.2.6 All reagent containers should be covered with Parafilm after insertion of the transmission lines to prevent contamination from airborne ammonia.

14.2.7 Reagent recipes from other automated wet chemistry analyzers should not be substituted.

14.2.8 If sample tube or standard container materials other than polystyrene are used, standards and samples in these containers should be analyzed to investigate absorption or contamination.

14.2.9 The blank in this method should not give a peak. If the blank peak is negative, the carrier is contaminated. If the blank peak is positive, the blank is contaminated.

14.2.10 If the detection limit is greater than that specified in the method the following outline should be followed.

- A. Verify standards preparation procedures.
- B. Verify that a 630 nm filter is being used.

- C. Verify that the sample loop is completely filled by running dye.
- D. Verify that the reagents are being added in the correct order.
 - E. Verify that the reagents are flowing smoothly. If not, check for obstructions in the manifold and replace worn tubing.
 - F. Prepare fresh reagents. Take care to be sure that the EDTA is completely dissolved.

14.2.11 If the samples are colored or are suspected to show a background absorbance, this interference should be subtracted. This can be done by using the following procedure:

- A. Calibrate the system in the normal manner.
- B. Disable the check standards or DQM features and analyze the samples.
- C. Place reagent and carrier lines in DI water and allow the baseline to stabilize.
- D. Inject samples again without recalibrating.
- E. Subtract the “background” concentration from the original concentration to give the correct concentration.

$$\text{Corrected Concentration} = \text{Original Concentration} - \text{Background}$$

Concentration

14.2.12 It is critical that the peak be detected on the “flat top” of the standard peaks. If the window is not on the “flat top”, the peak start time should be adjusted.

14.2.13 To ascertain that the standards peak is properly positioned in the window, observe the peaks on the system unit screen. The baseline trace should be flat for the duration of the baseline window. If the peak is not properly positioned, change the peak detection parameters. Low nutrient or artificial seawater, as is and spiked at 1 or 2 low levels, can be used to set timing parameters.

14.2.14 For low level analysis it is recommended that samples be analyzed in replicate from each sample cup. This is done by entering Replicates = 2 when entering sample information.

14.2.15 Use consumer bleaches with caution. Proprietary additives may contribute to staining of tubing and data quality.

14.2.16 Add reagents in the order that they appear on the manifold to reduce staining.

Data from Lachat instrument is back up periodically.

15. CALCULATIONS

15.1 Calibration is done by injecting standards. The data system will then prepare a calibration curve by plotting response versus standard concentration. Sample concentration is calculated from the regression equation.

15.2. Report only those values that fall between the lowest and highest calibration standards. Samples exceeding the highest standard should be diluted and reanalyzed.

15.3 Report results in mg/L or µg/L.

16. METHOD PERFORMANCE

16.1. Laboratory Reagent Blank (LRB) – The laboratory must analyze at least one LRB with each batch of samples. Data produced are used to assess contamination from the laboratory environment. Values that exceed the MDL indicate laboratory or reagent contamination should be suspected and corrective actions must be taken before continuing the analysis.

16.2. Laboratory Fortified Blank (LFB) – The laboratory must analyze at least one LFB (also referred to as a Quality Control Sample, QCS, or Laboratory Control Sample, LCS, with each batch of samples. Calculate accuracy as percent recovery (Section 18). If the recovery of any analyte falls outside the required control limits of 90-110%, that analyte is judge out of control, and the source of the problem should be identified and resolved before continuing analyses.

16.3. The laboratory must use LFB analyses data to assess laboratory performance against the required control limits of 90-110%.

16.4. Instruments Performance Check Solution (IPC) – For all determinations the laboratory must analyze the IPC (a mid-range check standard) and a calibration blank immediately following daily calibration, after every tenth sample (or more frequently, if required) and at the end of the sample run. Analysis of the IPC solution and calibration blank immediately following calibration must verify that the instrument is within +/- 10% of calibration. Subsequent analyses of the IPC solution must verify the calibration is still within +/- 10%. If the calibration cannot be verified within the specific limits, reanalyze the IPC solution. If the second analysis of the IPC solution confirms calibration to be outside the limits, sample analysis must be discontinued, the cause determined and/or in the case of drift the instrument recalibrated. All samples following the last acceptable IPC solution must be reanalyzed. The analysis data of the calibration blank and IPC solution must be kept on file with sample analyses data.

A demonstration of capability (DOC) should be performed each time there is a significant change in the chemistry of the method, a major modification to an existing instrument, or a new instrument is installed. A DOC is performed by each analyst designated to analyze samples using this method. An annual check must subsequently be performed and documented for each analyst using this method.

17. POLLUTION PREVENTION

17.1 Pollution prevention encompasses any technique that reduces or eliminates the quantity or toxicity of waste at the point of generation. Numerous opportunities for pollution prevention exist in laboratory operation. Whenever feasible, laboratory personnel should use pollution

prevention techniques to address their waste generation. When wastes cannot be feasibly reduced at the source, the recycling as the next best option. The quantity of chemicals purchased should be based on expected usage during its shelf life and disposal cost of unused material. Actual reagent preparation volumes should reflect anticipated usage and reagent stability. For information about pollution prevention consult the IEC Laboratory's Health and Safety Manual.

18. DATA ASSESSMENT AND ACCEPTANCE CRITERIA FOR QUALITY CONTROL MEASURES

18.1 Laboratory Spike Sample Matrix (LSM) – The laboratory must spike, in duplicate, minimum of 10% of routine samples. In each case the LSM aliquots must be duplicate of the aliquot used for sample analyses. The spiking level shall be at 1 to 5 times higher than the background concentration of the sample.

Calculate the percent recovery for each analyte, corrected for concentrations measured in the unfortified sample, and compare these values to the designated LFM recovery range 90-110%. Percent recovery may be calculated using the following equation.

$$R = \frac{C_1 - C}{s} \times 100$$

Where R= percent recovery, C_1 = fortified sample concentration, C= sample background concentration, s= concentration equivalent of analyte added to sample.

18.2 If the recovery of any analyte falls outside the designated LSM recovery range and the laboratory performance for that analyte is shown to be in control (sect. 18.3), the recovery problem encountered with the LSM is judge to be either matrix or solution related, not system related.

18.3 Computer the relative percent difference (RPD) between the two LSM results and compare the value to the designated RPD recovery range of 10%. The RPD may be calculated using the following equation:

$$RPD = \frac{(D_1 - D_2)}{(D_1 + D_2) / 2} \times 100$$

Where D1= concentration of analyte in the sample, D2= concentration of analyte in the second (duplicate) sample.

18.4 if the RPD falls outside the designated recovery range and the laboratory performance for that analyte is shown to be in control (sect 18.3), the recovery problem encountered with the

LSM is judge to be either matrix or solution related, not system related.

18.5 Where reference materials are available, they should be analyzed to provide additional performance data. The analysis of reference samples is a valuable tool for demonstrating the ability to perform the method acceptably.

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions.

19. CORRECTIVE ACTIONS FOR OUT-OF-CONTROL DATA

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions.

20. CONTINGENCIES FOR HANDLING OUT-OF-CONTROL OR UNACCEPTABLE DATA

See Quality Control, Section 12 for a description of contingencies for handling out-of-control or unacceptable data (corrective actions).

21. WASTE MANAGEMENT

The USEPA requires that laboratory waste management practice be conducted consistent with all applicable rules and regulations. Excess reagents, samples and method process wastes should be characterized and disposed of in an acceptable manner. IEC urges staff to protect the air, water and land by minimizing and controlling all releases from hoods and bench operations, complying with the letter and spirit of any water discharge permit and regulations, and by complying with all solid and hazardous waste regulations, particularly the hazardous waste identification rules and land disposal restrictions. For further information on waste management consult IEC's Health and Safety Manual and the College of Staten Island's Chemical Hygiene Plan.

22. REFERENCES

22.1. Lachat Instruments, September 2008. Determination of Ammonia (Phenolate) By Flow Injection Analysis Colorimetry. QuikChem® Method 10-107-06-1-B

22.2. U.S. Environmental Protection Agency, methods for Chemical Analysis of Water and Wastewater, EPA-600/4-79-020, Rev 2; August 1993, Method 350.1.

23. TABLES, DIAGRAMS, FLOWCHARTS, AND VALIDATION DATA:

NONE.

See Lachat Manual QuickChem Method 10-107-06-1-B (above).

Appendix C-3

Revision No. 3
December 2018
Page 1 of 15

Nitrate , Nitrite, Nitrate+Nitrite and Total Nitrogen

1) Test Method

This SOP is based on EPA method 353.2, Revision 2.0 for Nitrate and Nitrite and Lachat method 31-107-04-4-A for Total Nitrogen.

2) Applicable Matrix or Matrices

Non-potable waters.

3) Method Detection Limit

The MDL and reporting limit for Nitrate (NO₃ as N) is 0.095 mg/L (MDL) and 0.238 mg/L (RL)

The MDL and reporting limit for Nitrite (NO₂ as N) is 0.004 mg/L (MDL) and 0.01mg/L (RL)

The MDL and reporting limit for Nitrate-Nitrite (as N) is 0.098 mg/L (MDL) and 0.245 mg/L (RL)

4) Scope and Application

This method is applicable to surface and saline waters, and domestic and industrial waste waters. This analysis are performed by Flow Injection analysis (FIA).

5) Summary of Method

Nitrate is quantitatively reduced to nitrite by passage of the sample through a copperized cadmium column. The nitrite (reduced nitrate plus original nitrite) is then determined by diazotizing with sulfanilamide followed by coupling with N-(1-naphthyl) ethylenediamine dihydrochloride. The resulting water soluble dye has a magenta color which is read at 520 nm. Nitrite alone can also be determined by removing the cadmium column. Total Nitrogen can be determined by adding a persulfate digestion prior to analysis through the cadmium column.

6) Definitions

In waters and wastewaters the forms of nitrogen of greatest interest are, in order of decreasing oxidation state: nitrate, nitrite, ammonia and organic nitrogen. All these forms of nitrogen are biochemically interconvertible and are components of the nitrogen cycle. Nitrate generally occurs in trace quantities in surface water. It is an essential nutrient for many photosynthetic autotrophs

and in some cases has been identified as the growth-limiting nutrient. Nitrite is an intermediate oxidation state of nitrogen, both in the oxidation state of ammonia to nitrate and in the reduction of nitrate.

7) Interferences

7.1 Residual chlorine can produce a negative interference by limiting reduction efficiency. Before analysis, samples should be checked and if required, dechlorinated with sodium thiosulfate. Low results would be obtained for samples that contain high concentrations of iron, copper or other metals. In this method, EDTA is added to the buffer to reduce this interference.

7.2 Samples that contain large concentrations of oil and grease will coat the surface of the cadmium. This interference is eliminated by pre-extracting the sample with an organic solvent.

7.3 Sample turbidity may interfere. Remove turbidity first by filtration with 0.45 um pore diameter membrane filter prior to analysis

8) Safety

8.1 The toxicity and carcinogenicity of each reagent used in this method has not been fully established. Each chemical should be regarded as a potential health hazard and exposure to these compounds should be minimized by good laboratory practices, e.g. wear proper protective equipment, safety glasses, gloves, lab coat and working inside hoods whenever possible.

Refer to the Interstate Environmental Commission Laboratory Health and Safety Manual and CSI Chemical Hygiene Plan for specific guidelines.

For detailed explanations consult the Material Safety Data Sheets (MSDS), available in the Laboratory.

The following chemicals have the potential to be highly toxic or hazardous, for detailed explanations consult the MSDS:

Cadmium
Hydrochloric acid
Phosphoric acid
Sulfuric acid

9) Equipment and Supplies

9.1 Lachat 8500 QuikChem® Autoanalyzer , manifold reaction unit, colorimetric detector, 520 nm

wavelength filter and Omnion software data system.

9.2 Cadmium Reduction Column

9.3 Autoclave

9.4 Miscellaneous laboratory apparatus such as volumetric pipettes, flasks, etc.

10) Reagents and Standards

10.1 Digestion Reagent (Total Nitrogen only):

In a 1L volumetric flask dissolve 10 g potassium persulfate, 6.0 g boric acid, 3.0 g sodium hydroxide in approximately 800 mL DI water. Dilute to the mark and mix. Prepare fresh monthly and store in plastic.

10.1.1 Sodium persulfate Digestion Reagent recommended by Lachat for brackish and sea waters. (Total Nitrogen only):

In 1 volumetric flask, dissolve 50 g sodium persulfate in about 800 mL DI water. Stir until dissolved. Dilute to the mark, and invert to mix. Prepare fresh monthly. Store in plastic.

Analysis Reagents:

10.2 Sodium Hydroxide, 15N

Slowly add 150 g NaOH to 250 mL DI water with constant stirring. This solution will get very hot. Cool and store in plastic bottle.

10.3 Ammonium Chloride Buffer, pH 8.5

Dissolve 85.0 g ammonium chloride and 1.0 g disodium ethylenediamine tetraacetic acid dihydrate in approximately 800 mL DI water in a 1 L flask. Adjust the pH to 8.5 with 15 N sodium hydroxide. Dilute to the line and mix. Or preferably use Lachat Ammonia Chloride Buffer, pH 8.5 Reagent CAT#52003.

10.4 Sulfanilamide color reagent

Add approximately 600 mL DI water to a 1 L flask. Then add 100 mL of 85% phosphoric acid, 40.0 g sulfanilamide and 1.0 g N-(1-naphthyl) ethylenediamine dihydrochloride. Stir until dissolved, approximately 30 minutes. Dilute to the line and mix. Store in a dark bottle. This solution is stable for 1 month. Or preferably use Lachat Sulfanilamide Reagent 1 CAT#52022 and Sulfanilamide Reagent 2 CAT#52023. Mix together.

10.5 Dechlorinating reagent

Dissolve 0.35g sodium thiosulfate ($\text{Na}_2\text{S}_2\text{O}_3 \cdot 5\text{H}_2\text{O}$) in 100 ml DI water.

Standard Preparation: Prepare fresh weekly or purchase equivalent from approved vendors (obtain and retain certificates of analysis on file in the laboratory)

10.6 Stock Nitrate Standard, 100 mg/L

Dissolve 0.722 g potassium nitrate FISHER CAT# P383-500 in approximately 600 mL DI water in a 1 L flask. Add 2 mL chloroform, dilute to the line and mix. This solution is stable for 6 months.

10.7 Stock Nitrite Standard, 100 mg/L

Dissolve 0.493 g sodium nitrite FISHER CAT# S347-500 in approximately 800 mL DI water in a 1 L flask. Add 2 mL chloroform, dilute to the line and mix. Alternately, use a commercially prepared stock solution.

10.8 Working Nitrate Standard, 10 mg/L

Dilute 10.0 mL stock nitrate standard up to 100 mL with DI water.

10.9 Cadmium Efficiency Check Standard, Nitrate, 1.0 mg/L. Dilute 10 ml of 10 mg/L NO₃ standard up to 100 mL with DI water.

10.10 Working Nitrite Standard, 10 mg/L

Dilute 10.0 mL stock nitrite standard up to 100 mL with DI water.

10.11 Cadmium Efficiency Check Standard, Nitrite, 1.0 mg/L. Dilute 10 ml of 10 mg/L NO₂ standard up to 100 mL with DI water.

10.12 Working Standards:

Dilute the following amount of working stock standard to 100 mL. The upper and lower value of the standard curve should not be modified. If a project requires a different range the upper and lower standards may be modified but the reporting limit must also be adjusted accordingly. Samples above the high standard (+10%) must be diluted.

The mid-range standards listed below are recommended but may be modified by the analyst. If preparing standards for total nitrogen only, there is no need to use both the nitrate and nitrite working standard. Either one can be used but the total amount of standard used must remain the same as in the table below.

<u>mL Working Nitrate Stock</u>	<u>mL Working Nitrite Stock</u>	<u>mg/L NO₃+NO₂</u>
1.00	1.00	0.10
2.00	2.00	0.20
5.00	5.00	0.50
10.0	10.0	1.00
15.0	15.0	1.50

BS/BSD Solution:

Obtain solutions from ERA or other reliable sources. Prepare according to instructions supplied by the manufacturer.

All purchased and prepared standards and reagents are recorded in the chemical inventory logbook. All containers must be labeled with the Name, ID#, concentration, preparation date or date received, and expiration date (where applicable).

11) Sample Collection, Preservation, Shipment and Storage

For the Unified Water Study, samples are field-filtered according to UWS-approved QAPP and sampling SOP. Filters are then packaged, frozen and forwarded to IEC laboratory as per established UWS procedure and approved QAPP. See UWS QAPP and sampling SOP for details.

11.1 Samples are to be collected in plastic or glass containers and must be refrigerated to 4°C. The holding time for Nitrate and Nitrite as separate analytes is 48 hours.

11.2 Samples for Nitrate + Nitrite and Total Nitrogen analysis are to be preserved to pH <2 with sulfuric acid. The holding time for Nitrate + Nitrite and Total Nitrogen is 28 days. Samples collected from Western LIS monitoring surveys may be frozen for up to one year (after filtration) as per QAPP.

12) Quality Control

12.1 Calibration Curve

Acceptance Criteria - A minimum of 5 standards and a blank must be used to

For the Unified Water Study, samples are field-filtered according to UWS-approved QAPP and sampling SOP. Filters are then packaged, frozen and forwarded to IEC laboratory as per established UWS procedure and approved QAPP. See UWS QAPP and sampling SOP for details.

11.1 Samples are to be collected in plastic or glass containers and must be refrigerated to 4°C. The holding time for Nitrate and Nitrite as separate analytes is 48 hours.

11.2 Samples for Nitrate + Nitrite and Total Nitrogen analysis are to be preserved to pH <2 with sulfuric acid. The holding time for Nitrate + Nitrite and Total Nitrogen is 28 days. Samples collected from Western LIS monitoring surveys may be frozen for up to one year (after filtration) as per QAPP.

12) Quality Control

12.1 Calibration Curve

Acceptance Criteria - A minimum of 5 standards and a blank must be used to generate the calibration curve. The correlation coefficient must be >0.995. The % residual for each standard should be $0 \pm 10\%$ with the exception of the lowest standard which is at the reporting limit. The % residual for this standard may be $\pm 30\%$. A calibration curve should be generated at least every six months or following any major change or repair of the instrument. The calibration curve must be verified on each working day by the analysis of an IPC standard and blank.

Corrective Action - If the correlation coefficient of the calibration curve, consisting of at least five standards and a blank, is <0.995, the calibration is disallowed. The analysis must be terminated, and repeated after correcting the problem. % residuals for the standards are monitored so that any possible instrument or dilutor troubleshooting may be performed. If the IPC check does not pass, the instrument should be recalibrated.

12.2 Instrument Performance Check (IPC) Standard or Initial Calibration Verification (ICV) and Continuing Calibration Verification (CCV) Standard.

Acceptance Criteria - Analyze the IPC solution for all determinations immediately following each calibration, after every tenth sample (or more frequently, if required)

and at the end of the sample run. Analysis of the IPC solution immediately following calibration must verify that the instrument is within $\pm 10\%$ of the true value. Subsequent analyses of the continuing IPC solution must be within $\pm 10\%$ of the true value.

Corrective Action - If the calibration cannot be verified within the specified limits, analysis must be discontinued, the cause determined and/or in the case of drift the instrument re-calibrated. All samples following the last acceptable IPC solution must be reanalyzed.

12.3 Laboratory Reagent Blank (LRB), Prep Blank (PB), Initial Calibration Blank (ICB) and Continuing Calibration Blank (CCB)

Acceptance Criteria - Analyze a blank along with each batch of 20 or fewer samples. All LRB/PB/ICB/CCB results must be $<$ the Reporting Limit.

Corrective Action - If the results of the LRB/PB/ICB/CCB are $>$ the Reporting Limit then all associated samples with a concentration of $\leq 10\times$ the amount found in the LRB/PB/ICB/CCB should be reprepared and reanalyzed (sample results $\geq 10\times$ the amount found in the LRB/PB/ICB/CCB are not considered to be affected by the blank contamination or drift).

If the samples cannot be reprepared, then all affected sample results must be either 1) qualified accordingly, or 2) the reporting limit is raised to the amount found in the blank. Check with the team leader/section chief to determine which option should be used.

12.4 Laboratory Fortified Blank (LFB), Blank Spike/Blank Spike Duplicate (BS/BSD) or Quality Control Samples (QCS). (Also referred to as Laboratory Control Samples-LCS)

Acceptance Criteria - Analyze two LFB/BS/BSD/QCS samples with each batch of 20 or fewer samples. Calculate accuracy as percent recovery using the following equation:

$$\% \text{ Recovery} = \frac{\text{LFB/BS/BSD/QCS}}{s} \times 100$$

where:

LFB/BS/BSD/QCS = control sample results determined by laboratory

s = concentration equivalent of analyte added to fortify the LFB/BS/BSD/QCS solution.

The % recovery of the LFB/BS/BSD/QCS for samples analyzed under other programs should be within 85-115%. The relative percent difference (RPD) of the duplicates should not exceed 20% for aqueous standards.

Corrective Action - If the % recovery or RPD results are outside the required control limits, the affected samples should be reprepared and reanalyzed. If the samples cannot be reprepared, then all affected sample results must be qualified accordingly.

12.5 Laboratory Fortified Matrix (LFM) or Matrix Spike (MS) Recovery Acceptance Criteria. For samples analyzed under other programs, e.g. Ambient Water, prepare one LFM/MS per matrix for an analytical batch of 20 samples or less regardless of the number of different projects that comprise the analytical batch. In each case the LFM aliquot must be a duplicate of the aliquot used for sample analysis. Calculate the percent recovery, corrected for background concentration measured in the unfortified sample aliquot as per the equation below. The recovery should be 80- 120%. Calculate percent recovery using the following equation:

$$R = \frac{C_s - C}{s} \times 100$$

where: R = percent recovery,
Cs = fortified sample concentration,
C = sample background concentration, and
s = conc. equivalent of spike added to sample.

Corrective Action - If the % recovery of the LFM/MS is outside the required

control limits, and the laboratory performance is shown to be in control, the recovery problem encountered is judged to be matrix related, not system related. The native sample result of the sample used to produce the LFM/MS must be qualified accordingly.

Note: The % recovery of the LFM/MS is not evaluated if the result of the unfortified sample concentration is $>1X$ the level used to fortify the sample.

13) Calibration and Standardization

See Quality Control Section 12.1 for the procedure for preparing the calibration curve.

14) Procedure

14.1 Sample Preparation

Sample preparation is documented in the Sample Preparation Log Book

Screen all samples received under the NPDES program for chlorine prior to analysis. If chlorine is present, dechlorinate with sodium thiosulfate.

For samples requiring Total Nitrogen, pipet 10mL of sample into a culture tube. Add 5 mL of digestion solution, cap and vortex. Digest the samples in the autoclave for 30 minutes at 15 psi.

Note: If the final value of a sample or BS/BSD is higher than 2.0 mg/L, make an initial dilution prior to digestion since above this level there will be insufficient persulfate for complete oxidation.

14.2 Instrument Set-up

Turn on the Lachat instrument and install the proper manifold for the analysis to be performed. If nitrate and/or nitrite are to be reported separately, install both the nitrate + nitrite manifold and the nitrite only manifold. If only $\text{NO}_3 + \text{NO}_2$ or Total Nitrogen is to be reported, use only the $\text{NO}_3 + \text{NO}_2$ manifold. Consult the Lachat methods manual for installation details.

Make sure the cadmium column on the $\text{NO}_3 + \text{NO}_2$ manifold is off line, then pump Milli-Q water through the system and check for leaks and smooth flow. Perform any routine maintenance if necessary. Once instrument is stable, remove reagent lines from the water and place in the proper reagent receptacle. Turn on the cadmium reduction column once all reagents are flowing. **Do not run water thru the cadmium column!**

Instrument Maintenance and Repair records are entered into the Lachat Instrument Maintenance Log.

14.3 Sample Analysis

Click on the Omnion icon. From file, open the method file for the analysis being performed. Update the run worksheet with the correct sample IDs. Make any changes to the method that may be necessary, i.e. changes in standard concentrations, etc.

All Lachat methods have been configured to alert the user if the QC criteria Method Performance section 16 has been met. If the criteria is not met, follow the corrective action in the appropriate section.

Enter all standards and samples into the run worksheet protocol. Enter any necessary dilutions into the worksheet such as those done during the distillation step or those due to high level standards or samples.

When ready to start analysis, click on the start button at the top of the Windows screen.

Once the run is finished, click on tools, custom report and open format. Choose standard report and click OK. Preview the report and make any required changes before printing. Print the report and exit from the reporting area.

Data from the Lachat instrument is backed up periodically.

Close the flow to the cadmium column, then remove all reagent lines and place into a DI water receptacle. Flush system with water for at least 15 minutes. Remove lines from water and allow air to pump through system. Release tension on pump tubing by lowering arms on pump.

15) Calculations

15.1 Integration

Since the peak expectation window may shift within a method if any of the instrument conditions change, an analyst may need to reintegrate data after the run is finished. Some of the instrument conditions that affect timing are ones that will either speed up or slow down the flow of sample between the autosampler and detector such as changing pump tubing or removing a clog from a reagent line. If the peak expectation

window is adjusted for any sample in a run, the adjustment is made to each sample and standard in the entire run. When setting the peak expectation window, the entire area of the subject peak and only the area of the subject peak should be integrated. If any analytical run is altered after analysis such as modifying the peak expectation window, save the data file using the original date/time stamp and add the word reprocessed to the end so that the original and modified data file are saved. Refer to the Lachat User Manual for specific instructions on operating this software.

Sample results are calculated by the Omnion 3.0 software supplied with the Lachat autoanalyzer. The calculations are performed by determining the area of each sample or standard peak that falls within a set peak expectation window. The peak expectation window is determined separately for each method based upon the time needed for a sample to travel from the autosampler to detector and also the width of each peak.

The efficiency of the cadmium column must be calculated each time Nitrate, Nitrite and Nitrogen are analyzed. The efficiency should be between 90-110%. If the efficiency is outside of this range, stop the analysis and perform required system or column maintenance before continuing.

$$\% \text{ cadmium efficiency: } \frac{1.0 \text{ mg/L NO}_3 \text{ result}}{1.0 \text{ mg/L NO}_2 \text{ result}} \times 100$$

16) Method Performance

A demonstration of capability (DOC) should be performed each time there is a significant change in the chemistry of the method, a major modification to an existing instrument, or a new instrument is installed. A DOC is performed by each analyst designated to analyze samples using this method. An annual check must subsequently be performed and documented for each analyst using this method.

16.1 Demonstration of Capability

A demonstration of capability study must be performed and documented for each analyst using this method. The study should consist of the analysis of four standards which are from a source independent of the standard curve. The results of the standards must be within the acceptance criteria for BS/BSD samples in section 14. If samples are analyzed under more than one program, the acceptance criteria used should be the tighter limits. The % RSD should be within 20%. The results of the accuracy and precision study (true value, % recovery, standard deviation and % RSD) are maintained by the Quality Assurance Officer for each analyst and are located

in the IEC Laboratory Active Employee Training Manual.

Continuing Demonstration of Capability

An annual continuing demonstration of capability study must be performed and documented. It may consist of either successfully analyzing a PT sample or analyzing 2 sets of BS/BSD standards to within control limits as stated in section 12. The results of the continuing accuracy and precision study (true value, % recovery, standard deviation and % RSD or final report from the PT provider) are maintained by the Quality Assurance Officer for each analyst and are located in the IEC Laboratory Active Employee Training Manual.

16.2 Method Detection Limit (MDL)

An MDL Study was conducted for this method. The study is based on the requirements listed in 40 CFR Part 136 Appendix B. The MDL Study comprised the analysis of seven reagent grade water samples fortified at a level between 2-3x the detection limit. The results of the MDL determination (true value, average concentration, standard deviation and calculated MDL) are maintained by the Quality Assurance Officer for each method and are located in the Lachat analysis Log Book.

16.3 Limit of Quantitation (LOQ)

The Laboratory performs a Limit of Quantitation (LOQ) study on an annual basis for analytes associated with chemistry methods. The validity of LOQ is confirmed by successful analysis of a Laboratory Fortified Blank (LFB) at approximately 2X the reporting limit. The acceptance criteria for each analyte is $\pm 30\%$ of the true value. After this study is completed, it is reviewed and approved by the Laboratory Management. A summary of all LOQ study performance is maintained.

17) Pollution Prevention

Pollution prevention encompasses any technique that reduces or eliminates the quantity or toxicity of waste at the point of generation. Numerous opportunities for pollution prevention exist in laboratory operation. Whenever feasible, laboratory personnel should use pollution prevention techniques to address their waste generation. When wastes cannot be feasibly reduced at the source, the recycling as the next best option.

The quantity of chemicals purchased should be based on expected usage during its shelf life and disposal cost of unused material. Actual reagent preparation volumes should reflect anticipated usage and reagent stability.

For information about pollution prevention consult the IEC Laboratory's Health and Safety Manual

18) Data Assessment and Acceptance Criteria for Quality Control Measures

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions.

19) Corrective Actions for Out-of-Control Data

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions.

20) Contingencies for Handling Out-of-Control or Unacceptable Data

See Quality Control, Section 12 for a description of contingencies for handling out-of-control or unacceptable data (corrective actions)

21) Waste Management

The IEC requires that laboratory waste management practice be conducted consistent with all applicable rules and regulations. Excess reagents, samples and method process wastes should be characterized and disposed of in an acceptable manner. IEC urges staff to protect the air, water and land by minimizing and controlling all releases from hoods and bench operations, complying with the letter and spirit of any water discharge permit and regulations, and by complying with all solid and hazardous waste regulations, particularly the hazardous waste identification rules and land disposal restrictions.

IEC SOP Section No XXXXI
Revision No. 3
December 2018
Page **14** of **14**

22) References

EPA Methods for the Determination of Inorganic Substances in Environmental Samples, Rev. 2.0 August 1993, Method 353.2

Lachat QuikChem Automated Ion Analyzer Methods Manual, Method 10-107-04- 1-C. July 2008.

Standard Methods for the Examination of Water and Wastewater, Method 4500- $\text{NO}_3^- \text{I}$ Cadmium Reduction Flow Injection Method, 21st Edition, 2000

23) Tables, Diagrams, Flowcharts and Validation Data

None.

See Lachat Manual QuickChem Method 31-107-04-1-C (above).

Appendix C-4

SOP XXXXII
Revision No. 3
December 4, 2018
Page 1 of 15

TOTALPHOSPHOROUS/ORTHOPHOSPHATE

1. **Test method**

1.1 This SOP is based on EPA Method 365.1, Revision 2.0. Determination of phosphorous by semi-automated colorimetry.

2. **Applicable Matrix or Matrices**

This method is applicable to surface and saline waters, and domestic and industrial waste waters.

3. **Method Detection Limits**

MDL Orthophosphate is 0.049 mg/L Reporting Limit 0.121 mg/L
MDL Total Phosphorus is 0.074 mg/L Reporting Limit 0.186 mg/L

4. **Scope And Application**

4.1 This analysis is performed by Flow Injection analysis (FIA).

4.2 The orthophosphate method determines total orthophosphate or if the sample is filtered through a 0.45 micron pore size filter, the result is termed dissolved orthophosphate. The difference between the result of a sample determined directly and filtered is termed insoluble (particulate) orthophosphate

4.3 The Total Phosphorus method determines total phosphorus, or if the sample is filtered through a 0.45micron pore size filter, the result is termed total dissolved phosphorus. The method are based on reactions that are specific for the orthophosphate (PO_4^{3-}) ion.

5. **Summary of Method**

The orthophosphate ion (PO_4^{3-}) reacts with ammonium molybdate and antimony potassium tartrate under acidic conditions to form a complex. This complex is reduced with ascorbic acid to form a blue complex which absorbs light at 880 nm. The absorbance is proportional to the concentration of orthophosphate in the sample. Polyphosphates may be converted to the orthophosphate form by sulfuric acid digestion and organic phosphorous may be converted to orthophosphate by persulfate digestion. The determination of Total phosphorus (or total dissolved phosphorus) is performed by

performing this method on digested samples, and utilization of a Total Phosphorus manifold.

6. **Definitions**

Phosphorus is a nutrient that occurs in natural waters and in wastewaters almost solely as phosphates. Small amounts of orthophosphate or certain condensed phosphates are added to some water supplies during treatment. Larger quantities are added through use of detergents and fertilizers. Phosphorus is essential to the growth of organisms and can be the limiting nutrient for primary productivity of an organism.

7. **Interferences**

- 7.1 Orthophosphate: Only orthophosphate forms a blue color in this test. Polyphosphates and organic phosphorus compounds are not recovered. The sulfuric acid in the molybdate reagent does not have enough contact time with polyphosphates to hydrolyze them. The PO_4^{3-} ion reacts with ammonium molybdate and antimony potassium tartrate under acidic conditions to form an antimony-phosphomolybdate complex. This complex is reduced with ascorbic acid to form a blue complex which absorbs light at 880nm. The absorbance is proportional to the concentration of PO_4^{3-} in the digested sample. Sample turbidity must be removed by filtration immediately after sampling for orthophosphate samples
- 7.2 Total Phosphorus: Samples for dissolved total phosphorus should be filtered only after digestion. Silica forms a pale blue complex, which also absorbs at 880nm, this interference is generally insignificant as a silicate concentration of approximately 30 mg SiO_2/L would be required to produce a 0.005 mg P/L positive error in orthophosphate. Glassware contamination is a problem in low level phosphorus determinations. Glassware should be washed with 1:1 HCl and rinsed with deionized water. Commercial detergents should rarely be needed but, if they are used, use special phosphate-free preparations for lab glass.
- 7.3 High concentrations of iron may interfere by consuming some of the reducing agent, resulting in low recoveries. This interference is eliminated by the addition of sodium bisulfite. Sample color that absorbs in the photometric range used for analysis will also interfere
- 7.4 Arsenate is a positive interference. If it is known to be present it can also be eliminated by the addition of sodium bisulfite.
- 7.5 This method is very sensitive to contamination. Detergent can contain extremely high concentrations of phosphorus. All glassware must be scrupulously cleaned

and acid washed before use.

8. **Safety**

- 8.1 The toxicity and carcinogenicity of each reagent used in this method has not been fully established. Each chemical should be regarded as a potential health hazard and exposure to these compounds should be minimized by good laboratory practices, e.g. wear proper protective equipment, safety glasses, gloves, lab coat and working inside hoods whenever possible.
- 8.2 Refer to the IEC Laboratory Health and Safety Manual and College of Staten Island Chemical Hygiene Plan for specific guidelines.
- 8.3 For detailed explanations consult the Material Safety Data Sheets (MSDS), available in the Laboratory.
- 8.4 The following chemicals have the potential to be highly toxic or hazardous, for detailed explanations consult the MSDS.
 - 8.4.1 Sulfuric Acid
 - 8.4.2 Dodecyl Sulfate

9. **Equipment and Supplies**

- 9.1 Balance
- 9.2 Glassware-Class A volumetric flasks, pipettes or plastic containers as required. Samples may be stored in plastic or glass. All glassware used in the determination of phosphate should be washed with hot 1:1 HCl and rinsed with distilled water. Preferably, this glassware should be used only for the determination of phosphate and after use it should be rinsed with distilled water and kept covered until needed again. If this is done, the treatment with 1:1 HCl is only required occasionally.
- 9.3 Lachat Autoanalyzer, manifold reaction unit (heating unit), colorimetric detector, 880 nm wavelength filter and Omnion software data system.
- 9.4 Autoclave
- 9.5 Miscellaneous laboratory apparatus such as volumetric pipettes, flasks, etc.

10. **Reagents And Solutions**

- 10.1 Reagents
 - 10.1.1 Stock Ammonium Molybdate Solution

SOP XXXXII
Revision No. 3
December 4, 2018
Page 4 of 15

Dissolve 40.0 g ammonium molybdate tetrahydrate in approximately 800 mL of water in a 1 liter flask. Dilute to the line and mix. Store in plastic and refrigerate.

10.1.2 Stock Antimony Potassium Tartrate Solution

Dissolve 3.0g antimony potassium tartrate (potassium antimonyl tartrate hemihydrate) or 3.22 g antimony potassium tartrate (potassium antimonyl tartrate trihydrate) in approximately 800 mL of water in a 1 liter flask.

Dilute to the line and mix. Store in a dark bottle and refrigerate.

10.1.3 Molybdate Color Reagent

Add approximately 500 mL water to a 1 liter flask. Slowly add 21.0 mL of concentrated sulfuric acid while mixing. When the solution can be comfortably handled, add 72.0 mL of Stock Antimony Potassium Tartrate Solution and 213 mL of Ammonium Molybdate Solution. Dilute to 1 liter and mix. Use HACH LACHAT Color reagent CAT# 52002 as alternative.

10.1.4 Ascorbic Acid Reducing Solution, 0.33 M

Dissolve 60.0 g granular ascorbic acid in approximately 700 mL of water in a 1-liter flask. Dilute to the mark and mix. Add 1.0 g sodium dodecyl sulfate and mix. Prepare fresh weekly and discard if the solution becomes yellow. Use HACH LACHAT Ascorbic Acid reagent CAT# 52020 as alternative.

10.1.5 Carrier, Sulfuric Acid, 0.13M

Add approximately 500 mL water to a 1-liter flask. Add 7.2 mL of concentrated sulfuric acid. Dilute to the mark and mix.

10.1.6 Carrier, Dodecyl Sulfate (Recommended by Lachat for Brackish and Seawater analysis).

To 1 Liter volumetric flask containing about 900 mL of DI water, add 1.0g of dodecyl sulfate. Dilute to the mark. Stir to mix. Do not degas this reagent.

10.1.7 Ammonium Persulfate Digestion Solution

Dissolve 8.0 g of ammonium persulfate in 50 mL DI water. While mixing, slowly add 6.2 mL of concentrated sulfuric acid and dilute to 100 mL. Make fresh weekly.

10.1.8 Sodium Persulfate Digestion Solution (Recommended by Lachat for Brackish and Seawater Analysis).

SOP XXXXII
Revision No. 3
December 4, 2018
Page 5 of 15

In 1 Liter volumetric flask, dissolve 50 g sodium persulfate in about 800 mL DI water. Stir until dissolved. Dilute to the mark, and invert to mix. Prepare fresh monthly. Store in plastic.

If necessary, prevent bubble formation by degassing all prepared solutions except the standards with helium. Use He at 140kPa (20 lb/in²). Preferably, purchase equivalent solutions which should not require degassing.

10.2 Standard Preparation

10.2.1 Stock Phosphorous Standard, 100 mg/L

In a 1-liter volumetric flask, dissolve 0.4396g anhydrous potassium phosphate monobasic (KH₂PO₄) FISHER CAT# P382-500, that has been dried for one hour at 105°C.

10.2.2 Working stock standard, 10 mg/L

Dilute 10.0 mL of Stock Phosphorous Standard to 100 mL with DI water.

10.2.3 Working Standards

Dilute the following mL of working stock standard to 100 mL. The upper and lower value of the standard curve should not be modified. If a project requires a different range the upper and lower standards may be modified but the reporting limit must also be adjusted accordingly. Samples above the high standard (+10%) must be diluted. The mid-range standards listed below are recommended but may be modified by the analyst.

<u>mL Working Stock</u>	<u>mg/L Phosphorous</u>
50.0	5.00
40.0	4.00
20.0	2.00
10.0	1.00
5.0	0.500
1.0	0.100

Note: Select and prepare the required standards according to the analysis and linear range being used.

10.2.4 BS/BSD Solution

Obtain solutions from ERA, EMSL or other reliable sources. Prepare according to instructions provided by the supplier.

- 10.3 All purchased and prepared standards and reagents are recorded in the chemical inventory logbook (purchased) or the standards preparation logbook (prepared) which assigns a unique ID# to each. All containers must be labeled with the Name, ID#, concentration, preparation date and expiration date (where applicable).

11. Sample Collection, Preservation, Storage and Holding Time

For the Unified Water Study, samples are field-filtered according to UWS-approved QAPP and sampling SOP. Filters are then packaged, frozen and forwarded to IEC laboratory as per established UWS procedure and approved QAPP. See UWS QAPP and sampling SOP for details.

- 11.1 Sample containers may be of plastic or glass. All bottles must be thoroughly cleaned and acid rinsed. The volume collected should be sufficient to insure a representative sample and allow for replicate analysis if required. Samples for orthophosphate should be filtered immediately upon collection, with a maximum holding time of 48 hours.
- 11.2 Samples must be refrigerated to 4°C.
- 11.3 Samples for total phosphorous are preserved by the addition of sulfuric acid to a pH of <2 and analyzed within 28 days of collection. This is accomplished by adding no more than 2 mL concentrated H₂SO₄ per liter and verifying that the pH is less than 2. If the pH is still greater than 2, more sulfuric acid is added until the pH is <2.
- 11.4 Samples for orthophosphate are not preserved and must be analyzed within 48 hours of collection.
- 11.5 Digested samples may be store up to one month in screw cap tubes.
- 11.6 Samples collected from Western LIS monitoring survey may be frozen for up to one year (after filtration) as per QAPP.

12. Quality Control

a. Calibration Curve

Acceptance Criteria - A minimum of 5 standards and a blank must be used to generate the calibration curve. The correlation coefficient must be >0.995 . The % residual for each standard should be $0 \pm 10\%$ with the exception of the lowest standard which is at the reporting limit. The % residual for this standard may be $\pm 30\%$. A calibration curve should be generated at least every run or following any major change or repair of the instrument. The calibration curve must be verified on each working day by the analysis of an IPC standard and blank. Analyze at least five standards and a blank using a first order calibration curve. If any verification data exceeds the initial values by $\pm 10\%$, reanalyze the standards. If any portion of the range is shown to be nonlinear, sufficient standards must be used to clearly define the nonlinear portion. The LCR must be verified every six months or whenever a significant change in instrument response is observed.

Corrective Action - If the correlation coefficient of the calibration curve, consisting of at least five standards and a blank, is <0.995 , the calibration is disallowed. The analysis must be terminated, and repeated after correcting the problem. % residuals for the standards are monitored so that any possible instrument or dilutor troubleshooting may be performed. If the IPC check does not pass, the instrument should be recalibrated.

b. Instrument Performance Check (IPC) Standard or Initial Calibration Verification (ICV) and Continuing Calibration Verification (CCV) Standard.

Acceptance Criteria - Analyze the IPC solution for all determinations immediately following each calibration, after every tenth sample (or more frequently, if required) and at the end of the sample run. Analysis of the IPC solution immediately following calibration must verify that the instrument is within $\pm 10\%$ of the true value. Subsequent analyses of the continuing IPC solution must be within $\pm 10\%$ of the true value.

Corrective Action - If the calibration cannot be verified within the specified limits, analysis must be discontinued, the cause determined and/or in the case of drift the instrument re-calibrated. All samples following the last acceptable IPC solution must be reanalyzed.

c. Laboratory Reagent Blank (LRB), Prep Blank (PB), Initial Calibration Blank (ICB) and Continuing Calibration Blank (CCB)

Acceptance Criteria - Analyze a blank along with each batch of 20 or fewer samples. All LRB/PB/ICB/CCB results must be $<$ the Reporting Limit.

SOP XXXXII
Revision No. 3
December 4, 2018
Page 8 of 15

Corrective Action - If the results of the LRB/PB/ICB/CCB are > the Reporting Limit then all associated samples with a concentration of <10x the amount found in the LRB/PB/ICB/CCB should be reprepared and reanalyzed (sample results $\geq 10x$ the amount found in the LRB/PB/ICB/CCB are not considered to be affected by the blank contamination or drift).

If the samples cannot be reprepared, then all affected sample results must be either 1) qualified accordingly, or 2) the reporting limit is raised to the amount found in the blank. Check with the team leader/section chief to determine which option should be used.

- d. Laboratory Fortified Blank (LFB), Blank Spike/Blank Spike Duplicate (BS/BSD) or Quality Control Samples (QCS)

Acceptance Criteria - Analyze two LFB/BS/BSD/QCS samples with each batch of 20 or fewer samples. Calculate accuracy as percent recovery using the following equation:

$$\% \text{ Recovery} = \frac{\text{LFB/BS/BSD/QCS}}{s} \times 100$$

where:

LFB/BS/BSD/QCS = control sample results determined by laboratory

s = concentration equivalent of analyte added to fortify the LFB/BS/BSD/QCS solution.

The % recovery of the LFB/BS/BSD/QCS for samples analyzed under NPDES program should be within 10% of true value. The relative percent difference (RPD) of the duplicates should not exceed 20% for aqueous standards.

Corrective Action - If the % recovery or RPD results are outside the required control limits, the affected samples should be reprepared and reanalyzed. If the samples cannot be reprepared, then all affected sample results must be qualified accordingly.

- e. Laboratory Fortified Matrix (LFM) or Matrix Spike (MS) Recovery Acceptance

Criteria:

For samples analyzed under NPDES program, prepare one LFM/MS per matrix for an analytical batch of 10 samples or less regardless of the number of different projects that comprise the analytical batch. In each case the LFM aliquot must be a duplicate of the aliquot used for sample analysis. Calculate the percent recovery, corrected for background concentration measured in the unfortified sample aliquot as per the

SOP XXXXII
Revision No. 3
December 4, 2018
Page 9 of 15

equation below. The % recovery for NPDES or DW samples should be 90-110%. The recovery for all other programs should be 80- 120%.

Corrective Action - If the % recovery of the LFM/MS is outside the required

$$R = \frac{C_s - C}{s} \times 100$$

where: R = percent recovery,
Cs = fortified sample concentration,
C = sample background concentration, and
s = conc. equivalent of spike added to sample.

control limits, and the laboratory performance is shown to be in control, the recovery problem encountered is judged to be matrix related, not system related. The native sample result of the sample used to produce the LFM/MS must be qualified accordingly.

Note: The % recovery of the LFM/MS is not evaluated if the result of the unfortified sample concentration is >1X the level used to fortify the sample.

13. Calibration

See Section 12 Quality Control for Calibration Procedures.

14. Procedure

14.1 Sample Preparation

14.1.1 Sample preparation is documented in the Sample Digestion Log Book.

Digestion procedure for total phosphorous

14.1.2 Check sample pH before digestion and record on digestion bench sheet. Samples should be preserved to a pH of <2.

14.1.3 Measure 10 mLs of each standard and sample into a digestion tube and add 1.0 mL of persulfate digestion solution. Vortex.

14.1.4 Heat for 30 minutes in an autoclave at 121oC (15-20 psi).

14.1.5 Cool and remix on vortex mixer. If the samples are turbid after digestion, filter prior to analysis

14.2 Instrument Operating Conditions

14.2.1 Turn on the Lachat instrument and install the proper manifold for the analysis to be performed. Use the shorter heater coil and set the temperature to 37o C. Consult the Lachat methods manual for installation details if necessary.

14.2.2 Pump Milli-Q water through the system and check for leaks and smooth flow. Perform any routine maintenance if necessary. Once instrument is stable, remove reagent lines from the water and place in the proper reagent receptacle.

14.2.3 Instrument Maintenance and Repair records are entered into the Instrument Maintenance Log.

14.3 Sample Analysis

14.3.1 Click on the Omnion icon. From file, open the method file for the analysis being performed. Update the run worksheet with the correct sample IDs. Make any changes to the method that may be necessary, i.e. changes in standard concentrations, etc.

14.3.2 Place standards and samples as per the run worksheet protocol. Enter any necessary dilutions into the worksheet such as those done during the distillation step or those due to high level standards or samples. When diluting samples, use the appropriate volume of digestion solution so that the matrix matches that of the standard curve.

14.3.3 When ready to start analysis, click on the start button at the top of the Windows screen.

14.3.4 All Lachat methods have been configured to alert the user if the QC criteria in sections 14.1-14.3.5 has been met. If the criteria are not met, follow the corrective action in the appropriate section.

14.3.5 Once the run is finished, click on tools, custom report and open format. Choose standard report and click OK. Preview the report and make any required changes before printing. Print the report and exit from the reporting area.

14.3.6 Data from the Lachat instrument is backed up periodically.

14.3.7 Remove all reagent lines and place into a DI water receptacle. Flush system with water for at least 15 minutes. Remove lines from water and allow air to pump through system. Release tension on pump tubing by lowering arms on pump.

15. Calculations

15.1 Integration

Since the peak expectation window may shift within a method if any of the instrument conditions change, an analyst may need to reintegrate data after the run is finished. Some of the instrument conditions that affect timing are ones that will either speed up or slow down the flow of sample between the sampler and detector such as changing pump tubing or removing a clog from a reagent line. If the peak expectation window is adjusted for any sample in a run, the adjustment is made to each sample and standard in the entire run. When setting the peak expectation window, the entire area of the subject peak and only the area of the subject peak should be integrated. If any analytical run is altered after analysis such as modifying the peak expectation window, save the datafile using the original date/time stamp and add the word reprocessed to the end so that the original and modified datafile are saved. Refer to the Lachat User Manual for specific instructions on operating this software.

15.2 Calculations

Sample results are calculated by the Omnion 3.0 software supplied with the Lachat autoanalyzer. The calculations are performed by determining the area of each sample or standard peak that falls within a set peak expectation window. The peak expectation window is determined separately for each method based upon the time needed for a sample to travel from the sampler to detector and also the width of each peak.

16. Method Performance

A demonstration of capability (DOC) should be performed each time there is a significant change in the chemistry of the method, a major modification to an existing instrument, or a new instrument is installed. A DOC is performed by each analyst designated to analyze samples using this method. An annual check must subsequently be performed and documented for each analyst using this method.

a. Accuracy and Precision

i. Demonstration of Capability

A demonstration of capability study must be performed and documented for each analyst using this method. The study should consist of the analysis of four standards which are from a source independent of the standard curve. The results of the standards must be within the acceptance criteria for BS/BSD samples in section 14. If samples are analyzed under more than one program, the acceptance criteria used should be the tighter limits. The % RSD should be within 20%. The results of the accuracy and precision study (true value, % recovery, standard deviation and % RSD)

SOP XXXXII
Revision No. 3
December 4, 2018
Page 12 of 15

are maintained by the Quality Assurance Officer for each analyst and are located in the IEC Laboratory Active Employee Training Manual.

ii. Continuing Demonstration of Capability

An annual continuing demonstration of capability study must be performed and documented. It may consist of either successfully analyzing a PT sample or analyzing 2 sets of BS/BSD standards to within control limits as stated in section

13.1.1. The results of the continuing accuracy and precision study (true value, % recovery, standard deviation and % RSD or final report from the PT provider) are maintained by the Quality Assurance Officer for each analyst and are located in the IEC Laboratory Active Employee Training Manual.

b. Method Detection Limit (MDL)

An MDL Study was conducted for this method. The study is based on the requirements listed in 40 CFR Part 136 Appendix B. Specific procedures for conducting an MDL study can be found in SOP # G-8. The MDL Study comprised the analysis of seven reagent grade water samples fortified at a level between 2-3x the detection limit. The results of the MDL determination (true value, average concentration, standard deviation and calculated MDL) are maintained by the Quality Assurance Officer for each method and are located in the Laboratory.

c. Limit of Quantitation (LOQ)

The Laboratory performs a Limit of Quantitation (LOQ) study on an annual basis for analytes associated with chemistry methods. The validity of LOQ is confirmed by successful analysis of a Laboratory Fortified Blank (LFB) at approximately 2X the reporting limit. The acceptance criteria for each analyte is $\pm 30\%$ of the true value. After this study is completed, it is reviewed and approved by the Laboratory Management. A summary of all LOQ study performance is maintained in the Laboratory.

17. Pollution Prevention

- 17.1 Pollution prevention encompasses any technique that reduces or eliminates the quantity or toxicity of waste at the point of generation. Numerous opportunities for pollution prevention exist in laboratory operation. The IEC has established a preferred hierarchy of environmental management techniques that places pollution prevention as the management option of first choice. Whenever feasible, laboratory personnel should use pollution prevention techniques to address their waste generation.

SOP XXXXII
Revision No. 3
December 4, 2018
Page 13 of 15

- 17.2 The quantity of chemicals purchased should be based on expected usage during its shelf life and disposal cost of unused material. Actual reagent preparation volumes should reflect anticipated usage and reagent stability.

18. Data Assessment and Acceptance Criteria for Quality Control Measures

18.1 Laboratory Spike Sample Matrix (LSM) – The laboratory must spike, in duplicate, minimum of 10% of routine samples. In each case the LSM aliquots must be duplicate of the aliquot used for sample analyses. The spiking level shall be at 1 to 5 times higher than the background concentration of the sample.

Calculate the percent recovery for each analyte, corrected for concentrations measured in the unfortified sample, and compare these values to the designated LFM recovery range 90-110%. Percent recovery may be calculated using the following equation.

$$R = \frac{C_s - C}{s} \times 100$$

Where R= percent recovery, C_s = fortified sample concentration, C= sample background concentration, s= concentration equivalent of analyte added to sample.

18.2 If the recovery of any analyte falls outside the designated LSM recovery range and the laboratory performance for that analyte is shown to be in control (sect. 18.3), the recovery problem encountered with the LSM is judge to be either matrix or solution related, not system related.

18.3 Computer the relative percent difference (RPD) between the two LSM results and compare the value to the designated RPD recovery range of 10%. The RPD may be calculated using the following equation:

$$RPD = \frac{(D_1 - D_2)}{(D_1 + D_2) / 2} \times 100$$

Where D1= concentration of analyte in the sample, D2= concentration of analyte in the second (duplicate) sample.

18.4 If the RPD falls outside the designated recovery range and the laboratory performance for that analyte is shown to be in control (sect 18.3), the recovery problem encountered with the LSM is judge to be either matrix or solution related, not system related.

18.5 Where reference materials are available, they should be analyzed to provide additional performance data. The analysis of reference samples is a valuable tool for demonstrating the

ability to perform the method acceptably.

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions.

19. Corrective Actions for Out-of-Control Data

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions

20. Contingencies for Handling Out-Of-Control Data

See Quality Control, Section 12, for a discussion of QC acceptance criteria and corrective actions

21. Waste Management

The IEC requires that laboratory waste management practice be conducted consistent with all applicable rules and regulations. Excess reagents, samples and method process wastes should be characterized and disposed of in an acceptable manner. The agency urges laboratories to protect the air, water and land by minimizing and controlling all releases from hoods and bench operations, complying with the letter and spirit of any water discharge permit and regulations, and by complying with all solid and hazardous waste regulations, particularly the hazardous waste identification rules and land disposal restrictions.

22. References

Lachat QuikChem Automated Ion Analyzer Methods Manual, Method 10-115-01-1-E (2001) Determination of Total Phosphorus by Flow Injection Analysis (Acid Persulfate Digestion Method)

Lachat QuikChem Automated Ion Analyzer Methods Manual, Method 10-115-01-1-A (2007) Determination of Ortho Phosphate in Waters by Flow Injection Analysis Colorimetry

Standard Methods for the Examination of Water and Wastewater, Method 4500-P G. Flow Injection Analysis for Orthophosphate, 21st Edition, 1999.

Standard Methods for the Examination of Water and Wastewater, Method 4500-PH. Flow Injection Analysis for Total Phosphorous (Proposed), 20th Edition, 1998

SOP XXXXII
Revision No. 3
December 4, 2018
Page 15 of 15

23. Tables, Diagrams, Flowcharts, And Validation Data

See Lachat Methods 10-115-01-1-E (2001) and Lachat Method 10-115-01-1-A (2007) in Lachat methods manual